

© Copyright 2007, 2010 Applied Biosystems. All rights reserved.

Information in this document is subject to change without notice. Applied Biosystems assumes no responsibility for any errors that may appear in this document. This document is believed to be complete and accurate at the time of publication. In no event shall Applied Biosystems be liable for incidental, special, multiple, or consequential damages in connection with or arising from the use of this document.

NOTICE TO PURCHASER:

This instrument is Authorized for use in DNA sequencing and fragment analysis. This authorization is included in the purchase price of this instrument and corresponds to the up-front fee component of a license under process claims of U.S. patents and under all process claims for DNA sequence and fragment analysis of U.S. patents now or hereafter owned or licensable by Applied Biosystems for which an Authorization is required, and under corresponding process claims in foreign counterparts of the foregoing for which an Authorization is required. The running royalty component of licenses may be purchased from Applied Biosystems or obtained by using Authorized reagents purchased from Authorized suppliers in accordance with the label rights accompanying such reagents. Purchase of this instrument does not itself convey to the purchaser a complete license or right to perform the above processes. This instrument is also licensed under U.S. patents and apparatus and system claims in foreign counterparts thereof. No rights are granted expressly, by implication or by estoppel under composition claims or under other process or system claims owned or licensable by Applied Biosystems. For more information regarding licenses, please contact the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

Applied Biosystems 3130 Genetic Analyzer:

Covered by one or more of U.S. Pats. Nos. RE37,606; RE37,941.

Applied Biosystems 3130xl Genetic Analyzer:

Covered by one or more of U.S. Pats. Nos. RE37,606; RE37,941.

NOTICE TO PURCHASER

The purchase price of this Applied Biosystems 3130/3130xl Genetic Analyzer includes a grant of a limited, non-transferable license under U.S. patent claims and method claims of its foreign counterparts, and element claims of its foreign counterparts, to use this particular instrument for electrophoresis methods employing fluorescence as a means of detection. No other licenses or rights are hereby conveyed either expressly, by implication, or estoppel including, but not limited to, any claims to a composition.

The Applied Biosystems 3130 and 3130xl Genetic Analyzers includes patented technology licensed from Hitachi, Ltd. as part of a strategic partnership between Applied Biosystems and Hitachi, Ltd., as well as patented technology of Applied Biosystems.

Trademarks

Applied Biosystems, AmpFlSTR, BigDye, Coffler, GeneMapper, Identifiler, Profiler, Profiler Plus ID, SeqScape, SGM Plus, SNaPshot, and StockMarks are registered trademarks and AB (Design), Applera, Hi-Di, KB, POP, POP-4, POP-6, POP-7, SEfiler, and Yfiler are trademarks of Applied Biosystems or its subsidiaries in the U.S. and/or certain other countries.


Dell is a registered trademark of Dell Computer Corporation.

Microsoft, Windows, and Windows XP are registered trademarks of the Microsoft Corporation in the United States and other countries.

Oracle is a registered trademark of the Oracle Corporation.

All other trademarks are the sole property of their respective owners.

Part Number 4352715 Rev. D 06/2010

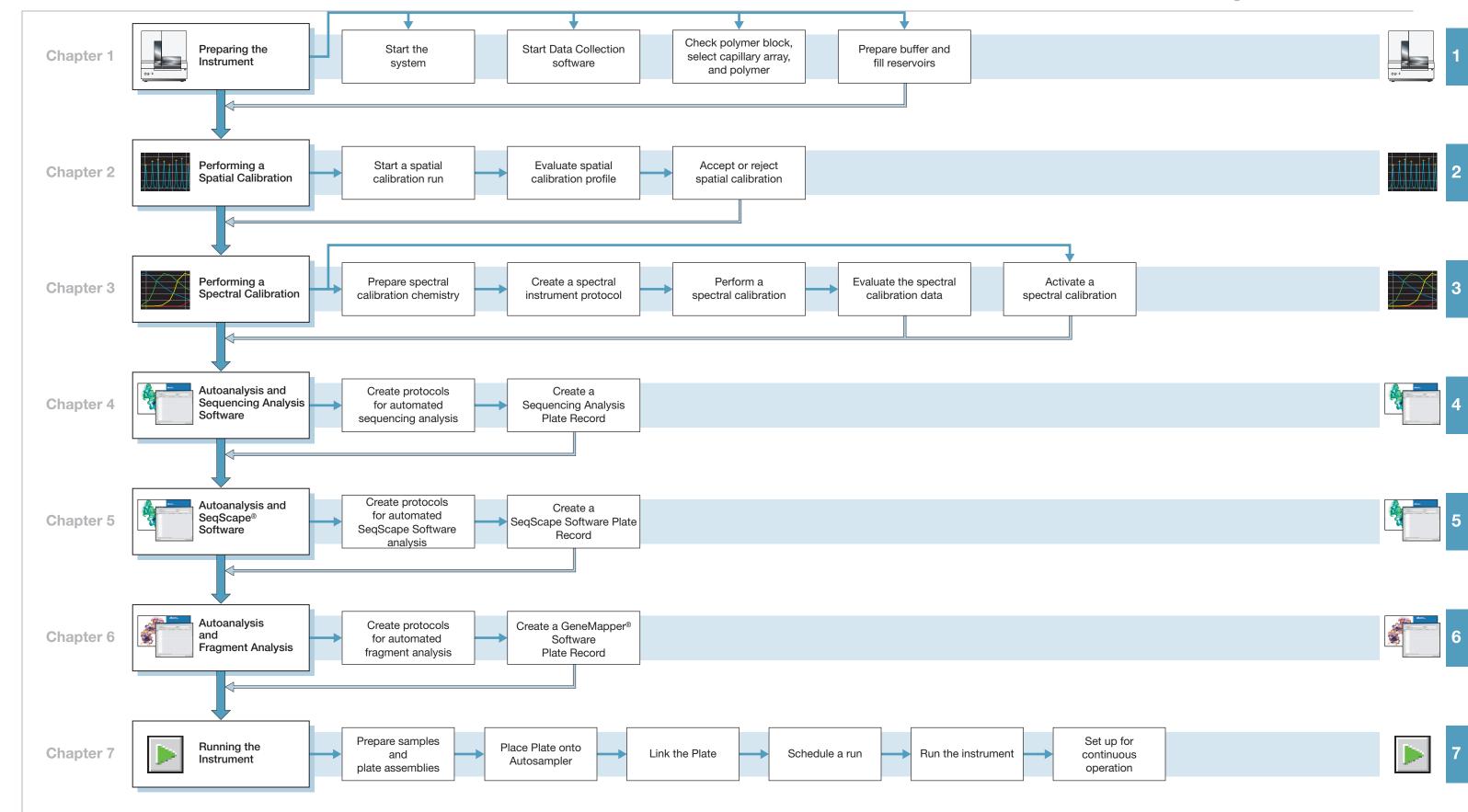

Contents

	Preface	IX
	How to Use This Guide	іх
	How to Obtain More Information	
	How to Obtain Support	x
	Safety	xiii
	Safety Conventions Used in This Document	xiv
	Symbols on Instruments	
	Safety Labels on Instruments	xv
	General Instrument Safety	xvi
	Chemical Safety	xvii
	Chemical Waste Safety	
	Electrical Safety	xx
	Physical Hazard Safety	xx
	Laser Safety	xx
	Workstation Safety	xxi
	Safety and Electromagnetic Compatibility (EMC) Standards	xxii
Chapter 1	Preparing the Instrument	1
	Instrument Description	2
	Application Summary Tables	3
	Barcode Readers	6
	Starting the 3130/3130x/ Genetic Analyzer System	7
	Data Collection Software	9
	Preparing the Instrument	11
	Preparing Buffer and Filling Reservoirs	16
Chapter 2	Performing a Spatial Calibration	21
	Spatial Calibration	22
	Creating a Spatial Calibration File	
	Evaluating a Spatial Calibration File	
	Examples of Spatial Profiles	
	Examples of Spatial Follows	

Chapter 3	Performing a Spectral Calibration	27
	Spectral Calibration	28
	Preparing the Spectral Calibration Chemistry	31
	Loading Samples	
	Creating a Spectral Instrument Protocol	
	Performing a Spectral Calibration	
	Evaluating the Spectral Calibration Data	
	Activating a Spectral Calibration	
	Examples of Passing Sequencing Spectral Calibrations	
	Examples of Fassing Fragment Analysis opecital Gallbrations	
Chapter 4	Autoanalysis and	
	Sequencing Analysis Software	53
	Sequencing Analysis Software	
	About Plate Records and Sequencing Analysis	
	Creating Protocols for Automated Sequencing Analysis	
	Analysis Protocol for Sequencing Analysis	
	Results Group for Sequencing Analysis	
	Tilling Out a dequencing Analysis Flate Necold	
Chapter 5	Autoanalysis and SeqScape Software	81
	SeqScape Software	82
	About Plate Records and Seqscape Software	83
	Creating Protocols for Automated SeqScape Software Analysis	
	Filling Out a SeqScape Software Plate Record	101
Chapter 6	Autoanalysis and Fragment Analysis	105
σπαρισι σ		
	GeneMapper Software	
	Creating Protocols for Automated Fragment Analysis	
	Creating a GeneMapper Software Plate Record	
Chapter 7	Running the Instrument	127
	Preparing Samples and Plate Assemblies	128
	Linking and Unlinking a Plate	128
	Running the Instrument	
	Controlling the Run	
	Set Up for Continuous Operation	
	Working with Data in The Run History View	
	Viewing Data from a Completed Run in the Data Collection Software	140

	Index	3
	384-Well Plate Mapping	
	Capillary Array Map	
	Injection Scheduling	
Appendix A	Plate Mapping	161
	Viewing Analyzed Data	160
	Using Autoanalysis Manager Software	155
	Viewing Autoextraction Results	147

Preface

How to Use This Guide

Purpose of This Guide

The Applied Biosystems 3130/3130xl Genetic Analyzers Getting Started Guide provides step-by-step instructions for preparing and analyzing a sample. It is designed to help you quickly learn how to use the Applied Biosystems 3130/3130xl Genetic Analyzers.

Audience

This manual is written for principle investigators and laboratory staff who are planning to operate and maintain the Applied Biosystems 3130/3130xl Genetic Analyzers.

Assumptions

This guide assumes you have the following background:

- Familiarity with Microsoft® Windows® XP operating system.
- Knowledge of general techniques for handling DNA samples and preparing them for electrophoresis.
- A general understanding of hard drives, data storage, file transfers, and copying and pasting.

If you want to integrate the 3130/3130xl Genetic Analyzers into your existing laboratory data flow system, you need networking experience.

Text Conventions

This guide uses the following conventions:

- **Bold** indicates user action. For example:
 - Type **0**, then press **Enter** for each of the remaining fields.
- *Italic* text indicates new or important words and is also used for emphasis. For example:
 - Before analyzing, *always* prepare fresh matrix.
- A right arrow bracket (>) separates successive commands you select from a dropdown or shortcut menu. For example:

Select File > Open > Spot Set.

Right-click the sample row, then select View Filter > View All Runs.

User Attention Words

Two user attention words appear in Applied Biosystems user documentation. Each word implies a particular level of observation or action as described below:

- **Note** Provides information that may be of interest or help but is not critical to the use of the product.
- **IMPORTANT!** Provides information that is necessary for proper instrument operation, accurate chemistry kit use, or safe use of a chemical.

Examples of the user attention words appear below:

Note: The size of the column affects the run time.

IMPORTANT! To verify your client connection to the database, you need a valid Oracle user ID and password.

Safety Alert Words

Safety alert words also appear in user documentation. For more information, see "Safety Alert Words" on page xiv.

How to Obtain More Information

Related Documentation

The following related document is shipped with the system:

- Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide (P/N 4352716) - Contains information on maintaining the instrument and computer system, troubleshooting tables to diagnose performance failures and advance reference content.
- Applied Biosystems 3130/3130xl Genetic Analyzers Site Preparation Guide (P/N 4352717) - Contains the space, environmental, and electrical requirements to support the 3130/3130xl Genetic Analyzer system.
- Applied Biosystems 3730/3730xl DNA Analyzers and Applied Biosystems 3130/3130xl Genetic Analyzers AB Navigator Software Administrator Guide (P/N 4359472) - Contains information and procedures for the administrator maintaining the computer system and software files of the Applied Biosystems 3130/3130xl Genetic Analyzers.
- Applied Biosystems 3130/3130xl Genetic Analyzers Quick Reference Card (P/N 4362825) Contains a flowchart on how to run your samples and instrument, a table of maintenance tasks, and a Data Collection software reference guide.

Note: For additional documentation, see "How to Obtain Support" on page xi.

Send Us Your Comments

Applied Biosystems welcomes your comments and suggestions for improving its user documents. You can e-mail your comments to:

techpubs@appliedbiosystems.com

How to Obtain Support

For the latest services and support information for all locations, go to http://www.appliedbiosystems.com, then click the link for **Support**.

At the Support page, you can:

- Search through frequently asked questions (FAQs)
- Submit a question directly to Technical Support
- Order Applied Biosystems user documents, MSDSs, certificates of analysis, and other related documents
- Download PDF documents
- Obtain information about customer training
- Download software updates and patches

In addition, the Support page provides access to worldwide telephone and fax numbers to contact Applied Biosystems Technical Support and Sales facilities.

Preface

How to Obtain Support

Safety

This safety section describes the following:
Safety Conventions Used in This Document xiv
Symbols on Instruments
Safety Labels on Instrumentsxv
General Instrument Safetyxvi
Chemical Safetyxvii
Chemical Waste Safety
Electrical Safety
Physical Hazard Safetyxx
Laser Safety
Workstation Safetyxxi
Safety and Electromagnetic Compatibility (FMC) Standards xxii

Safety Conventions Used in This Document

Safety Alert Words

Four safety alert words appear in Applied Biosystems user documentation at points in the document where you need to be aware of relevant hazards. Each alert word—**IMPORTANT, CAUTION, WARNING, DANGER**—implies a particular level of observation or action, as defined below:

Definitions

IMPORTANT! – Indicates information that is necessary for proper instrument operation, accurate chemistry kit use, or safe use of a chemical.

CAUTION – Indicates a potentially hazardous situation that, if not avoided, may result in minor or moderate injury. It may also be used to alert against unsafe practices.

WARNING – Indicates a potentially hazardous situation that, if not avoided, could result in death or serious injury.

DANGER – Indicates an imminently hazardous situation that, if not avoided, will result in death or serious injury. This signal word is to be limited to the most extreme situations.

Except for IMPORTANTs, each safety alert word in an Applied Biosystems document appears with an open triangle figure that contains a hazard symbol. *These hazard symbols are identical to the hazard icons that are affixed to Applied Biosystems instruments* (see "Safety Symbols" on page xvi).

Examples

The following examples show the use of safety alert words:

IMPORTANT! You must create a separate a Sample Entry Spreadsheet for each 96-well microtiter plate.

CAUTION The lamp is extremely hot. Do not touch the lamp until it has cooled to room temperature.

WARNING CHEMICAL HAZARD. Formamide. Exposure causes eye, skin, and respiratory tract irritation. It is a possible developmental and birth defect hazard. Read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

DANGER ELECTRICAL HAZARD. Failure to ground the instrument properly can lead to an electrical shock. Ground the instrument according to the provided instructions.

Symbols on Instruments

Electrical Symbols on Instruments

The following table describes the electrical symbols that may be displayed on Applied Biosystems instruments.

Symbol	Description
	Indicates the On position of the main power switch.
0	Indicates the Off position of the main power switch.
Φ	Indicates the On/Off position of a push-push main power switch.
Ť	Indicates a terminal that may be connected to the signal ground reference of another instrument. This is not a protected ground terminal.
(Indicates a protective grounding terminal that must be connected to earth ground before any other electrical connections are made to the instrument.
~	Indicates a terminal that can receive or supply alternating current or voltage.
=	Indicates a terminal that can receive or supply alternating or direct current or voltage.

Safety Symbols

The following table describes the safety symbols that may be displayed on Applied Biosystems instruments. Each symbol may appear by itself or in combination with text that explains the relevant hazard (see "Safety Labels on Instruments" on page xvi). These safety symbols may also appear next to DANGERS, WARNINGS, and CAUTIONS that occur in the text of this and other product-support documents.

Symbol	Description
<u> </u>	Indicates that you should consult the manual for further information and to proceed with appropriate caution.
<u></u>	Indicates the presence of an electrical shock hazard and to proceed with appropriate caution.
<u></u>	Indicates the presence of a hot surface or other high-temperature hazard and to proceed with appropriate caution.
*	Indicates the presence of a laser inside the instrument and to proceed with appropriate caution.
	Indicates the presence of moving parts and to proceed with appropriate caution.

Safety Labels on Instruments

The following CAUTION, WARNING, and DANGER statements may be displayed on Applied Biosystems instruments in combination with the safety symbols described in the preceding section.

English	Francais
CAUTION Hazardous chemicals. Read the Material Safety Data Sheets (MSDSs) before handling.	ATTENTION Produits chimiques dangeureux. Lire les fiches techniques de sûreté de matériels avant la manipulation des produits.
CAUTION Hazardous waste. Read the waste profile (if any) in the site preparation guide for this instrument before handling or disposal.	ATTENTION Déchets dangereux. Lire les renseignements sur les déchets avant de les manipuler ou de les éliminer.
CAUTION Hazardous waste. Refer to MSDS(s) and local regulations for handling and disposal.	ATTENTION Déchets dangereux. Lire les fiches techniques de sûreté de matériels et la régulation locale associées à la manipulation et l'élimination des déchets.
WARNING Hot lamp.	AVERTISSEMENT Lampe brûlante.
WARNING Hot. Replace lamp with an Applied Biosystems lamp.	AVERTISSEMENT Composants brûlants. Remplacer la lampe par une lampe Applied Biosystems.
CAUTION Hot surface.	ATTENTION Surface brûlante.
DANGER High voltage.	DANGER Haute tension.

English	Français
WARNING To reduce the chance of electrical shock, do not remove covers that require tool access. No user-serviceable parts are inside. Refer servicing to Applied Biosystems qualified service personnel.	AVERTISSEMENT Pour éviter les risques d'électrocution, ne pas retirer les capots dont l'ouverture nécessite l'utilisation d'outils. L'instrument ne contient aucune pièce réparable par l'utilisateur. Toute intervention doit être effectuée par le personnel de service qualifié de Applied Biosystems.
DANGER Class 3B laser radiation present when open and interlock defeated. Avoid direct exposure to laser beam.	DANGER Class 3B rayonnement laser en cas d'ouverture et d'une neutralisation des dispositifs de sécurité. Eviter toute exposition directe avec le faisceau.
DANGER Class 3B laser radiation when open. Avoid direct exposure to laser beam.	DANGER Class 3B rayonnement laser en cas d'ouverture. Eviter toute exposition directe avec le faisceau.
DANGER Class 2(II) laser radiation present when open and interlock defeated. Do not stare directly into the beam	DANGER de Class 2(II) rayonnement laser en cas d'ouverture et d'une neutralisation des dispositifs de securite. Eviter toute exposition directe avec le faisceau.
DANGER Class 2(II) laser radiation present when open. Do not stare directly into the beam.	DANGER de Class 2(II) rayonnement laser en cas d'ouverture. Eviter toute exposition directe avec le faisceau.
DANGER Class 2(II) LED when open and interlock defeated. Do not stare directly into the beam.	DANGER de Class 2(II) LED en cas d'ouverture et d'une neutralisation des dispositifs de securite. Eviter toute exposition directe avec le faisceau.
DANGER Class 2(II) LED when open. Do not stare directly into the beam.	DANGER de Class 2(II) LED en cas d'ouverture. Eviter toute exposition directe avec le faisceau.
CAUTION Moving parts.	ATTENTION Parties mobiles.

General Instrument Safety

WARNING PHYSICAL INJURY HAZARD. Use this product only as specified in this document. Using this instrument in a manner not specified by Applied Biosystems may result in personal injury or damage to the instrument.

Moving and Lifting the Instrument caution PHYSICAL INJURY HAZARD. The instrument is to be moved and positioned only by the personnel or vendor specified in the applicable site preparation guide. If you decide to lift or move the instrument after it has been installed, do not attempt to lift or move the instrument without the assistance of others, the use of appropriate moving equipment, and proper lifting techniques. Improper lifting can cause painful and permanent back injury. Depending on the weight, moving or lifting an instrument may require two or more persons.

Moving and Lifting Stand-Alone Computers and Monitors

WARNING Do not attempt to lift or move the computer or the monitor without the assistance of others. Depending on the weight of the computer and/or the monitor, moving them may require two or more people.

Things to consider before lifting the computer and/or the monitor:

- Make sure that you have a secure, comfortable grip on the computer or the monitor when lifting.
- Make sure that the path from where the object is to where it is being moved is clear of obstructions.
- Do not lift an object and twist your torso at the same time.
- Keep your spine in a good neutral position while lifting with your legs.
- Participants should coordinate lift and move intentions with each other before lifting and carrying.
- Instead of lifting the object from the packing box, carefully tilt the box on its side and hold it stationary while someone slides the contents out of the box.

Operating the Instrument

Ensure that anyone who operates the instrument has:

- Received instructions in both general safety practices for laboratories and specific safety practices for the instrument.
- Read and understood all applicable Material Safety Data Sheets (MSDSs). See "About MSDSs" on page xviii.

Chemical Safety

Chemical Hazard Warning

WARNING CHEMICAL HAZARD. Before handling any chemicals, refer to the Material Safety Data Sheet (MSDS) provided by the manufacturer, and observe all relevant precautions.

WARNING CHEMICAL HAZARD. All chemicals in the instrument, including liquid in the lines, are potentially hazardous. Always determine what chemicals have been used in the instrument before changing reagents or instrument components. Wear appropriate eyewear, protective clothing, and gloves when working on the instrument.

About MSDSs

Chemical manufacturers supply current Material Safety Data Sheets (MSDSs) with shipments of hazardous chemicals to *new* customers. They also provide MSDSs with the first shipment of a hazardous chemical to a customer after an MSDS has been updated. MSDSs provide the safety information you need to store, handle, transport, and dispose of the chemicals safely.

Each time you receive a new MSDS packaged with a hazardous chemical, be sure to replace the appropriate MSDS in your files.

Obtaining MSDSs

You can obtain from Applied Biosystems the MSDS for any chemical supplied by Applied Biosystems. This service is free and available 24 hours a day.

To obtain MSDSs:

- 1. Go to https://docs.appliedbiosystems.com/msdssearch.html
- 2. In the Search field, type in the chemical name, part number, or other information that appears in the MSDS of interest. Select the language of your choice, then click Search.
- **3.** Find the document of interest, right-click the document title, then select any of the following:
 - **Open** To view the document
 - **Print Target** To print the document
 - Save Target As To download a PDF version of the document to a destination that you choose
- **4.** To have a copy of a document sent by fax or e-mail, select **Fax** or **Email** to the left of the document title in the Search Results page, then click **RETRIEVE DOCUMENTS** at the end of the document list.
- **5.** After you enter the required information, click **View/Deliver Selected Documents Now**.

Chemical Safety Guidelines

To minimize the hazards of chemicals:

- Read and understand the Material Safety Data Sheets (MSDS) provided by the chemical manufacturer before you store, handle, or work with any chemicals or hazardous materials. (See "About MSDSs" on page xviii.)
- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing). For additional safety guidelines, consult the MSDS.
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use only with adequate ventilation (for example, fume hood). For additional safety guidelines, consult the MSDS.
- Check regularly for chemical leaks or spills. If a leak or spill occurs, follow the manufacturer's cleanup procedures as recommended on the MSDS.
- Comply with all local, state/provincial, or national laws and regulations related to chemical storage, handling, and disposal.

Chemical Waste Safety

Chemical Waste Hazard

CAUTION HAZARDOUS WASTE. Refer to Material Safety Data Sheets and local regulations for handling and disposal.

WARNING CHEMICAL WASTE HAZARD. Wastes produced by Applied Biosystems instruments are potentially hazardous and can cause injury, illness, or death.

Chemical Waste Safety Guidelines

To minimize the hazards of chemical waste:

- Read and understand the Material Safety Data Sheets (MSDSs) provided by the manufacturers of the chemicals in the waste container before you store, handle, or dispose of chemical waste.
- Provide primary and secondary waste containers. (A primary waste container holds the immediate waste. A secondary container contains spills or leaks from the primary container. Both containers must be compatible with the waste material and meet federal, state, and local requirements for container storage.)
- Minimize contact with chemicals. Wear appropriate personal protective equipment when handling chemicals (for example, safety glasses, gloves, or protective clothing). For additional safety guidelines, consult the MSDS.
- Minimize the inhalation of chemicals. Do not leave chemical containers open. Use
 only with adequate ventilation (for example, fume hood). For additional safety
 guidelines, consult the MSDS.
- Handle chemical wastes in a fume hood.
- After emptying the waste container, seal it with the cap provided.
- Dispose of the contents of the waste tray and waste bottle in accordance with good laboratory practices and local, state/provincial, or national environmental and health regulations.

Waste Disposal

If potentially hazardous waste is generated when you operate the instrument, you must:

- Characterize (by analysis if necessary) the waste generated by the particular applications, reagents, and substrates used in your laboratory.
- Ensure the health and safety of all personnel in your laboratory.
- Ensure that the instrument waste is stored, transferred, transported, and disposed of according to all local, state/provincial, and/or national regulations.

IMPORTANT! Radioactive or biohazardous materials may require special handling, and disposal limitations may apply.

Electrical Safety

DANGER ELECTRICAL SHOCK HAZARD. Severe electrical shock can result from operating the Applied Biosystems 3130/3130xl Genetic Analyzers without its instrument panels in place. Do not remove instrument panels. High-voltage contacts are exposed when instrument panels are removed from the instrument.

Power

DANGER ELECTRICAL HAZARD. Grounding circuit continuity is vital for the safe operation of equipment. Never operate equipment with the grounding conductor disconnected.

DANGER ELECTRICAL HAZARD. Use properly configured and approved line cords for the voltage supply in your facility.

DANGER ELECTRICAL HAZARD. Plug the system into a properly grounded receptacle with adequate current capacity.

Overvoltage Rating

The Applied Biosystems 3130/3130xl Genetic Analyzers have an installation (overvoltage) category of II, and is classified as portable equipment.

Physical Hazard Safety

Moving Parts

WARNING PHYSICAL INJURY HAZARD. Moving parts can crush and cut. Keep hands clear of moving parts while operating the instrument. Disconnect power before servicing the instrument.

Laser Safety

Laser Classification

The Applied Biosystems 3130/3130xl Genetic Analyzers use an Argon laser. Under normal operating conditions, the instrument laser is categorized as a Class I laser. When safety interlocks are disabled during certain servicing procedures, the laser can cause permanent eye damage, and, therefore, is classified under those conditions as a Class 3B laser.

The Applied Biosystems 3130/3130xl Genetic Analyzers has been tested to and complies with 21 CFR, 1040.10 and 1040.11, as applicable."

The Applied Biosystems 3130/3130xl Genetic Analyzers have been tested to and complies with standard EN60825-1, "Radiation Safety of Laser Products, Equipment Classification, Requirements, and User's Guide."

Laser Safety Requirements

To ensure safe laser operation:

- The system must be installed and maintained by an Applied Biosystems Technical Representative.
- All instrument panels must be in place on the instrument while the instrument is operating. When all panels are installed, there is no detectable radiation present. If any panel is removed when the laser is operating (during service with safety interlocks disabled), you may be exposed to laser emissions in excess of the Class 3B rating.
- Do not remove safety labels or disable safety interlocks.

Additional Laser Safety Information

Refer to the user documentation provided with the laser for additional information on government and industry safety regulations.

WARNING LASER HAZARD. Lasers can burn the retina causing permanent blind spots. Never look directly into the laser beam. Remove jewelry and other items that can reflect the beam into your eyes. Do not remove the instrument top or front panels. Wear proper eye protection and post a laser warning sign at the entrance to the laboratory if the top or front panels are removed for service.

WARNING LASER BURN HAZARD. An overheated laser can cause severe burns if it comes in contact with the skin. DO NOT operate the laser when it cannot be cooled by its cooling fan. Always wear appropriate laser safety goggles.

Workstation Safety

Correct ergonomic configuration of your workstation can reduce or prevent effects such as fatigue, pain, and strain. Minimize or eliminate these effects by configuring your workstation to promote neutral or relaxed working positions.

CAUTION MUSCULOSKELETAL AND REPETITIVE MOTION

HAZARD. These hazards are caused by potential risk factors that include but are not limited to repetitive motion, awkward posture, forceful exertion, holding static unhealthy positions, contact pressure, and other workstation environmental factors.

To minimize musculoskeletal and repetitive motion risks:

- Use equipment that comfortably supports you in neutral working positions and allows adequate accessibility to the keyboard, monitor, and mouse.
- Position the keyboard, mouse, and monitor to promote relaxed body and head postures.

Safety and Electromagnetic Compatibility (EMC) Standards

This section provides information on:

- U.S. and Canadian Safety Standards
- · Canadian EMC Standard
- European Safety and EMC Standards
- Australian EMC Standards

U.S. and Canadian Safety Standards

This instrument has been tested to and complies with standard UL 3101-1, "Safety Requirements for Electrical Equipment for Laboratory Use, Part 1: General Requirements."



This instrument has been tested to and complies with standard CSA 1010.1, "Safety Requirements for Electrical Equipment for Measurement, Control, and Laboratory Use, Part 1: General Requirements."

Canadian EMC Standard

This instrument has been tested to and complies with ICES-001, Issue 3: Industrial, Scientific, and Medical Radio Frequency Generators.

European Safety and EMC Standards

Safety



This instrument meets European requirements for safety (Low Voltage Directive 73/23/EEC). This instrument has been tested to and complies with standards EN 61010-1:2001, "Safety Requirements for Electrical Equipment for Measurement, Control and Laboratory Use, Part 1: General Requirements" and EN 61010-2-010, "Particular Requirements for Laboratory Equipment for the Heating of Materials."

EMC

This instrument meets European requirements for emission and immunity (EMC Directive 89/336/EEC). This instrument has been tested to and complies with standard EN 61326 (Class B), "Electrical Equipment for Measurement, Control and Laboratory Use – EMC Requirements."

Australian EMC Standards

This instrument has been tested to and complies with standard AS/NZS 2064, "Limits and Methods Measurement of Electromagnetic Disturbance Characteristics of Industrial, Scientific, and Medical (ISM) Radio-frequency Equipment."




Safety

Safety and Electromagnetic Compatibility (EMC) Standards




Preparing the Instrument

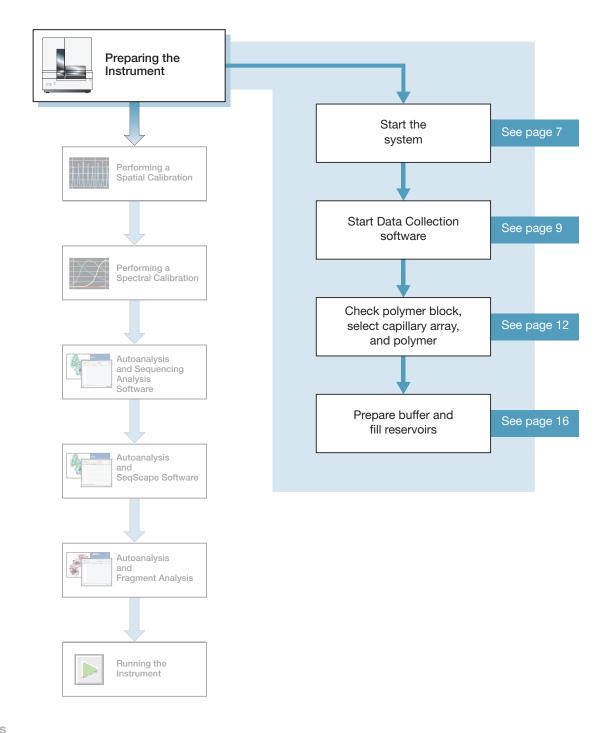

Instrument Description

Instrument Layout A typical laboratory layout is shown below.

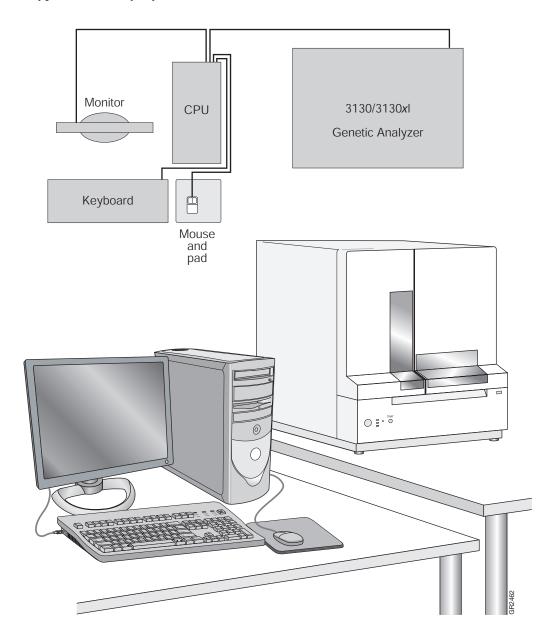

Notes			



Application Summary Tables

Sequencing Resolution Performance and **Specifications**

Decide what combination of capillary array and polymer from the table below matches your resolution and performance specifications.

Time of Divis	Capillary	Polymer		Run Sample Sample (number sample Sample Sample (min) Sample Sample (min) Sample Sample Sample (min) Sample	24 hr Throughput (number of samples)		KB Basecaller QV ₂₀ LOR ^{a b}
Type of Run	Length (cm)	Туре	Module		3130 <i>xl</i> Genetic Analyzer		
Ultra rapid	36	POP-4 [™]	UltraSeq36_POP4	40	144	576	400
		POP-7 [™]	UltraSeq36_POP7	35	164	656	500
Rapid	36	POP-6 [™]	RapidSeq36_POP6	60	96	384	500
		POP-7	RapidSeq36_POP7		96	384	600
Fast	50	POP-7	FastSeq50_POP7	60	96	384	700
Standard	50	POP-4	StdSeq50_POP4	100	56	224	600
		POP-6	StdSeq50_POP6	150	36	144	600
		POP-7	StdSeq50_POP7	120	48	192	850
Long read	80	POP-4	LongSeq80_POP4	210	24	96	700
		POP-7	LongSeq80_POP7	170	32	128	950

a Length of Read (LOR) is the usable range of high-quality or high-accuracy bases determined by Quality Values (QV) generated by KB[™] Basecaller v1.2. The LOR is determined by using a sliding window of 20 bases, which has an average QV > 20.

- N	ь	$\overline{}$	4.	~
1	A (U	U	E:

b 98.5% basecalling accuracy, less than 2% Ns.

Chapter 1 Preparing the Instrument Application Summary Tables

Fragment Analysis Resolution Performance and Specifications

Decide what combination of capillary array and polymer from the table below matches your resolution and performance specifications. See "Fragment Analysis Kits and Run Modules" on page 5 for information on capillary, polymer, and run module.

	Capillary	Polymer Type	Run Time (min)	24 hr Throughput (GT ^a)		Resolution	Specification
Run Modules	Length (cm)			3130 Analyzer	3130 <i>xl</i> Analyzer	(bp)	(SD) ^b
High Throughput, Sma	II Size Fragme	ent Analysis		-	-		+
FragmentAnalysis22 _POP4	22	POP-4	20	5,760	23,040	250	0.15
SNP22_POP4	22		20	5,760	23,040	120	0.50
Standard Fragment An	alysis						
FragmentAnalysis36 _POP4	36	POP-4	45	2,560	10,240	500	0.15
HIDFragmentAnalysis 36_POP4			45	2,560	10,240	500	0.15
SNP36_POP4			30	3,840	15,360	120	0.50
FragmentAnalysis36 _POP7		POP-7	35	3,290	13,170	500	0.15
FragmentAnalysis50 _POP4	50	POP-4	65	1,760	7,040	500	0.15
FragmentAnalysis50 _POP6		POP-6	90	1,200	4,800	500	0.15
FragmentAnalysis50 _POP7		POP-7	50	2,300	9,220	500	0.15

a 20 GT (Genotypes)/capillary/injection.

Notes			

b 1bp resolution at 99.99% accuracy.

Fragment Analysis Kits and Run Modules

The table below lists the Applied Biosystem kit types, with the available run module(s) and dye sets.

		Module								
Application/Kit	SNP22_POP4	SNP36_POP4	HTSNP36_POP7	FragmentAnalysis22_POP4	FragmentAnalysis36_POP4	FragmentAnalysis36_POP7	FragmentAnalysis50_POP4	FragmentAnalysis50_POP6	FragmentAnalysis50_POP7	HIDFragmentAnalysis36_POP4
SNaPshot® Multiplex System	E5	E5		ΐ	ь	ь	ь́	ь́	ம்	H
Custom oligos		-		D, F, G5						
Linkage Mapping Set (human) v2.5				D, G5						
ABI PRISM® Mouse Mapping Set v1.0				D, G5						
4-Dye Stockmarks® kits (bovine and canine)					F					
5-Dye Stockmarks® kit (equine)					G5					
AFLP® kits					F					
4-Dye AmpF <i>l</i> STR [®] kits										F
5-Dye AmpFℓSTR kits										G5

AmpF&STR Kit Table

Kits	HIDFragmentAnalysis 36_POP4
AmpF&STR COfiler® Kit	F
AmpF&STR Profiler Plus® Kit	
AmpF <i>t</i> STR Profiler Plus <i>ID</i> Kit	
AmpF&STR SGM Plus® Kit	
Other 4-Dye AmpFlSTR Kits	
AmpFℓSTR SEfiler [™] Kit	G5
AmpFtSTR Identifiler® Kit	
AmpFℓSTR Yfiler [™] Kit	
Other 5-Dye AmpFlSTR Kits	

Notes			

Barcode Readers

DANGER ELECTRICAL HAZARD. Power off the instrument and the computer before connecting an external barcode reader to the instrument.

External Barcode Readers

KEYENCE BL-80VE



An external barcode reader can be used with the 3130/3130xl instrument. With the KEYENCE BL-80VE (see photo above), which connects to the instrument computer keyboard, you can scan barcodes into any text box in the Data Collection software. The KEYENCE BL-80VE barcode reader uses an LED as a light source.


Notes			

Starting the 3130/3130xl Genetic Analyzer System


Starting the Computer Workstation

IMPORTANT! You must start the computer workstation before starting the instrument.

1. Power on the monitor.

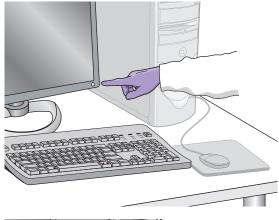

2. Power on the computer.

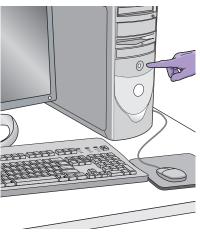

- **3.** In the Log On to Windows dialog box:
 - a. Enter the user name.
 - **b.** If applicable, enter a password.

Note: If the computer is connected to a network, you do not need to log on to the network before starting the instrument.

c. Click OK .



Starting the 3130/3130x/ Genetic Analyzer

- **1.** Ensure that the:
 - · Oven door is closed and locked
 - Instrument doors are closed

Note: If the doors are open during power on, the yellow light will continue to flash until you close the doors.

2. Ensure that the computer is powered on.


Note: The computer must be on and running to allow the instrument to copy the firmware from the computer.

3. Power on the instrument by pressing the on/off button on the front of the instrument.

Note: While the instrument is booting up and performing self-checks, the yellow status light blinks.

4. Ensure the green status light is on and not flashing before proceeding.

Note: If the green status light does not come on, start the Data Collection software and view the log. The pathway to the log is: E:\AppliedBiosystems\UDC\DataCollection\ Log\Instrument Name

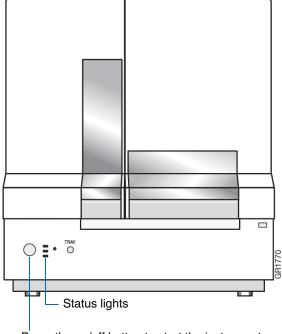

Press the on/off button to start the instrument




Data Collection Software

IMPORTANT! Do not rename the computer. The instrument computer was assigned a unique name before the 3130/3130xl genetic analyzer Data Collection software was installed. Do not rename the computer once the Data Collection software has been installed. Doing so may cause the Data Collection software to malfunction.

Starting the 3130/3130xl Data Collection Software

1. Select Start > All Programs > Applied
Biosystems > Data Collection > Run 3130
Data Collection v3.0 or Run 3130xl Data
Collection v3.0 to display the Service Console.
By default, all applications are off, indicated by
the red circles. They launch automatically with
the 3130/3130xl Data Collection software.

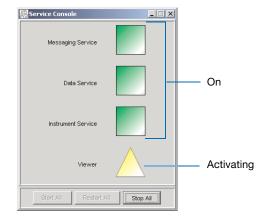

As each application activates, the red circles (off) change to yellow triangles (activating), and then to green squares (on) when they are fully functional.


When all the applications are running (displaying all green squares—this could take several minutes), the Foundation Data Collection window displays.






2. Click + to expand subfolders in the left tree pane. All application folders are now visible.

Select items from this tree pane to open manager, viewer, and other program windows.

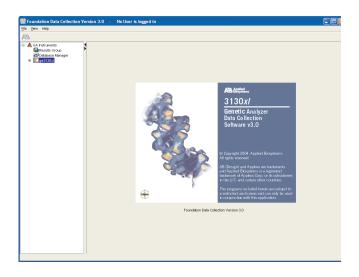

3. Click + to expand subfolders in the left tree pane. All application folders are now visible.


Select items from this tree pane to open manager, viewer, and other program windows.

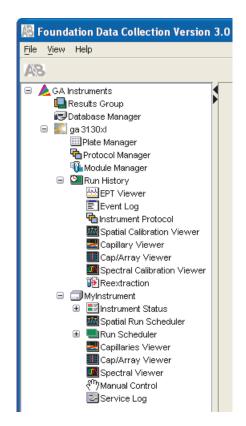



Preparing the Instrument

Instrument Doors and Interior

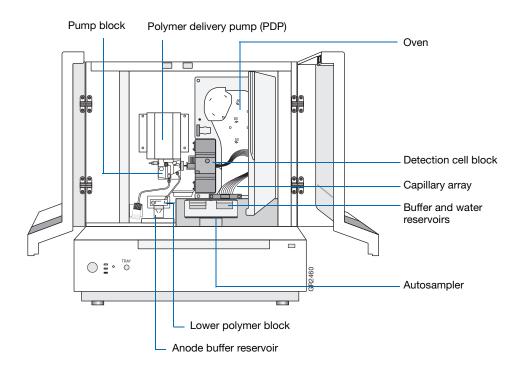


Part	Function
Anode buffer reservoir	Contains 16 mL of 1X running buffer.
Buffer and water reservoirs (four)	Each contain 16 mL of 1X running buffer or water.
Autosampler	Holds the sample plates and reservoirs and moves to align the samples, water, or buffer with the capillaries.
Capillary array	Enables the separation of the fluorescent-labeled DNA fragments by electrophoresis. It is a replaceable unit composed of 4 or 16 silica capillaries.
Detection cell block and heater	Holds the capillaries in place for laser detection.
Lower polymer block	Contains the buffer valve, anode electrode, and anode buffer reservoir.
Oven	Maintains uniform capillary array temperature.
Polymer delivery pump (PDP)	Pumps polymer into the array and performs maintenance procedures.
Pump block	Includes the displacement pump chamber, piston water seal, array attachment point (array port), and connection to the lower polymer block through the interconnect tube.

Inspecting the Instrument

- **1.** Open the instrument doors.
- **2.** Perform the daily maintenance tasks -see Chapter 1 in the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide.*
- **3.** Install clean drip trays.

Polymer Delivery Pump

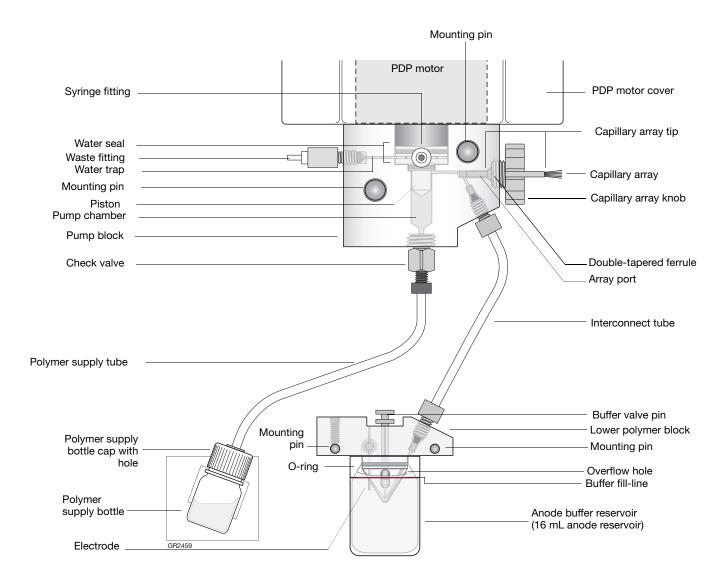

Notes			



Inspecting the Instrument

- **1.** Open the instrument doors.
- **2.** Perform the daily maintenance tasks -see the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide* for more detail on maintenance.
- **3.** Install clean drip trays.

Installing or Replacing the Capillary Array

IMPORTANT! The capillary array length defined in the wizard must match the array length you are using for correct autoanalysis results.

IMPORTANT! Wear gloves while handling polymer, the capillary array, septa, or buffer reservoirs.



CAUTION CHEMICAL HAZARD. POP

polymer may cause eye, skin, and respiratory tract irritation. Please read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Use for research and development purposes only.

 Install a capillary array by clicking <instrument name>, selecting Install Array Wizard and following the prompts.

IMPORTANT! To install or replace an array that is a different length than the one you were using, reset the active spectral calibration (see page 47) or create a new spectral calibration for your dye set and array length combination (see page 37).

2. Optional: Select **Update Cap Array Info** wizard to correct any errors or update capillary array and serial number information.

Wizards Help

Install Array Wizard
Change Polymer Type Wizard
Replenish Polymer Wizard
Bubble Remove Wizard
Water Wash Wizard
Instrument Shutdown Wizard
Autosampler Calibration Wizard
Update Cap Array Info

When to Replenish or Change Polymer

IMPORTANT! Always replace polymer that has been on the instrument longer than one week.



If polymer on the instrument	Then	
has been on less than one week and is in sufficient quantity to complete your runs ^a	Remove all bubbles, and then proceed with instrument preparation.	
has been on less than one week, and insufficient in quantity to complete your runs	Add fresh polymer to the polymer supply by following the Replenish Polymer Wizard.	
has been on longer than one week		
is the wrong type (a change between POP-4, POP-6, and/or POP-7 polymers is required)	Replace the installed polymer type with a different type by following the Change Polymer Type Wizard.	

a A 3130x/ genetic analyzer run uses 50 to 80 μ L of polymer and a 3130 genetic analyzer run uses ~25 to 40μ L of polymer.

Before Using the Polymer

- **1.** Remove the polymer from 4 °C storage.
- **2.** Loosen the cap and bring the polymer to room temperature.
- **3.** To dissolve deposits, tighten the cap and gently swirl the polymer.

Replenishing the Polymer

IMPORTANT! Wear gloves while handling polymer, the capillary array, septa, or buffer reservoirs.



CAUTION CHEMICAL HAZARD. POP **polymer** may cause eye, skin, and respiratory tract irritation. Please read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Use for research and development purposes only.

1. Click < **Instrument Name**> in the tree pane.

Notes		

Preparing the Instrument



2. Select Wizards > Replenish Polymer Wizard to replenish polymer.

IMPORTANT! The polymer type defined in the wizard must match the polymer type that you are using.

Changing Polymer Type

IMPORTANT! Wear gloves while handling polymer, the capillary array, septa, or buffer reservoirs.



CAUTION CHEMICAL HAZARD. POP **polymer** may cause eye, skin, and respiratory tract irritation. Please read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves. Use for research and development purposes only.

- 1. Click < Instrument Name > in the tree pane.
- 2. Select Wizards > Change Polymer Type Wizard to change to a different polymer.

Wizards Help

Install Array Wizard Change Polymer Type Wizard Replenish Polymer Wizard Bubble Remove Wizard Water Wash Wizard Instrument Shutdown Wizard Autosampler Calibration Wizard

Update Cap Array Info

Wizards Help

Install Array Wizard Change Polymer Type Wizard Replenish Polymer Wizard Bubble Remove Wizard Water Wash Wizard Instrument Shutdown Wizard Autosampler Calibration Wizard Update Cap Array Info

Preparing Buffer and Filling Reservoirs

Required Materials


Genetic Analyzer Buffer with EDTA may cause eye, skin, and respiratory tract irritation. Read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

Required materials to prepare 1X running buffer:

- 10X Genetic Analyzer Buffer (P/N 402824)
- Purified (distilled or deionized) water
- 50 mL graduated cylinder

Preparing Buffer for a Single Run


- **1.** Add 5 mL of 10X Genetic Analyzer buffer into a graduated cylinder.
- **2.** Add purified water to bring the total volume up to 50 mL.
- 3. Mix well.

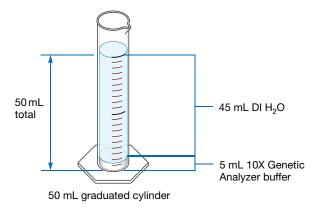

Storing Buffer

Store 1X running buffer at:

- 2 to 8 °C for up to 1 month
- Room temperature for 1 week

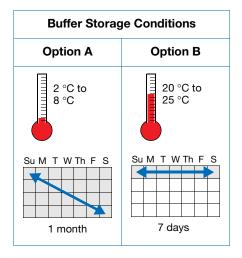

N	Л	$\overline{}$	+	$\overline{}$	0
ı,	И	U	U	U	J



Replacing Buffer and Water

Replace the 1X running buffer in the anode and cathode buffer reservoirs daily, or before each batch of runs.

IMPORTANT! Failing to replace buffer may lead to loss of resolution and data quality.

IMPORTANT! To clean and refill the reservoirs, the autosampler is brought to the forward position, thereby leaving the capillary tips exposed to the air. Do not leave the autosampler in this position for an extended time because the capillaries can dry out and the array may be damaged.

Filling the Water and Cathode Buffer Reservoirs


IMPORTANT! Wear gloves while performing the following procedure, and any other time you handle the capillary array, septa, or buffer reservoirs.



CAUTION CHEMICAL HAZARD. 1X

Genetic Analyzer Buffer with EDTA may cause eye, skin, and respiratory tract irritation. Read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

- **1.** Verify the oven and instrument doors are closed.
- **2.** Press the Tray button on the outside of the instrument to bring the autosampler to the forward position
- **3.** Wait until the autosampler stops at the forward position, then open the instrument door.

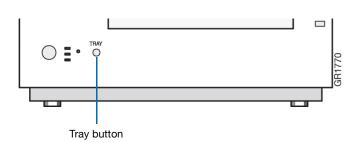



Chapter 1 Preparing the Instrument Preparing Buffer and Filling Reservoirs

- **4.** Remove the cathode buffer and water reservoirs from the instrument.
- **5.** Dispose of remaining fluids and rinse out the reservoirs with deionized water.


Note: Follow your company's waste disposal practices for appropriate disposal procedures.


- **6.** Rinse the cathode reservoir with 1X running buffer, and then fill to the line with 1X running buffer (about 16 mL).
- **7.** Fill the two water reservoirs to the line with quality deionized water (about 16 mL).


CAUTION Be sure that the septa fit securely and flush on the tops of the reservoirs to prevent damaging the capillary tips.

- **8.** Place a clean reservoir septa on each reservoir, and dry the outside of the reservoirs using a lint-free tissue wipe.
- **9.** Place the reservoirs into position on the autosampler as shown.
- **10.** Close the instrument doors.

Note: Closing the doors returns the autosampler to the last known position, placing the tips of the capillaries in water or buffer.

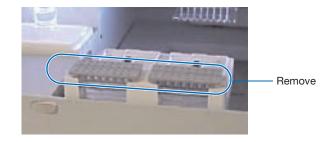

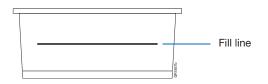

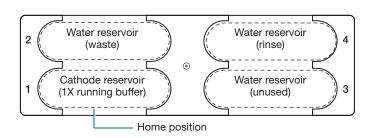





Filling the Anode Buffer Reservoir

Change the anode buffer:


- Every 24 hours
- Before each run or batch of runs
- Every time you replenish the polymer or change polymer type

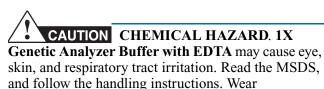

appropriate protective eyewear, clothing, and gloves.

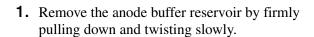


- **2.** Discard the used buffer appropriately.
- **3.** Clean and rinse the reservoir with deionized water, and then rinse with buffer.

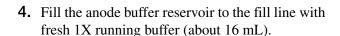


Note: The meniscus should line up with the fill line.

- **5.** Put the anode buffer reservoir on the instrument.
- **6.** If the reservoir fills with fluid, repeat this procedure to discard and replace the running buffer.


Note: The reservoir could fill during bubble removal.

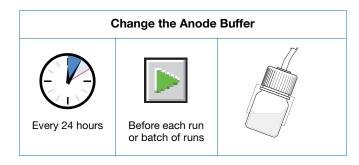



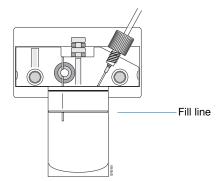




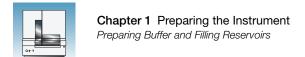

Notes	
Notes	
20	A 1' 1 B'
20	Applied Biosystems 3130/3130x/ Genetic Analyzers Getting Started Guide



Performing a Spatial Calibration

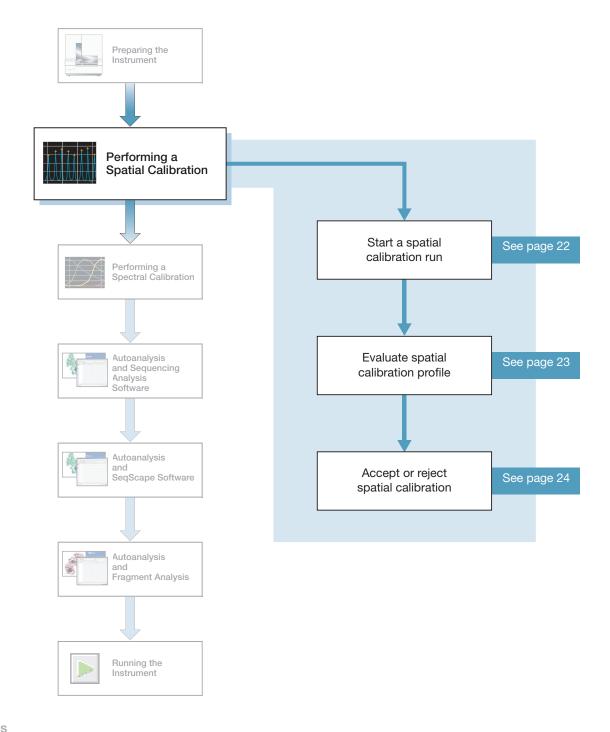

Spatial Calibration

The 3130/3130xl genetic analyzer Data Collection software uses images collected during the spatial calibration to establish a relationship between the signal emitted by each capillary and the position where that signal falls on and is detected by the CCD camera.

When to Perform the Calibration

You are required to perform a spatial calibration when you:


- Install or replace a capillary array
- Temporarily remove the capillary array from the detection block
- Move the instrument

Creating a Spatial Calibration File

1. In the tree pane of the Data Collection software, click

▲ GA Instruments > ■ ga3130 or ga3130xl > □ instrument

name > ■ Spatial Run Scheduler.

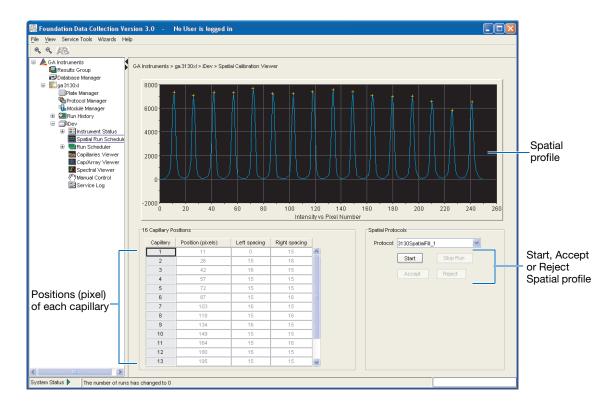



- **2.** In the Spatial Protocols section, select one of the following:
 - If the capillaries contain fresh polymer, select Protocol > 3130SpatialNoFill 1
 - Otherwise, select **Protocol** > 3130SpatialFill_1


Note: You do not need to fill the capillaries each time you perform a spatial calibration.

3. Click Start.

The calibration run lasts approximately:

- 2 min. without filling the capillaries
- 6 min. when filling the capillaries

Note: The spatial profile window turns black when you start a spatial calibration.

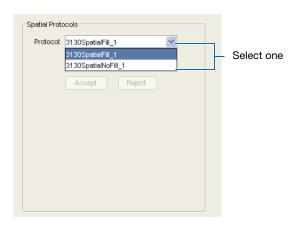




Evaluating a Spatial Calibration File


1. Evaluate the spatial calibration profile using the following criteria:

Peak Attribute	Acceptable Criteria
Height	Similar heights for all peaks.
Orange crosses	One orange cross marking the top of every peak. No misplaced crosses.
Shape	Single sharp peak for each capillary. Small shoulders are acceptable.
Spacing	The difference between adjacent positions is 13 to 16 pixels. Theoretical spacing between capillaries is 15.

Spatial calibration profile for 3130xl system

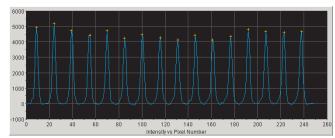

Spatial calibration profile for 3130 system

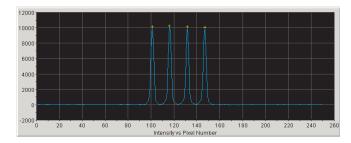





2. Examine each row in the 16 or 4 Capillary Positions table and verify that the values in both the Left spacing and Right spacing columns range between 13 to 16 pixels.

To move the cross:

- **a.** Type a new value in the Positions (pixels) box for the capillary of interest.
- **b.** Click outside of that box or press **Enter**.


Capillary	Position (pixels)	Left spacing	Right spacing	
1	10	0	15	1
2	25	15	16	
3	41	16	15	
4	56	15	16	
5	72	16	15	
6	87	15	15	1
7	102	15	15	
8	117	15	15	
9	132	15	16	
10	148	16	15	
11	163	15	15	1
12	178	15	16	
13	194	16	15	

	ry Position (pixels)	Left spacing	Right spacing
96	96	0	15
111	111	15	15
126	126	15	15
141	141	15	.0
ht spacing	eft spacing and ght spacing blumns		

Accepting or Rejecting a Spatial Calibration

If the calibration:


- Passed, click Accept to write the calibration data to the database and .ini file.
- Failed, click Reject, then see the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide.

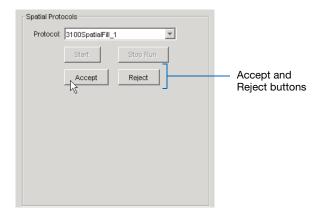




Examples of Spatial Profiles


Passing Profiles 3130 Genetic Analyzer

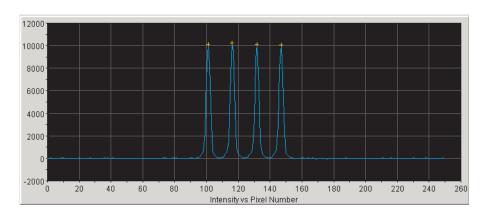

3130xl Genetic Analyzer

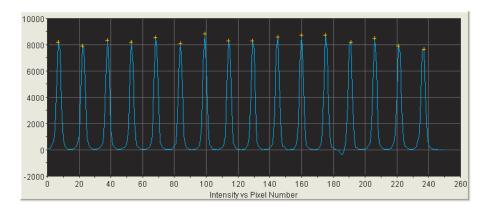

Failing Profile 3130xl Genetic Analyzer

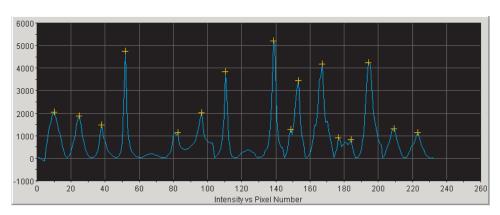




Notes____



Performing a Spectral Calibration

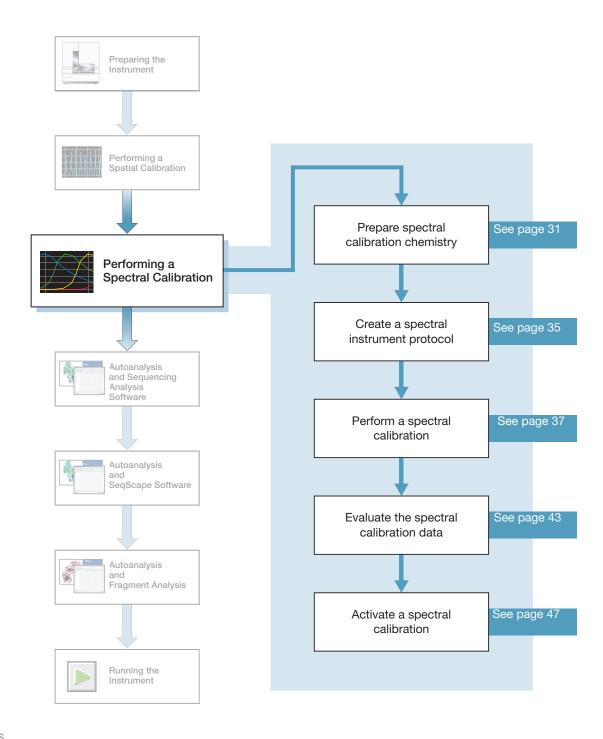

Spectral Calibration

A spectral calibration creates a matrix that is used during a run to reduce raw data from the instrument to the 4-dye or 5-dye data stored in sample files. Performing a spectral calibration is similar to performing a sample run except that calibration standards are run in place of samples and a spectral calibration module is used in place of a run module.

When to Perform the Calibration

Perform a spectral calibration:

- When you use a new dye set on the instrument
- When you change capillary array length (or polymer type for fragment analysis)
- After the laser or CCD camera has been realigned/replaced by a service engineer
- If you begin to see a decrease in spectral separation (pull-up and/or pull-down peaks) in the raw or analyzed data

What Happens?

Run the spectral standards in all 16 or 4 capillaries. The Data Collection software then:

- Collects the data and stores it in 16 or 4 separate temporary files
- Analyzes the data and generates a matrix for each capillary
- Stores the spectral calibration data for the dye set run

Changing Capillary Array Lengths and Polymer Type

For each dye set, a unique spectral calibration is used for each different capillary array length.

- For every sequencing dye set, you must create a separate spectral calibration for each capillary array length you use.
- For every fragment analysis dye set, you must create a separate spectral calibration for each capillary array length and polymer type combination you use.

See "Activating a Spectral Calibration" on page 47, on how to switch calibrations once calibrations are performed for each dye set on each capillary length.

Supported Sequencing Chemistries

New basecaller, mobility files, and run modules will be developed for the 3130/3130xl series instruments solely in support of the ABI PRISM® BigDye® Terminator v1.1 and v3.1chemistry kits. For more information, please contact Applied Biosystems Technical Support.

ABI PRISM® BigDye® Terminator v1.0, v2.0 and v3.0 chemistry kits have been discontinued.

Notes			



Types of Calibration **Standards**

There are two types of spectral calibration standards:

- Matrix standards for fragment analysis or sequencing A tube that contains four or five fragments each labeled with a different single dye.
- BigDye® v3.1 or BigDye® v1.1 Terminator Sequencing Standard A tube of a standard chemistry reaction that contains multiple labeled fragments in each of the four dyes

Dye Set Tables

Determine the correct dye set and matrix standard set for the application you are using by using the tables below.

Sequencing Chemistry Dye Sets, Calibration Standards, and Chemistry File


Sequencing Chemistry	Dye Set	Spectral Calibration Standard	Chemistry File
BigDye® Terminator v3.1 Cycle	Z_BigDyeV3	BigDye® v3.1 Matrix Standards	Matrix Standard
Sequencing Kit • ABI PRISM® dGTP BigDye® Terminator v 3.0 Cycle Sequencing Ready Reaction Kita		BigDye® v3.1 Terminator Sequencing Standard	Sequence Standard
BigDye® Terminator v1.1 Cycle	E_BigDyeV1	DS-01 Matrix Standards	Matrix Standard
Sequencing Kit • ABI PRISM® BigDye® Primer Cycle Sequencing Kits		BigDye® v1.1 Terminator Sequencing Standard	Sequence Standard
ABI PRISM® dGTP BigDye® Terminator Cycle Sequencing Ready Reaction Kita			
ABI PRISM [®] dRhodamine Terminator Cycle Sequencing Kit		dRhodamine Matrix Standards Kit	Matrix Standard

a dGTP kits are not supported on capillary electrophoresis instruments due to compressions on certain sequence context regions; you can run the kits if compression is not an issue.

Fragment Analysis Dye Sets, Calibration Standards, and Chemistry File

Fragment Analysis Chemistry	Dye Set	Spectral Calibration Standard	Chemistry File
Custom oligos	D	DS-30 Matrix Standards	Matrix Standard
ABI PRISM® Mouse Mapping Set v1.0	D	DS-31 Matrix Standards	
Custom oligos			
AFLP® kits	F	DS-32 Matrix Standards	
Stockmarks® Kits 4-dye (bovine and canine)			
AmpFtSTR® COfiler® Kit			
AmpFtSTR® Profiler Plus® Kit			
AmpFtSTR® Profiler Plus® ID Kit			
AmpFtSTR® SGM Plus® Kit			
Other 4-Dye AmpFlSTR Kits			
ABI PRISM® SNaPshot® Multiplex System	E5	DS-02 Matrix Standards	
ABI PRISM® Linkage Mapping Set v2.5	G5	DS-33 Matrix Standards	
Stockmarks® Kit 5-dye (equine)			
Custom Oligos			
AmpFtSTR® Identifiler® Kit			
 AmpFℓSTR® SEfiler[™] Kit 			
 AmpFℓSTR® Yfiler[™] Kit 			
Other 5-Dye AmpFlSTR Kits			

Notes			

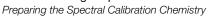




Preparing the Spectral Calibration Chemistry

Preparing the Calibration Standard

1. Prepare one of the following:

WARNING CHEMICAL HAZARD. **Formamide** causes eye, skin, and respiratory tract irritation. It is a possible reproductive and birth defect hazard. Read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

- ABI PRISM BigDye Terminator v1.1 or v3.1 Sequencing Standard:
 - Remove a tube of the Sequencing Standard from the freezer.
 - Add 170 μL of Hi-Di[™] formamide to resuspend the BigDye Terminator v1.1 or v3.1 Sequencing Standard.
- Sequencing or fragment analysis matrix standards:
 - Remove a tube of the matrix standard from the refrigerator.
 - Mix thoroughly, then spin briefly in a microcentrifuge.
 - Follow the matrix standard insert for matrix standard and Hi-Di formamide ratios.
- **2.** Vortex thoroughly. Briefly centrifuge the mixture.
- **3.** Heat the standard tube at 95 °C for 5 minutes to denature the DNA.
- **4.** Cool the tubes on ice for 2 minutes.

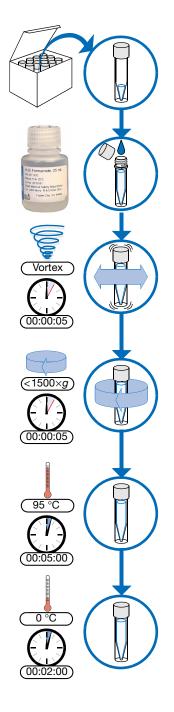

Loading Samples

chemicals on the instrument, including liquid in the lines, are potentially hazardous. Please read the MSDS, and follow the handling instructions. Wear

appropriate eyewear, protective clothing, and gloves when working on the instrument.



Do not use warped



Note: The efficient way to lay out samples in a plate is illustrated here. See Appendix A, "Plate Mapping," for the relationship between sample position and priority in scheduled injections.

To load samples:

- **1.** Add the denatured samples to the wells of a 96-or 384-well reaction plate:
 - If using a 3130xl genetic analyzer:
 - 96-well plate Add 10 μL of denatured standard to wells A1 through H2.
 - **384-well plate** Add 5 μ L of denatured standard into alternating wells of the plate:


Row 1: A1, C1, E1, ...K1, M1, O1

Row 2: Empty

Row 3: A3, C3, E3, ...K3, M3, O3

- If using 3130 genetic analyzer:
 - **96-well plate** Add 10 μ L of denatured standard to wells A1, B1, C1 and D1.
 - **384-well plate** Add 5 μ L of denatured standard into alternating wells of the plate:

Row 1: A1, C1, E1 and G1

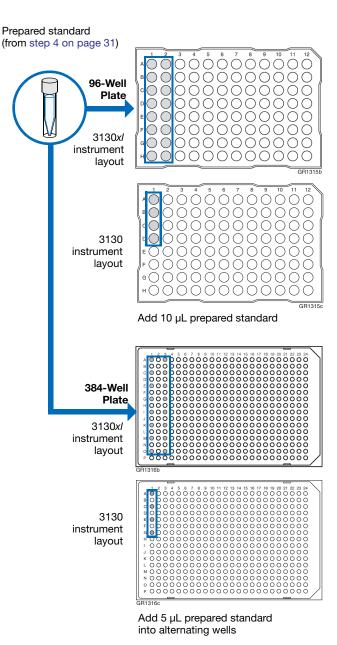





- **2.** Seal the plate:
 - **a.** Place the plate on a clean, level surface.
 - **b.** Lay the septa flat on the plate.
 - c. Align the holes in the septa strip with the wells of the plate, then firmly press downward onto the plate.

IMPORTANT!: Do not heat plates that are sealed with septa.

3. To prevent damage to the capillary array, inspect the plate and septa to verify the septa fits snugly and flush on the plate.

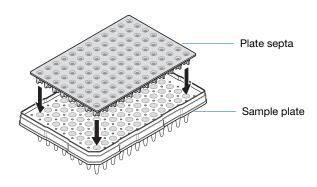

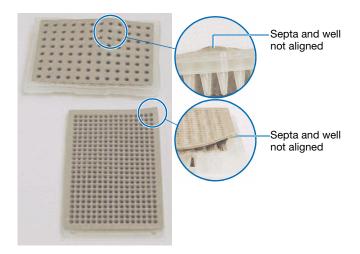


Prepared standard

- **4.** Briefly centrifuge the plate.
- **5.** Remove the plate from the centrifuge and verify that each sample is positioned correctly in the bottom of its well.

If the reagents of any well contain bubbles or are not located at the bottom of the well, repeat steps 4 and 5.


6. Leave the plate on ice until you are ready to prepare the plate assembly and place the assembly on the autosampler.

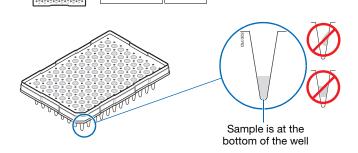

<1500 × g


Chapter 3 Performing a Spectral Calibration *Loading Samples*


- **7.** Assemble the plate assembly:
 - **a.** Place the sample plate into the plate base.
 - **b.** Snap the plate retainer onto the plate and plate base.

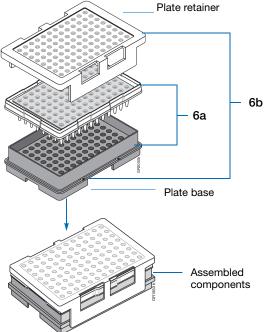

8. Verify that the holes of the plate retainer and the septa strip are aligned. If not, re-assemble the plate assembly (see step 7).

IMPORTANT! Damage to the array tips will occur if the plate retainer and septa strip holes do not align correctly.

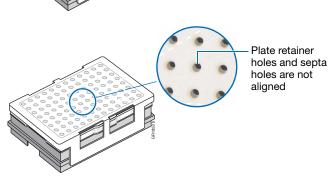


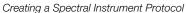





Creating a Spectral Instrument Protocol


In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or ga3130 > ➡ Protocol Manager to open the Protocol Manager window.

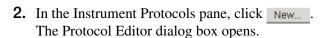

- **3.** Complete the Protocol Editor dialog box.
 - **a.** Type a name for the protocol.
 - **b.** Type a description for the protocol (optional).
 - c. Select Spectral in the Type drop-down list.
 - **d.** Select the correct dye set for your run. See "Dye Set Tables" on page 29.

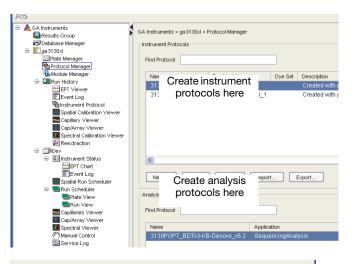

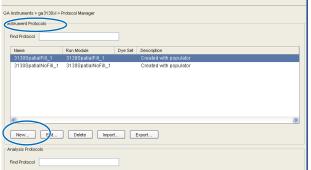

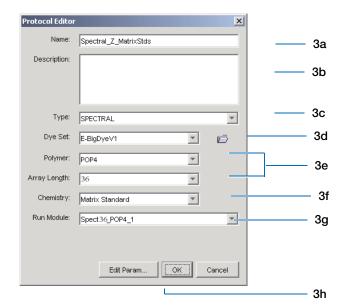





- e. Select the Polymer and Array Length from the appropriate drop-down list. See table, "Polymer, Array Length and Spectral Run Modules."
- f. Select your chemistry file in the Chemistry drop-down list. Failure to select the correct chemistry file for your spectral calibration samples results in a failing spectral run. See "Dye Set Tables" on page 29 for matching chemistry files.

Note: The chemistry file for fragment analysis dye sets defaults to the Matrix Standard.

g. In the Run Module drop-down list, select the run module. See the table "Polymer, Array Length and Spectral Run Modules."

Note: The modules list is filtered based on the polymer type, then the array length you selected in step e. You may have only one run module option available.

h. Click ok .

Polymer, Array Length and Spectral Run Modules


Polymer Type	Array Length (cm)	Run Module
POP-4	22	Spect22_POP4
	36	Spect36_POP4
		SpectSQ36_POP4
	50	Spect50_POP4
	80	Spect80_POP4
POP-6	36	Spect36_POP6
	50	Spect50_POP6
POP-7	36	Spect36_POP7
	50	Spect50_POP7
	80	Spect80_POP7


Performing a Spectral Calibration

Creating the Plate Record

- **1.** In the tree pane of the Data Collection software, click A GA Instruments > ga3130xl or $3130 > \square$ instrument name $> \square$ Plate Manager.
- **2.** Click New... to open the New Plate dialog box.
- Results Group ■ Database Manage N= 34 30√I Plate Manager Plate Manager Protocol Manag Module Manager Run History

- **3.** Complete the New Plate dialog box:
 - a. Enter a name for the plate.
 - **b.** Optional: Enter a description for the plate record.
 - **c.** In the Application drop-down list, select Spectral Calibration.
 - d. In the Plate Type drop-down list, select 96-Well or 384-Well
 - e. Enter a name for the owner.
 - f. Enter a name for the operator.
 - g. Click OK .
- 4. In the Spectral Calibration Plate Editor dialog
 - **a.** In the Sample Name column, enter a sample name, then click the next cell. The value 100 automatically displays in the Priority column.
 - b. Optional: In the Comments column, enter any additional comments or notations for the sample at the corresponding position of the plate.
 - c. In the **Instrument Protocol 1** column, select a protocol from the list or create a new protocol (see step 2 on page 35)
- **5.** Highlight the entire row.

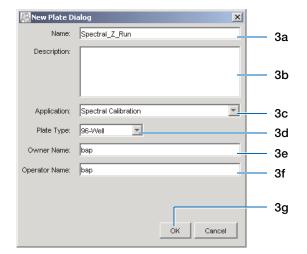

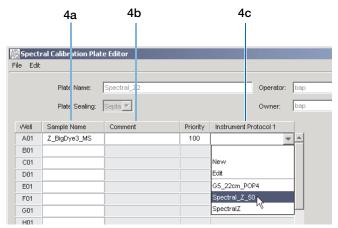




Chapter 3 Performing a Spectral Calibration Performing a Spectral Calibration

6. Select Edit > Fill Down Special.

Based on your plate type (96- or 384-well) and capillary array (16 or 4 capillaries), the software automatically fills in the appropriate well numbers for a single run.

7. Click OK .

You have successfully created the plate record for the spectral calibration plate.

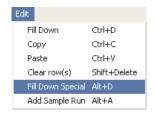





Placing the Plate Assembly into the Instrument

- **1.** Verify the oven and front doors are closed.
- **2.** Press the Tray button and wait for the autosampler to stop at the forward position.
- **3.** Open the front doors.
- **4.** Place the plate assembly on the autosampler in position A or B for the 3130*xl* genetic analyzer and position B for the 3130 genetic analyzer.

Note: There is only one orientation for the plate, with the notched end of the plate base away from you.

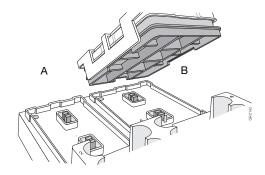

3130xl genetic analyzer

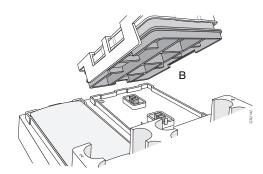


3130 genetic analyzer

- **5.** Ensure the plate assembly fits flat in the autosampler. Failure to do so may allow the capillary tips to lift the plate assembly off of the autosampler.
- **6.** Close the instrument doors.

Note: Closing the doors returns the autosampler to the home position, placing the tips of the capillaries in buffer.

Running the Spectral Calibration Plate

In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or 3130 > ■ instrument name > ■ Run Scheduler > ■ Plate View.

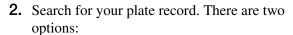


- Select Barcode in the Type of Search dropdown list.
 - In the Scan or Type Plate ID text box, type in the plate name and click Search to find the plate.
 - Or if you have a limited number of plates in the database, click Find All .

All plates in the database display in plate record section.


- Perform an advanced search by selecting Advanced in the Type of Search drop-down list.
 - Use the drop-down list to define search conditions for a category or multiple categories (Run Name, Results Group Name, Plate Name, etc.)

Note: Your entry for Plate Name will also be set as the Plate ID.

- For each category with a condition selected, type a value (primary search string) in the Value 1 column.
- Click search . All plates in the database that match the search criteria display in the plate record section.

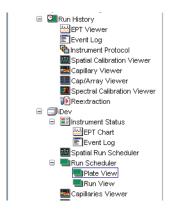



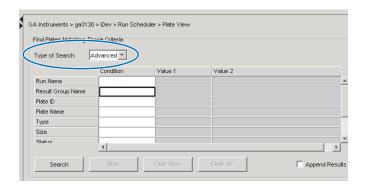




- **3.** Link the plate.
 - **a.** Select the plate record you want to run.
 - b. Click the plate position indicator that matches the plate you want linked.
 The plate map color will change from yellow to green when it is successfully linked.

Note: The 3130 genetic analyzer has only one plate position (Bay B) to link a plate record.

- **4.** In the toolbar of the Data Collection software window, click to begin the run.
- **5.** The Processing Plates dialog box opens. Click OK |

Note: The instrument may pause before running the plate to raise the oven temperature.

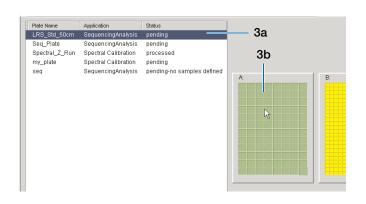



Spectral Calibration Approximate Run Times

Capillary Length (cm)	Run Type	Ramp Time (min)	Approximate Total Run Time (min)
22	Spec22_POP4	3	21
36	Spec36_POP4	10	35
	SpecSQ36_POP4	10	48
	Spec36_POP6	10	54
	Spec36_POP7	5	30
50	Spec50_POP4	7.5	110
	Spec50_POP6	10	95
	Spec50_POP7	7.5	52
80	Spec80_POP4	10	140
	Spec80_POP7	10	120

Notes____



Viewing the Pass/Fail Status After the Run

After the spectral calibration run, the pass or fail status of each capillary is recorded in the Event Log section of the Instrument Status window.

- **1.** In the tree pane of the Data Collection software, click ▲ GA Instruments > **Second ga3130xl** or ga3130 > instrument name > **■** Instrument Status > **■** Event Log.
- 2. In the Event Messages section of the window, view the status of each capillary.

Note: The window below shows the dye set G5 status results.



				Cap # Pass/fail status Q-value Condition Number
vent Messages				
Туре	Date	Time	Publisher	Description
w into	07707703	10.41.51	IDev	Finished Saving Spectral campiation data
Info	07/07/03	15:41:30	iDev	Saving spectral calibration data
🚳 Info	07/07/03	15:41:30	iDev	Capillary 16 successfully calibrated : q=0.988 c=9.12
🚳 Info	07/07/03	15:41:30	iDev	Capillary 15 successfully calibrated : q=0.986 c=9.15
Info	07/07/03	15:41:30		Run completed
Info	07/07/03	15:41:30	iDev	Capillary 14 successfully calibrated : q=0.986 c=9.01
Info	07/07/03	15:41:30	iDev	Capillary 13 successfully calibrated : q=0.988 c=8.99
Info	07/07/03	15:41:29	iDev	Capillary 12 successfully calibrated : q=0.989 c=8.87

Each capillary should have a Q-value above 0.95 and a Condition number range shown in the table.

Note: If the entire spectral calibration failed, see the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide for troubleshooting information.

Dye Set	Acceptable Condition Number Range	Q-Value
Sequencin	0.95	
Z_BigDyeV3	3 to 5	
E_BigDyeV1		
Fragment		
D	4 to 8.5	
F	6 to 12	
E5	2.5 to 4	
G5	8.5 to 14.5	
Any 4- or 5-Dye set	1 to 20	0.80




Evaluating the Spectral Calibration Data

IMPORTANT! Evaluate the spectral calibration profile for each capillary, even if the Spectral Calibration Results box indicates that they passed.

Note: Pages 49 to 50 for examples of passing sequencing spectral calibration profiles. See pages 51 to 52 for examples of passing fragment analysis spectral calibration profiles.

Evaluating the Spectral Profile and Raw Data

In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or ga3130 > ■ instrument name > ■ Spectral Viewer.

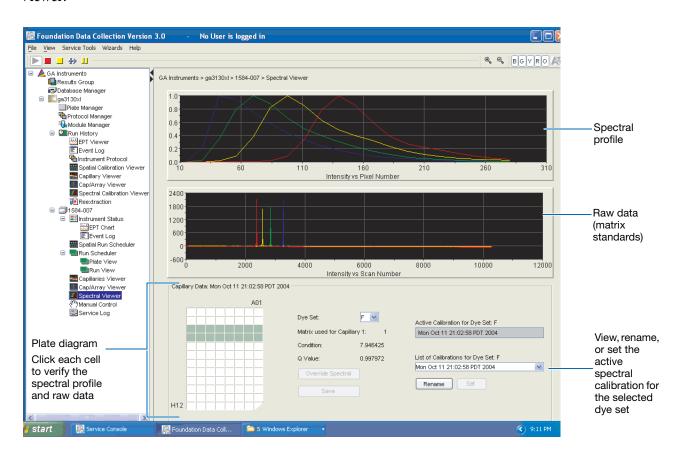





2. In the Dye Set drop-down list, select the dye set you just created.

Note: If the spectral calibration failed (no spectral profiles are created), see the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide* for more information.

3. In the plate diagram, select a well on the plate diagram to view the capillary spectral results.


Note: A failing capillary is automatically assigned the spectral profile of its nearest passing capillary.

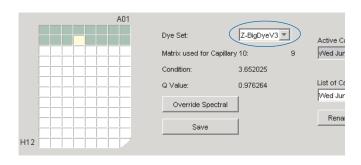

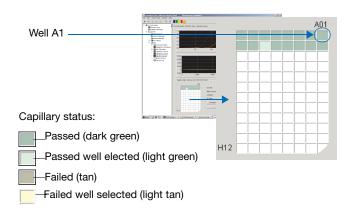


- **4.** Evaluate the spectral profile and raw data for the selected capillary:
 - **a.** Verify that the order of the peaks in the spectral profile from left to right are:
 - 4-dye: blue-green-yellow-red
 - 5-dye: blue-green-yellow-red-orange

Do the peaks in the profile appear in the correct order?	Then
Yes	Go to step b.
No	The calibration run has failed. See the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide.

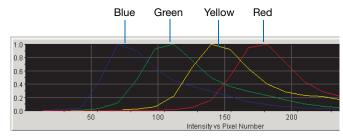

Example of a 4-dye spectral profile

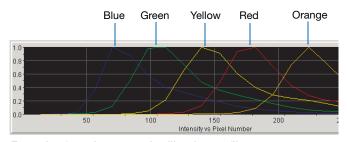

Example of a 5-dye spectral calibration profile



b. Verify that the order of the peaks in the raw data profile from left to right are:

Sequencing


- 4-dye: red-yellow-blue-green

Fragment Analysis


- 4-dye: red-yellow-green-blue

- 5-dye: orange-red-yellow-green-blue


Are peaks in the wrong order or extraneous peaks adversely affecting the spectral profile?	Then
Yes	The calibration run has failed. See the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide.
No	Go to step c.

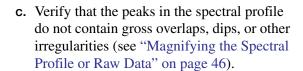

Are the peaks separate and distinct?	Then
Yes	The capillary has passed. Go to step 5.
No	The calibration run has failed. See the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide.

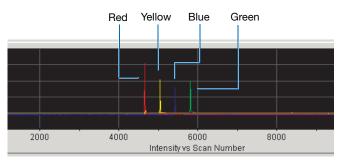


Example of a 4-dye sequencing raw data profile

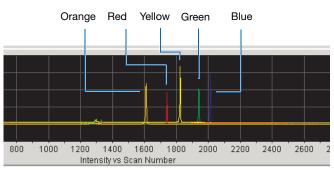


Example of a 5-dye fragment analysis raw data profile

Peaks are distinct, regular and in the proper order - pass

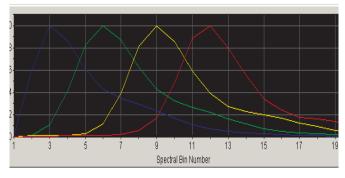


Red peak is not distinct, regular or in the proper order - fail

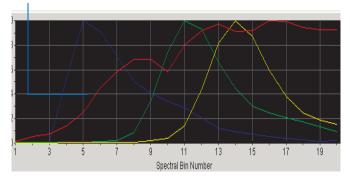



- **5.** Repeat steps 3 and 4 for each capillary in the array.
- **6.** Optional: Rename the spectral run. The spectral file default name is the day, date, and time of the run.
 - a. Click Rename
 - b. In the Rename Calibration dialog box, enter a descriptive name for the spectral calibration including the dye set, array length and polymer type (optional).
 - c. Click OK .



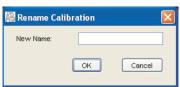

Magnifying the Spectral Profile or Raw Data

1. In the tree pane of the Data Collection software, click

▲ GA Instruments > ■ ga3130xI or ga3130 > ■ instrument name > ■ Spectral Viewer.


name > **||||| Spectral viewer**.

2. In the spectral profile or raw data display, click-drag the cursor to create a box around the area of interest.


3. Release the mouse button.

The Data Collection software displays the selected region.

4. Press **r** to reset the view.

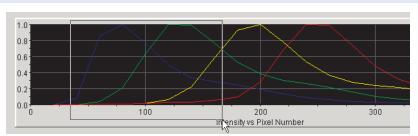

Selecting an area to magnify in a spectral profile

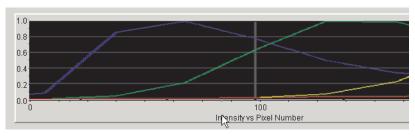

Magnified area of that spectral profile



Activating a Spectral Calibration

IMPORTANT! A run cannot start unless a calibration file that matches the dye set and capillary array length combination to be used for the run, is active.

IMPORTANT! While the software does not force you to recalibrate spectrally after changing polymer types, it is strongly recommended that you perform a spectral calibration after a polymer type change.

IMPORTANT! Whenever you install or replace an array or switch dye sets, you must either set the active spectral calibration for that dye set and array length combination or perform a new spectral calibration and then activate it.

IMPORTANT! Every new spectral calibration is automatically the active one for that dye set.

You may activate a previously created spectral calibration for a run as long as the calibration matches the dye set and array length combination (and, for fragment analysis, the polymer type) that you wish to use. Use an active spectral calibration for:

- Sequencing analysis applications that require a separate spectral calibration (for the same dye set) for different capillary array lengths and polymer type
- Fragment analysis applications that require a separate spectral calibration (for the same dye set) for different capillary array lengths and polymer type
- Repeat spectral calibrations where the original calibration is better than the second one

Setting an Active Spectral Calibration

1. In the tree pane of the Data Collection software, click

▲ GA Instruments > \blacksquare ga3130xl or ga3130 > \blacksquare instrument name > \blacksquare Spectral Viewer.


IMPORTANT! If the Spectral Viewer window is blank and deactivated, then either: —The spectral calibration for that dye set is not in the database or,

- -You changed the array length and you do not have a spectral calibration file activated for that dye set and array length combination.
- 2. In the Dye Set drop-down list, select a dye set.


Notes		



Chapter 3 Performing a Spectral Calibration *Activating a Spectral Calibration*

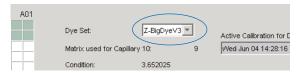

3. In the List of Calibrations for Dye Set drop-down list, select the spectral calibration you want to use. The spectral profile and raw data is displayed.

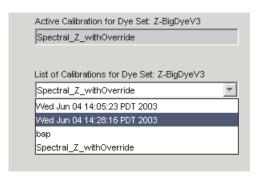


4. If the spectral calibration is acceptable, then click set . Otherwise, run a new spectral calibration.




Examples of Passing Sequencing Spectral Calibrations

Dye Set Z Created from Matrix Standard

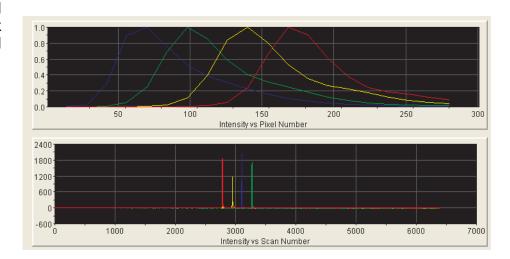


Dye Set Z Created from a Sequencing Standard

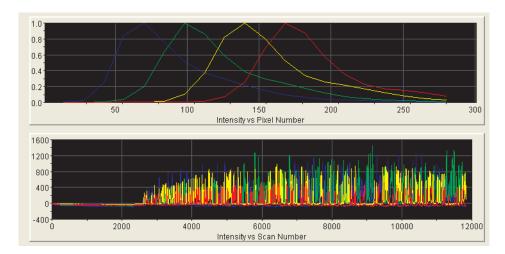




Dye Set E Created from Matrix Standard Set **DS-01**

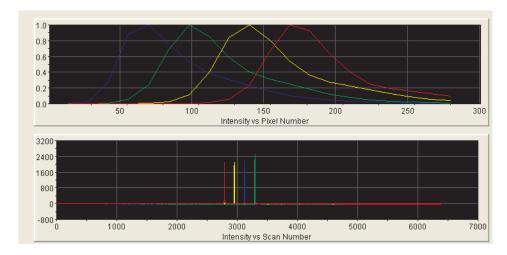


Dye Set E Created from a Sequencing Standard

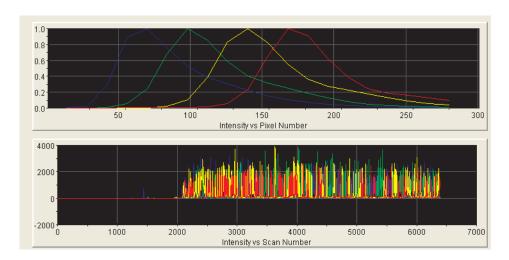




Examples of Passing Fragment Analysis Spectral Calibrations

Dye Set G5 Created from Matrix Standard Set DS-33

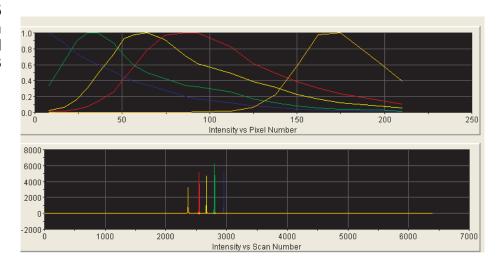


Dye Set F Created from Matrix Standard Set DS-32

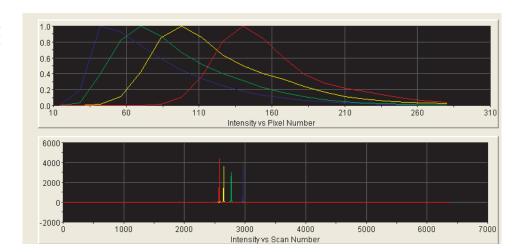




Dye Set D Created from Matrix Standard Set DS-30

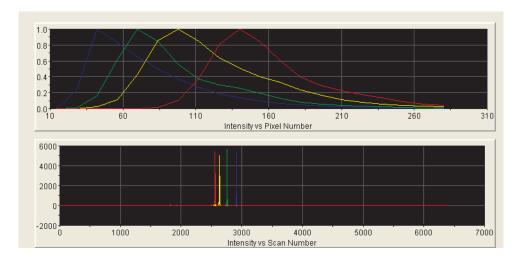


Dye Set E5 Created from Matrix Standard Set DS-02

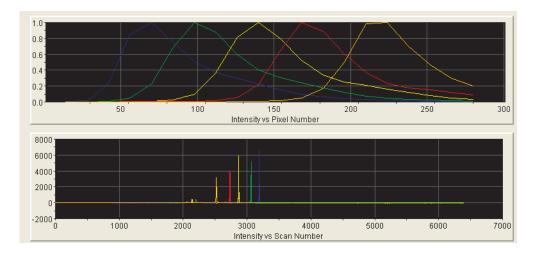



Autoanalysis and Sequencing Analysis Software

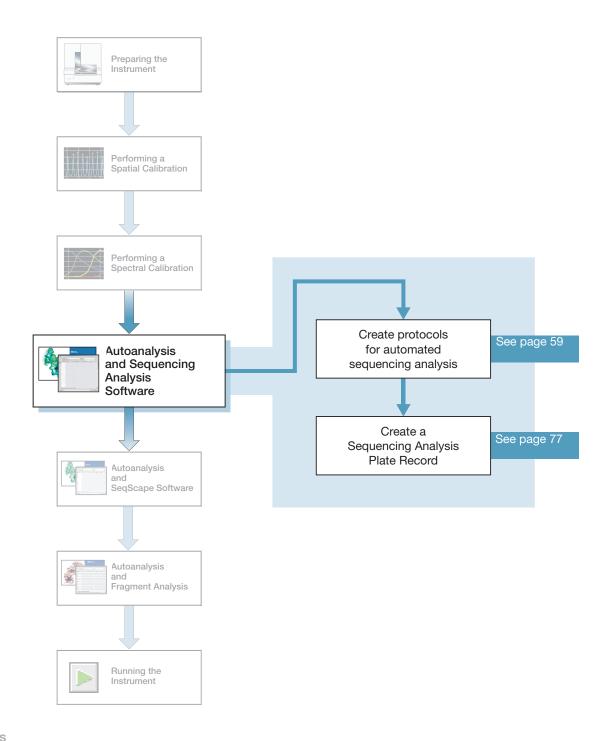

Sequencing Analysis Software

See the *Applied Biosystems DNA Sequencing Analysis Software v5.1 User Guide* (P/N 4346366) for details on sequencing analysis or the *Sequencing Analysis Software v5.2 User Bulletin* titled *New Features* (P/N 4358355) for an update on Sequencing Analysis Software v5.2.

Analyze your sequencing samples automatically using autoanalysis or manually.

Autoanalysis

Perform autoanalysis of sequencing samples using features of the 3130/3130xl genetic analyzer Data Collection software and Sequencing Analysis Software v5.2 or higher.

The Sequencing Analysis Software must be installed and registered with the 3130/3130xl genetic analyzer Data Collection software before you can create files required for autoanalysis.

Autoanalysis can only be performed on the same computer that collected the sample files. If you perform autoanalysis on samples, but wish to edit/review results on another computer, transfer the analysis protocol to the Sequencing Analysis Software database. To analyze samples on another computer, transfer the files to that location.

Manual Analysis

If the run is not set up for autoanalysis, see the *Applied Biosystems DNA Sequencing Analysis Software v5.1 User Guide* and the *Sequencing Analysis Software v5.2 User Bulletin - New Features*.

File-Naming Convention


Some alphanumeric characters are not valid for user names or file names. Do not use the characters below:

spaces

\/:*?"<>|

IMPORTANT! An error message is displayed if you use any of these characters. You must remove the invalid character to continue.

Notes		

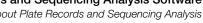



About Plate Records and Sequencing Analysis

Overview

Plate records are data tables in the instrument database that contain the following information:

- Plate name, type, and owner
- Position of the sample on the plate (well number)
- Sample Name
- Mobility file (in Analysis Protocol)
- Comments about the plate and about individual samples
- Name of the run module and Dye set information (run modules specify the conditions to run the samples)
- Name of the Analysis Protocol

Note: A plate record is similar to a sample sheet or an injection list that you may have used with other ABI PRISM® instruments.

When to Create a Plate Record

Create a plate record for each plate of samples for the following types of runs:

- Spectral calibrations
- · Sequencing Analysis Software
- SeqScape Software analysis
- GeneMapper® software
- Mixed (sequencing and fragment analysis samples see the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide for creating a mixed plate record)

For runs to begin, you must create a plate record for and link it to a plate loaded on the instrument. However, you can create plate records for new plates while a run is in progress.

About a Sequencing **Analysis Plate** Record

The Plate Editor displays an empty plate record for the application selected in the New Plate dialog box. The data fields within a given plate record vary depending on the selected application. This section describes the sequencing analysis plate record's data fields.


The table below describes required plate editor fields:

Parameters	Description	See Page
Instrument Protocol	Contains the run module and dye set needed to run the instrument.	59
Analysis Protocol	Contains everything needed to analyze sequencing data.	62
Results Group	Defines the file type, the file name, file save locations, analysis software and autoanalysis.	70



About Plate Records and Sequencing Analysis

Elements of a Sequencing Analysis Plate Record

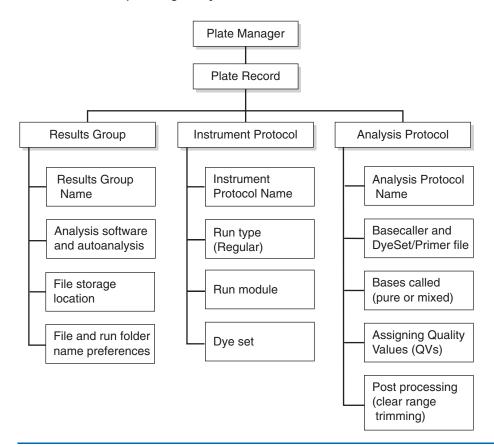


IMPORTANT! For data collection and autoanalysis to succeed, each run of samples must have an Instrument Protocol, an Analysis Protocol, and a Results Group assigned within a plate record.

Notes_____



A Blank Sequencing Analysis Plate Record

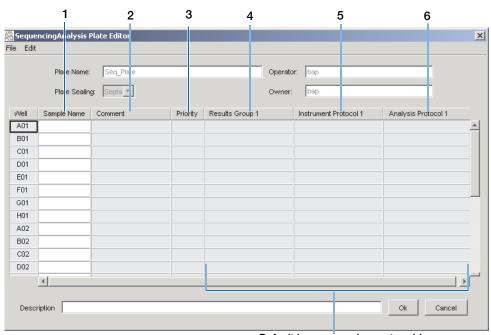

Default is one sample run, to add additional runs see page 79

Columns inserted in a plate record for a sequencing analysis run

Number and Column	Description
1. Sample Name	Name of the sample
2. Comment	Comments about the sample (optional)
3. Priority	A default value of 100 to each sample. Changing the value to a smaller number increases the priority of that set of 16 or 4 samples to run before the others in the injection list.
4. Results Group	Some options:
	New: Opens the Results Group Editor dialog box
	 Edit: Opens the Results Group Editor dialog box for the Results Group listed in the cell
	None: Sets the cell to have no selected Results Group
	Select one of the available Results groups from the list
	Note: You must have a Results Group selected for each sample entered in the Sample Name column.
	See "Results Group for Sequencing Analysis" on page 70.

About Plate Records and Sequencing Analysis

Number and Column	Description
5. Instrument	New: Opens the Protocol Editor dialog box.
Protocol	Edit: Opens the Protocol Editor dialog box for the Instrument Protocol listed in the cell.
	None: Sets the cell to have no selected protocol.
	List of Instrument Protocols: In alpha-numeric order.
	Note: You must have an Instrument Protocol selected for each sample entered in the Sample Name column.
	See "Instrument Protocol for Sequencing Analysis" on page 59.
6. Analysis Protocol	New: Opens the Analysis Protocol Editor dialog box.
	Edit: Opens the Analysis Protocol Editor dialog box for the Instrument Protocol listed in the cell.
	None: Sets the cell to have no selected protocol.
	List of Analysis Protocols: In alpha-numeric order
	Note: You must have an Analysis Protocol selected for each sample entered in the Sample Name column.
	See "Analysis Protocol for Sequencing Analysis" on page 62.

	0		



Creating Protocols for Automated Sequencing Analysis

If the appropriate instrument protocol, analysis protocol, and results group have been created, proceed to "Filling Out a Sequencing Analysis Plate Record" on page 77.


Instrument Protocol for Sequencing Analysis

An instrument protocol contains all the settings necessary to run the instrument: the protocol name, type of run, run module, and dye set.

WARNING Do not edit the instrument protocol currently being used, while the instrument is running.

Creating an Instrument Protocol

In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or ga3130 > ♠ Protocol Manager.

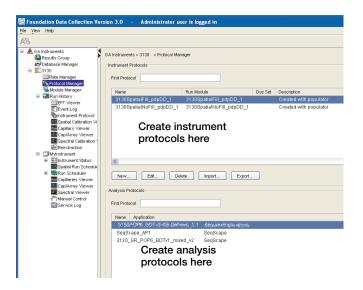

Chapter 4 Autoanalysis and Sequencing Analysis Software




Creating Protocols for Automated Sequencing Analysis


2. In the Instrument Protocol section, click New...


The Protocol Editor opens.

- **3.** Complete the Protocol Editor:
 - **a.** Type a name for the protocol.
 - **b.** Type a description for the protocol (optional).
 - **c.** Select **Regular** in the Type drop-down list.

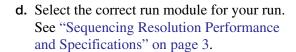

Note: To customize a run module, see the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide.

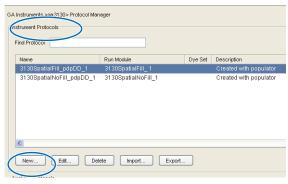

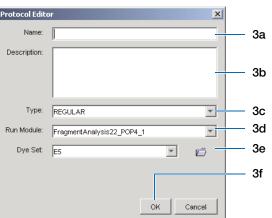



e. Select the correct Dye Set for your run by using the table below.

			POP-Polym	-		P-6 ymer		POP	-7 Pol	ymer	
Chemistry	Dye Set	UltraSeq36	StdSeq50	LongSeq80	RapidSeq36	StdSeq50	UltraSeq36	RapidSeq36	FastSeq50	StdSeq50	LongSeq80
BigDye® Terminator v3.1 Cycle Sequencing Kit	Z_BigDye	1	1	1	1	1	1	1	1	1	1
ABI PRISM® dGTP BigDye® Terminator v3.0 Cycle Sequencing Ready Reaction Kit	V3	_	_	_	_	_	_	-	-	-	_
BigDye® Terminator v1.1 Cycle Sequencing Kit	E_BigDye	1	1	1	1	1	1	1	1	1	1
ABI PRISM® dGTP BigDye® Terminator Cycle Sequencing Kit*	V1	_	_	_	_	_	_	-	-	-	_
ABI PRISM® dRhodamine Dye Terminator Cycle Sequencing Ready Reaction Kit		1	1	1	1	1	_	_	_	_	_
ABI PRISM® BigDye® Primer Cycle Sequencing Kits		_	_	_	1	1	_	_	_	_	_

^{*}dGTP kits are not supported on capillary electrophoresis instruments due to compressions on certain sequence context regions; you can run the kits if you do not care about the compression issues.

f. Click ok .

Importing an Instrument Protocol

- 1. Click import in the Instrument Protocols pane of the Protocol Editor window to display the standard File Import dialog box.
- **2.** Navigate to the location of the .xml file you want to import.
- **3.** Select the .xml file and click **Open**.

Exporting an Instrument Protocol

- **1.** In the Instrument Protocols pane, highlight the protocol you want to export.
- **2.** Click Export to display the standard File Export dialog box.
- **3.** Browse to the desired folder location.

4.	Click	Save	
••	CHCK	Save	

-Instrument Proto	cols					
Find						
Name		Run Module			Dye Set	
	1				1	
New	Edit	Delete	Import	Export		

Analysis Protocol for Sequencing Analysis

An analysis protocol contains all the settings necessary for analysis and post processing:

- Protocol name The name, description of the analysis protocol, and the sequence file formats to be used
- Basecalling settings The basecaller,
 DyeSet/Primer file, and analysis stop point to be used
- Mixed Bases (Optional) When two bases are found at the same position. Define the percent value of the second highest to the highest peak
- Clear Range The high quality sequence remaining after trimming low quality sequences typically found at the beginning and ends of the sequence. The clear range can be based on base positions, sample quality values, and/or number of ambiguities (Ns) present

Note: If you created an appropriate analysis protocol in the Sequencing Analysis software, you can use it in Data Collection software.

IMPORTANT! Do not delete an analysis protocol during a run while it is being used for that run. Autoanalysis will not be performed if you do so.


Creating an Analysis Protocol

See the *Applied Biosystems DNA Sequencing Analysis Software v5.2 User Guide* (P/N 4346366) for more information on analysis protocols.

1. In the Analysis Protocol section of the Protocol Manager, click New...


If more than one analysis application is installed on the data collection computer, the Analysis Applications dialog box opens.

Note: When you name your analysis protocol, assign a version name to help you distinguish your analysis protocols.

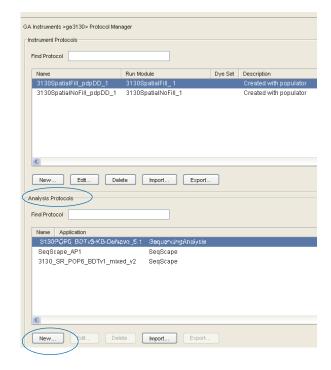




2. Select Sequencing Analysis, then click OK to open the Analysis Protocol Editor dialog box.

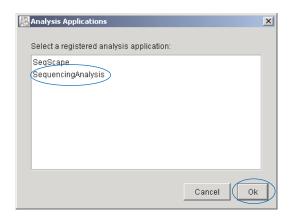

3. In the **General** tab:

- **a.** Enter a unique name and description for the new protocol.
- **b.** Select the appropriate Sequence File formats settings.

Option	Creates
Write .Seq File check box	A .seq file for printing the sequence as text file or for use in other software.
	 Use ABI format with Applied Biosystems software. Use FASTA format with other software.
Write Standard Chromatogram Format file (.scf)	A .scf file to use with other software. When created, the .scf extension is not appended to the file name.
Write Phred (.phd.1) File	A .phd.1 file to use with other software if you used the KB [™] basecaller.

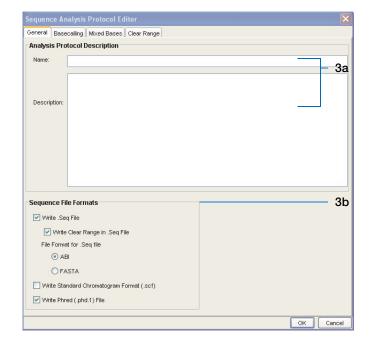




Analysis Protocol for Sequencing Analysis

4. Select the **Basecalling** tab.

a. See "3130/3130xl Genetic Analyzer Basecaller and DyeSet/Primer Files" on page 66 to select your basecaller and DyeSet primer file.

Note: Sequencing Analysis Software and 3130/3130*xl* genetic analyzer Data Collection software filter .mob file choices to match the chosen .bcp file.

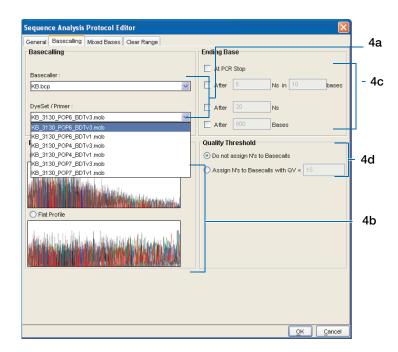

b. In the Processed data pane, select True or Flat Profile.

Option	Used to display data as processed traces scaled
True Profile	Uniformly - the average height of peaks in the region of strongest signal is about equal to a fixed value. The profile of the processed traces will be very similar to that of the raw traces.
Flat Profile	Semi-locally - the average height of peaks in any region is about equal to a fixed value. The profile of the processed traces will be flat on an intermediate scale (> about 40 bases).
	Note: This option is applied to data that is analyzed with the KB [™] basecaller only. If you use the ABI basecaller, the profile option reverts to True Profile.

- **c.** If desired, select one or more stop points for data analysis.
- d. Select your Threshold Quality option.

Option	Function
© Call all bases and assign QV	With the KB [™] basecaller, use this setting to assign a base to every position, as well as the QV.
Assign 'N' for bases with QV < 15	With the KB basecaller, use this setting to assign Ns to base with QVs less than the set point. The QV is still displayed.

Notes_____




5. Select the **Mixed Bases** tab.

Note: This function is active with the KB Basecaller only.

- a. For mixed bases only, select **Use Mixed Base Identification**.
- b. Use the default setting of 25% or change the detection level by entering a new value or dragging the % line up or down.

Note: Do not use less than 15% as your detection limit.

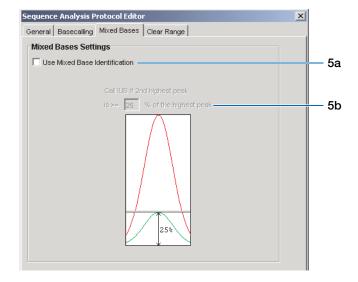


6. Select the Clear Range tab.

Note: The clear range is the region of sequence that remains after excluding the low-quality or error prone sequence at both the 5′ and 3′ ends.

Select one or more Clear Range methods. If you apply multiple methods, the smallest clear range results.

7. Click ok to save the protocol and close the Sequence Analysis Protocol Editor dialog box.

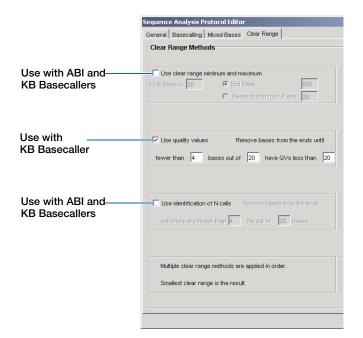

3130/3130xl Genetic Analyzer Basecaller and DyeSet/Primer Files

Basecaller and DyeSet/Primer Files Used with BigDye® Terminator Chemistry and KB Basecalling

BigDye® Terminator v1.1 Cycle Sequencing Kit POP-4™ UltraSeq36_POP4 StdSeq50_POP4 LongSeq80_POP4 KB_3130_POP4_BDTv1.mob	KB.bcp	
LongSeq80_POP4		
<u> </u>		
The state of the s		
POP-6 [™] RapidSeq36_POP6 KB_3130_POP6_BDTv1.mob		
StdSeq50_POP6		
POP-7 [™] UltraSeq36_POP7 KB_3130_POP7_BDTv1.mob		
RapidSeq36_POP7		
FastSeq50_POP7		
StdSeq50_POP7		
LongSeq80_POP7		
BigDye® Terminator v3.1 POP-4 UltraSeq36_POP4 KB_3130_POP4_BDTv3mob		
Cycle Sequencing Kit StdSeq50_POP4		
LongSeq80_POP4		
POP-6 RapidSeq36_POP6 KB_3130_POP6_BDTv3.mob		
StdSeq50_POP6		
POP-7 UltraSeq36_POP7 KB_3130_POP7_BDTv3.mob		
RapidSeq36_POP7		
FastSeq50_POP7		
StdSeq50_POP7		
LongSeq80_POP7		

Basecaller and DyeSet/Primer Files Used with BigDye® Terminator Chemistry and ABI Basecalling

DNA Sequencing Chemistry	Polymer	ABI Basecalling Run Module	Basecaller	DyeSet/Primer
BigDye [®] Terminator v1.1	POP-4 [™]	UltraSeq36_POP4	Basecaller- 3130POP4UR.bcp	DT3130POP4LR{BD}v1. mob
Kit		LongSeq80_POP4	Basecaller- 3130POP4_80cmv3.bcp	
	POP-6 [™]	RapidSeq36_POP6	Basecaller- 3130POP6RRv2.bcp	DT3130POP6{BD}v2.mob
		StdSeq50_POP6	Basecaller- 3130POP6SR.bcp	

Notes			



Basecaller and DyeSet/Primer Files Used with BigDye® Terminator Chemistry and ABI Basecalling (continued)

BigDye Terminator v3.1	POP-4	UltraSeq36_POP4	Basecaller- 3130POP4UR.bcp	DT3130POP4{BDv3}v1.m ob	
Cycle Sequencing Kit		LongSeq80_POP4	Basecaller- 3130POP4_80cmv3.bcp		
	POP-6	RapidSeq36_POP6	Basecaller- 3130POP6RRv2.bcp	DT3130POP6{BDv3}v1.m ob	
		StdSeq50_POP6	Basecaller- 3130POP6SRv2.bcp		
ABI PRISM® dRhodamine	mine minator LongSeq80_POP4 cing	UltraSeq36_POP4	Basecaller- 3130APOP4UR.bcp	DT3130POP4{dRhod}v2. mob	
Dye Terminator Cycle Sequencing		LongSeq80_POP4	Basecaller- 3130POP4_80cmv3.bcp		
Ready Reaction Kit		Basecaller- 3130POP6RRv2.bcp	DT3130POP6{dRhod}v2. mob		
		StdSeq50_POP6	Basecaller- 3130POP6SR.bcp		

Dye Primer Chemistry

Basecaller and DyeSet/Primer Files Used for Dye Primer Chemistry

DNA Sequencing Chemistry	Polymer	ABI Basecalling Run Module	Basecaller	DyeSet/Primer
ABI PRISM® BigDye® Primer Cycle Sequencing Kit	POP-6	RapidSeq36_POP6	Basecaller- 3130POP6RRv2.bcp	DP3130POP6{BD -21M13}v1.mob
		StdSad		Basecaller- 3130POP6SR.bcp

Notes		



Editing and Deleting Analysis Protocols


Editing an Analysis Protocol


- 1. In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to edit.
- 2. Click Edit....
- **3.** Make changes in the General, Basecalling, Mixed Bases and Clear Range tabs, if appropriate.
- 4. Click to save the protocol and close the Analysis Protocol Editor dialog box.

Deleting an Analysis Protocol

IMPORTANT! Do not delete an analysis protocol during a run while it is being used for that run. Autoanalysis will not be performed if you do so. You must first delete any plate records using the analysis protocol before you can delete or modify the analysis protocol for these plate records.

- 1. In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to delete.
- **2.** Click Delete to display the Deletion Confirmation dialog box.
- 3. Click Yes .

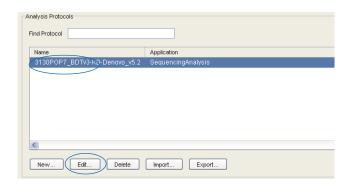

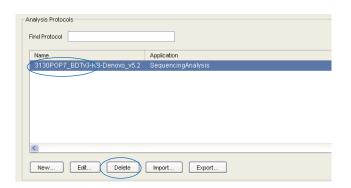




Exporting an Analysis Protocol


- **1.** In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to export.
- **2.** Click Export to display the standard File Export dialog box.
- **3.** Browse to the desired folder location.
- 4. Click Save.

Importing an Analysis Protocol

- 1. Click mort to display the standard File Import dialog box.
- **2.** Browse for the .xml file to import and click **Open**.

Note: For multiple applications, select the appropriate application to associate with the analysis protocol.

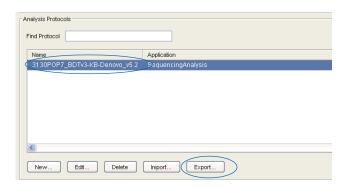

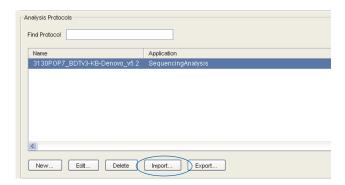



Results Group for Sequencing Analysis

A Results Group is a component within Data Collection that organizes samples and certain user settings under a single name. It is called a Results Group because it is used to name, sort, and deliver samples that result from a run.

Creating a Results Group

- **1.** In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ Results Group.
- 2. Click New... to display the Results Group Editor window.




GA Instruments > Results Group


Default_Results_Group GM_Results_Group MJD_Results_Group

Find Results Group

Name

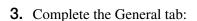

- a. Type a unique Results Group Name. The name can be used in grouping sample files.
- **b.** Type a Results Group Owner (optional). The owner name can be used in naming and sorting sample files.
- **c.** Type a Results Group Comment (optional).

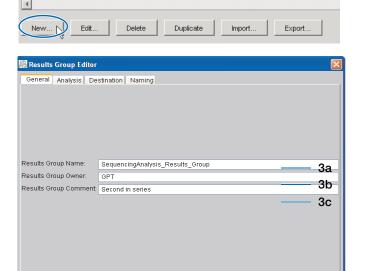

OK Cancel

Owner Comment



- **4.** Select the **Analysis** tab, then:
 - a. Select Sequencing Analysis from the Analysis Type drop-down list.
 - b. In the Analysis Actions section, select **Do** Autoanalysis, if you want your data automatically analyzed after a run.


Note: Login ID and password are not required for Sequencing Analysis software.


5. Select the **Destination** tab, then use the default destination or define a new location for data storage.

To use	Then
default location*	skip to step 6.
custom location	complete steps a-b.

- a. Click Use Custom Location, then click Browse... to navigate to a different save location.
- **b.** Click Test to test the Location path name connection:

If it	Then a message box displays
Passes	Test succeeded: <"path">.
Fails	Test failed:<"path">.

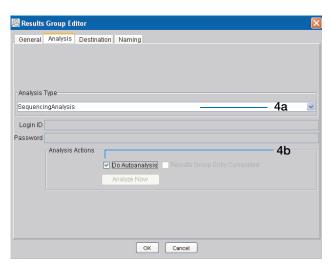

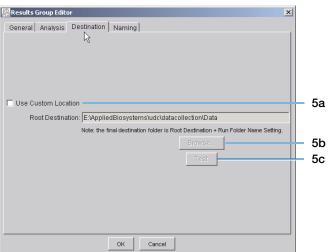

*Sample File Destinations

Locations where sample files are placed during extraction:

- Default Destination, default folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\<instrument type>\ <instrument name>\run folder
- Default Destination, custom folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\top custom folder\ subfolders, etc.
- Custom Destination, default folder naming: Destination\<instrument type>\<instrument name>\run folder
- Custom Destination, custom folder naming: Destination\top customer folder\subfolders, etc.

Chapter 4 Autoanalysis and Sequencing Analysis Software




Results Group for Sequencing Analysis

6. Select the Naming tab.

Use the Naming tab to customize sample file and run folder names. You have two options:

- Use the default names for the folder and sample file.
- Use custom names for the folder and sample file. See "Optional: Completing Sample File Name Format Pane" on page 73 for elements of the Naming tab.

Note: Sample name, run folder name, and path name, *combined*, can total no more than 250 characters.

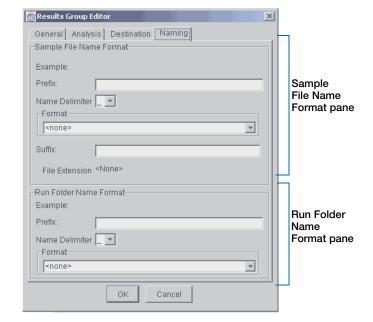

Sample File Name Format

MJD_ .ab1

Example:




Optional: Completing Sample File Name Format Pane


Note: The default sample file naming for sequencing analysis is:

seq_<capillary number>_<wellposition>.ab1


1. Select the Naming tab.

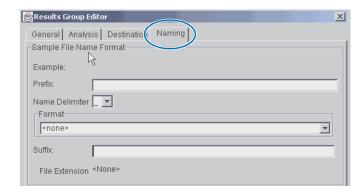


- **2.** Type a prefix for the file name in the Prefix box. Your entry is shown in the Example line.
- Select the symbol that separates the Format elements in the file name from the Name Delimiter drop-down list. Only one delimiter symbol may be chosen.

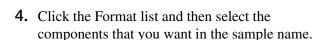

Note: All the samples from a single run can be placed in the same run or results folder, so the name of every sample from a single run should be different. Most of the Format options are not different between samples, so select at least one of the options that make the sample names unique within a run.

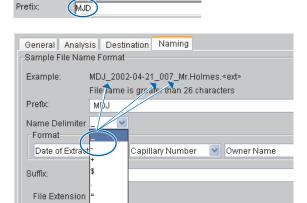

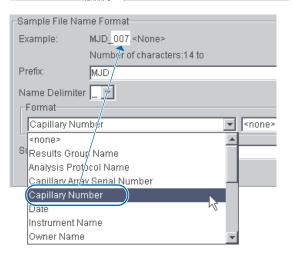



Results Group for Sequencing Analysis


If a unique identifier is not included in the name, a warning message displays. Select from the elements to make a unique sample file name. Examples are:

- Well position
- Capillary number
- Run sequence number
- Plate Quadrant

Note: some identifiers, such as well position, capillary number and plate quadrant are not unique in multiple sample instances. Plate Quadrant is not a unique identifier for a 96-well plate.


As you select the elements for the file name, they are placed in the Example line. An additional element drop-down menu displays allowing you the option of selecting an additional element.

	Warning messa	age
Sample File Nar	ne Format	
Example:	BasecallerProtocol.saz.ab1	
	INVALID NAME: Filename does not have a unique identifier	in it.
Prefix:	\(\frac{1}{2}\)	
Name Delimiter		
Analysis Prote	ocol Name	
Suffix:		
File Extension	ab1	

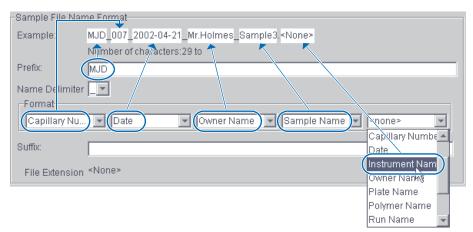

The names of the Format elements eventually truncate, but the Example field remains visible (up to 72 characters).

Note: Resize the Results Group window horizontally to view the element options.

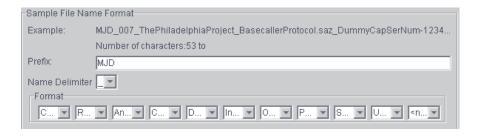



5. Type the suffix for the file name in the Suffix box.

Note: The File Extension field displays the file extension generated from the Analysis Type specified on the Analysis tab (see page 71). For example, Sequencing Analysis produces sample files with an .ab1 extension.

Optional: Run Folder/Sub-Folder Name Format Pane

Note: The default run folder naming for sequencing analysis is Run_<Instrumentname>_<Dateand TimeofRun>_<RunSeq#>.

Follow the same steps described for the Sample File Name Format pane (see page 73) to change the subfolder name within the run folder. Make the subfolder name unique, such as using the Run Folder name.

Saving a Results Group

Click OK from any tab once all the elements within the Results Group have been chosen.


Note: Even if you create a custom run folder location, the system generates a separate default run folder that contains the log file.

Importing and Exporting a Results Group

Results Groups can be imported from, or exported to, XML files allowing sharing of identical Results Groups between instruments.

Importing a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- 2. Click mport to display the standard File Import dialog box.

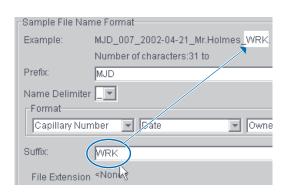

Chapter 4 Autoanalysis and Sequencing Analysis Software



Results Group for Sequencing Analysis

3. Navigate to the file you want to import.

Note: Import file type is .xml (XML file).

4. Click Open .

Exporting a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- **2.** Click the Results Group name to select it.
- 3. Click Export .

A standard file export dialog box displays with the chosen Results Group name.

- **4.** Navigate to the location where you want to save the exported file.
- 5. Click Save

Note: If there is a name conflict with a Results Group that already exists at the save location, then duplicate the Results group. Duplication copies the settings into a similar Results Group without the risk of user error when copying it manually (see procedure below).

Duplicating a Results Group

- **1.** Click the Results Group name.
- 2. Click Duplicate .

Note: When you duplicate a Results Group, you are asked to type a name for the new Results Group and for the analysis application type.


Notes		



Filling Out a Sequencing Analysis Plate Record

Creating a Sequencing Analysis Plate Record

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > ga3130xl or ga3130 > Plate Manager.
- **2.** Click New... to display the New Plate Dialog dialog box.
- **3.** Complete the information in the New Plate Dialog:
 - **a.** Type a name for the plate.
 - **b.** Type a description for the plate (optional).
 - **c.** Select your sequencing application in the Application drop-down list.
 - d. Select **96-well** or **384-well** in the Plate Type drop-down list.
 - **e.** Type a name for the owner and operator.
 - f. Click open the Sequencing Analysis Plate Editor.

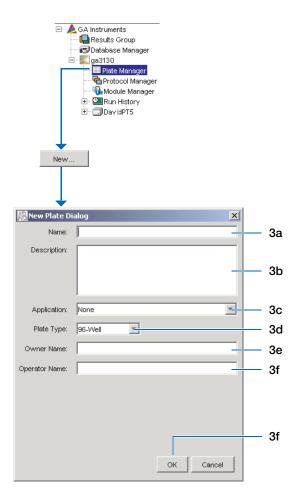






Filling Out a Sequencing Analysis Plate Record

Completing a Sequencing Analysis Plate Record

- 1. In the **Sample Name** column of a row, enter a sample name, then click the next cell. The value 100 automatically displays in the Priority column.
- **2.** In the **Comments** column, enter any comments or notations for the sample.
- **3.** In the **Priority** column, change the priority value, if desired. A lower number for each 4 or 16 set, receives a higher run priority.
- **4.** In the **Results Group 1** column, select a group from the drop-down list (see page 70) or create a new Results Group.
- **5.** In the **Instrument Protocol 1** column, select a protocol from the drop-down list (see page 59) or create a new instrument protocol.
- **6.** In the **Analysis Protocol 1** column, select a protocol from the drop-down list (see page 62) or create a new analysis protocol.
- **7.** Complete the plate record based on the samples loaded in your plate:
 - For the single runs that use the same samples and protocols Highlight the entire row, then select **Edit** > **Fill Down Special**. Based on your plate type (96- or 384-well) and capillary array (16 or 4 capillaries), the software automatically fills in the appropriate well numbers for a single run or
 - For entire plates that use the same samples and protocols – Highlight the entire row, then select Edit > Fill Down to fill down the entire plate editor.
 - For plates with different samples and protocols complete the entries manually.

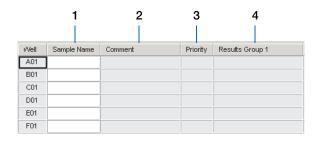

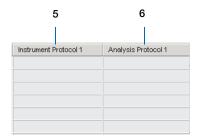

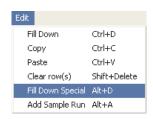


	ſΊ		

8. To do more than one injection of the same sample, select **Edit > Add Sample Run**.

Results Group, Instrument Protocol and Analysis Protocol columns are added to the right end of the plate record.

- **9.** Complete the columns for the additional runs.
- **10.** Click OK .


Note: After clicking OK in the Plate Editor, the completed plate record is stored in the Plate Manager database. The plate record can be searched for, edited, duplicated, exported, or deleted in the Plate Manager.

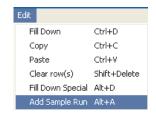

Chapter 4 Autoanalysis and Sequencing Analysis Software Filling Out a Sequencing Analysis Plate Record



Autoanalysis and SeqScape Software

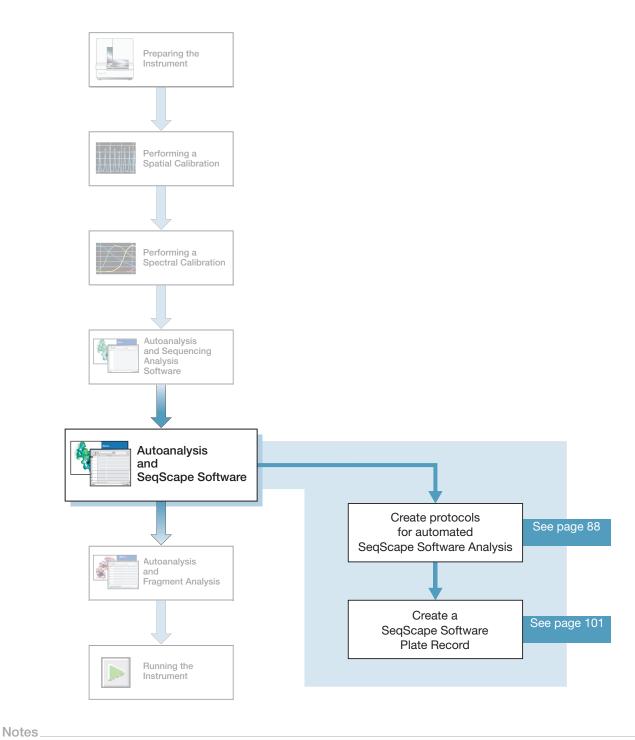

SeqScape Software

Analyzing your sequencing samples may be done by autoanalysis or manual analysis.

Autoanalysis

Use the ABI PRISM® SeqScape® software to automatically analyze Applied Biosystems 3130/3130xl Genetic Analyzers sequencing data. Perform autoanalysis only on the same instrument computer that collected the sample files. You can configure the software packages to perform data collection and then data analysis without requiring user interaction.

Install and register the SeqScape software with the 3130/3130xl genetic analyzer Data Collection software before creating required files for autoanalysis.

See Appendix B in the *SeqScape Software User Guide v2.5* (P/N 4359442), for more details on setting up autoanalysis.

Required Software Packages

Autoanalysis requires three software packages:

• 3130/3130xl genetic analyzer Data Collection software

The Data Collection software runs the instrument and collects fluorescent data from samples. For autoanalysis to occur, the software must communicate with downstream software.

Analysis parameters used for data collection as well as that created in SeqScape software are accessible in Data Collection software.

Autoanalysis Manager

The Autoanalysis Manager software is part of the integration between the data collection, SeqScape, and GeneMapper[®] software. It can queue messages and track their processing status. Each message is considered a batch job, whether it contains a single sample, samples from a result group, or an entire run of samples.

Autoanalysis Manager is installed by Seqscape or GeneMapper software when loaded on a system with Data Collection software.

• A version of SeqScape software without a user interface

This SeqScape software version is identical to the regular software version except that there is no user interface. The Autoanalysis Manager opens and uses this version of software to analyze the data in the projects.

The automated processing version and the standard version of SeqScape software are installed from the SeqScape Software installation CD.

IMPORTANT! When installing SeqScape software on a computer that is connected to a 3130/3130xl genetic analyzer, the Data Collection software must be running. Otherwise, the SeqScape software does not register with the Data Service. See Chapter 2 in the SeqScape Software User Guide v2.5, for information on installing the software.

Notes		



Importing Reports and Projects After Autoanalysis

In the SeqScape Software v2.5, select **Tools > Options** tab. You can set the options to automatically import:

- Reports after autoanalysis
- Projects after autoanalysis

See the SegScape Software User Guide v2.5 (P/N 4359442) for more details.

File-Naming Convention

Some alphanumeric characters are not valid for user names or file names. Do not use the invalid characters below:

spaces

\/:*?"<>|

IMPORTANT! An error message is displayed if you use any of these characters. You must remove the invalid character to continue.

Manual Analysis

See the SegScape Software User Guide v2.5 (P/N 4359442) for information on performing manual analysis.

About Plate Records and Segscape Software

Requirements

Successful automatic analysis requires that:

- SeqScape software is installed properly
- SeqScape software is registered and the appropriate user IDs have been created
- Autoanalysis Manager software is running
- The 3130/3130xl genetic analyzer is set up to run, and samples are prepared
- Files for a Data Collection software plate record are available:
 - An instrument protocol
 - An analysis protocol
 - A results group assigned within a plate record

When to Create a Plate Record

Create a plate record for each plate of samples for the following types of runs:

- Spectral calibrations
- Sequencing Analysis Software
- SeqScape Software analysis
- GeneMapper® software
- Mixed (sequencing and fragment analysis samples see the *Applied Biosystems* 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide for creating a mixed plate record)

Notes		

Chapter 5 Autoanalysis and SeqScape Software



About Plate Records and Segscape Software

For runs to begin, you must create a plate record for and link it to a plate loaded on the instrument. However, you can create plate records for new plates while a run is in progress.

SeqScape Software Plate Record Overview

The Plate Editor displays an empty plate record for the application selected in the New Plate dialog box. The data fields within a given plate record vary depending on the application. This section describes the SeqScape Plate Record's data fields.


IMPORTANT! For data collection and SeqScape software autoanalysis to succeed, each run of samples must have an instrument protocol, an analysis protocol, and a results group assigned within a plate record. You are also required to have a project, project template, and specimen information.

The table below describes required plate editor fields.

Parameters	Description	See Page	
Instrument Protocol	nstrument Protocol Contains the run module and dye set needed to run the instrument.		
	Created in Data Collection software.		
Analysis Protocol	Alysis Protocol Contains everything needed to analyze sequencing data.		
	Created in Data Collection software or SeqScape software. If you prefer to use analysis protocols created in SeqScape software, then select "Always use this protocol" - in the SeqScape software analysis defaults settings.		
Results Group	Defines the file type, the file name, file save locations, default analysis protocols linked to sample injections, and user name and password.	98	
	Created in Data Collection software.		
Project	A group of related sequences that share the same reference or for which there is no explicit reference.	See the SeqScape Software User Guide v2.5 for more information	
Project Template	Contains a reference data group (RDG), analysis defaults, and display settings.		
Specimen	The container that holds all the sample data as assembled contigs from a biological source or PCR product		

Notes		

Elements of a SeqScape Plate Record

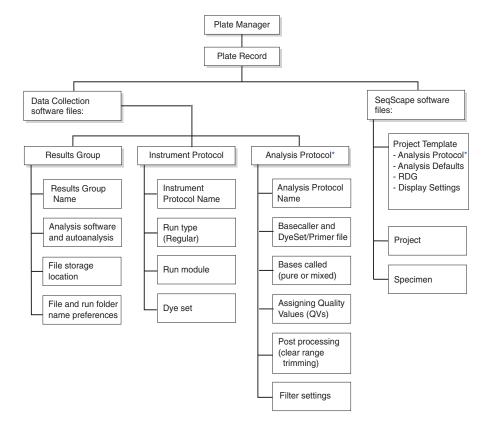


*You can create Analysis Protocols in either SeqScape or Data Colletion software



About Plate Records and Seqscape Software

SeqScape Software Plate Record

A Seqscape software plate record displaying the columns and descriptions is shown below.

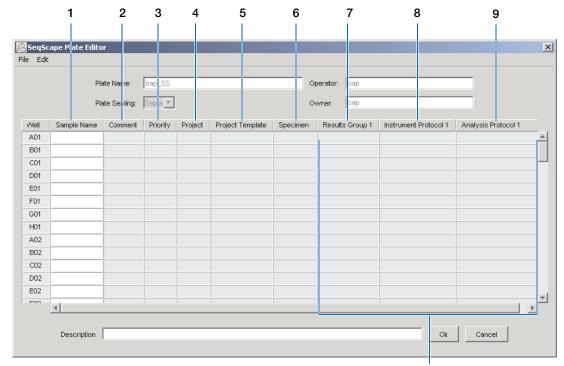

Default is one sample run, to add runs see page 103

	Number and Column	Description
1.	Sample Name	Name of the sample.
2.	Comment	Comments about the sample (optional).
3.	Priority	A default value of 100 to each sample. Changing the value to a smaller number causes that set of 16 or 4 samples to run to before the others in the injection list.
4.	Project	Select one of the available Project from the list that was created in Data Collection or SeqScape software.
5.	Project Template	Completed automatically based on the Project selected.
6.	Specimen	Select one of the available Specimen from the list that was created in Data Collection or SeqScape software.

Notes		



	Number and Column	Description
7.	Results Group	Some options:
		New: Opens the Results Group Editor dialog box
		Edit: Opens the Results Group Editor dialog box for the Results Group listed in the cell
		None: Sets the cell to have no selected Results Group
		Select one of the available Results groups from the list
		Note: You must have a Results Group selected for each sample entered in the Sample Name column.
		See "Creating a Results Group" on page 98.
8.	Instrument Protocol	New: Opens the Protocol Editor dialog box.
		Edit: Opens the Protocol Editor dialog box for the Instrument Protocol listed in the cell.
		None: Sets the cell to have no selected protocol.
		List of Instrument Protocols: In alpha-numeric order.
		Note: You must have an Instrument Protocol selected for each sample entered in the Sample Name column.
		See "Creating an Instrument Protocol" on page 88.
9.	Analysis Protocol	New: Opens the Analysis Protocol Editor dialog box.
		Edit: Opens the Analysis Protocol Editor dialog box for the Instrument Protocol listed in the cell.
		None: Sets the cell to have no selected protocol.
		List of Analysis Protocols: In alpha-numeric order
		Note: You must have an Analysis Protocol selected for each sample entered in the Sample Name column.
		See "Creating an Analysis Protocol" on page 93.

-11	v	()	110	н	3
	-	_	-	_	4

Creating Protocols for Automated SeqScape Software Analysis

If the appropriate SeqScape software protocols and results group have been created, proceed to "Filling Out a SeqScape Software Plate Record" on page 101.


Instrument Protocol for SeqScape Software

An instrument protocol contains all the necessary settings to run the instrument:


- · Protocol name
- Type of run
- Run module
- Dye set

Creating an Instrument Protocol


1. In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130 or ga3130xl > ♠ Protocol Manager.

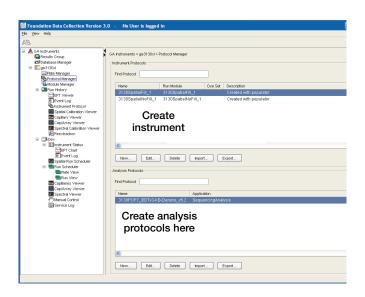

2. In the Instrument Protocols section, click New... to open the Protocol Editor.

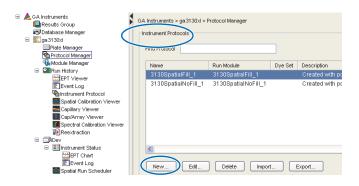

- **3.** Complete the Protocol Editor:
 - **a.** Type a name for the protocol.
 - **b.** Type a description for the protocol (optional).
 - c. Select Regular in the Type drop-down list.

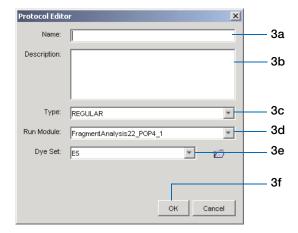

d. Select the correct run module for your run. See "Sequencing Resolution Performance and Specifications" on page 3 to see a list of modules or the Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide to modify a default module.

Chapter 5 Autoanalysis and SeqScape Software

Creating Protocols for Automated SeqScape Software Analysis

e. Select the correct Dye Set for your run using the table below.


			POP-4	-		P-6 mer		POP	-7 Pol	ymer	
Chemistry	Dye Set	UltraSeq36	StdSeq50	LongSeq80	RapidSeq36	StdSeq50	UltraSeq36	RapidSeq36	FastSeq50	StdSeq50	LongSeq80
BigDye® Terminator v3.1 Cycle Sequencing Kit	Z_BigDye	1	1	1	1	1	1	1	1	1	1
ABI PRISM® dGTP BigDye® Terminator v3.0 Cycle Sequencing Ready Reaction Kit	V3	_	_	_	_	_	_	_	_	_	_
BigDye® Terminator v1.1 Cycle Sequencing Kit	E_BigDye	1	1	1	1	1	1	1	1	1	1
ABI PRISM® dGTP BigDye® Terminator Cycle Sequencing Kit*	V1	_	_	_	_	_	_	_	_	_	_
ABI PRISM® dRhodamine Dye Terminator Cycle Sequencing Ready Reaction Kit		1	1	1	1	1	_	_	_	_	_
ABI PRISM® BigDye® Primer Cycle Sequencing Kits		_	_	_	1	1	_	_	_	_	_

^{*}dGTP kits are not supported on capillary electrophoresis instruments due to compressions on certain sequence context regions; you can run the kits if you do not care about the compression issues.

_	~		4
f.	Click	0K	ı.

Importing an Instrument Protocol

1. Click import in the Instrument Protocols pane of the Protocol Editor window to display the standard File Import dialog box.

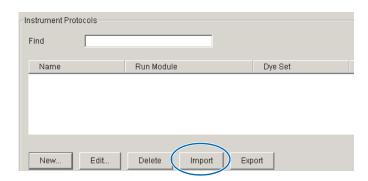

- **2.** Navigate to the location of the .xml file you want to import.
- **3.** Select the .xml file and click **Open**.

Exporting an Instrument Protocol

- **1.** In the Instrument Protocols pane, highlight the protocol you want to export.
- **2.** Click Export to display the standard File Export dialog box.
- **3.** Browse to the desired folder location.
- 4. Click Save

5

Analysis Protocol for SeqScape Software

About Analysis Protocols

An analysis protocol contains all the settings necessary for analysis and post processing:

- **Protocol name** The name, description of the analysis protocol, and the sequence file formats to be used
- **Basecalling settings** The basecaller, DyeSet/Primer file, and analysis stop point to be used
- **Mixed Bases** (Optional) When two bases are found at the same position. Define the percent value of the second highest to the highest peak
- Clear Range The high quality sequence remaining after trimming low quality sequences typically found at the beginning and ends of the sequence. The clear range can be based on base positions, sample quality values, and/or number of ambiguities (Ns) present
- **Filter** The settings that define which samples will be used in the assembly

Note: If you created an appropriate analysis protocol in the SeqScape software, you can use it in the Data Collection software. See Appendix B in the SeqScape Software User Guide v2.5 (P/N 4359442) for more information on autoanalysis and analysis protocols.

IMPORTANT! Do not delete the current analysis protocol being used, during a run. Autoanalysis will not be performed if you do so.

IMPORTANT! You are required to select the Always use this Analysis Protocol in Analysis Defaults > Analysis Protocol in the Seqscape software if you want to use it for SeqScape analysis. See the *SeqScape Software User Guide v2.5* (P/N 4359442) for more information.

Notes		

Creating Protocols for Automated SeqSc


Creating an Analysis Protocol


IMPORTANT! If you created an appropriate analysis protocol in SeqScape software, you can use it in the Data Collection software. You can also create an analysis protocol in the SeqScape software.

1. In the Analysis Protocol section of the Protocol Manager, click New... .


If more than one analysis application is installed on the data collection computer, the Analysis Applications dialog box opens.

2. Select **SeqScape**, then click OK to open the Analysis Protocol Editor dialog box.

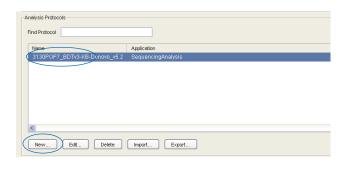

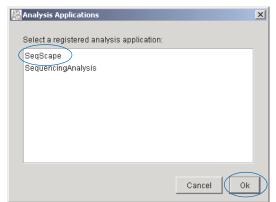

3. In the **General** tab, enter a unique name and description (optional) for the new protocol.

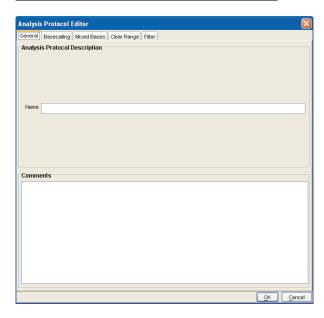

Notes

5




4. Select the **Basecalling** tab, then:

a. Select the appropriate basecaller and DyeSet primer based on the chemistry, capillary array length and polymer type you are using. The dye set primer files displayed are based on the basecaller selected. See "3130/3130xl Genetic Analyzer Basecaller and DyeSet/Primer Files" on page 66 for a table on basecallers and dyeset/primer files.

Note: SeqScape Software and the Data Collection software filter .mob file choices to match the chosen .bcp file.

b. In the Processed data pane, select True or Flat Profile.

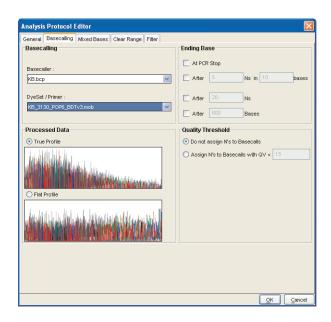

Option	Function
© True Profile	Displays data as processed traces scaled uniformly. The average peak height in the region of strongest signal is about equal to a fixed value. The profile of the processed traces is similar to that of the raw traces.
⊙ Flat Profile	Displays data as processed traces scaled semi-locally. The average peak height in any region is about equal to a fixed value. The profile of the processed traces is flat on an intermediate scale (> about 40 bases).
	Note: This option is applied to data analyzed with the KB [™] basecaller only. If you use the ABI basecaller, the profile option reverts to True Profile.

- **c.** If desired, select one or more stop points for data analysis.
- d. Select your Threshold Quality option.

Option setting	Function using KB basecaller
Call all bases and assign QV	Assigns a base to every position, as well as the QV.
Assign 'N' for bases with QV < 15	Assigns Ns to bases with QVs less than the set point. The QV will still be displayed.

Notes		

5. Select and fill in the **Mixed Bases** tab:


Note: This function is active with the KB basecaller only.


- a. For mixed bases only, select Use Mixed **Base Identification.**
- **b.** Use the default setting of 25% or change the detection level by entering a new value or dragging the % line up or down.

Note: Do not use less than 15% as your detection limit.

- **6.** Select the Clear Range tab, then, if desired, select one or more stop points for data analysis. Select the recommended settings, which are the default settings:
 - a. Use Quality values, fewer than 4 bases out of 20 have QVs less than 20.
 - **b.** Use reference trimming.

Note: The clear range is the region of the sequence that remains after excluding the lowquality or error-prone sequence at both the 5' and 3' ends.

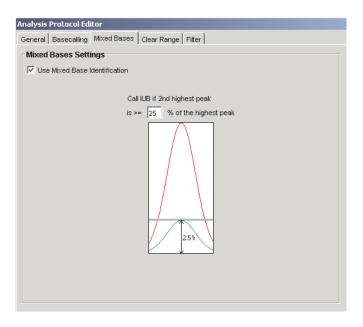

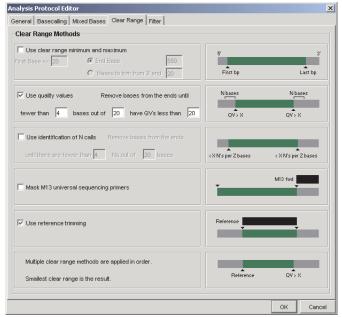

Notes

5

Chapter 5 Autoanalysis and SeqScape Software Creating Protocols for Automated SeqScape Software Analysis




7. Select the **Filter** tab, then, if desired, change one or more of the settings. Select the recommended settings, which are the default settings:

a. Maximum Mixed Bases (%): 20

b. Maximum Ns (%): 10.0

c. Minimum Clear Length (bp): 50


d. Minimum Sample Score: 20

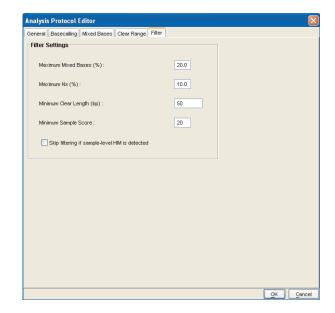

8. Click ok to save the protocol and close the Analysis Protocol Editor dialog box.

Editing an Analysis Protocol


- 1. In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to edit.
- 2. Click Edit...
- **3.** Make changes in the General, Basecalling, Mixed Bases and Clear Range tabs, as appropriate.
- **4.** Click or to save the protocol and close the Analysis Protocol Editor dialog box.

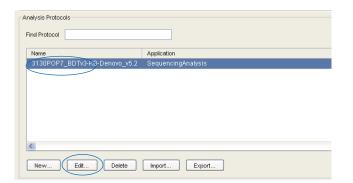

Deleting an Analysis Protocol

IMPORTANT! Do not delete an analysis protocol during a run while it is being used for that run or Autoanalysis will not be performed. Also, you must first delete any plate records using the analysis protocol before you can delete or modify the analysis protocol for these plate records.

- **1.** In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to delete.
- Click Delete .The Deletion Confirmation dialog box displays.
- 3. Click Yes .

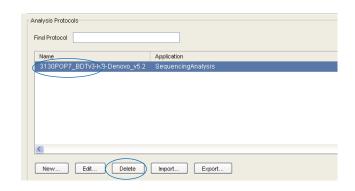

Exporting an Analysis Protocol

- 1. In the Analysis Protocols pane in the Analysis Protocol Manager, highlight the protocol you want to export.
- **2.** Click Export to display the standard File Export dialog box.
- **3.** Browse to the desired folder location.
- 4. Click Save.

Importing an Analysis Protocol

- 1. Click mort to display the standard File Import dialog box.
- **2.** Browse to the .xml file to import and click **Open**.

Note: For multiple applications, select the appropriate application to associate with the analysis protocol.

5

Results Group for SeqScape Software

A Results Group is a component within Data Collection that organizes samples and certain user settings under a single name. It is called a Results Group because it is used to name, sort, and deliver samples that result from a run.

Creating a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- 2. Click New....

The Results Group Editor window displays.


3. Complete the General tab:

- a. Type a Results Group Name. The name can be used in naming and sorting sample files. It must be unique (see page for a list of accepted characters).
- b. Optional: Type a Results Group Owner. The owner name can be used in naming and sorting sample files.
- **c.** Optional: Type a Results Group Comment.


4. Select the **Analysis** tab, then:

- a. Select SeqScape_computer name in the Analysis Type drop-down list.
- **b.** Select **Do Autoanalysis** in the Analysis Actions section.
- **c.** Type a valid SeqScape Software Login ID and Password in the text boxes.

Note: Failure to use the proper login and password causes your samples not to be analyzed automatically.

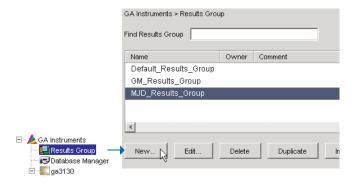



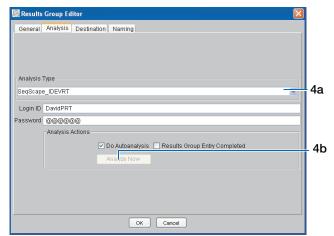

	ſΊ			


5. Select the **Destination** tab, then use the default destination or define a new location for data storage.


To use	Then
default location*	skip to step 6
custom location	complete steps a-c

- a. Click Use Custom Location, then click
 Browse... to navigate to a different save location.
- b. Click Test to test the Location path name connection:

If it	Then a message box displays
Passes	Test succeeded: <"path">.
Fails	Test failed:<"path">.

- 6. Select the Naming tab, then define custom names for the sample file and the run folder name. See "Optional: Completing Sample File Name Format Pane" on page 73 for more detailed information.
- 7. Click or to save and close the Results Group Editor.

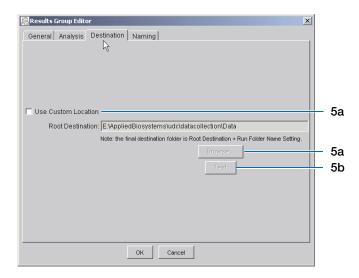

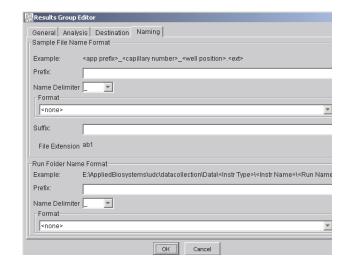

*Sample File Destinations

Locations where sample files are placed during extraction:

- Default Destination, default folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\<instrument type>\<instrument name>\run folder
- Default Destination, custom folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\top custom folder\ subfolders, etc.
- Custom Destination, default folder naming: Destination\<instrument type>\<instrument name>\run folder
- Custom Destination, custom folder naming: Destination\top customer folder\subfolders, etc.

Chapter 5 Autoanalysis and SeqScape Software Creating Protocols for Automated SeqScape Software Analysis

Importing and Exporting a Results Group

Note: Importing or exporting of SeqScape software Results Groups are not supported for transfer between separate computers.

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- **2.** Click Import to open a File Import dialog box.
- **3.** Navigate to the file you want to import.

Note: Import file type is .xml (XML file).

4. Click Open .

Note: When you import or duplicate a Results Group, you are asked to type a name for the new Results Group and for the analysis application type.

Exporting a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- **2.** Click the Results Group name to select it.
- 3. Click Export .

A standard File Export dialog box displays with the chosen Results Group name.

- **4.** Navigate to the location where you want to save the exported file.
- 5. Click Save .

Note: If there is a name conflict with a Results Group that already exists at the save location, the Results groups can be duplicated to copy settings into a similar Results Group without the risk of user error when copying it manually (see "Duplicating a Results Group" on page 100).

Duplicating a Results Group

t it.

Notes		



2. Click Duplicate .

Note: When you import or duplicate a Results Group, you are asked to type a name for the new Results Group and for the analysis application type.


Filling Out a SeqScape Software Plate Record

Creating a SeqScape Software Plate Record

- 1. Click the **Plate Manager** icon in the navigation pane.
- 2. Click New... .

The New Plate Dialog dialog box opens.

- **3.** Complete the New Plate dialog box:
 - **a.** Type a name for the plate.
 - **b.** Type a description for the plate (optional).
 - c. Select SeqScape_computer name in the Application drop-down list.
 - d. Select 96-well or 384-well in the Plate Type drop-down list.
 - **e.** Type a name for the owner.
 - **f.** Type a name for the operator.
 - g. Click OK . The SeqScape Plate Editor opens.

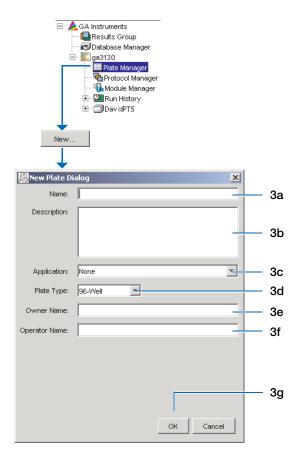




Completing a SeqScape Software Plate Record


- 1. In the **Sample Name** column of a row, enter a sample name, then click the next cell. The value 100 automatically display in the Priority column.
- **2.** In the **Comments** column, enter any additional comments or notations for the sample.
- **3.** In the **Priority** column, change the priority value, if desired. A lower number for each 4 or 16 set, receives a higher run priority. (See the *Applied Biosystems 3130/3130xl Maintenance, Troubleshooting, and Reference Guide* for changing priority values).
- 4. In the Project column, select a project from the drop-down list or create a new project with the corresponding project template.
 Based on the project you select, the project
- **5.** In the **Specimen** column, select or create a specimen.


template is filled in automatically.

- **6.** In the **Results Group 1** column, select a group from the drop-down list (see page 98).
- **7.** In the **Instrument Protocol 1** column of the row, select a protocol from the drop-down list (see page 88).
- **8.** In the **Analysis Protocol 1** column of the row, select a protocol from the drop-down list (see page 93).
- **9.** To complete the rest of the plate record based on the samples loaded in your plate, do one of the following:
 - For the same samples and protocols –
 Highlight the entire row, then select Edit >
 Fill Down Special.

Based your plate type (96- or 384-well) and capillary array (16 or 4 capillaries), the software automatically fills in the appropriate well numbers for a single run.

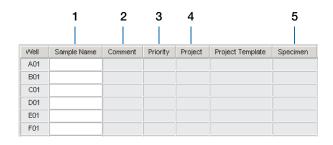

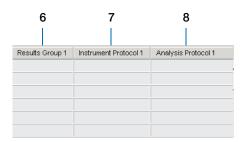

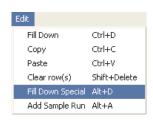



- For the same samples and protocols Highlight the entire row, then select **Edit** > Fill Down.
- For the different samples and protocols Complete the entries manually.
- **10.** If you want to do more than one run, then select Edit > Add Sample Run.


Additional Results Group, Instrument Protocol and Analysis Protocol columns are added to the right end of the plate record.

Add additional runs by selecting **Edit > Add** Sample Run again.

- **11.** Complete the columns for the additional runs.
- **12.** Click ok to save, then close the plate record.

Note: After clicking OK within the Plate Editor, the completed plate record is stored in the Plate Manager database. The plate record can be searched for, edited, exported, or deleted in the Plate Manager.


For sample preparation and running the instrument, see Chapter 7, "Running the Instrument."

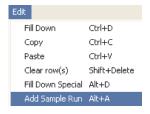

Chapter 5 Autoanalysis and SeqScape Software Filling Out a SeqScape Software Plate Record



Autoanalysis and Fragment Analysis

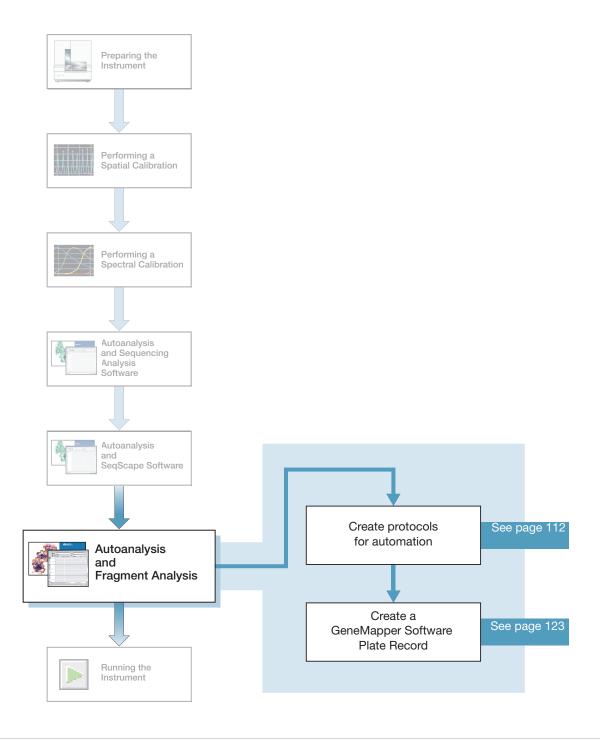

GeneMapper Software

Analyze your fragment analysis samples automatically using autoanalysis or manually.

Note: This chapter is written for both GeneMapper[®] Software v3.7 or higher and GeneMapper[®] *ID* Software v3.2. Graphic examples are from GeneMapper software v3.7.

GeneMapper Software v3.7 and GeneMapper ID Software v3.2

Perform autoanalysis of fragment analysis samples with features of the 3130/3130xl Data Collection and GeneMapper software using the same instrument computer that collected the sample files or on a remote computer.

- Autoanalysis can be performed on the same instrument computer that collected the sample files or on another computer using the remote analysis feature of GeneMapper software.
- After autoanalysis, if you wish to edit/review results on another computer, then
 transfer the GeneMapper software project, analysis methods, size standards, panel,
 and bin set information to the other GeneMapper software database. All
 components need to be exported and imported individually.
- When completing the plate record, fill in the instrument protocol and other information for Data Collection software to complete the run.
- When creating a new Results Group for a set of samples to be autoanalyzed, check the Do Autoanalysis check box and enter the appropriate GeneMapper Software user name and password. For remote analysis, define a default location for sample file storage.

File-Naming Convention

Some alphanumeric characters are not valid for user names or file names. Do not use the invalid characters below:

spaces

\/:*?"<>|

IMPORTANT! An error message is displayed if you use any of these characters. You must remove the invalid character to continue.

Manual Analysis

For information on manual analysis, see the *GeneMapper Software v3.7 User Guide* (PN 4359413) or *GeneMapper ID Software V3.1 User Guide* (PN 4338775), or the User Bulletin titled *GeneMapper ID Software v3.2*, Subject: *New Features and Installation Procedures for GeneMapper ID Software v3.2* (PN 4352543).

Notes		



About Fragment Analysis and Data Collection

When GeneMapper software is installed on a computer that has 3130/3130xl genetic analyzer Data Collection software, two applications are available through the Results Group Editor (see page 115):

- GeneMapper-Generic
- GeneMapper-<Computer Name>

GeneMapper-Generic

Use GeneMapper-Generic to generate .fsa files, without performing autoanalysis. When completing the Sample Sheet, fill in basic information for the Data Collection software to complete the run; all other GeneMapper software related fields are text entries. Text entries are useful if you are using other software analysis applications or choose to analyze your samples in GeneMapper software on another computer, but do not have the same entries in the GeneMapper software database stored on the Data Collection computer. For example, if you have a customized size standard definition on another GeneMapper software computer, you can type in that size standard name in the size standard text field and it will populate that column in your GeneMapper software project.

GeneMapper-<Computer Name>

Use GeneMapper-<Computer Name> for autoanalysis. The Size Standard, Analysis Method, and Panel columns in the Sample Sheet window read directly from the GeneMapper software database. Create these components in GeneMapper software prior to setting up the plate record for a run. New entries cannot be created for these columns once you are in the plate editor dialog box. If you create a new GeneMapper software component while the plate record dialog box is open, the columns will not update. The plate record must be closed and reopened to update the GeneMapper software components. For more information see, "Creating Protocols for Automated Fragment Analysis" on page 112.

Notes		

About GeneMapper Software Plate Records

A plate record is similar to a sample sheet or an injection list that you may have used with other ABI PRISM® instruments.

Plate records are data tables in the instrument database that store information about the plates and the samples such as:

- Plate name, type, and owner
- Position of the sample on the plate (well number)
- Comments about the plate and about individual samples
- Dye set information (in Instrument protocol)
- Name of the run module where run modules specify information about how samples are run (in Instrument protocol)

When to Create a Plate Record

Create a plate record for each plate of samples for the following types of runs:

- Spectral calibrations
- Sequencing Analysis Software
- SeqScape Software analysis
- GeneMapper® software
- Mixed (sequencing and fragment analysis samples see the *Applied Biosystems* 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide for creating a mixed plate record)

For runs to begin, you must create a plate record for and link it to a plate loaded on the instrument. However, you can create plate records for new plates while a run is in progress.

Files needed to create a GeneMapper software plate record:


Parameters	Description	
Instrument Protocol	Contains everything needed to run the instrument.	112
Results Group	Defines the file type, the file name, autoanalysis, and file save locations that are linked to sample injections.	115

IMPORTANT! For data collection and auto-analysis to be successful, each run of samples must have an Instrument Protocol, a Results Group and files created in GeneMapper software assigned within a plate record.


Notes		



Elements of a GeneMapper software plate record

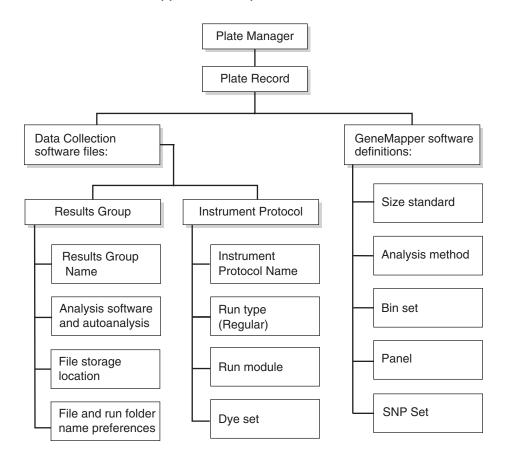



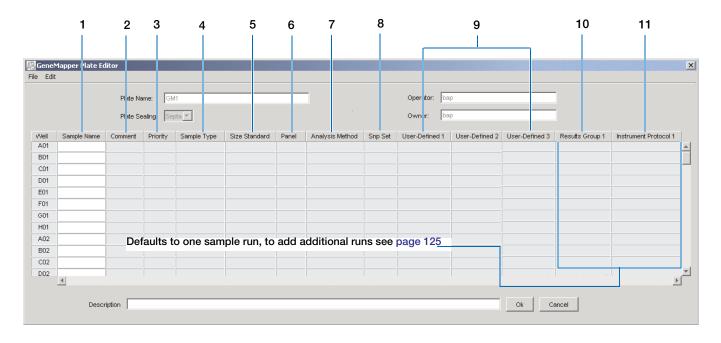

Blank GeneMapper software plate record

Columns inserted in a plate record for a fragment analysis run

	Number and Column	Description
1.	Sample Name	Name of the sample
2.	Comment	Comments about the sample (optional)
3.	Priority	A default value of 100 to each sample. Changing the value to a smaller number causes that set of 16 or 4 samples to run to before the others in the injection list.
4.	Sample Type	Use to identify the sample as Sample, Positive Control, Allelic Ladder or Negative Control.
5.	Size Standard	GeneMapper-Generic (optional):
	IMPORTANT! For GeneMapper- <computer< td=""><td>Manually enter size standards in the text field</td></computer<>	Manually enter size standards in the text field
	Name> ONLY:	GeneMapper- <computer name="">:</computer>
	Create the Size Standard, Panel, and Analysis Method in GeneMapper software before creating a new plate	Select a saved size standard from the drop-down list
6.	Panel	GeneMapper-Generic (optional):
	IMPORTANT! For GeneMapper- <computer< td=""><td>Manually enter panels in the text field</td></computer<>	Manually enter panels in the text field
	Name> ONLY:	GeneMapper- <computer name="">:</computer>
	Create the Size Standard, Panel, and Analysis Method in GeneMapper software before creating a new plate	Select a saved panel from the drop-down list

Notes			

Chapter 6 Autoanalysis and Fragment Analysis About GeneMapper Software Plate Records



	Number and Column	Description
7.	Analysis Method IMPORTANT! For GeneMapper <computer name=""> ONLY: Create the Size Standard, Panel, and Analysis Method in GeneMapper software before creating a new plate</computer>	 GeneMapper-Generic (optional): Manually enter analysis methods in the text field GeneMapper-<computer name="">: Select a saved analysis method from the drop-down list</computer>
8.	SNP Set IMPORTANT! For GeneMapper <computer name=""> ONLY: The SNP set, created in the GeneMapper software, links a SNP name to a marker name.</computer>	 GeneMapper-Generic (optional): Manually enter SNP set in the text field GeneMapper-<computer name=""></computer>
9.	Three User-defined columns	Optional text entries
10.	Results Group	 Options: New: Opens the Results Group Editor dialog box Edit: Opens the Results Group Editor dialog box for the Results Group listed in the cell None: Sets the cell to have no selected Results Group Select one of the available Results groups from the list Note: You must select a Results Group for each sample entered in the Sample Name column. See "Results Group for Fragment Analysis" on page 115.
11.	Instrument Protocol	 Options: New: Opens the Protocol Editor dialog box. Edit: Opens the Protocol Editor dialog box for the Instrument Protocol listed in the cell. None: Sets the cell to have no selected protocol. List of Instrument Protocols: In alpha-numeric order. Note: You must select an Instrument Protocol for each sample entered in the Sample Name column. See "Results Group for Fragment Analysis" on page 115.

- N	л.	$\overline{}$	4.	$\overline{}$	è
1	V	U	U	U	S

Creating Protocols for Automated Fragment Analysis

If the appropriate fragment analysis protocols and results group have been created, proceed to "Creating a GeneMapper Software Plate Record" on page 123.


Instrument Protocol for Fragment Analysis

An instrument protocol contains all the settings necessary to run the instrument:


- · Protocol name
- Type of run
- Run module
- Dye set

Creating an Instrument Protocol

1. In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or ga3130 > ➡ Protocol Manager.

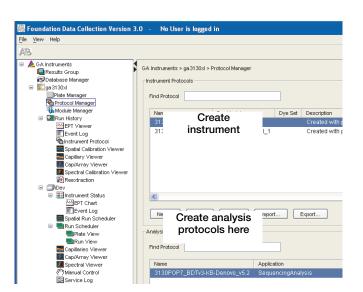


2. In the Instrument Protocols section, click New... to open the Protocol Editor.

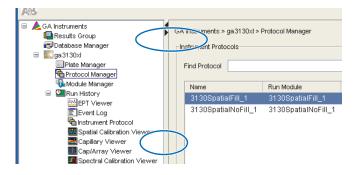



- **3.** Complete the Protocol Editor:
 - **a.** Type a name for the protocol.
 - **b.** Type a description for the protocol (optional).
 - c. Select Regular in the Type drop-down list.

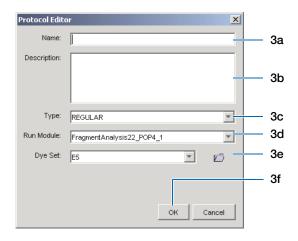

d. Using the table below, select the run module for your run. To customize a run module, see the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide* on modifying a module.

Application or Kit	Capillary Array Length (cm)	Run Module
SNaPshot® Multiplex System	22	SNP22_POP4_1
	36	SNP36_POP4_1
• LMS v2.5	22	FragmentAnalysis22_POP4_1
ABI PRISM® Mouse Mapping Set v1.0	36	FragmentAnalysis36_POP4_1
Custom oligos		FragmentAnalysis36_POP7_1
Stockmarks AFLP®	50	FragmentAnalysis50_POP4_1
		FragmentAnalysis50_POP6_1
		FragmentAnalysis50_POP7_1
 AmpFℓSTR® COfiler® AmpFℓSTR Profiler Plus® AmpFℓSTR SGM Plus® AmpFℓSTR Profiler Plus ID AmpFℓSTR SEfiler™ AmpFℓSTR Yfiler™ Other 4-Dye AmpFℓSTR AmpFℓSTR Identifiler Other 5-Dye AmpFℓSTR 	36	HIDFragmentAnalysis36_POP4_1

Chapter 6 Autoanalysis and Fragment Analysis Creating Protocols for Automated Fragment Analysis


e. Using the table below, select the correct Dye Set for your run.

Application or Kit	Dye Set	Matrix Standard Set
Custom oligos	D	DS-30
ABI PRISM Mouse Mapping Set v1.0 Custom oligos	D	DS-31
 AmpFlSTR COfiler AmpFlSTR Profiler Plus AmpFlSTR SGM Plus AmpFlSTR Profiler Plus ID Other 4-Dye AmpFlSTR AFLP Stockmarks -Canine and Bovine 	F	DS-32
ABI PRISM SNaPshot Multiplex System	E5	DS-02
 ABI PRISM Linkage Mapping Set v2.5 Custom Oligos AmpFℓSTR Identifiler™ AmpFℓSTR SEfiler AmpFℓSTR Yfiler Other 5-Dye AmpFℓSTR Stockmarks-Equine 	G5	DS-33

f. Click OK.

Importing an Instrument Protocol

1. Click in the Instrument Protocols pane of the Protocol Editor window to display the standard File Import dialog box.

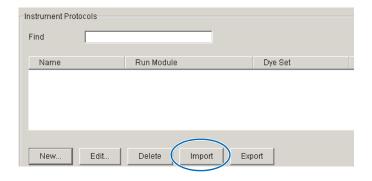

- **2.** Navigate to the location of the .xml file you want to import.
- **3.** Select the .xml file and click **Open**.



Exporting an Instrument Protocol


- **1.** In the Instrument Protocols pane, highlight the protocol you want to export.
- **2.** Click Export to display the standard File Export dialog box.
- **3.** Browse to the desired folder location.
- 4. Click Save .

Results Group for Fragment Analysis

A Results Group is a component within Data Collection that organizes samples and certain user settings under a single name. It is called a Results Group because it is used to name, sort, and deliver samples that result from a run.


Creating a Results Group for Autoanalysis

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- **2.** Click New... to open the Results Group Editor window.

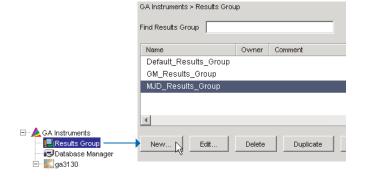




- **3.** Complete the General tab:
 - **a.** Type a unique Results Group Name. The name can be used in naming and sorting sample files.
 - **b.** Optional: Type a Results Group Owner. The owner name can be used in naming and sorting sample files.
 - **c.** Optional: Type a Results Group Comment.

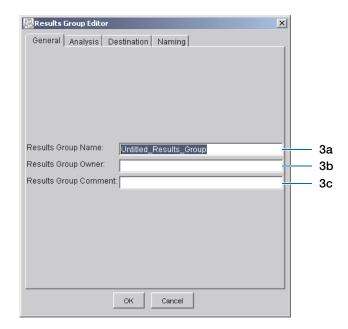

- 4. Select the Analysis tab, then:
 - **a.** Click **Analysis Type** and then select one of the following:

If You Select	Then
None	Only raw data files are generated.
GeneMapper- Generic	Autoanalysis is not enabled and only .fsa files are generated.
GeneMapper- <computer Name></computer 	Autoanalysis of completed runs is enabled. Steps b, c, and d below apply only to GeneMapper- <computer name=""> (not GeneMapper-Generic).</computer>

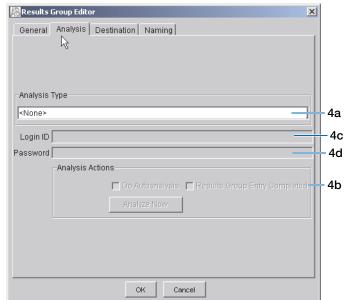



b. Select an option in the Analysis Actions section using the table below.


If You Select	Then	Use with Setting from Automated Processing Tab (page 118)
Do Autoanalysis	Samples are analyzed after each run of 16 or 4 samples.	When every run completes
Do Autoanalysis and Results Entry Group Complete	Samples are analyzed after all pending samples using the same results group have been run.	Only when the result group is complete

- c. Type the Login ID.
- d. Type the login password.

Note: The login ID and password relate to the GeneMapper software UserName and Password. These items can only be created through the GeneMapper software Options Users tab. Autoanalysis will fail if the incorrect username and password are entered.

5. Select the **Destination** tab, then use the default destination or define a new location for data storage.

To use a	Then
default location*	skip to step 6.
custom location	complete steps a-b.
Use for remote analysis using GeneMapper v3.7	

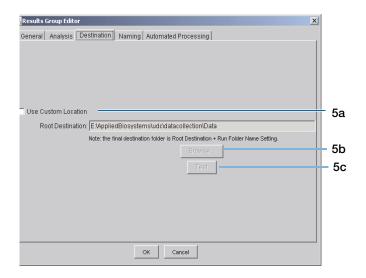

*Sample File Destinations

Locations where sample files are placed during extraction:

- Default Destination, default folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\<instrument type>\<instrument name>\run folder
- Default Destination, custom folder naming: E:\AppliedBiosystemsUDC\DataCollection\data\top custom folder\ subfolders, etc.
- Custom Destination, default folder naming: Destination\<instrument type>\<instrument name>\run folder
- · Custom Destination, custom folder naming: Destination\top customer folder\subfolders, etc.

Notes		



- a. Click Use Custom Location, then click Browse... to navigate to a different save location.
- **b.** Click Test to test the Location path name connection:

If it	Then a message box displays
Passes	Test succeeded: <"path">.
Fails	Test failed:<"path">.


6. Select the **Naming** tab to customize sample file and run folder names.

Note: The default sample file naming for fragment analysis is:

frag_<capillary number>_<wellposition>.fsa

Note: Sample name, run folder name, and path name, combined, can total no more than 250 characters.

For defining the elements of the Naming tab, see page 119.

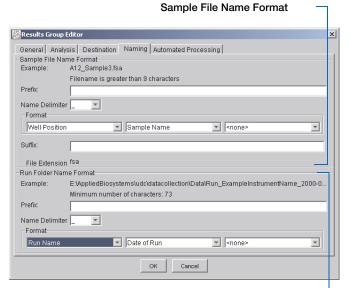

Run Folder Name Format pane

7. Select the Automated Processing tab.

In "Autoanalysis is performed" select when you want your samples autoanalyzed using the table below.



autoanalysis option

Select an

If You Select	Then	Use with Settings from Analysis Tab (page 117)
Only when the result group is complete	Samples are analyzed after all pending samples using the same results group have been run.	Do Autoanalysis and Results Entry Group Complete
When every run completes	Samples are analyzed after each run of 16 or 4 samples.	Do Autoanalysis


8. Click OK to save the Results Group.

Notes		



Completing the Sample File Name Format Pane

- 1. Type a prefix for the file name in the **Prefix** box. Anything that you type here is shown in the Example line.
- 2. Click the Name Delimiter list choose the symbol that will separate the Format elements in the file name (see step 3 below). Only one delimiter symbol may be chosen.

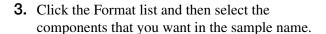


Note: All the samples from a single run can be placed in the same run or results folder; the name of every sample from a single run should be different. Most of the Format options will not be different between samples; select at least one of the options to make the sample names unique within a run.


If a unique identifier is not included in the name, a warning message displays. Select from the elements to make a unique sample file name. Examples are:


- Well position
- Capillary number
- Run sequence number
- Plate Quadrant

Note: Some identifiers, such as well position, capillary number and plate quadrant are not unique in multiple sample instances. Plate Quadrant is not a unique identifier for a 96-well plate.



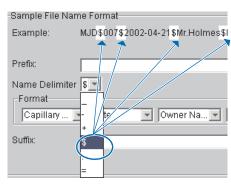

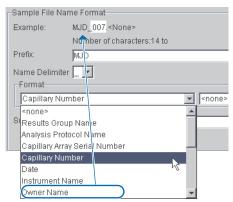

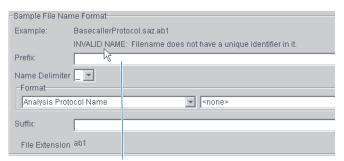


Warning message


Notes			



As you select the elements for the file name, they are placed in the Example line. An additional element drop-down menu displays allowing you the option of selecting an additional element.

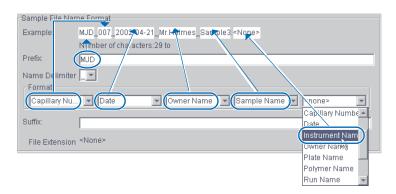

The names of the Format elements eventually truncate, but the Example field remains visible (up to 72 characters).

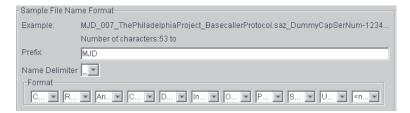


4. Click the Suffix box (optional) and type the suffix for the file name.

The File Extension field displays the file extension generated from the Analysis Type specified on the Analysis tab (see page 116). For example, Sequencing Analysis produces sample files with an .ab1 extension.

Run Folder/Sub-Folder Name Format Pane

Follow the same steps described above for the Sample File Name Format pane (see page 119) to change the sub-folder name within the run folder.

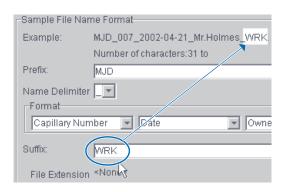



Importing and Exporting a Results Group

Importing or exporting of GeneMapper software Results Groups are not supported for transfer between separate computers.

Importing a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- 2. Click Import .

A standard File Import dialog box displays.

3. Navigate to the file you want to import.

Note: Import file type is .xml (XML file).

4. Click Open

Note: When you import or duplicate a Results Group, you are asked to type a name for the new Results Group and for the analysis application type.

Exporting a Results Group

- 1. In the tree pane of the Data Collection software, click ▲ GA Instruments > Results Group.
- **2.** Click the Results Group name to select it.
- 3. Click Export .

A standard file export dialog box displays with the chosen Results Group name.

- **4.** Navigate to the location where you want to save the exported file.
- 5. Click Save

Note: If there is a name conflict with a Results Group that already exists at the save location, the Results groups can be duplicated in order to copy settings into a similar Results Group without the risk of user error when copying it manually (see "Duplicating a Results Group" on page 122).



Duplicating a Results Group

- **1.** Click the Results Group.
- 2. Click Duplicate .

Note: When you import or duplicate a Results Group, you are asked to type a name for the new Results Group and for the analysis application type.

Notes			




Creating a GeneMapper Software Plate Record

Creating the Plate Record for Autoanalysis


- In the tree pane of the Data Collection software, click ▲ GA Instruments > ga3130xl or ga3130 > Plate Manager.
- **2.** Click New... to open New Plate Dialog.
- **3.** Complete the information in the New Plate Dialog:
 - **a.** Type a name for the plate.
 - **b.** Type a description for the plate (optional).
 - **c.** Select your GeneMapper application in the Application drop-down list.
 - d. Select **96-well** or **384-well** in the Plate Type drop-down list.
 - e. Type a name for the owner.
 - **f.** Type a name for the operator.
 - g. Click OK.


 The GeneMapper software Plate Editor opens.

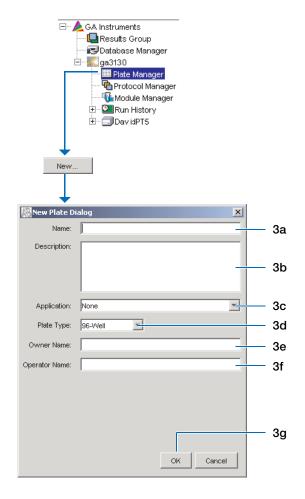


Completing a GeneMapper Software Plate Record for Autoanalysis

- **1.** In the Sample Name column of a row, enter a sample name, then click the next cell. The value 100 automatically display in the Priority column.
- **2.** In the Comment column, enter any additional comments or notations for the sample.
- **3.** In the Priority column, change the priority value, if desired.
- **4.** In the Sample Type column, select a sample type from the drop-down list.
- **5.** In the Size Standard column, select a size standard from the drop-down list.
- **6.** In the Panel column, select a panel from the drop-down list.
- **7.** In the Analysis Method column, select a method from the drop-down list.
- **8.** In the Snp Set column, select a SNP set from the drop-down lis if applicable, otherwise select None.
- **9.** Enter text for User-Defined columns 1 to 3.
- **10.** In the Results Group 1 column, create a new Results Group or select a group from the drop-down list (see page 115).
- **11.** In the Instrument Protocol 1 column, create a new instrument protocol or select a protocol from the drop-down list (see page 112).

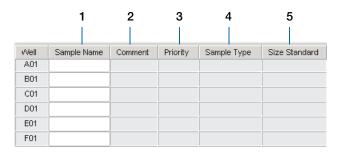

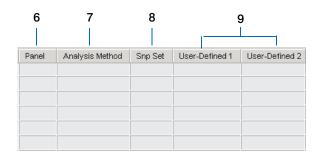

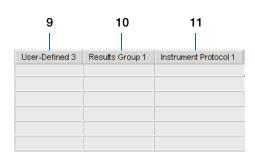



12. To complete the plate record:

For the same samples and protocols –
 Highlight the entire row, then select Edit >
 Fill Down Special.

Based on your plate type (96- or 384-well) and capillary array (16 or 4 capillaries), the software automatically fills in the appropriate well numbers for a single run.


- For the same samples and protocols –
 Highlight the entire row, then select Edit >
 Fill Down.
- For the different samples and protocols Complete the entries manually.
- **13.** To do more than one run, select **Edit > Add Sample Run**.


Additional Results Group, Instrument Protocol and Analysis Protocol columns are added to the right end of the plate record.

Add additional runs by selecting **Edit > Add Sample Run** again.

- **14.** Complete the columns for the additional runs.
- **15.** Click ok to save, then close the plate record.

IMPORTANT! After clicking OK in the Plate Editor, the plate record is stored in the Plate Manager database and can now be located, edited, exported, or deleted.

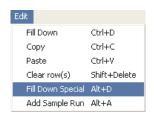

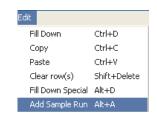

- N	ь	$\overline{}$	4.	~
1	A (U	U	E:




Chapter 6 Autoanalysis and Fragment Analysis

Creating a GeneMapper Software Plate Record

Notes			



Running the Instrument

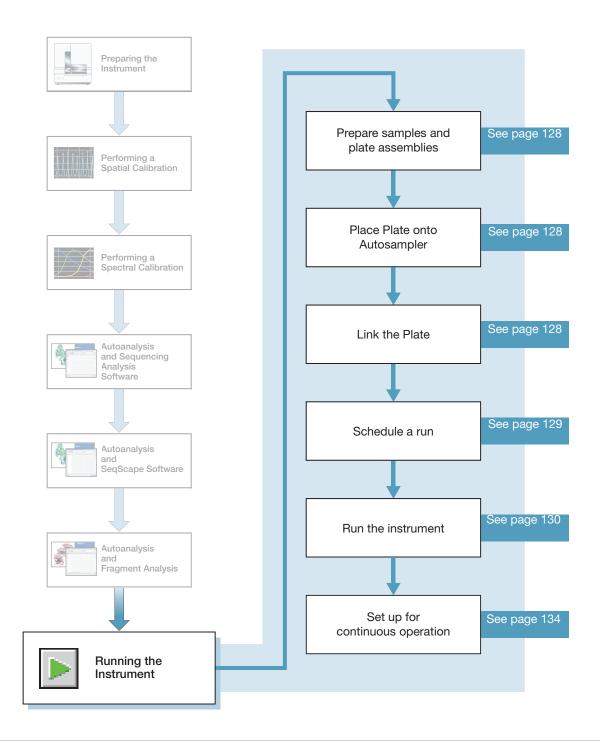

Preparing Samples and Plate Assemblies

Review Table

Generalized sample preparation and plate assembly procedures are discussed in Chapter 3, "Performing a Spectral Calibration," on page 27.

Topic	See
Sample Preparation	"Preparing the Calibration Standard" on page 31 and "Loading Samples" on page 32 to review handling samples.
	Follow your individual kit protocol to prepare sample load volume:
	10 to 30 µL for a 96-well plate
	5 to 15 μL for a 384-well plate
Sealing plates and plate assembly	"Seal the plate:" on page 33.
Placing plate assemblies in the instrument	"Placing the Plate Assembly into the Instrument" on page 39.
Searching for plate records to link to a plate	"Running the Spectral Calibration Plate" on page 40.


Linking and Unlinking a Plate

Link a plate on the autosampler to your newly created plate record before running the plate.

Select the desired plate record, then click the plate position indicator corresponding to the plate position in the instrument. The plate position (A or B) displays in the link column.

Note: The 3130 genetic analyzer has only one plate position (B) to link a plate record.


The plate position indicator changes from yellow to green when linked and the green run button becomes active.

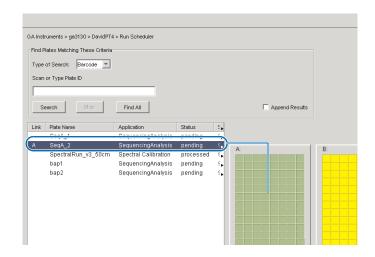

Notes____

Unlinking a Plate Record

- **1.** Click the plate record that you want to unlink.
- 2. Click Unlink .

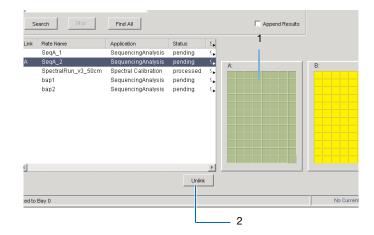


Viewing the Run Schedule

For more information on run scheduling and sample injection order, see "Plate Mapping" on page 161.

To verify that runs are scheduled correctly:

- In the tree pane of the Data Collection software, click ▲ GA Instruments > ga3130xl or ga3130 > instrument name > Run Scheduler > Run View.
- **2.** Select a row for any run. The corresponding wells to be injected for that run are highlighted in the plate diagram.

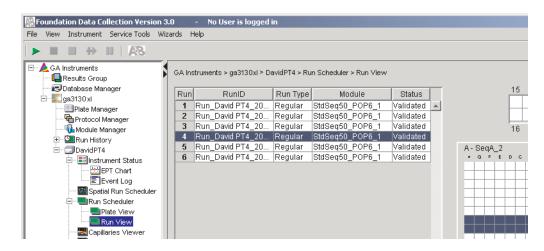

Running the Instrument

Starting the Run

1. Verify that the active spectral calibration matches your dye set and capillary array length for all scheduled runs.

To change the active spectral calibration, see "Activating a Spectral Calibration" on page 47.

To create a new spectral calibration, see "Creating a Spectral Instrument Protocol" on page 35.

- **2.** Click the green button in the toolbar.
- **3.** The Processing Plates dialog box opens, then click OK .

The software automatically performs a run validation:

- If the validation passes, the run starts
- If any of the validation test fails, the run does not start. Check the event log for information.




Starting the Autoanalysis Manager for SeqScape and/or GeneMapper Software

Use the Autoanalysis Manager (AAM) software with SeqScape® and/or GeneMapper® software to automatically analyze the data.

Select Start > Programs > Applied Biosystems > Autoanalysis Manager > Autoanalysis Manager 3.0.

Note: Autoanalysis Manager does not start automatically. Open Autoanalysis Manager to receive messages from the data collection software for autoanalysis in SeqScape and/or GeneMapper software.

The Autoanalysis Manager window opens.

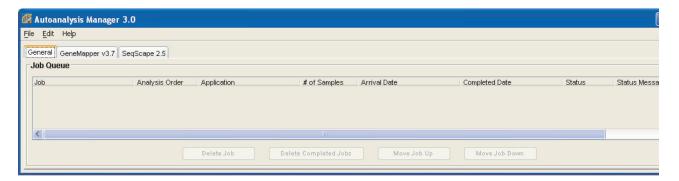

2. Quit the SeqScape and/or GeneMapper software.

No other interaction with the AAM software is needed until the runs finish. See "Using Autoanalysis Manager Software" on page 155 on how to use the Autoanalysis Manager.

Basic Run Module Steps

Automated module steps when the run starts

Module Steps	Approximate Time
Turn Oven On	N/A
Wait for oven to equilibrate Initialize autosampler	1 min 40 sec
Fill Array	3-4 min
PreRun	3 min
Inject samples	10 to 30 sec
Start separation Ramp voltage	10 min
Collect Data	Variable
Run ends: Leave oven on Laser to idle	Until next run starts
Total time prior to separation:	1

Total time prior to separation:

- Oven unheated: ~25 min
- Oven at set temperature: ~6.5 min

To customize a run module, see the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting and Reference Guide.*

Note: The PostBatch Utility, which runs automatically, turns off both the oven and the laser at end of a batch of runs.

DNA Sequencing Run Times

Approximate run times of common DNA sequencing analysis runs

Type of Run	Run Module	Run Time (min)
Ultra rapid	UltraSeq36_POP4	40
	UltraSeq36_POP7	35
Rapid	RapidSeq36_POP6	60
	RapidSeq36_POP7	
Fast	FastSeq50_POP7	60
Standard	StdSeq50_POP4	100
	StdSeq50_POP6	150
	StdSeq50_POP7	120
Long read	LongSeq80_POP4	210
	LongSeq80_POP7	170

Notes			

Fragment Analysis Run Times

Approximate run times of common fragment analysis runs

Type of Run	Run Modules	Run Time (min)
High Throughput, Small Size	FragmentAnalysis22_POP4	20
Fragment Analysis	SNP22_POP4	20
Standard Fragment Analysis	FragmentAnalysis36_POP4	45
	HIDFragmentAnalysis36_POP4	45
	SNP36_POP4	30
	FragmentAnalysis36_POP7	35
	FragmentAnalysis50_POP4	65
	FragmentAnalysis50_POP6	90
	FragmentAnalysis50_POP7	50

For the latest size standard modules (for example, run modules for the LIZ 1200 system), go to http://www.appliedbiosystems.com, then click the link for Support, then Software Downloads, then GeneScanTM Size Standards, then Run Modules.

Controlling the Run

Using the Toolbar

Use the toolbar at the top of the data collection software window to control the run.



Click	Description
•	Starts the run
Start Run	
	Stops the current run, and all other scheduled runs
Stop	
	Completes the current run, then stops all other scheduled runs
Stop After Current Run	

Click	Description
43-	Stops the current run, then starts the other scheduled runs
Skip to Next Run	
11	Pauses the current run ^a
Pause Run	

a Pausing the instrument for too long, especially after sample injection, will adversely affect data quality.

Set Up for Continuous Operation

Overview Use the continuous run feature to create and link a plate during a run. You can:

- Run one or more plates
- Remove the plate(s) once samples have run
- Link and run additional plates

When using the continuous run feature:

- Only mount or unmount plates when the instrument is paused.
- Create the plate record before pausing the instrument to make linking more efficient, reducing long pauses.
- Link new plates after a run has resumed.

Adding or Replacing a Plate During a Run

During a run, you can mount and unmount plates while the instrument is paused. The plate record can be created, then linked after the run has been resumed.

If the plate is in use, see "Adding, Replacing, or Removing a Plate During a Run" on page 135.

IMPORTANT! Pausing the instrument for too long, especially after sample injection, will affect data quality.

Notes		

Adding, Replacing, or Removing a Plate During a Run

Replace a completed plate or add a new plate to an unused plate bay (3130xl genetic analyzer only).

- **1.** Prepare your plate and create the plate record.
- **2.** Click [1] (Pause).

In the Pause dialog box, read the pause warning, then click ok to pause the run.

The Resume run dialog box opens when the run is paused.

IMPORTANT! Do not click OK to resume the run. Temporarily ignore the dialog box.

- **3.** Remove the old plate, if applicable.
 - a. Press the Tray button to bring the autosampler forward.
 - **b.** Open the instrument door.
 - **c.** Remove the old plate.
- **4.** Mount the new plate.
- **5.** Close the door.
- **6.** Click **OK** in the Resume Run dialog box.
- 7. In the Completed Run dialog box, click ok to continue if the samples have been injected, or click | Cancel | to abort the run and return the instrument to an idle state.

IMPORTANT! If you click OK, the instrument will continue running the current run regardless of whether the samples have actually been injected or not. If the samples have not been injected, the samples will be injected from the new plate.









8. Search for the plate record, then link the new plate.

The new plate runs after the current plate completes all scheduled injections.

Note: If you unmount the currently running plate prior to the first frame of data being collected but after sample injection (clicked OK to continue), the plate status changes to processed even though the run is actually continuing.

Notes		




Viewing Data During a Run

Run Scheduler > Plate View

In the tree pane of the Data Collection software, click

▲ GA Instruments > \blacksquare ga3130xl or ga3130 > \blacksquare instrument name > \blacksquare Run Scheduler > \blacksquare Plate View.


Note: The **note** Run Scheduler and **note** Plate View windows display the same information.

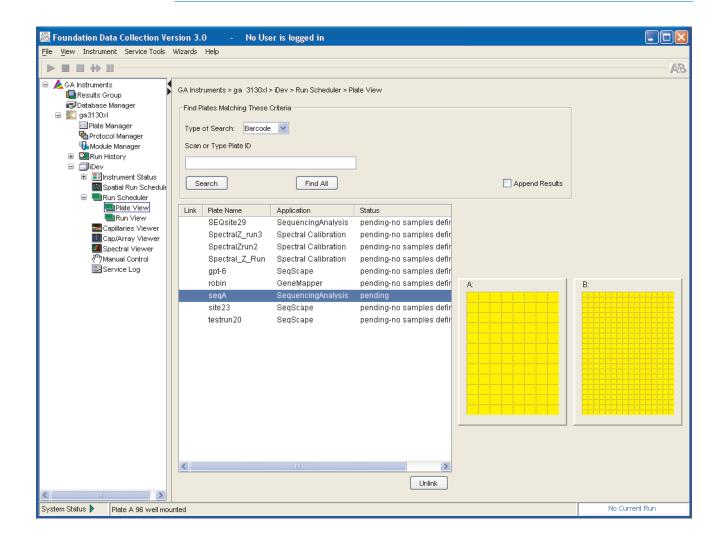

Run Scheduler > Run View

In the tree pane of the Data Collection software, click

△ GA Instruments > ■ ga3130xl or ga3130 > ■ instrument name > ■ Run Scheduler > ■ Run View to monitor the status of the scheduled runs.

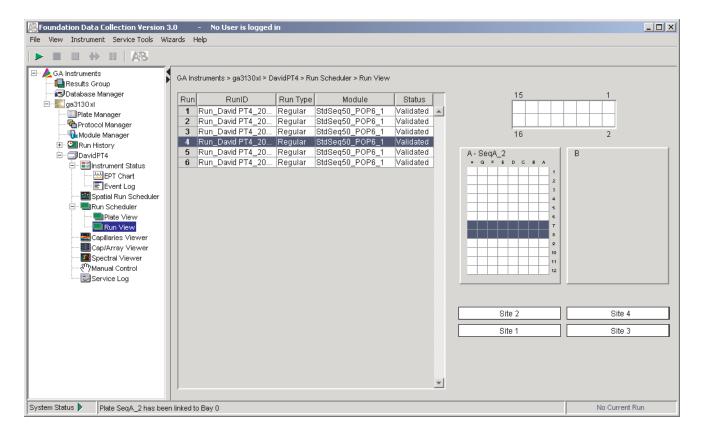


Note: For default load maps, see Appendix A, "Plate Mapping."



Instrument Status

In the tree pane of the Data Collection software, click

🛕 GA Instruments > 彲 ga3130xl or ga3130 > 🗊 instrument name > 📰 Instrument **Status** to monitor the status of the instrument or the current run.

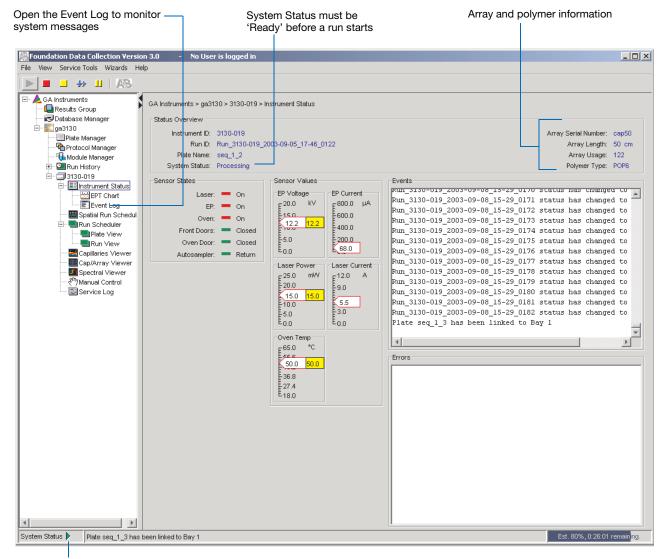

System Status changes from green to flashing red when errors occur, see Event Log.

Instrument Condition Sensor States Pane

The color of the indicator provides a quick way to check the status of the item to the left. See the table below for a definition of each color.

For	A green box indicates	A red box indicates	A yellow box indicates
Laser	Laser is off	Laser is on	Laser is idle
EP	Electrophoresis is off	Electrophoresis is on	_
Oven	Oven is off	Oven is on	_
Front Doors	Doors are closed	Doors are open	_
Oven Door	Door is closed	Door is open	_
Autosampler	Autosampler is homed	Autosampler is forward	_

Events Box

The Events box lists the:

- Instrument's recent actions
- Status of each capillary as passed or failed at the end of a spectral calibration
- Calibration data at the end of a spatial calibration

Some of the events listed in the Events box provide information for service engineers.

Errors Box

The Errors box lists errors that have occurred during the current run.

Some of the error messages provide information for service engineers. A "fatal" error usually requires that you restart the data collection software.


Notes		

Instrument Status > EPT Chart

In the tree pane of the Data Collection software, click ▲ GA Instruments >

ga3130xl or ga3130 > □ instrument name > □ Instrument Status > □ EPT

Chart. The EPT chart displays real-time electrophoresis (EP) data during a run.

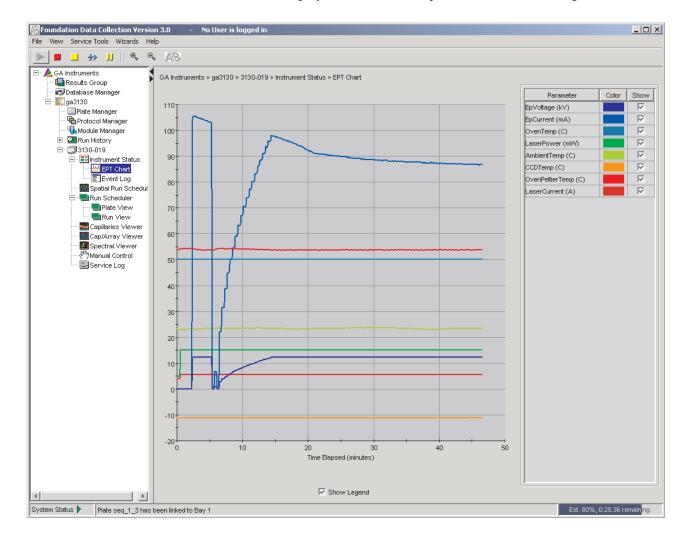


Notes____

Instrument Status > Event Log


In the tree pane of the Data Collection software, click ▲ GA Instruments > ■ ga3130xl or ga3130 > ■ instrument name > ■ Instrument Status > ■ Event Log. The Event log itemizes events such as errors and general information for all data collection steps.

Clear error messages by clicking **Clear Errors**. The System Status light flashes red until all errors are cleared. Take corrective action based on error message.

Note: This view can also be used to monitor spectral calibration results in real time to verify the capillary-by-capillary processing status.

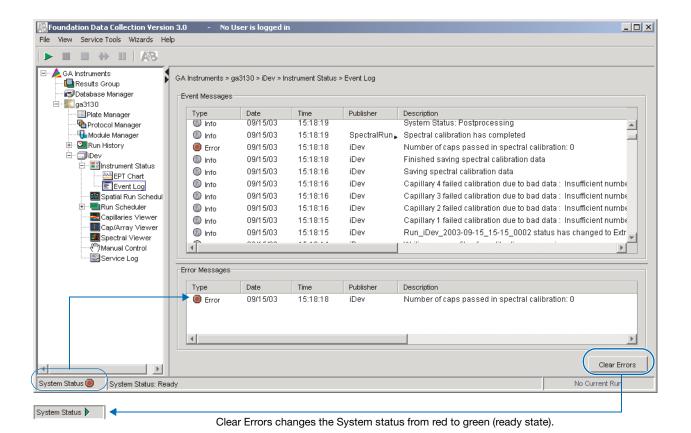

In the tree pane of the Data Collection software, click GA Instruments > ga3130 or ga3130xl > instrument name > Capillaries Viewer. Use the Capillary Viewer to examine the quality of the raw data during a run for several capillaries at once.

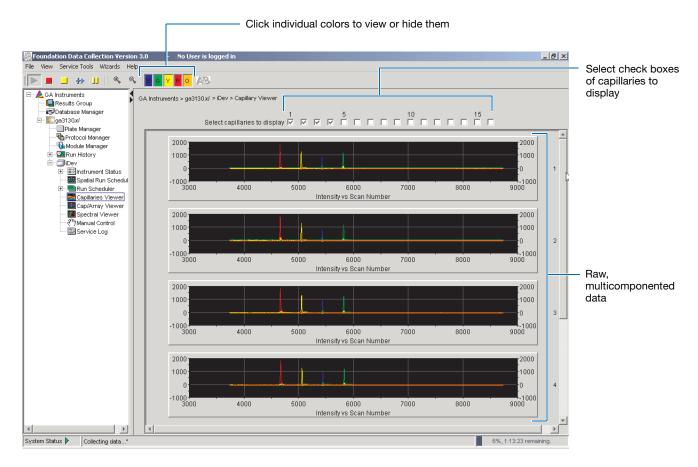

Check Boxes

Select the check boxes of the capillaries to view their electropherograms. The capillaries are displayed in the order in which the boxes are checked. The more boxes that are selected, the slower the refresh window rate.

Raw Data

An electropherogram is a graph of relative dye concentration against time, plotted for each dye. The displayed raw data has been corrected for spectral overlap (multicomponented).

How to Zoom

To zoom in and out:


- 1. Select a rectangular area over the area of interest by holding down the mouse button. Release the mouse button to zoom in.
- **2.** Click to return to full view.

Notes

143

Cap/Array Viewer

In the tree pane of the Data Collection software, click \triangle GA Instruments $> \blacksquare$ ga3130 or ga3130x $l > \blacksquare$ instrument name $> \blacksquare$ Cap/Array Viewer. Use the window during a run to examine the quality of your data, which is displayed as color data for the entire capillary array. You can view all the capillaries (vertical axis) as a function of time/scan numbers (horizontal axis).

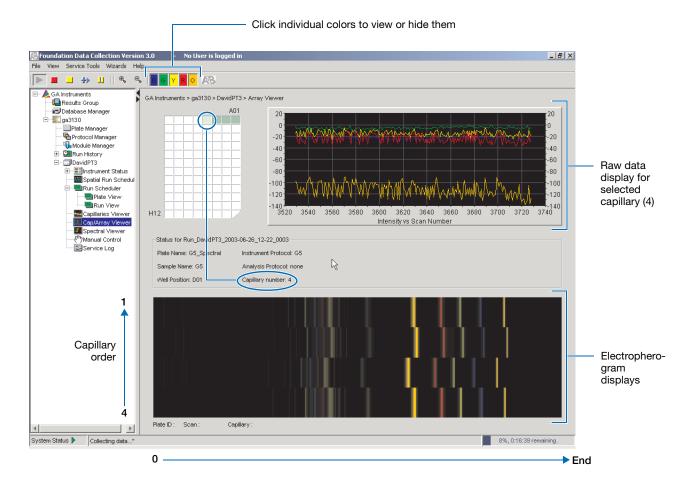

How to Zoom

To zoom in and out:

- **1.** Select a rectangular area over the area of interest by holding down the mouse button. Release the mouse button to zoom in.
- **2.** Click to return to full view.



Working with Data in The Run History View

Run History Components

Elements of the Run History Utility

Use the Run History utility only with completed runs stored in the local 3130/3130xl genetic analyzer Data Collection database. It does not provide real-time viewing of collecting runs.

In the left tree pane, click the icon next to the function to launch it.

Elements Within the Run History Utility

Icon	Element	Displays the
<u>uų</u>	EPT Viewer	Electrophoresis data of the run, such as voltage, current, and oven temperature profiles.
	Event Log	Event messages associated with the run.
4	Instrument Protocol	Instrument protocol and run module settings used for the run.
747	Spatial Calibration Viewer	Spatial calibration associated with the run.
<u> </u>	Capillaries Viewer	Individual electropherogram for each sample in the run.
****	Cap/Array Viewer	Array view of the run.
I	Spectral Viewer	Spectral calibration used for the run.
1	Reextraction	Extraction and analysis status for all samples in the run. See "Viewing Autoextraction Results" on page 147 for more details.

Note: If Cleanup Database has been used, you cannot view processed data in Run History.

Notes		


Viewing Data from a Completed Run in the Data Collection Software

View data using two formats within the 3130/3130xl genetic analyzer Data Collection software under the Run History icon:

- In the Cap/Array Viewer window (in much the same way that you might view the gel file output from an ABI PRISM® slab gel instrument).
- In the Capillary Viewer window, capillary-bycapillary.

Viewing Data from a Completed Run

- In the tree pane of the Data Collection software, click ▲ GA Instruments > ga3130xl or ga3130 > Run History to select the run you want to view.
- **2.** Search for your run by either Barcode or Advanced search.
- **3.** After choosing the run, click the run history elements, such as Cap/Array Viewer or the Capillary Viewer from the left tree pane to look at the run data.

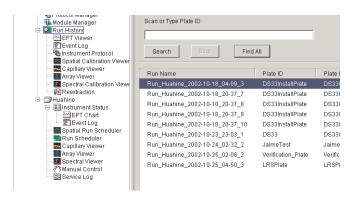

Viewing Autoextraction Results

Overview

After a run is completed, extraction and analysis is performed automatically, using the settings in the Plate Editor and the Results Group. View the results of extraction and analysis in the Reextraction Panel. Reextract samples with the same settings, or with different Analysis Protocols or Results Groups. This process can be useful for many reasons:

- The destination location may not have been available during extraction.
- Some samples may have failed analysis and a different Analysis Protocol might be more successful.
- Samples might be saved in different locations, or with no analysis at all to save space.
- Sample names may need to be edited.

Sample File Destinations

Locate the Results Group's Naming Tab and check the Destination Tab and Run Folder Name Format to find the sample file destination.


Effects of Changes Made in the Reextraction Panel Changes made in the Reextraction Panel to a Results Group, Analysis protocol, sample names, or Comments, also change in the original plate record. The original plate information is overwritten.

Notes		

Selecting and Queuing Samples for Extraction

Queue individual samples for reextraction to experiment with different Analysis Protocols for samples that have failed initial extraction.

- **1.** Click [(Run History).
- **2.** Enter the plate name for a plate that has been completed, or click **Search**. Plates that have runs still pending cannot be reextracted. All the runs from that plate appear in the window.
- **3.** Select a run from the list.

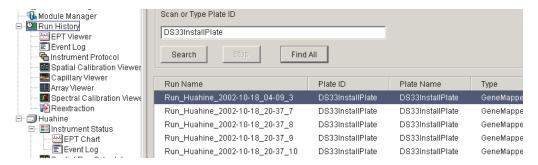

- **4.** Click (Reextraction) in the left tree pane to display the Reextraction window.
- **5.** Click the check boxes in the Extract column to select the samples to be reextracted.
- **6.** Click **Extract** to start the reextraction.

Note: Reextracted sample files are saved in the original folder that data was extracted to unless this destination was changed through the Results Group.




Elements of the Reextraction Window

All the samples are displayed with the results of extraction and analysis.

Note: Sort the columns of the re-extraction panel by holding the shift key and then clicking on a column header.


Reextraction Window for Sequencing Analysis

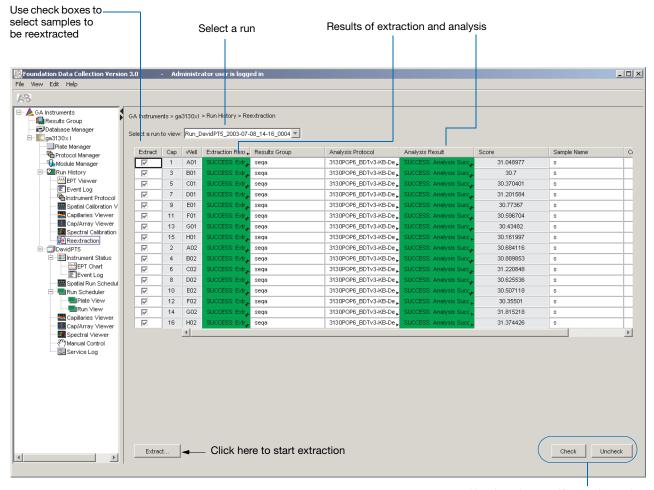

Use these buttons if several samples are highlighted

B. II		

Reextraction Window for Fragment Analysis

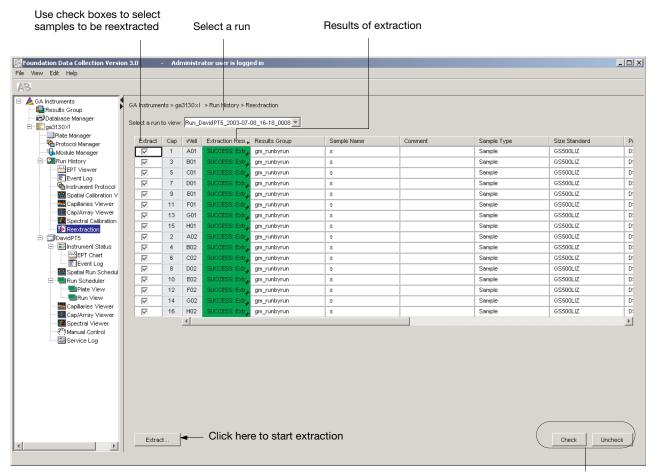

Use these buttons if several samples are highlighted

Notes_____


Results Column

The results of extraction and analysis are color coded in the Results column. The following table lists the colors and their values for Sequencing Analysis.


Color	Value
Red	Extraction or analysis failed
	Note: Descriptive messages can be viewed by resizing the Results column to view all text.
Yellow	Warnings for extraction or analysis
	Note: Descriptive messages can be viewed by resizing the Results column to view all text.
Green	Successful extraction (with no analysis intended), or successful extraction and analysis.

The Results column, by default, shows only the beginning of any processing message. The entire message and the sample file location can be viewed by expanding the cell. There is a tooltip view for each sample results message.

Tooltip view. Access by placing the cursor over the sample of interest

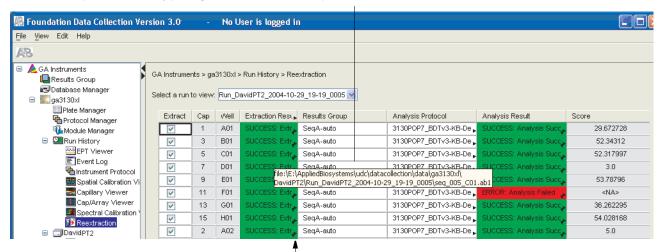

Drag the cell's edge to expand the column

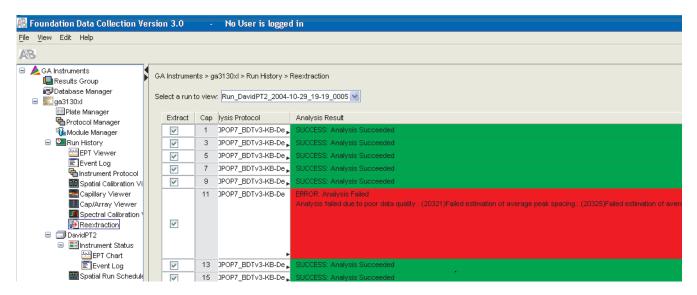

Expanded column

Quality Column

The Quality column represents the quality values for an entire sequence. Quality Values are only assigned to analyzed samples when using the KB^{TM} Basecaller.

Colors displayed and their associated value range

Color	Quality Value Range		
Red	< 15		
Orange	≥ 15 and < 20		
Yellow	≥ 20 and < 30		
Green	> 30		
Note: For more information on KB Basecaller and Quality Values, see the Applied			

The column is empty (white) if:

- Analysis was not performed
- · Analysis failed
- ABI Basecaller was used for analysis. This basecaller does not assign Quality Values.

Biosystems DNA Sequencing Analysis Software v5.1 User Guide, PN 4346366.



Results Group and Analysis Protocol Columns

You can edit and make changes for reextraction to the Results Group and the Analysis Protocol (Analysis Method in the GeneMapper® software).

Note: Select an entire column in the Reextraction window by clicking on the column header. For example, clicking on the Extract column header selects all samples. Clicking the Uncheck or Check buttons at the bottom of the window, enables or disables the check boxes for each sample. The filldown command (Ctrl+D) works the same here as in the Plate Editor for easier information input.

Sorting the Samples

- **1.** Hold down the shift key while clicking on the column header to sort samples according to the column properties.
- **2.** Shift-click again to sort them in reverse order. Sort by:
 - Capillary number
 - Well position
 - Results
 - Quality
 - Extract column

Note: For example, bring all of the failed analysis or extraction samples to the top of the column. Now they can be viewed without having to scroll down to each sample individually.

Notes			

Reextracting Selected Samples

- **1.** Expand the Results column cells for any yellow or red results to see a description of the warning or failure.
- 2. If desired, select a new Results Group, or edit the current one. Selecting a new Results Group allows you to turn off autoanalysis, change the samples and folder naming options, the location where they are placed, and the owner of the Results Group.
- **3.** If desired, change the Analysis Protocol to experiment with different ways of analyzing the sample, using a different basecaller for example.
- **4.** Check the check box in the Extract column for the samples you wish to extract again.
- 5. Click Extract.

Note: Reextraction creates an entirely new sample file and does not replace the previously saved sample file. The presence of a previous sample file has no effect on the creation of a new sample file. If the same naming options that are used for reextraction are identical to those used previously, a number is appended to the filename. For example, if the first sample is, "sample 01.ab1" then the second sample would be, "sample 01 (1).ab1."

Notes			

Using Autoanalysis Manager Software

Use Autoanalysis Manager software with SeqScape and/or GeneMapper software. The Autoanalysis Manager software is installed by the SeqScape or GeneMapper software installation CD.

Autoanalysis occurs in the following sequence:

- **1.** When data collection software finishes a run, the Message Service sends the message "Run Completed."
- **2.** The Autoanalysis Manager receives the message, and the job is submitted. The job appears in the General tab.
- **3.** The Autoanalysis Manager polls for jobs every 2 minutes and opens the automated processing version of SeqScape and/or GeneMapper software to analyze the data.
- **4.** At the end of analysis, the status in the Autoanalysis Manager is updated.

Files Created

The data collection software stores the sample files in the location specified in the results group. The Autoanalysis Manager copies the files into the Data Store for SeqScape or GeneMapper processing.

To maintain sufficient storage space on your hard drive, delete or transfer unnecessary sample files to a different network or CD.

Autoanalyzed Samples

Note: If your GeneMapper or SeqScape software was opened during autoanalysis, then close your analysis application and relaunch the Autoanalysis Manager.


Once an internal message from the instrument is received by the Autoanalysis Manager, it opens the automated processing version of SeqScape and/or GeneMapper software to autoanalyze the samples. The standard user version of SeqScape and/or GeneMapper software must be closed in order for autoanalysis to begin.

Components

The Autoanalysis Manager has two or three tabs:

- General tab
- GeneMapper tab, if GeneMapper v3.7 or GeneMapper ID v3.2 software is installed
- SeqScape tab, if SeqScape software is installed

General Tab The General tab shows the jobs that have been submitted and their status.

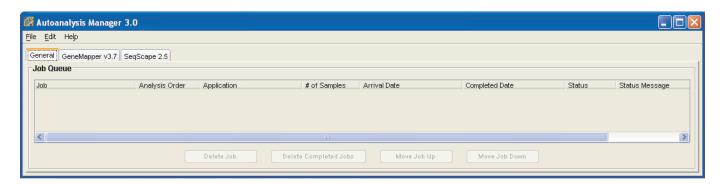

Command Buttons in the General Tab




Command Button Functions in the General Tab

Button Name	Function
Delete Job	Deletes an individual job/project from the Autoanalysis Manager list.
	Does not delete sample files, SeqScape or GeneMapper software project.
Delete Completed Jobs	Deletes all completed jobs/projects from the Autoanalysis Manager list. Only successful jobs are deleted.
	Does not delete sample files, SeqScape or GeneMapper software projects.
Move Job Up	The active job/project is always given a queue number of 1. Once
Move Job Down	job 1 is finished analyzing, job 2 becomes job 1 and all other numbers are changed accordingly. Use the Move Up/Down buttons if you want to rearrange the analysis order.

Notes____

GeneMapper Software Tab

The GeneMapper software tab shows the jobs that have been submitted and their status.

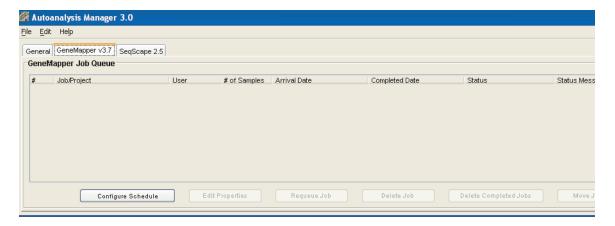

Command Buttons in the GeneMapper Software Tab



Command Button Functions in the GeneMapper Software Tab

Button Name	Function
Configure Schedule	Next Analysis Time:
	Enables you to set a start time for autoanalysis. Before this time arrives, autoanalysis of projects will not occur.
	Periods restricting automated analysis:
	Enables you to set times during which autoanalysis will not occur. Useful if you know that you are going to be reviewing data during a certain time period and don't want to be bothered by the "Runs ready for processing" dialog box. Runs build up in the queue until the restricting time period is over at which time the runs will be autoanalyzed.
	Automatic Deletion of Completed Jobs:
	Enables you to set the software to automatically delete successfully completed jobs. Jobs that failed or have not been analyzed will not be deleted. Only the Autoanalysis Manager job is deleted, sample files and GeneMapper software projects are not.
Edit Properties	Enables you to change the following settings:
	Job/Project Name
	UserName: GeneMapper software UserName
	 Password: Matching password for GeneMapper software UserName
	 Queue position: Enter a new queue position number for the project


Notes			

Button Name	Function
Requeue Job	Samples that need to be autoanalyzed have queue numbers listed in the # column.
	 Samples that are already analyzed or failed, have a blank cell in the # column.
	To resubmit a job for autoanalysis, use the Requeue Job button to assign a queue number to that job.
Delete Job	Deletes an individual job/project from the AutoAnalysis Manager list. Does not delete sample files or GeneMapper software project.
Delete Completed Jobs	Deletes all completed jobs/projects from the AutoAnalysis Manager list. Only successful jobs are deleted. Does not delete sample files or GeneMapper software projects.

ΝΙ	_	4	_	_
ΙVΙ	\cap	Т	е	S
	_	•	_	_

SeqScape Software Tab

The SeqScape software tab shows the jobs, project, and status information.

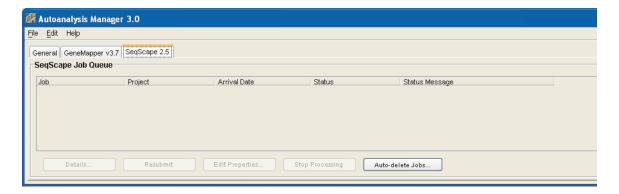

Command Button Functions in the SeqScape Software Tab

Button Name	Function	
Details	Displays the project in the navigation pane	
	Details for Job "Run_SSIntegration_2002-08-23_16-59_1 - SeqScape_Results_Gr □ □ HLA-C-3100 □ □ 360.2 □ seq_005_F01_10301 □ seq_007_E01_10301 □ seq_011_C01_10301 □ seq_013_B01_10301 □ seq_013_B01_10301	
Resubmit	Submits a job for analysis	
Edit Properties	Edits the name and password (active only if analysis failed)	
Delete	Deletes a job from the Autoanalysis Manager	


Viewing Analyzed Data

Locating Sample Files

Analyzed sample files are extracted into a run folder defined in the Naming tab of your Results Group, and placed in a location defined in the Destination tab.

The default location is:

E:\AppliedBiosystems\UDC\Data Collection\Data\ <instrument type>\<instrument name>\<Run Name>

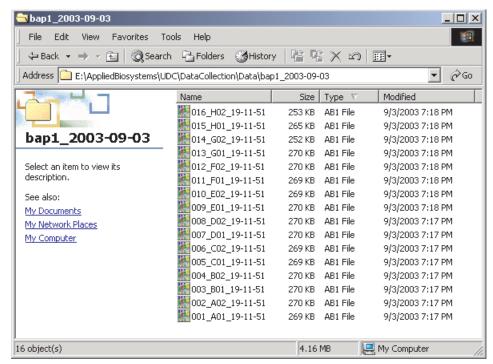

Locating Sample Files

Locate the reextracted data in the location defined by the Results Group or the default destination location.

Viewing Sample Files

Use the Sequencing Analysis, SeqScape, or the GeneMapper Software to view reextracted sample electropherogram data, both raw, and analyzed.

Sample file type	File extension
Sequencing, SeqScape	.ab1
Fragment analysis	.fsa

Plate Mapping

Injection Scheduling

Samples are scheduled for injection based on their position in a 96-or 384-well plate and the capillary number. Study the appropriate plate mapping patterns described in "96-Well Plate Mapping" or "384-Well Plate Mapping" to arrange your samples according to your preferred injection priority.

The injection schedule system is based on the:


- Order of the linked plates (3130xl instrument only). See page 128 for linking plate information.
- Sample priority value in the plate record.

If all priorities are set to 100 (default), then runs are scheduled as outlined below. See the *Applied Biosystems 3130/3130xl Genetic Analyzers Maintenance, Troubleshooting, and Reference Guide* on "Run Priority Scheduling".

Capillary Array Map

The capillary numbers in the capillary array are shown in the layout below. The capillary array layout is the same for both 96- and 384-well plates.

Capillary Number Layout

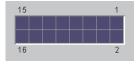




96-Well Plate Mapping

3130xl Genetic Analyzer


For a 96-well plate, injections are made from every well in two consecutive rows, starting with an odd row. A full 96-well plate requires six runs to inject all samples. See the examples below where samples (grey) and capillary number (number) positions can affect your number of runs.

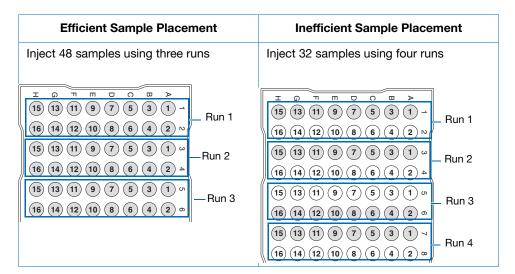

3130 Genetic Analyzer

For a 96-well plate, injections are made from four consecutive wells in a row. A full plate of 96 sample requires 24 runs to inject all samples once.

In the following example of a 96-well plate, the gray circles represent samples and the number in the well indicates capillary number. It takes four runs to inject 16 samples.

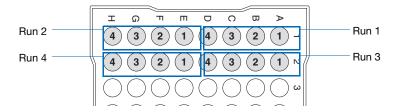


384-Well Plate Mapping

3130xl Genetic Analyzer

For a 384-well plate, injections are made from every other well and every other row. A full plate of 384 samples requires 24 runs to inject all samples once.


The example below is the injection pattern for the first four injections, starting with well A01. The light gray circles represent samples and the dark gray circles indicate the injection pattern.

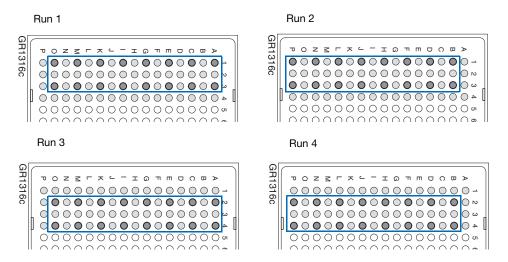

3130 Genetic Analyzer

For a 384-well plate, injections are made from every other well. A full plate of 384 sample requires 96 runs to inject all samples once.

The example below is the injection pattern for the first six injections, starting with well A01. The light gray circles represent samples and the dark gray circles indicate the injection pattern.

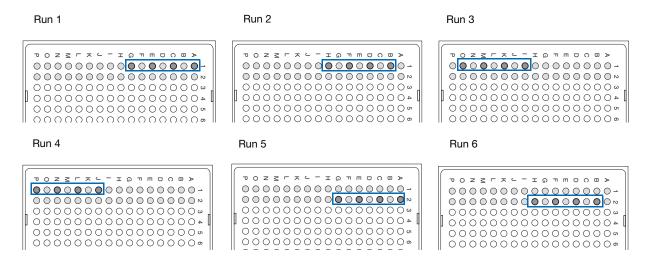

384-Well Plate Mapping		
Notes		

Appendix A

Index

Numerics	GeneMapper Software, computer name 107
3130 basecaller and DyeSet/Primer files 66, 67	manually, fragment analysis documentation 106 manually, Seqscape software documentation 83
	manually, sequencing documentation 54
A	sequencing 54
ABI PRISM® BigDye® Primer Cycle Sequencing Kits	Autoanalysis Manager
table, dye sets and standards 29	components 155
ABI PRISM® dGTP BigDye® Terminator Cycle	explained 82 general tab 156
Sequencing Ready Reaction Kit, table of dye	starting 131
set and standards 29	using 155
ABI PRISM® dGTP BigDye® Terminator v 3.0 Cycle	autoextraction
Sequencing Ready Reaction Kit	overview 147
table of dye set and standards 29	Automated Processing tab 118
ABI PRISM® dRhodamine Terminator Cycle Sequencing	automated SeqScape analysis 82
Kit	automatic analysis, before you start 83
table of information 29	autosampler 18
Activating spectral calibration 47	function 11
AmpFISTR kits matrix standard 30	
table of dye sets, filter, modules 5	В
analysis protocol	barcode reader 6
deleting 68, 97	barcode, type of search 40
editing 68, 96	basecaller and DyeSet/Primer files 66
export 69, 97	basecaller files 66, 67
general description 55 import 69, 97	BigDye® Terminator chemistry, information tables 66
options 58	BigDye® Terminator v3.1 Cycle Sequencing Kit
options, SeqScape software 87	table, dye set and standard 29
analysis, manual 54, 83, 106	biohazardous waste, handling xx
analyzing GeneMapper Software data 160	bold text, when to use ix
anode buffer reservoir	buffer
filling 19	hazards 32
function 11	making 16
Applied Biosystems	buffer fill-line 12
contacting xi	buffer reservoir, anode, illustration of 12
customer feedback on documentation x Information Development department x	buffer valve pin, illustration of 12
Technical Support xi	• ,
array length, polymer type, table 36	C
array port, illustration of 12	calibration standards
assembling the plate 34	types, fragment analysis 30
assumptions for using this guide ix	types, sequencing 29
Australian EMC standards xxiii	Canadian safety standards xxiii
autoanalysis	cap/array viewer 144
	capillaries viewer 143

capillary array function 11	viewing analyzed sequencing data 160 viewing complete run 146
illustration of 12 installing or replacing 13	Data Collection software starting 9
capillary array knob and tip, illustration of 12	data collection software, starting 9
capillary array length	destinations 147
polymer type, table 36	detection cell block and heater
cathode buffer reservoir, filling 17	function 11
CAUTION	documentation
description xiv example xv	GeneMapper Software User Guide (PN) 106 related x
Change Polymer Type Wizard, using 15	double-tapered ferrule, illustration of 12
characters	dye primer chemistry, information table 67
acceptable 54, 83, 106	dye set
check valve diagram of 12	fragment analysis 30, 114 table, sequencing 29
chemical safety xviii	Dye Terminator chemistry, files 66
chemical safety guidelines xix	DyeSet/Primer files 66, 67
chemical waste	list of 64, 94
hazards xx safety xx	E
safety guidelines xx	E dye set, sequencing kit table 61
chemical waste disposal, guidelines xx	electrical hazard symbol xvi
computer	electrical safety xxi
assigning a unique name 9 starting 7	electrical shock hazards xxi
condition number, spectral calibration 42	
conventions 42	electrical symbols, on instruments xv
bold text ix	electrode, illustration of 12
IMPORTANTS! ix	electromagnetic compatibility standards. See EMC
in this guide ix	standards
italic text ix	electropherogram, definition 143
menu commands ix Notes ix	EMC standards xxiii
user attention words ix	Australian xxiii Canadian xxiii
conventions, safety xiv	European xxiii
creating	EPT chart 141
analysis protocol for autoanalysis 62, 93	ergonomic safety xxii
GeneMapper Software plate record for	errors box 140
autoanalysis 123	European EMC standards xxiii
SeqScape Software plate record 101	European safety standards xxiii
Sequencing Analysis plate record for autoanalysis 77	event log 142
customer feedback, on Applied Biosystems	icon, displays 145
documents x	events, box 140
D	F
DANGER	file extension 160
description xiv	file naming, invalid characters 54
example xv	fill down special, function 38
data	fill line, anode buffer reservoir 19
how to zoom 143 viewing 137	filter set

fragment analysis 5 find all plate search 40	physical xxi repetitive motion xxii
•	repetitive motion - xxii
fragment analysis	1
data analysis 160	1
filter set 5	illustration of 12
kit types 5 passing profiles, spectral calibration 51	IMPORTANT
run modules 113	description xiv
run, average time 133	example xv
function 11	IMPORTANTS!, description ix
Tunction 11	Information Development department, contacting x
G	Install Array Wizard
GeneMapper Software	using 13
computer name 107	installation category xxi
elements 109	instrument
generic 107	condition, status 140
instrument protocol 108	diagram of interior 11
plate records 108	documentation x
results group 108	EPT status 141
GeneMapper Software tab 157	layout 2 operation 130
GeneMapper TM Software and GeneMapper TM ID	setup 12
Software 106	starting the system 8
general hazard symbol xvi	instrument operational safety, instructions for xviii
guidelines	instrument protocol
chemical safety xix	creating 35
chemical waste disposal xx	creating for SeqScape software 88
chemical waste safety xx	creating, for fragment analysis 109, 112
waste disposal xx	general description 55
	icon, displays 145
H	options, SeqScape software 87
hard drive space, outcometic checking, 122	options,s equencing 58
hard drive space, automatic checking 132	instrument status 139
hazard icons	interconnect tube 12
accompanying safety alert words xiv	invalid characters in names 54, 83, 106
components xiv	
described xiv	italic text, when to use ix
in documents xiv on instruments xiv, xvi	
See also hazard symbols	K
See also safety symbols	KB basecalling run modules 66
hazard symbols	C
electrical xvi	1
general xvi	<u> </u>
hot surface xvi	labels, instrument safety xvi
in documents xv	laser hazard xxi
laser hazard xvi	symbol xvi
moving parts xvi	warnings xxii
on instruments xvi	laser safety xxi
See also hazard icons	laser classification xxi
See also safety symbols	requirements xxii
hazards	Launching Data Collection software 9
chemical waste xx	layout, instrument 2
electrical shock xxi	length of read, definition 3
laser xxi	
moving parts xxi moving/lifting instrument xvii	linking a plate 128

LOR definition 3	P
lower polymer block 11	pausing a run 133
lower polymer block, illustration of 12	PDP motor cover, illustration of 12
Luer fitting, illustration of 12	physical hazard safety xxi
	physical hazards xxi
M	piston, illustration of 12
manual analysis, fragment analysis documentation 106	-
manual analysis, Seqscape software documentation 83	plate assembly diagram 34
manual analysis, sequencing documentation 54	placing in instrument 39
matrix standard	plate editor, filling in 37
fragment analysis 30	plate manager, access 37
matrix standards	plate record
fragment analysis 114	creating 37
preparing for sequencing 31 sequencing 29	creating for GeneMapper Software autoanalysis 123
menu commands, conventions for describing ix	creating for SeqScape Software 101
mixed plate record 55	creating for Sequencing Analysis autoanalysis 77
modules	discussed 55, 83, 108 elements, sequencing analysis 56
fragment analysis 4, 5	linking and unlinking 128
KB basecalling 66	mixed 55
sequencing 3	search 40
spectral 36	plate view 137
mounting pin, illustration of 12	plates
moving and lifting	linking and unlinking from plate records 128
computer xviii monitor xviii	sealing 33
moving and lifting instrument, safety xvii	polymer
moving parts	array length, table 36 changing to new type 15
hazard xxi	replenishing or changing 14
hazard symbol xvi	when to add or change 14
safety xxi	polymer delivery pump
MSDSs	diagram 12
description xviii	polymer supply bottle cap with hole, illustration of 12
obtaining xix referring to xix, xx	polymer supply bottle, illustration of 12
when to review xviii	polymer supply tube illustration of 12
	protocol manager, access 35
N	pump block 12 function 11
Naming	pump chamber, illustration of 12
tab, fragment analysis 118	
Notes, description ix	Q
0	quality column 152
O-ring, illustration of 12	quality value range 152
oven	Q-value, spectral calibration 42
function 11	
overflow hole, illustration of 12	R
overvoltage category (rating) xxi	radioactive waste, handling xx
overvoltage rating xxi	reader, barcode 6

reextraction	conventions xiv		
editing Results Group, Analysis Protocol, Analysis	electrical xxi		
Method 153	ergonomic xxii		
results column 151	instrument xvii		
selected samples 154	instrument operation xviii		
selecting 148	laser xxi		
window elements 149	moving and lifting instrument xvii		
reextraction panel	moving parts xxi physical hazard xxi		
effects of changes made in 147	standards xxiii		
repetitive motion hazard xxii	workstation xxii		
Replenish Polymer Wizard			
using 15	safety alert words		
reservoirs	accompanying hazard icons xiv CAUTIONS xiv		
filling 17	DANGERS xiv		
positions on the autosampler 18	description xiv		
resolution	IMPORTANTS xiv		
fragment analysis 4	WARNINGS xiv		
	safety labels, on instruments xvi		
results column, reextraction 151	-		
Results Group	safety requirements, laser xxii		
creating for sequencing analysis 70	safety standards		
creating, fragment analysis 115	Canadian xxiii		
creating, SeqScape software 98	European xxiii		
editor, GeneMapper 107	U.S. xxiii		
general description 55	safety symbols		
importing and exporting 100 options 57	on instruments xvi		
options, SeqScape software 165	See also hazard symbols		
	sample file 147		
run	default location 160		
fragment analysis 133	type 160		
requires plate record 55, 84, 108 sensor states 140	samples		
sequencing, average time 132	autoanalyzing 155		
starting 130	loading 32		
starting 150 starting, stopping, skipping, pausing 133	sensor states 140		
viewing data 137	SeqScape software tab 159		
run history utility 145	SeqScape Software v2.5		
run modules	autoanalysis 82		
	plate editor 84		
automated steps, description 132 selecting for fragment analysis 113	plate record elements 85		
selecting for fragment analysis 113	user guide P/N 82		
run schedule, viewing 129	sequencing		
_	automated, creating protocols 59		
run scheduler	automation, required files 55		
spatial 22	data, automating the analysis 82		
verifying schedule 129	matrix standards 31		
run time	passing profiles, spectral calibration 49		
fragment analysis 4	plate editor 55		
spectral calibration 41	run time 132		
running buffer, making and storing 16	specifications 3		
	spectral run modules 36		
S	viewing analyzed data 160		
	sequencing chemistry		
safety	supported kits 28		
before operating the instrument xvii	service console, using 9		
chemical xviii chemical waste xx	set (activate) spectral calibration 47		
CHCHICAL WASIC AA	or (see and operation of		

software, starting 9	U
spatial calibration	unlinking a plate 129
accepting or rejecting 24	
creating 22	Update Cap Array Info
evaluating profiles 23	using 13
examples of profiles 25	US safety standards xxiii
viewer 145	user attention words, described ix
when required 22	utility, run history 145
specifications	
fragment analysis 4 sequencing 3	V
	viewer
spectral calibration	cap/array 144
creating plate record 37 evaluating results 43	capillaries 143
pass/fail status 42	icons and display information 145
passing profiles, fragment analysis 51	spectral 145
passing profiles, sequencing 49	•
performing a 37	W
preparing standards for 31	
run times 41	WARNING, description xiv
set spectral calibration 47	waste disposal, guidelines xx
when to perform 28	water seal, illustration of 12
spectral calibration standard, types 29	water trap
spectral viewer 145	illustration of 12
spectral viewer, blank or deactivated 47	wizard
standards	Replenish Polymer Wizard, using 15
EMC xxiii	Update Cap Array Info, using 13
safety xxiii	workstation safety xxii
starting	
run 133	Z
spectral calibration run 40	7 due set seguencing kit tehle 61
status	Z dye set, sequencing kit table 61
instrument 139	
service console 9	
status lights	
on instrument startup 8	
stopping a run 133	
supported sequencing chemistry kits 28	
symbols	
hazard xvi	
hot surface xvi	
symbols on instruments	
electrical xv	
safety xvi	
-	
T	
tables	
summary 3	
Technical Support, contacting xi	
text conventions ix	
toolbar 133	
training, information on xi	

Headquarters

850 Lincoln Centre Drive Foster City, CA 94404 USA Phone: +1 650.638.5800

Toll Free (In North America): +1 800.345.5224

Fax: +1 650.638.5884

Worldwide Sales and Support

Applied Biosystems vast distribution and service network, composed of highly trained support and applications personnel, reaches 150 countries on six continents. For sales office locations and technical support, please call our local office or refer to our Web site at www.appliedbiosystems.com.

www.appliedbiosystems.com



Applied Biosystems is committed to providing the world's leading technology and information for life scientists.

6/2010 Part Number 4352715 Rev. D