

TSQ Series

Hardware Manual

70111-97163 Revision E June 2014



Thermo
SCIENTIFIC

© 2014 Thermo Fisher Scientific Inc. All rights reserved.

Foundation, Ion Max, TSQ, TSQ Quantum Access, TSQ Quantum Access MAX, TSQ Vantage, TSQ Vantage EMR, TSQ Vantage AM, TSQ Quantum Ultra, TSQ Quantum Ultra EMR, and TSQ Quantum Ultra AM are trademarks, and Accela, TSQ Quantum, and Thermo Scientific are registered trademarks of Thermo Fisher Scientific Inc.

The following are registered trademarks in the United States and possibly other countries: Swagelok is a registered trademark of the Crawford Fitting Company. Tygon is a registered trademark of Saint-Gobain Performance Plastics Company. Restek is a registered trademark of Restek Corporation. Microsoft and Windows are registered trademarks of Microsoft Corporation. Convectron and Granville-Phillips are registered trademarks of Brooks Automation, Inc. Intel and Pentium are registered trademarks of Intel Corporation. KEL-F is a registered trademark of 3M Company. Rheodyne and Upchurch Scientific are registered trademarks of IDEX Health Science, LLC. Kalrez and Viton are registered trademarks of DuPont Performance Elastomers LLC. Delrin, Teflon, and Tefzel are registered trademarks of E.I. du Pont de Nemours & Co. Alconox is a registered trademark of Alconox, Inc. Gastight is a registered trademark of Hamilton Company. Kimwipe is a registered trademark of Kimberly-Clark Corporation.

PEEK is a trademark of Victrex plc. Edwards is a trademark of Edwards Limited.

All other trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.

Thermo Fisher Scientific Inc. provides this document to its customers with a product purchase to use in the product operation. This document is copyright protected and any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of Thermo Fisher Scientific Inc.

The contents of this document are subject to change without notice. All technical information in this document is for reference purposes only. System configurations and specifications in this document supersede all previous information received by the purchaser.

This document is not part of any sales contract between Thermo Fisher Scientific Inc. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale, which Terms and Conditions of Sale shall govern all conflicting information between the two documents.

Release history: Revision A, February 2008; Revision B, June 2008; Revision C, March 2009; Revision D, October 2010; Revision E, June 2014.

Software version: (Thermo) Foundation 2.0 and later, TSQ 2.5 and later

For Research Use Only. Not for use in diagnostic procedures.



Regulatory Compliance

Thermo Fisher Scientific performs complete testing and evaluation of its products to ensure full compliance with applicable domestic and international regulations. When the system is delivered to you, it meets all pertinent electromagnetic compatibility (EMC) and safety standards as described below.

Changes that you make to your system might void compliance with one or more of these EMC and safety standards. Changes to your system include replacing a part or adding components, options, or peripherals not specifically authorized and qualified by Thermo Fisher Scientific. To ensure continued compliance with EMC and safety standards, replacement parts and additional components, options, and peripherals must be ordered from Thermo Fisher Scientific or one of its authorized representatives.

- [TSQ Quantum Access MAX](#)
- [TSQ Quantum Ultra](#)
- [TSQ Vantage](#)
- [TSQ Quantum Access](#)

TSQ Quantum Access MAX

EMC Directive 2004/108/EEC

EMC compliance has been evaluated by TUV Rheinland of North America, Inc.

EN 55011: 2007, A2: 2007	EN 61000-4-4: 2004
EN 61000-3-2: 2006	EN 61000-4-5: 2005
EN 61000-3-3: 1995, A1: 2001, A2: 2005	EN 61000-4-6: 2007
EN 61000-4-2: 1995, A1: 1999, A2: 2001	EN 61000-4-11: 2004
EN 61000-4-3: 2006	EN 61326-1: 2006

FCC Class A: CFR 47 Part 15: 2007

Low Voltage Safety Compliance

This device complies with European Union Directive 2006/95/EC implemented by 61010-1: 2001.



TSQ Quantum Ultra

EMC Directive 89/336/EEC as amended by 92/31/EEC and 93/68/EEC

EMC compliance has been evaluated by Underwriters Laboratories Inc.

EN 55011: 1998	EN 61000-4-4: 1995, A1: 2001, A2: 2001
EN 61000-3-2: 1995, A1: 1998, A2: 1998, A14: 2000	EN 61000-4-5: 1995, A1: 2001
EN 61000-3-3: 1998	EN 61000-4-6: 2001
EN 61000-4-2: 2000	EN 61000-4-11: 1994, A1: 2001
EN 61000-4-3: 2002	EN 61326-1: 1998

FCC Class A, CFR 47 Part 15: 2005

CISPR 11: 1999, A1: 1999, A2: 2002

Low Voltage Safety Compliance

This device complies with Low Voltage Directive 73/23/EEC and harmonized standard EN 61010-1: 2001.

TSQ Vantage

EMC Directive 2004/108/EC

EMC compliance has been evaluated by TUV Rheinland of North America, Inc.

EN 55011: 1998, A1: 1999, A2: 2002	EN 61000-4-4: 2004
EN 61000-3-2: 2006	EN 61000-4-5: 2006
EN 61000-3-3: 1995, A1: 2001, A2: 2005	EN 61000-4-6: 2001
EN 61000-4-2: 2001	EN 61000-4-11: 2004
EN 61000-4-3: 2006	EN 61326-1: 2006

FCC Class A, CFR 47 Part 15: 2007

CISPR 11: 1999, A1: 1999, A2: 2002

Low Voltage Safety Compliance

This device complies with the Low Voltage Directive 2006/95/EC and harmonized standard EN 61010-1.

TSQ Quantum Access

EMC Directive 89/336/EEC, 92/31/EEC, 93/68/EEC

EMC compliance has been evaluated by TUV Rheinland of North America, Inc.

EN 55011: 1998, A1: 1999, A2: 2002

EN 61000-3-2: 2000

EN 61000-3-3: 1995, A1: 2001

EN 61000-4-2: 2001

EN 61000-4-3: 2002

EN 61000-4-4: 1995, A1: 2000, A2: 2001

EN 61000-4-5: 2001

EN 61000-4-6: 2003

EN 61000-4-11: 2001

EN 61326: 1997, A1: 1998, A2: 2001, A3: 2003

FCC Class A, CFR 47 Part 15: 2005

CISPR 11: 1999, A1: 1999, A2: 2002

Low Voltage Safety Compliance

This device complies with the Low Voltage Directive EN 61010-1:2001 and harmonized standard EN 61010-1: 2001.

FCC Compliance Statement

THIS DEVICE COMPLIES WITH PART 15 OF THE FCC RULES. OPERATION IS SUBJECT TO THE FOLLOWING TWO CONDITIONS: (1) THIS DEVICE MAY NOT CAUSE HARMFUL INTERFERENCE, AND (2) THIS DEVICE MUST ACCEPT ANY INTERFERENCE RECEIVED, INCLUDING INTERFERENCE THAT MAY CAUSE UNDESIRE OPERATION.

Notice on Lifting and Handling of Thermo Scientific Instruments



CAUTION Read and understand the various precautionary notes, signs, and symbols contained inside this manual pertaining to the safe use and operation of this product before using the device.

For your safety, and in compliance with international regulations, the physical handling of this Thermo Fisher Scientific instrument **requires a team effort** to lift and/or move the instrument. This instrument is too heavy and/or bulky for one person alone to handle safely.



Notice on the Proper Use of Thermo Scientific Instruments

In compliance with international regulations: Use of this instrument in a manner not specified by Thermo Fisher Scientific could impair any protection provided by the instrument.

Notice on the Susceptibility to Electromagnetic Transmissions

Your instrument is designed to work in a controlled electromagnetic environment. Do not use radio frequency transmitters, such as mobile phones, in close proximity to the instrument.

For manufacturing location, see the label on the instrument.

WEEE Compliance

This product complies with the European Union's Waste Electrical & Electronic Equipment (WEEE) Directive 2002/96/EC. It is marked with the following symbol:



Thermo Fisher Scientific is registered with B2B Compliance ([B2Bcompliance.org.uk](https://www.b2bcompliance.org.uk)) in the UK and with the European Recycling Platform ([ERP-recycling.org](https://www.eur-ecycling.org)) in all other countries of the European Union and in Norway.

If this product is located in Europe and you want to participate in the Thermo Fisher Scientific Business-to-Business (B2B) Recycling Program, send an email request to weee.recycle@thermofisher.com with the following information:

- WEEE product class
- Name of the manufacturer or distributor (where you purchased the product)
- Number of product pieces, and the estimated total weight and volume
- Pick-up address and contact person (include contact information)
- Appropriate pick-up time
- Declaration of decontamination, stating that all hazardous fluids or material have been removed from the product

For additional information about the Restriction on Hazardous Substances (RoHS) Directive for the European Union, search for RoHS on the Thermo Fisher Scientific European language websites.

IMPORTANT This recycling program is **not** for biological hazard products or for products that have been medically contaminated. You must treat these types of products as biohazard waste and dispose of them in accordance with your local regulations.

Conformité DEEE

Ce produit est conforme avec la directive européenne (2002/96/EC) des Déchets d'Equipements Electriques et Electroniques (DEEE). Il est marqué par le symbole suivant:









Thermo Fisher Scientific s'est associé avec une ou plusieurs sociétés de recyclage dans chaque état membre de l'Union Européenne et ce produit devrait être collecté ou recyclé par celle(s)-ci. Pour davantage d'informations, rendez-vous sur la page www.thermoscientific.fr/rohs.






WEEE Konformität

Dieses Produkt entspricht der EU Waste Electrical & Electronic Equipment (WEEE) Richtlinie 2002/96/EC. Es ist mit dem folgenden Symbol gekennzeichnet:



Thermo Fisher Scientific hat Vereinbarungen mit Verwertungs-/Entsorgungsfirmen in allen EU-Mitgliedsstaaten getroffen, damit dieses Produkt durch diese Firmen wiederverwertet oder entsorgt werden kann. Weitere Informationen finden Sie unter www.thermoscientific.de/rohs.

CAUTION Symbol	CAUTION	VORSICHT	PRECAUCIÓN	MISE EN GARDE
	Risk electric shock: This instrument uses voltages that can cause electric shock and/or personal injury. Before servicing, shut down the instrument and disconnect it from line power. While operating the instrument, keep covers on. Do not remove the protective covers from the printed circuit board assemblies (PCBAs).	Stromschlaggefahr: Dieses Gerät arbeitet mit Spannungen, die Stromschläge und/oder Personenverletzungen verursachen können. Vor Wartungsarbeiten muss das Gerät abgeschaltet und vom Netz getrennt werden. Betreiben Sie das Gerät nicht mit abgenommenen Abdeckungen. Nehmen Sie die Schutzabdeckungen von Leiterplatten nicht ab.	Riesgo de descargas eléctricas: Este instrumento utiliza voltajes que pueden causar descargas eléctricas y/o lesiones personales. Antes de revisar o reparar el instrumento, apáguelo y desconéctelo de la red eléctrica. Mantenga colocadas las cubiertas mientras se utiliza el instrumento. No retire las cubiertas protectoras del circuito impreso completo (PCBA).	Risque de choc électrique : l'instrument utilise des tensions susceptibles de provoquer une électrocution et/ou des blessures corporelles. Il doit être arrêté et débranché de la source de courant avant toute intervention. Ne pas utiliser l'instrument sans ses couvercles. Ne pas enlever les capots de protection des cartes à circuit imprimé (PCBA).
	Chemical hazard: Wear gloves and other protective equipment, as appropriate, when handling toxic, carcinogenic, mutagenic, corrosive, or irritant chemicals. Use approved containers and proper procedures to dispose of waste oil and when handling wetted parts of the instrument.	Gefahr durch Chemikalien: Tragen Sie beim Umgang mit toxischen, karzinogenen, mutagenen, ätzenden oder reizenden Chemikalien Schutzhandschuhe und weitere geeignete Schutzausrüstung. Verwenden Sie bei der Entsorgung von verbrauchtem Öl und beim Umgang mit medienberührenden Komponenten die vorgeschriebenen Behälter, und wenden Sie ordnungsgemäße Verfahren an.	Peligro por sustancias químicas: Cuando manipule sustancias químicas, tóxicas, carcinogénicas, mutágenas, corrosivas o irritantes, utilice guantes y otro equipo de protección. Utilice siempre recipientes homologados y siga los procedimientos adecuados cuando deseche aceite residual o manipule partes mojadas del instrumento.	Danger lié aux produits chimiques : porter des gants et d'autres équipements de protection appropriés pour manipuler les produits chimiques toxiques, cancérogènes, mutagènes, corrosifs ou irritants. Utiliser des récipients homologués et des procédures adéquates pour la mise au rebut des huiles usagées et lors de la manipulation des pièces de l'instrument en contact avec l'eau.
	Hot surface: Before touching, allow any heated components to cool.	Heiße Oberflächen: Lassen Sie heiße Komponenten vor der Berührung abkühlen.	Superficies calientes: Antes de tocar los componentes calientes, espere a que se enfrién.	Surface chaude : laisser refroidir les composants chauffés avant toute manipulation.
	Flammable substances hazard: Use care when operating the system in the presence of flammable substances.	Gefahr durch entzündbare Substanzen: Beachten Sie die einschlägigen Vorsichtsmaßnahmen, wenn Sie das System in Gegenwart von entzündbaren Substanzen betreiben.	Peligro por sustancias inflamables: Tenga mucho cuidado cuando utilice el sistema cerca de sustancias inflamables.	Danger lié aux substances inflammables : agir avec précaution lors de l'utilisation du système en présence de substances inflammables.
	Risk of eye injury: Eye injury could occur from splattered chemicals, airborne particles, or sharp objects. (Sharp objects that customers might install in the instrument include fused-silica tubing, the autosampler needle, and so on.) Wear safety glasses when handling chemicals or servicing the instrument.	Augenverletzungsrisiko: Verspritzte Chemikalien, Schwebstoffpartikel oder scharfe Objekte können Augenverletzungen verursachen. (Scharfe Objekte, die Kunden möglicherweise im Gerät installieren, sind z. B. Quarzglas-Kapillaren, die Nadel des Autosamplers, usw.) Tragen Sie beim Umgang mit Chemikalien oder bei der Wartung des Gerätes eine Schutzbrille.	Riesgo de lesiones oculares: Las salpicaduras de sustancias químicas, las partículas flotantes en el aire y los objetos afilados pueden causar lesiones oculares. (Entre los objetos afilados que los clientes pueden instalar en el instrumento se encuentran tubos de sílice fundida, agujas del muestreador automático, etc.). Para manipular sustancias químicas o realizar tareas de mantenimiento, utilice gafas de seguridad.	Risque de lésion oculaire : les projections chimiques, les particules en suspension dans l'air et les objets tranchants peuvent entraîner des lésions oculaires. (Les objets tranchants pouvant être installés par les clients dans l'instrument comprennent les tubes en silice fondue, les aiguilles du passeur automatique, etc.). Porter des lunettes de protection lors de toute manipulation de produit chimique ou intervention sur l'instrument.
	General hazard: A hazard is present that is not included in the other categories. This symbol also appears on the instrument. For details about the hazard, refer to the instrument manual. When the safety of a procedure is questionable, contact Technical Support for Thermo Scientific San Jose products.	Allgemeine Gefahr: Es besteht eine weitere Gefahr, die nicht in den vorstehenden Kategorien beschrieben ist. Dieses Symbol wird auch auf dem Gerät angebracht. Einzelheiten zu dieser Gefahr finden Sie in den Gerätehandbüchern. Wenn Sie sich über die Sicherheit eines Verfahrens im Unklaren sind, setzen Sie sich, bevor Sie fortfahren, mit dem technischen Support für Thermo Scientific San Jose Produkte in Verbindung.	Peligro general: Existen peligros que no se incluyen en las otras categorías. Este símbolo también aparece en el instrumento. Si desea obtener más información sobre estos peligros, consulte el manual del instrumento. En caso de duda sobre la seguridad de un procedimiento, póngase en contacto con el personal de servicio técnico de los productos Thermo Scientific San Jose.	Danger d'ordre général : indique la présence d'un risque n'appartenant pas aux catégories citées plus haut. Ce symbole figure également sur l'instrument. Pour plus de détails sur ce danger potentiel, se reporter au manuel de l'instrument. Si la sûreté d'une procédure est incertaine, contacter l'assistance technique pour les produits Thermo Scientific San Jose.

CAUTION Symbol	CAUTION	VORSICHT	PRECAUCIÓN	MISE EN GARDE
	Laser hazard: This instrument uses a laser that is capable of causing personal injury. This symbol also appears on the instrument. For details about the hazard, refer to the instrument manual.	Gefahr durch Laserstrahlen: Der in diesem Gerät verwendete Laser kann zu Verletzungen führen. Dieses Symbol wird auch auf dem Gerät angebracht. Einzelheiten zu dieser Gefahr finden Sie in den Gerätehandbüchern.	Peligro por láser: Este instrumento utiliza un láser que puede producir lesiones personales. Este símbolo también aparece en el instrumento. Si desea obtener más información sobre el peligro, consulte el manual del instrumento.	Danger lié au laser : l'instrument utilise un laser susceptible de provoquer des blessures corporelles. Ce symbole figure également sur l'instrument. Pour plus de détails sur ce danger potentiel, se reporter au manuel de l'instrument.
	Ultra violet light hazard: Do not look directly at the ultra-violet (UV) light or into the UV source. Exposure can cause eye damage. Wear UV eye protection.	Gefahr durch UV-Licht: Richten Sie Ihren Blick nicht direkt auf ultraviolettes Licht (UV-Licht) oder in die UV-Quelle. Dies kann zu Augenschäden führen. Tragen Sie eine UV-Schutzbrille.	Peligro por luz ultravioleta: No mire directamente a una luz ultravioleta (UV) ni a una fuente UV. La exposición puede causar daños oculares. Lleve protección ocular para UV.	Danger lié aux rayons ultraviolets : ne jamais regarder directement la lumière ultraviolette (UV) ou la source d'UV. Une exposition peut entraîner des lésions oculaires. Porter des protections oculaires anti-UV.
	Sharp object: Avoid physical contact with the object.	Scharfes Objekt: Vermeiden Sie den physischen Kontakt mit dem Objekt.	Objeto puntiagudo: Evite el contacto físico con el objeto.	Objet tranchant : éviter tout contact physique avec l'objet.
	Pinch point: Keep hands away from this area.	Quetschgefahr: Halten Sie Ihre Hände von diesem Bereich fern.	Puntos de pinzamiento: Mantenga las manos apartadas de esta área.	Risque de pincement : éloigner les mains de cette zone.
	Heavy objects: Never lift or move the instrument by yourself; you can suffer personal injury or damage the equipment. For specific lifting instructions, refer to the instrument manual.	Schweres Objekt: Bewegen und heben Sie das Gerät niemals allein an; dies kann zu Verletzungen oder zur Beschädigung des Geräts führen. Spezifische Anweisungen zum Anheben finden Sie im Gerätehandbuch.	Objeto pesado: Nunca levante ni mueva el instrumento por su cuenta, podría sufrir lesiones personales o dañar el equipo. Para obtener instrucciones específicas sobre levantamiento, consulte el manual del instrumento.	Objet lourd : ne jamais soulever ou déplacer l'instrument seul sous peine de blessure corporelle ou d'endommagement de l'instrument. Pour obtenir des instructions de levage spécifiques, se reporter au manuel de l'instrument.
	Trip obstacle: Be aware of cords, hoses, or other objects located on the floor.	Stolpergefahr: Achten Sie auf Kabel, Schläuche und andere Objekte auf dem Fußboden.	Tropiezo con obstáculos: Tenga en cuenta los cables, mangueras u otros objetos colocados en el suelo.	Risque de trébuchement : faire attention aux câbles, tuyaux et autres objets situés sur le sol.
	When the safety of a procedure is questionable, contact Technical Support for Thermo Scientific San Jose products.	Wenn Sie sich über die Sicherheit eines Verfahrens im unklaren sind, setzen Sie sich, bevor Sie fortfahren, mit Ihrer lokalen technischen Unterstützungsorganisation für Thermo Scientific San Jose Produkte in Verbindung.	En caso de duda sobre la seguridad de un procedimiento, póngase en contacto con el personal de servicio técnico de los productos Thermo Scientific San Jose.	Si la sûreté d'une procédure est incertaine, contacter l'assistance technique pour les produits Thermo Scientific San Jose.

CAUTION Symbol	CAUTION	警告	危険警告
	Risk electric shock: This instrument uses voltages that can cause electric shock and/or personal injury. Before servicing, shut down the instrument and disconnect it from line power. While operating the instrument, keep covers on. Do not remove the protective covers from the printed circuit board assemblies (PCBAs).	感電の危険性: この機器では、感電および/または身体傷害を引き起こすおそれのある電圧を使用しています。整備点検の前には、機器の電源を切り、電源コードを抜いてください。機器の作動中は、カバーを付けたままにしてください。プリント基板アセンブリ (PCBA) から保護カバーを取り外さないでください。	触电危险: 本仪器所用电压可能导致电击或人身伤害。进行维修服务前, 务必关闭仪器电源并断开其电源连接。操作此仪器时, 不要卸下顶盖。勿卸下印刷电路板组件 (PCBA) 的保护盖。
	Chemical hazard: Wear gloves and other protective equipment, as appropriate, when handling toxic, carcinogenic, mutagenic, corrosive, or irritant chemicals. Use approved containers and proper procedures to dispose of waste oil and when handling wetted parts of the instrument.	化学的危険性: 毒性、発癌性、変異原性、腐食性、または刺激性のある化学薬品を取り扱うときは、必要に応じて手袋などの保護具を着用します。廃油を処分したり、機器の接液部品を取り扱うときは、認可された容器を使用し、適切な手順に従います。	化学品危险: 当处理毒性、致癌性、致突变性、腐蚀性或者刺激性化学品时, 佩戴手套和其他保护性设备。当处理浸湿的仪器部件以及废油时, 使用认可的容器和合适的步骤。
	Hot surface: Before touching, allow any heated components to cool.	高温面: 触れる前に、加熱した部品を冷ましてください。	热表面: 待高温部件冷却之后再行维修。
	Flammable substances hazard: Use care when operating the system in the presence of flammable substances.	可燃性物質の危険性: 可燃性物質があるところでシステムを作動させる場合は十分注意してください。	易燃物危险: 在有易燃物质的场地操作该系统时, 务必小心谨慎。
	Risk of eye injury: Eye injury could occur from splattered chemicals, airborne particles, or sharp objects. (Sharp objects that customers might install in the instrument include fused-silica tubing, the autosampler needle, and so on.) Wear safety glasses when handling chemicals or servicing the instrument.	眼外傷の危険性: 飛散した化学薬品、浮遊粒子、または鋭利な物体によって眼外傷を負うおそれがあります (機器に取り付けられる可能性がある鋭利な物体は、ヒューズドシリカ、オートサンプラーニードルなどです)。化学薬品を取り扱ったり、機器を整備点検するときは、保護メガネを着用します。	眼睛伤害风险: 眼睛受伤可能源自飞溅的化学品、空气中的颗粒, 或者锋利的物体。(安装在仪器内的锋利物体包括熔融石英管、自动进样器的进样针等。) 处理化学品或对仪器进行维修服务时, 务必戴上防护眼镜。
	General hazard: A hazard is present that is not included in the other categories. This symbol also appears on the instrument. For details about the hazard, refer to the instrument manual. When the safety of a procedure is questionable, contact Technical Support for Thermo Scientific San Jose products.	一般的な危険性: それぞれのカテゴリに当てはまらない危険があります。この標識記号は機器にも表示されています。この危険の詳細については、機器のマニュアルを参照してください。手順の安全性にご不明な点がある場合は、Thermo Scientific San Jose 製品のテクニカルサポートまでお問い合わせください。	普通危険: 未归入其他类别的危险。此符号也会在仪器上出现。有关此危险的详细信息, 参阅适当的仪器手册。若对任何步骤的安全事项有疑问, 联系 Thermo Scientific San Jose 产品的技术支持中心。

CAUTION Symbol	CAUTION	警告	危険警告
	Laser hazard: This instrument uses a laser that is capable of causing personal injury. This symbol also appears on the instrument. For details about the hazard, refer to the instrument manual.	レーザー光線の危険性: この機器では、身体傷害を引き起こすおそれのあるレーザーを使用しています。この標識記号は機器にも表示されています。この危険の詳細については、機器のマニュアルを参照してください。	激光危险: 本仪器所用激光会导致人身伤害。此符号也会在仪器上出现。有关此危险的详细信息，参阅适当的仪器手册。
	Ultra violet light hazard: Do not look directly at the ultra-violet (UV) light or into the UV source. Exposure can cause eye damage. Wear UV eye protection.	紫外光の危険性: 紫外 (UV) 光または UV 光源を直接見ないでください。照射によって眼損傷を引き起こすおそれがあります。UV 保護メガネを着用します。	紫外光危険: 不要直视紫外 (UV) 光或者紫外光源。直视可能导致眼睛伤害。佩戴紫外线防护眼镜。
	Sharp object: Avoid physical contact with the object.	鋭利な物体: 物体との身体的接触を避けてください。	锋利物体: 避免直接接触锋利的物体。
	Pinch point: Keep hands away from this area.	ピンチポイント: この部分には手を挟まれないようにしてください。	夹点: 勿将手放在此部位。
	Heavy objects: Never lift or move the instrument by yourself; you can suffer personal injury or damage the equipment. For specific lifting instructions, refer to the instrument manual.	重量物: 1人で機器を持ち上げたり移動しないでください。身体傷害を負ったり、機器を損傷するおそれがあります。具体的な持ち上げ方法については、機器のマニュアルを参照してください。	重物: 切勿独自提起或移动本仪器；可能遭受人身伤害或损坏仪器。有关具体的提起说明，参阅仪器手册。
	Trip obstacle: Be aware of cords, hoses, or other objects located on the floor.	作業の障害物: 床にあるコード、ホース、その他の物体に注意してください。	绊倒危险: 注意地面上的线、管或其他物品。
	When the safety of a procedure is questionable, contact Technical Support for Thermo Scientific San Jose products.	手順の安全性にご不明な点がある場合は、Thermo Scientific San Jose 製品のテクニカルサポートまでお問い合わせください。	如对安全程序有疑问，联系 Thermo Scientific San Jose 产品的技术支持中心。

Contents

	Preface	xvii
	Related Documentation	xvii
	Safety and Special Notices	xviii
	Safety Precautions	xviii
	Solvent and Gas Purity Requirements	xix
	Service Philosophy	xix
	Level of Repair	xx
	Contacting Us	xxi
Chapter 1	Introduction	1
	Ionization Modes	3
	Ion Polarity Modes	3
	Scan Modes	3
	Q1MS and Q3MS Scan Modes	5
	Product Scan Mode	5
	Parent Scan Mode	6
	Neutral Loss Scan Modes	7
	Data Dependent Scan Mode	9
	Scan Types	10
	Full Scan	10
	Selected Ion Monitoring	10
	Selected Reaction Monitoring	11
	AutoSIM	11
	Data Types	11
	Profile Data Type	12
	Centroid Data Type	12
	Mass/Charge Range	12
Chapter 2	Functional Description	13
	Autosampler	14
	Liquid Chromatograph	15
	Syringe Pump	15
	Divert/Inject Valve	16

	Mass Spectrometer	18
	Controls and Indicators	19
	API Source	22
	Ion Source Interface	22
	Ion Optics	24
	Mass Analyzer	28
	Ion Detection System	35
	Vacuum System and Inlet Gases Hardware	37
	Electronic Assemblies	43
	Data System	44
	Computer Hardware	45
	Data System / Mass Spectrometer / LC Communication, Contact Closure, and Line Power Connections	45
	Data System / Local Area Network Interface	46
	Printer	46
Chapter 3	System Shutdown, Startup, and Reset	49
	Shutting Down the System in an Emergency	49
	Placing the System in Standby Condition	51
	Shutting Down the System Completely	52
	Starting Up the System after a Complete Shutdown	54
	Starting Up the LC	54
	Starting Up the Data System	54
	Starting Up the Mass Spectrometer	54
	Starting Up the Autosampler	55
	Setting Up Conditions for Operation	56
	Resetting the Mass Spectrometer	57
	Resetting the Data System	57
	Resetting the Data System by Using the Windows Shutdown and Restart Procedure	58
	Resetting the Data System by Turning the Personal Computer Off Then On	58
	Turning Off Selected Mass Spectrometer Components	59
Chapter 4	Daily Operation	63
	Things to Do Before Operating the TSQ System	63
	Checking the Argon and Nitrogen Supplies	64
	Checking the ESI Fused-Silica Sample Tube for Elongation	64
	Checking the System Vacuum Levels	64
	Things to Do After Operating the TSQ System	66
	Flushing the Sample Transfer Line, Sample Tube, and API Probe	66
	Placing the System in Standby Condition	67
	Cleaning the Ion Sweep Cone and Ion Transfer Tube	68
	Emptying the Solvent Waste Bottle	68

Chapter 5	Removing and Reinstalling the Ion Source Housing	69
	Removing the Ion Max or Ion Max-S Ion Source Housing	69
	Installing the Ion Max or Ion Max-S Ion Source Housing	71
	Ion Max and Ion Max-S Housing Maintenance	73
Chapter 6	Maintenance	75
	Frequency of Cleaning	77
	Maintaining the Ion Sweep Cone and Ion Transfer Tube	79
	Removing and Cleaning the Ion Transfer Tube	79
	Maintaining the Ion Source Interface Assembly	84
	Shutting Down and Venting the System	84
	Removing the Ion Max or Ion Max-S Ion Source Housing	84
	Removing the Ion Source Interface	85
	Removing the S-Lens and Exit Lens (TSQ Vantage)	86
	Cleaning the S-Lens and Exit Lens (TSQ Vantage)	87
	Removing the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)	87
	Cleaning the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)	88
	Reinstalling the S-Lens and Exit Lens (TSQ Vantage)	89
	Reinstalling the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)	89
	Reinstalling the Ion Source Interface Assembly	90
	Reinstalling the Ion Max or Ion Max-S Ion Source Housing	90
	Starting Up the System	90
	Cleaning the Q00 Ion Optics	91
	Shutting Down and Venting the System	91
	Removing the Ion Max or Ion Max-S Ion Source Housing	91
	Removing the Ion Source Interface Assembly	91
	Removing the Ion Optics Cage Assembly	92
	Disassembling the Ion Optics Cage Assembly	96
	Cleaning Q00 and L0	97
	Reassembling the Ion Optics Cage Assembly	98
	Reinstalling the Ion Optics Cage Assembly	99
	Reinstalling the Ion Source Interface Assembly	100
	Reinstalling the Ion Max or Ion Max-S Ion Source Housing	100
	Starting Up the System	100
	Maintaining the Forepump	101
	Cleaning the Fan Filters	101
Chapter 7	Diagnostics and PCB and Assembly Replacement	103
	Running the TSQ System Diagnostics	104
	Replacing a Fuse	105
	Replacing PCBs and Power Supplies	105

Chapter 8	Replaceable Parts.....	107
	MS Accessory Kit	107
	TSQ Vantage Instrument Specific Accessory Kit	108
	TSQ Quantum Ultra, Ultra AM, and Ultra EMR Instrument Specific Accessory Kit	108
	HESI-II Probe Kit	109
	TSQ Quantum Access Instrument Specific Accessory Kit	109
	TSQ Accurate Mass Calibration Compounds Kit	109
	Fittings, Ferrules, and Sample Loops	109
	Index	111

Preface

This *TSQ Series Hardware Manual* contains a description of the modes of operation and principle hardware components of your Thermo Scientific™ TSQ™ system. In addition, this manual provides step-by-step instructions for cleaning and maintaining your mass spectrometer.

Contents

- [To suggest changes to documentation or to Help](#)
- [Safety and Special Notices](#)
- [Safety Precautions](#)
- [Solvent and Gas Purity Requirements](#)
- [Service Philosophy](#)
- [Level of Repair](#)
- [Contacting Us](#)

❖ To suggest changes to the documentation or to the Help

Complete a brief survey about this document by clicking the button below.
Thank you in advance for your help.



Related Documentation

In addition to this manual, Thermo Fisher Scientific provides the following for TSQ Series mass spectrometers:

- *TSQ Series Preinstallation Requirements Guide*
- *TSQ Series Getting Connected Guide*
- *TSQ Series Getting Started Guide*

- *H-ESI Probe User Guide*
- *HESI-II Probe User Guide*
- *Ion Max and Ion Max-S API Source Hardware Manual*
- Help available from within the software

Safety and Special Notices

Make sure you follow the precautionary statements presented in this guide. The safety and other special notices appear in boxes.

Safety and special notices include the following:



CAUTION Highlights hazards to humans, property, or the environment. Each CAUTION notice is accompanied by an appropriate CAUTION symbol.

IMPORTANT Highlights information necessary to prevent damage to software, loss of data, or invalid test results; or might contain information that is critical for optimal performance of the system.

Note Highlights information of general interest.

Tip Highlights helpful information that can make a task easier.

Safety Precautions

Observe the following safety precautions when you operate or perform service on the mass spectrometer.



CAUTION Do not perform any servicing other than that contained in the *TSQ Series Hardware Manual*. To avoid personal injury or damage to the instrument, do not perform any servicing other than that contained in the *TSQ Series Hardware Manual* or related manuals unless you are authorized to do so.



CAUTION Shut down the mass spectrometer and disconnect it from line power before you service it. High voltages capable of causing personal injury are used in the instrument. Some maintenance procedures require that the mass spectrometer be shut down and disconnected from line power before service is performed. Do not operate the mass spectrometer with the top or side covers off. Do not remove protective covers from printed circuit boards (PCBs).



CAUTION Respect heated zones. Treat heated zones with respect. The ion transfer capillary, the H-ESI vaporizer, and the APCI vaporizer might be very hot and might cause severe burns if they are touched. Allow heated components to cool before you service them.



CAUTION Place the mass spectrometer in Standby (or Off) before you open the atmospheric pressure ionization (API) source. The presence of atmospheric oxygen in the API source when the mass spectrometer is On could be unsafe. (The TSQ system automatically turns the mass spectrometer Off when you open the API source; however, it is best to take this added precaution.)



CAUTION Provide an adequate fume exhaust system. It is your responsibility to provide an adequate fume exhaust system. Samples and solvents that are introduced into the TSQ mass spectrometer will eventually be exhausted from the forepump. Therefore, the forepump should be connected to a fume exhaust system. Consult local regulations for the proper method of exhausting the fumes from your system.



CAUTION Use care when changing the vacuum pump oil. Treat drained vacuum pump oil and pump oil reservoirs with care. Hazardous compounds introduced into the system might have become dissolved in the pump oil. Always use approved containers and procedures for disposing of waste oil. Whenever a pump has been operating on a system used for the analysis of toxic, carcinogenic, mutagenic, or corrosive/irritant chemicals, the pump must be decontaminated by the user and certified to be free of contamination before repairs or adjustments are made by a Thermo Fisher Scientific San Jose Customer Support Engineer or before it is sent back to the factory for service.

Solvent and Gas Purity Requirements

Use the highest purity solvents available. The TSQ mass spectrometer is extremely sensitive to solvent impurities. Some solvent impurities are transparent to UV/Visible detectors, but are easily detected by the TSQ mass spectrometer. Liquid chromatography grade is the minimum acceptable purity. Higher grade solvents are preferred. Distilled water is recommended. Deionized water contains chemicals and is not recommended.

You can order high quality solvents from Fisher Global Chemicals, telephone (800) 766-7000.

The TSQ mass spectrometer uses argon as a collision gas. The argon must be high purity (99.995%). The required gas pressure is 135 ± 70 kPa (20 ± 10 psig). Thermo Fisher Scientific has found that particulate filters are often contaminated and are therefore not recommended.

Service Philosophy

Servicing the TSQ system consists of performing procedures required to maintain system performance standards, prevent system failure, restore the system to an operating condition, or all of the above. Routine and preventive maintenance procedures are documented in this manual.

The user is responsible for routine and preventive maintenance during and after the warranty period. Regular maintenance increases the life of the system, maximizes the up-time of your system, and allows you to achieve optimum system performance.

Only a Thermo Fisher Scientific customer support engineer can perform services not described in this manual.

Level of Repair

Thermo Fisher Scientific's service philosophy for the TSQ system calls for troubleshooting to the lowest part, assembly, or module listed in [Chapter 8, "Replaceable Parts."](#)

For mechanical failures: A mechanical assembly typically is to be repaired to the level of the smallest item listed in [Chapter 8, "Replaceable Parts."](#)

Contacting Us

There are several ways to contact Thermo Fisher Scientific for the information you need.

For Thermo Scientific™ products	Access by phone, fax, email, or website
Technical Support	<p>(U.S.) Phone: 1 (800) 532-4752 Fax: 1 (561) 688-8736</p> <p>Email: us.techsupport.analyze@thermofisher.com</p> <p>Web—for product support, technical documentation, and knowledge bases: www.thermoscientific.com/support</p>
Customer Service (Sales and service)	<p>(U.S.) Phone: 1 (800) 532-4752 Fax: 1 (561) 688-8731</p> <p>Email: us.customer-support.analyze@thermofisher.com</p> <p>Web—for product information: www.thermoscientific.com/lc-ms</p> <p>Web—for customizing your service request:</p> <ol style="list-style-type: none"> 1. From any Products & Services web page, click Contact Us. 2. In the Contact Us box, complete the information requested, scroll to the bottom, and click Send.
User Documentation	<p>Web—for downloading documents: mssupport.thermo.com</p> <ol style="list-style-type: none"> 1. On the Terms and Conditions web page, click I Agree. 2. In the left pane, click Customer Manuals. 3. To locate the document, click Search and enter your search criteria. For Document Type, select Manual. <p>Email—to send feedback directly to Technical Publications: techpubs-lcms@thermofisher.com</p> <p>Web—to complete a survey about this Thermo Scientific document: www.surveymonkey.com/s/PQM6P62</p>

Introduction

TSQ Series mass spectrometers are members of the Thermo Scientific family of mass spectrometers. The TSQ Series mass spectrometer is an advanced analytical instrument that includes a syringe pump, a divert/inject valve, a mass spectrometer, and a data system.

Contents

- [Ionization Modes](#)
- [Ion Polarity Modes](#)
- [Scan Modes](#)
- [Scan Types](#)
- [Data Types](#)
- [Mass/Charge Range](#)

In a typical analysis, a sample can be introduced in the following ways:

- Using the syringe pump (direct infusion)
- Using the divert/inject valve fitted with a sample loop and liquid chromatograph (LC) (flow injection analysis)
- Using the divert/inject valve and high-performance liquid chromatograph (HPLC) fitted with a column (LC/MS)

In a typical analysis, an LC introduces a sample. The LC separates the sample into its various components. The components elute from the LC and pass into the mass spectrometer where they are analyzed.

Analysis by direct infusion or flow injection provides no chromatographic separation of components in the sample before they pass into the mass spectrometer. The data from the mass spectrometer is then stored and processed by the data system.

The TSQ mass spectrometer consists of an atmospheric pressure ionization (API) source, ion optics, triple-stage mass analyzer, and ion detection system. A vacuum manifold encloses the ion optics, mass analyzer, ion detection system, and part of the API source. Ionization of the sample takes place in the API source. The specific process used to ionize the sample is known

as the ionization mode. The ion optics transmit the ions produced in the ion source into the mass analyzer, where they are separated according to their mass-to-charge ratio. The polarity of the potentials applied to the lenses in the ion source and ion optics determines whether they transmit positively charged ions or negatively charged ions to the mass analyzer. You can configure the TSQ mass spectrometer to analyze positively or negatively charged ions (called the positive or negative ion polarity mode).

The TSQ instrument's triple-stage mass analyzer performs either one or two stages of mass analysis:

- The TSQ system is operated as a conventional mass spectrometer with one stage of mass analysis. The ion source ionizes the sample and the ion products are subjected to mass analysis in the first rod assembly. The second and third rod assemblies transmit the resulting mass-selected ions to the ion detection system.¹
- The TSQ system is operated as a tandem mass spectrometer with two stages of mass analysis. The ion source ionizes the sample and the ion products are mass analyzed by the first rod assembly. In this case, however, mass-selected ions exiting the first rod assembly collide with an inert gas in the second rod assembly and fragment to produce a set of ions known as product ions. (A chamber called the collision cell surrounds the second rod assembly. The collision cell can be pressurized with an inert gas.) The product ions undergo further mass analysis in the third rod assembly to detect selected ions. Two stages of mass analysis yield far greater chemical specificity than a single stage can achieve, because of the system's ability to select and determine two discrete but directly related sets of masses.

In a first stage of mass analysis the TSQ systems can be used to elucidate the structures of pure organic compounds and the structures of the components within mixtures. Furthermore, in a second stage of mass analysis, the mass spectrometer can fragment and separate each ionic fragment of a molecule formed in the ion source to build up an entire structure for the molecule, piece by piece. Thus, TSQ systems make investigating all pathways for the formation and fragmentation of each ion in the mass spectrum possible.

The two stages of mass analysis, with resultant reduction of chemical noise in the final mass spectrum, allow for very selective and sensitive analysis.

Each sequence of single- or triple-stage mass analysis of the ions is called a scan. The TSQ mass spectrometer uses several different scan modes and different scan types to filter, fragment, or transmit ions in the mass analyzer. Along with the ionization and ion polarity modes, the ability to vary the scan mode and scan type affords the user great flexibility in the instrumentation for solving complex analytical problems.

¹ The instrument can also be used as a single-stage mass spectrometer by transmitting the ions through the first and second rod assemblies followed by mass analysis in the third rod assembly.

Ionization Modes

The atmospheric pressure ionization (API) source forms gas phase sample ions from sample molecules that elute from the LC or are introduced by the syringe pump. You can operate the API source in the electrospray ionization (ESI) mode or, optionally, the heated electrospray ionization (H-ESI), nanospray ionization (NSI), atmospheric pressure photo ionization (APPI), or atmospheric pressure chemical ionization (APCI) mode. Refer to the *Ion MAX and Ion MAX-S API Source Hardware Manual*, *HESI-II Probe User Guide*, *H-ESI Probe User Guide*, *Ion MAX APPI Source Operator's Manual*, or *Nanospray Ion Source Operator's Manual* for more information regarding the ionization modes.

Ion Polarity Modes

You can operate the TSQ mass spectrometer in either of two ion polarity modes: positive or negative. Both positively charged and negatively charged ions form in the ion source of the mass spectrometer. The TSQ mass spectrometer can control whether positive ions or negative ions are transmitted to the mass analyzer for mass analysis by changing the polarity of the potentials applied to the ion source and ion optics. The ion optics deliver the ions produced in the ion source, in a collimated beam, to the mass analyzer.

The information obtained from a positive ion mass spectrum is different from and complementary to the information from a negative ion spectrum. Thus, the ability to obtain both positive ion and negative ion mass spectra aids you in the qualitative analysis of your sample. You can choose the ion polarity mode and ionization mode to obtain maximum sensitivity for the particular analyte of interest.

Scan Modes

You can operate the TSQ mass spectrometer in a variety of scan modes. The most commonly used scan modes can be divided into two categories: single mass spectrometry (MS) scan modes and MS/MS scan modes. The scan modes in each category are as follows:

- MS scan modes: Q1MS and Q3MS scan modes
- MS/MS scan modes product scan mode, parent scan mode, neutral loss scan mode
- Data-dependent scan mode

The scan modes that can be employed depend on the number and type of rod assemblies and the voltages applied to the rod assemblies.

The TSQ system mass analyzer has three rod assemblies.² The first and third rod assemblies, Q1 and Q3, are quadrupoles, and the second rod assembly, Q2, is a square-profile quadrupole.

Rod assemblies can operate in either of two capacities:

- As ion transmission devices
- As mass analyzers

If only RF voltage is applied, a rod assembly serves as an ion transmission device that passes all ions within a large range of mass-to-charge ratios (that is, virtually all ions present).

When you apply both RF and dc voltages to a rod assembly, the separation of ions of different mass-to-charge ratios occurs. This separation allows the rod assembly to serve as a mass analyzer.

On the TSQ mass spectrometer, the quadrupole rod assemblies can operate with RF and dc voltages or with only RF voltage. That is, Q1 and Q3 can act either as mass analyzers or ion transmission devices. The Q2 rod assembly operates exclusively with RF voltage. Thus, Q2 is always an ion transmission device. For a summary of how the rod assemblies function in several of the major scan modes, see [Table 1](#).

Table 1. Summary of scan modes

Scan mode	Q1 quadrupole	Q2 collision cell	Q3 quadrupole
Q1MS	Scan ^a	Pass all ions ^b	Pass all ions
Q3MS	Pass all ions	Pass all ions	Scan
Product	Set ^c	Fragment ions ^d , then pass all fragments	Scan
Parent	Scan	Fragment ions, then pass all fragments	Set
Neutral Loss	Scan	Fragment ions, then pass all fragments	Scan

^a Scan = full scan or transmission of selected ions

^b Pass all ions or fragments = pass ions or fragments within a wide range of mass-to-charge ratios

^c Set = set to pass ions of a single mass-to-charge ratio or a set of mass-to-charge ratios

^d Fragment ions = collisions with argon gas cause ions to fragment

² A rod assembly is a regular array of metal rods. See “Mass Analyzer” on [page 28](#) for a discussion of the rod assemblies used on the TSQ Series instrument.

Q1MS and Q3MS Scan Modes

The Q1MS and Q3MS scan modes perform only one stage of mass analysis. The mass spectrum obtained is equivalent to the mass spectrum obtained from an instrument with a single mass analyzer. In the one stage of analysis, ions formed in the ion source enter the analyzer assembly. One of the mass analyzers (Q1 or Q3) is scanned to obtain a complete mass spectrum. The other rod assemblies (Q2 and Q3, or Q1 and Q2, respectively) act as ion transmission devices. In the Q1MS scan mode, Q1 is used as the mass analyzer; in the Q3MS scan mode, Q3 is used as the mass analyzer.

Product Scan Mode

Product scan mode performs two stages of analysis. In the first stage, ions formed in the ion source enter Q1, which is set to transmit ions of one mass-to-charge ratio. Ions selected by this first stage of mass analysis are called parent ions. (As a result, Q1 is referred to as the parent mass analyzer, and the mass-to-charge ratio of ions transmitted by the parent mass analyzer is referred to as the parent set mass.) Parent ions selected by Q1 then enter Q2, which is surrounded by the collision cell.


Note For convenience, when referring to the first, second, and third rod assemblies as pieces of hardware, you can call them Q1, Q2, and Q3, respectively. However, for clarity when discussing their function in MS/MS scan modes, you can refer to them as the parent mass analyzer, collision cell (ion transmission device surrounded by the collision cell), and product mass analyzer, respectively.

In the second stage of analysis, ions in the collision cell can fragment further to produce product ions. Two processes produce product ions: by unimolecular decomposition of metastable ions or by interaction with argon collision gas present in the collision cell. This latter step is known as collision-induced dissociation (CID). Ions formed in the collision cell enter Q3 (the product mass analyzer) for the second stage of mass analysis. Q3 is scanned to obtain a mass spectrum that shows the product ions produced from the fragmentation of the selected parent ion.

A mass spectrum obtained in the product scan mode (product mass spectrum) is the mass spectrum of a selected parent ion.

Figure 1 illustrates the product scan mode.

Figure 1. Illustration of product scan mode

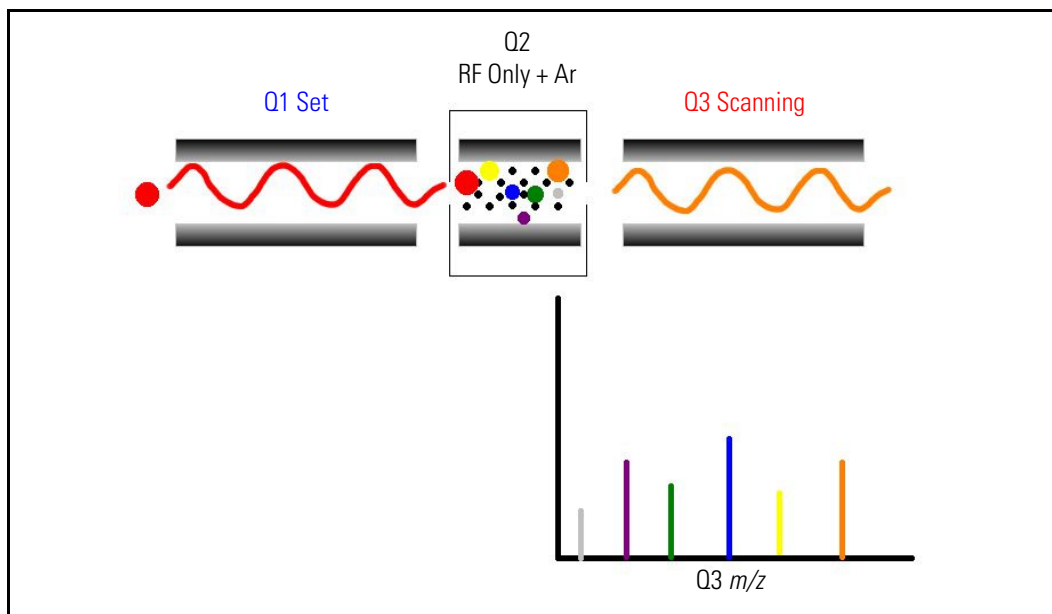


Parent Scan Mode

The parent scan mode also uses two stages of analysis. In the first stage, ions formed in the ion source are introduced into the parent mass analyzer, which is scanned to transmit parent ions sequentially into the collision cell.

In the second stage of analysis, in the collision cell, parent ions can fragment to produce product ions by unimolecular decomposition of metastable ions or by collision-induced dissociation. Ions formed in the collision cell enter the product mass analyzer, which transmits a selected product ion. (The product set mass is the mass-to-charge ratio of ions transmitted by the product mass analyzer.)

The resultant spectrum shows all the parent ions that fragment to produce the selected product ion. Note that for a mass spectrum obtained in the parent scan mode (parent mass spectrum), data for the mass-to-charge ratio axis is obtained from Q1 (the parent ions), whereas data for the ion intensity axis is obtained from Q3 (from monitoring the product ion).

Figure 2 illustrates the parent scan mode.

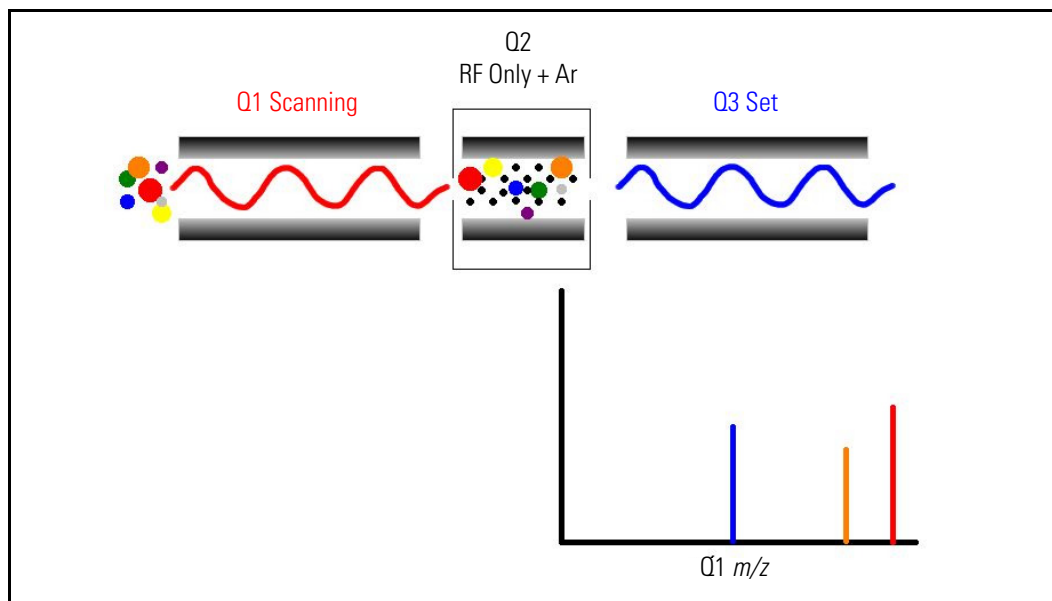Figure 2. Illustration of the parent scan mode

Experiments that employ the parent scan mode (parent experiments) can be used in structure and fragmentation studies as well as in survey analyses of mixtures. In general, parent experiments detect all compounds that decompose to a common fragment. The experiments are useful for the rapid detection of a series of structural homologs (for example, substituted aromatics, phthalates, steroids, or fatty acids) that have a common fragment ion (for example, m/z 149 for the phthalates).

Neutral Loss Scan Modes

In the neutral loss scan mode, the two mass analyzers (Q1 and Q3) are linked together so that they are scanned at the same rate over mass ranges of the same width. The respective mass ranges, however, are offset by a selected mass so that the product mass analyzer scans a selected number of mass units lower than the parent mass analyzer.

As a result, in the neutral loss scan mode, there are two stages of mass analysis. In the first stage, the parent mass analyzer separates ions formed in the ion source by mass-to-charge ratio. Then the ions are introduced sequentially into the collision cell.

In the second stage of analysis, ions admitted to the collision cell can fragment further by metastable ion decomposition or by CID to produce product ions. The product mass analyzer then separates these product ions by their mass-to-charge ratio. [Figure 3](#) illustrates the neutral loss scan mode. Examples of compounds with a common neutral loss fragment appear in [Figure 4](#).

For an ion to be detected, between the time the ion leaves Q1 and enters Q3, it must lose a neutral moiety whose mass (the neutral loss mass) is equal to the difference in the mass ranges being scanned by the two mass analyzers. Thus, a neutral loss mass spectrum is a spectrum that shows all the parent ions that lose a neutral species of a selected mass.

Note that a neutral gain (or association) experiment can also be performed in which the mass range scanned by Q3 is offset by a selected mass above the mass range scanned by Q1.

For a neutral loss (or neutral gain) mass spectrum, as for a parent mass spectrum, Q1 (the parent ion) provides data for the mass-to-charge ratio axis, whereas Q3 (the product ion being monitored) provides data for the ion intensity axis.

Experiments that use the neutral loss scan mode (neutral loss experiments) are useful when a large number of compounds are being surveyed for common functionality. Neutral moieties are frequently lost from substituent functional groups (for example, CO_2 from carboxylic acids, CO from aldehydes, HX from halides, and H_2O from alcohols).

Figure 3. Illustration of the neutral loss scan mode


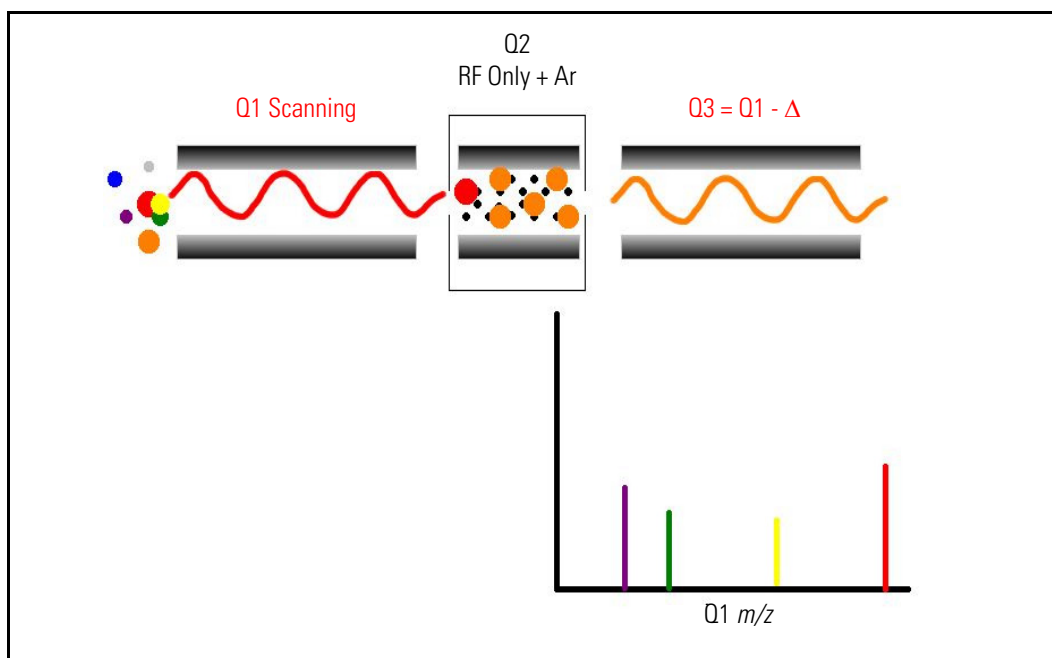


Figure 4. Examples of compounds with a common neutral-loss fragment

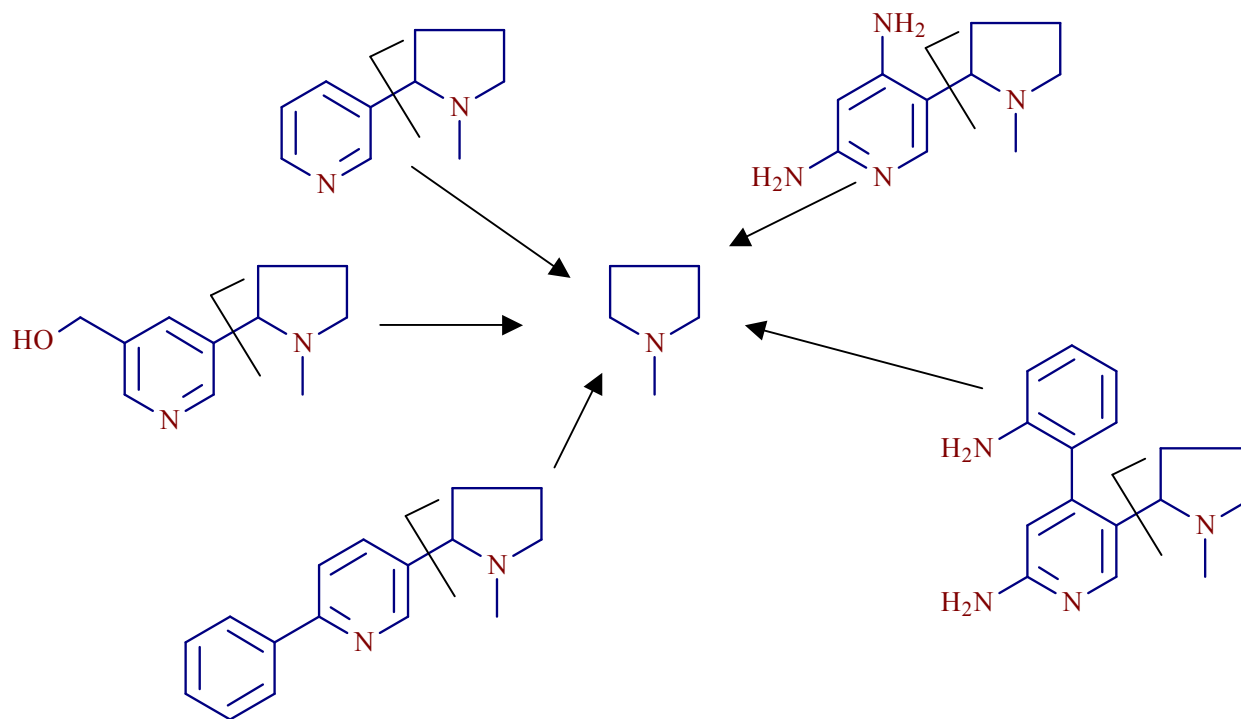

Data Dependent Scan Mode

The TSQ mass spectrometer uses the information in a data-dependent scan mode experiment to make automatic decisions about the next step of the experiment without input from a user. In data-dependent scan mode you specify criteria to select one or more ions of interest on which to perform subsequent scans, such as MS/MS. You can approach the setup of data-dependent experiments in either of two ways:

- If you have some idea of what the parent ion is, or if you expect a certain kind of parent, you can set up a list of possible parent ions. Then, when one of the parent ions you specified is detected, you can acquire product spectra and analyze the information. Conversely, you can also set up a list of ions that you do *not* want selected for fragmentation.
- If you have little information about your compound, you can set up the parameters of a data-dependent experiment so that if the intensity of the ion signal is above a specified threshold, the TSQ system generates product spectra. Later, you can decide if the information is useful.

Because a data-dependent scan needs to use a target ion from a previous scan, the first scan event cannot be a data-dependent scan.

Scan Types

TSQ systems can be operated with a variety of scan types. The most common scan types are as follows:

- [Full Scan](#)
- [Selected Ion Monitoring \(SIM\)](#)
- [Selected Reaction Monitoring \(SRM\)](#)
- [AutoSIM](#)

Full Scan

The full-scan scan type provides a full mass spectrum of each analyte. With full scan, the scanning mass analyzer is scanned from the first mass to the last mass, without interruption, in a given scan time.

Full-scan experiments are used to determine or confirm the identity of unknown compounds or the identity of each component in a mixture of unknown compounds. (Generally, a full mass spectrum is needed to determine the identity of an unknown compound.)

The full-scan scan type gives you more information about an analyte than does SIM, but a full scan does not yield the sensitivity that the other two scan types can achieve. With full scan, you spend less time monitoring the signal for each ion than you do in SIM or SRM. Full scan provides greater information but lower sensitivity than the other two scan types.

To use the SIM or SRM, you need to know what ions or reactions you are looking for before you can perform an experiment with these scan types. Thus, you might use a full scan for SIM to determine the identity of an analyte and to obtain its mass spectrum and a full scan for SRM to determine the mass spectrum and product mass spectra for parent ions of interest. Then, you might use SIM or SRM to do routine quantitative analysis of the compound.

Selected Ion Monitoring

Selected ion monitoring (SIM) monitors a particular ion or set of ions. SIM experiments are useful in detecting small quantities of a target compound in a complex mixture when you know the mass spectrum of the target compound. Thus, SIM is useful in trace analysis and in the rapid screening of a large number of samples for a target compound.

Because SIM monitors only a few ions, it can provide lower detection limits and greater speed than the full-scan modes. SIM achieves lower detection limits because more time is spent monitoring significant ions that are known to occur in the mass spectrum of the target analyte. SIM can achieve greater speed because it monitors only a few ions of interest; SIM does not monitor regions of the spectrum that are empty or have no ions of interest.

SIM can improve the detection limit and decrease analysis time, but it can also reduce specificity. Because SIM monitors only specific ions, any compound that fragments to produce those ions will appear to be the target compound. The result can be a false positive.

Selected Reaction Monitoring

Selected reaction monitoring (SRM) monitors a particular reaction or set of reactions, such as the fragmentation of an ion or the loss of a neutral moiety.

SRM monitors a limited number of parent or product ion pairs. In product-type experiments, a parent ion is selected as usual, but generally only one product ion is monitored. SRM experiments are normally conducted with the product scan mode.

As does SIM, SRM provides for the very rapid analysis of trace components in complex mixtures. However, because SRM selects two sets of ions, it obtains specificity that is much greater than what SIM can obtain. Any interfering compound would not only have to form an ion source product (parent ion) of the same mass-to-charge ratio as the selected parent ion from the target compound, but that parent ion would also have to fragment to form a product ion of the same mass-to-charge ratio as the selected product ion from the target compound.

AutoSIM

In the scan type known as AutoSIM, the mass spectrometer automatically selects the most intense masses (m/z values) in a survey scan, builds a SIM scan list for them, and then acquires and records ion current at only these selected masses. AutoSIM scans can be performed on any full scan in any scan mode but not on data-dependent scans.

There might be times when the scan ranges of two (or more) selected masses overlap. If this happens, both masses are placed in one SIM window. In EZ Tune, the SIM table in the Define Scan dialog box displays the center mass for this new scan window, not each selected mass.

Data Types

You can acquire and display mass spectral data (intensity versus mass-to-charge ratio) with the TSQ mass spectrometer in one of two data types:

- [Profile Data Type](#)
- [Centroid Data Type](#)

Profile Data Type

In the profile data type, you can see the shape of the peaks in the mass spectrum. Each atomic mass unit is divided into many sampling intervals. The intensity of the ion current is determined at each of the sampling intervals. The profile data type displays the intensity at each sampling interval with the intensities connected by a continuous line. In general, use the profile scan data type when you tune and calibrate the mass spectrometer so that you can easily see and measure mass resolution.

Centroid Data Type

The centroid data type displays the mass spectrum as a bar graph and sums the intensities of each set of multiple sampling intervals. This sum is displayed versus the integral center of mass of the sampling intervals. In general, use the centroid scan data type for data acquisition for faster scan speed. Data processing is also much faster for centroid data.

Mass/Charge Range

The TSQ mass spectrometer can operate in a mass/charge range of 10 to 3000 Da (TSQ Quantum[™] Access[™], TSQ Quantum Access MAX[™], TSQ Vantage EMR[™], and TSQ Quantum Ultra EMR[™] mass spectrometers), or 10 to 1500 Da (TSQ Vantage[™], TSQ Vantage AM[™], TSQ Quantum Ultra[™] and TSQ Quantum Ultra AM[™] mass spectrometers).

Functional Description

This chapter describes the principal components of the TSQ system and their respective functions.

Contents

- [Autosampler](#) (optional)
- [Liquid Chromatograph](#) (optional)
- [Mass Spectrometer](#)
- [Data System](#)


A functional block diagram of the TSQ mass spectrometer is shown in [Figure 5](#). A sample transfer line connects the LC to the mass spectrometer. The autosampler and LC are usually installed on the left of the mass spectrometer. The syringe pump and divert/inject valve are integrated into the mass spectrometer cabinet.

In a typical analysis, a sample can be introduced in any of the following ways:

- Using the syringe pump (direct infusion)
- Using the divert/inject valve fitted with a loop and an LC (flow injection analysis)
- Using a divert/inject valve and LC fitted with a column (LC/MS)

In analysis by LC/MS, a sample is injected onto an LC column. The sample then separates into its various components. The components elute from the LC column and pass into the mass spectrometer where they are analyzed.

The Electrospray (ESI), heated electrospray (H-ESI), nanospray (NSI), atmospheric pressure photo ionization (APPI), and atmospheric pressure chemical ionization (APCI) techniques each ionize sample molecules at atmospheric pressure. The ion optics focus and accelerate the resulting sample ions into the mass analyzer where they are analyzed according to their mass-to-charge ratios. An ion detection system then produces a signal proportional to the number of ions detected. The system electronics receive and amplify the ion current signal from the ion detection system. That signal is then passed on to the data system for further processing, storage, and display. The data system provides the primary TSQ mass spectrometer user interface.

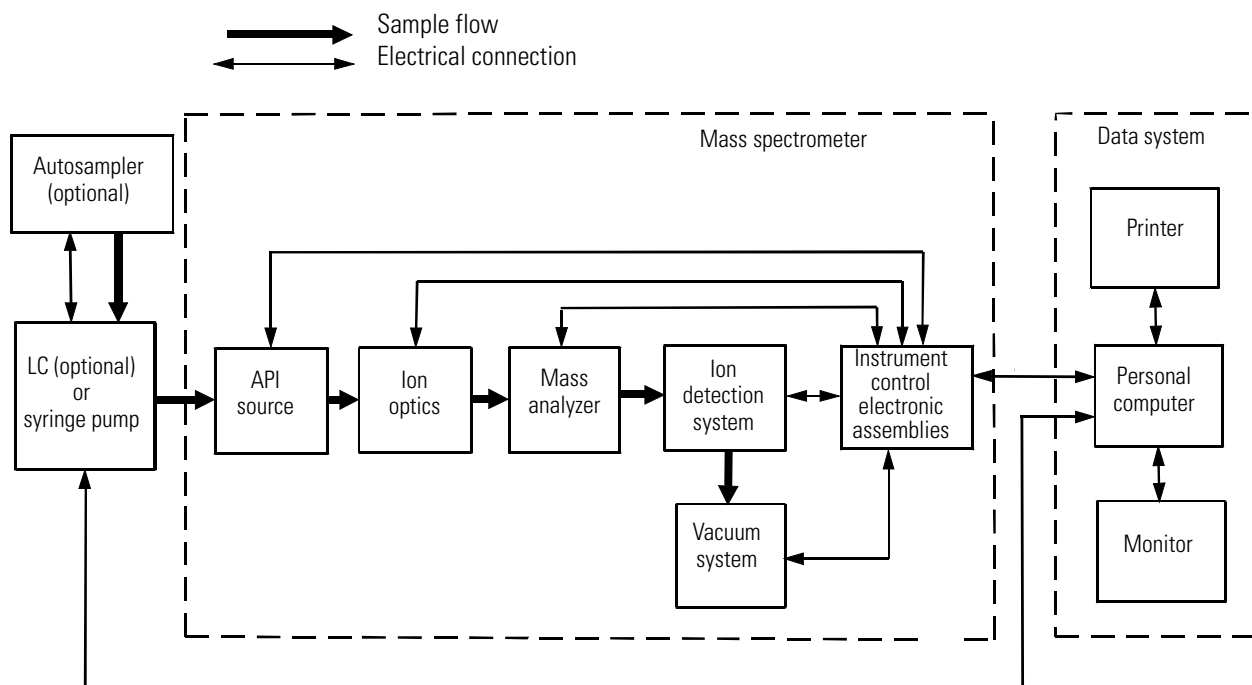Figure 5. Functional block diagram of the TSQ system

Autosampler

The (optional) autosampler injects samples automatically into the liquid chromatograph (LC) inlet. The TSQ data system computer can control most autosamplers. With an autosampler, you can automate your LC/MS analyses.

Autosampler Start/Stop signals with the TSQ mass spectrometer are provided by contact closure. Refer to the *TSQ Series Getting Connected Guide* for information on connecting an autosampler to the TSQ mass spectrometer by contact closure.

You configure the autosampler from the data system computer. Select the appropriate autosampler button in the Instrument Configuration window, which is available by choosing **Start > All Programs > Thermo Foundation > Instrument Configuration**. See Instrument Configuration Help for a description of autosampler configuration options.

You also use the data system to set up the autosampler to inject samples. Choose **Start > All Programs > Thermo Xcalibur > Xcalibur** and click **Instrument Setup** to open the Instrument Setup window. Then, click the appropriate autosampler icon to open the Autosampler page. For instructions on running your autosampler, refer to the Help.

If your autosampler is controlled from a Thermo Scientific data system, refer to the interconnect diagrams in [Figure 33 on page 47](#) and [Figure 34 on page 48](#), and the *Getting Connected Guide* for the autosampler, for information on connecting the autosampler to the data system computer, mass spectrometer, and LC.

For front-panel (keypad) operation (if any) and maintenance procedures, refer to the documentation provided with the autosampler.

Liquid Chromatograph

The (optional) liquid chromatograph (LC) separates a sample mixture into its chemical components by liquid chromatography. In liquid chromatography, the sample mixture partitions between a solid stationary phase of large surface area and a liquid mobile phase that percolates over the stationary phase. The molecular structure of each component of the mixture determines in which order each component elutes from the LC and enters the mass spectrometer.

You can control most LCs and the corresponding UV detectors directly from the TSQ data system computer. You can set the flow rate from 0 to 1000 $\mu\text{L}/\text{min}$ or more. Refer to the *TSQ Series Getting Connected Guide* for information on connecting an LC to the TSQ mass spectrometer.

You configure the LC from the data system computer. Select the appropriate LC button in the Instrument Configuration window, which is available by choosing **Start > All Programs > Thermo Foundation > Instrument Configuration**. See Instrument Configuration Help for a description of LC configuration options, refer to Xcalibur Help.

You also use the data system to set up the LC. Choose **Start > All Programs > Thermo Xcalibur > Xcalibur** and click **Instrument Setup** to open the Instrument Setup window. Then, click the appropriate LC icon to open the LC page. For instructions on running your LC, refer to the Help.

If your LC is controlled from a Thermo Scientific data system, refer to the interconnect diagrams in [Figure 33](#) on [page 47](#) and [Figure 34](#) on [page 48](#), and the *Getting Connected Guide* for the LC, for information on connecting the LC to the data system computer.

For front-panel (keypad) operation (if any) and maintenance procedures, refer to the documentation provided with the LC.


Syringe Pump

The TSQ mass spectrometer includes an electronically-controlled, integrated syringe pump. See [Figure 6](#). The syringe pump delivers sample solution from the syringe into the atmospheric pressure ionization (API) source. When the syringe pump is operating, a motor drives a pusher block that depresses the plunger of the syringe at a rate of one percent of the syringe volume per minute. Liquid flows out of the syringe needle and into the sample transfer line as the plunger is depressed. A syringe holder holds the syringe in place. Refer to the *TSQ Series Getting Started Guide* or the *TSQ Series Getting Connected Guide* for instructions on setting up the syringe pump.

You can start and stop the syringe pump from the EZ Tune window, which is available by choosing **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune**. Refer to EZ Tune Help for instructions on operating the syringe pump from the data system. You can also start and stop the syringe pump by pressing the syringe pump button. By holding down the button, you can run the syringe pump in Purge mode, where the flow rate is five percent of the syringe volume per second.

The syringe pump light-emitting diode (LED) (see [Figure 9](#) and [Figure 10](#)) illuminates green whenever the syringe pump is pumping. The LED illuminates yellow if the syringe pump is at the end of its travel.

Figure 6. Syringe pump

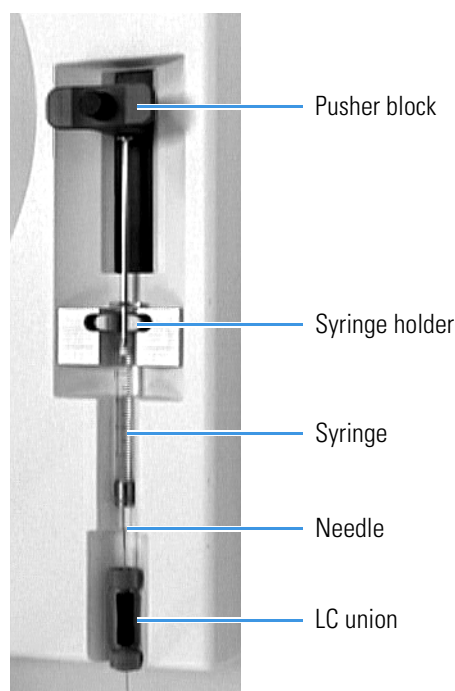

Divert/Inject Valve

The Rheodyne™ 7750E-185 divert/inject valve is a motorized, stainless steel, six-port valve that switches between two positions: Load and Inject. The divert/inject valve is located on the front of the TSQ mass spectrometer above the API source. See [Figure 7](#). You can configure (plumb) the divert/inject valve as a loop injector (for flow injection analysis) or as a divert valve. See [Figure 8](#). The divert valve enables switching the solvent front, gradient endpoint, or any portion of the LC run to waste. The *TSQ Series Getting Started Guide* provides procedures for plumbing the valve in the loop injector or divert valve configuration.

You can control the divert/inject valve from the data system. You specify the parameters of the divert/inject valve on the Divert Valve page, accessible from the Instrument Setup window. For instructions on operating the divert/inject valve from the data system, refer to the Help.

You can also use the divert/inject valve button to divert the LC flow between the mass spectrometer and waste when the valve is in the divert valve configuration, or switch between load and inject modes when the valve is in the loop injector configuration.

Figure 7. Divert/inject valve



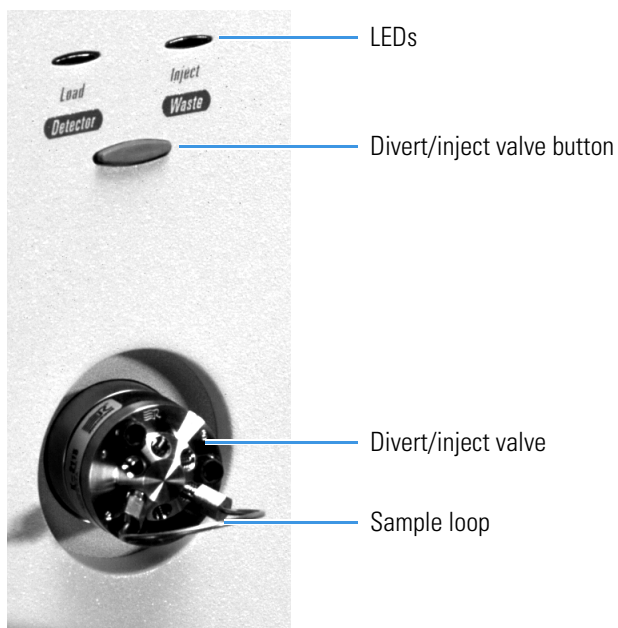


Figure 8. Divert/inject valve plumbed as a loop injector and as a divert valve

Plumbed as a loop injector

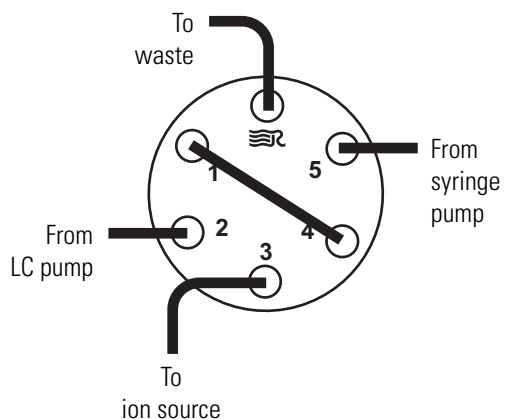

Plumbed as a divert valve

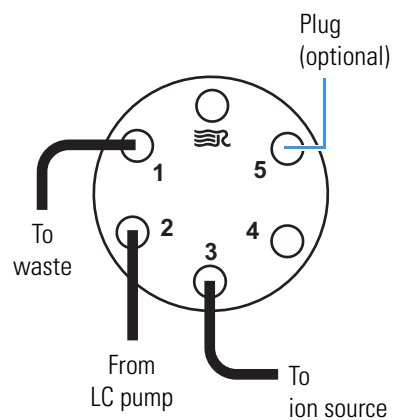

Mass Spectrometer

The mass spectrometer provides sample ionization and mass analysis of injected samples or samples eluted from a liquid chromatograph. The TSQ mass spectrometer uses a triple-quadrupole mass analyzer with an ion source external to the mass analyzer. Several important features of the TSQ mass spectrometer are as follows:

- High sensitivity and resolution
- m/z 10 to 3000 mass range for TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Vantage and TSQ Quantum Ultra mass spectrometers with extended mass range (EMR) electronic assemblies installed; and m/z 10 to 1500 mass range for TSQ Vantage, TSQ Vantage AM, TSQ Quantum Ultra, and TSQ Quantum Ultra AM mass spectrometers without extended mass range electronic assemblies
- ESI, H-ESI, NSI, APPI, and APCI ionization modes
- Positive and negative ion polarity modes
- MS and MS/MS scan modes
- Full-scan, SIM, SRM, AutoSIM, and data-dependent scan types

The mass spectrometer includes the following components:

- [Controls and Indicators](#)
- [API Source](#)
- [Ion Source Interface](#)
- [Ion Optics](#)
- [Mass Analyzer](#)
- [Ion Detection System](#)
- [Vacuum System and Inlet Gases Hardware](#)
- [Electronic Assemblies](#)

Controls and Indicators

Six LEDs are located at the upper right side of the front panel of the mass spectrometer. See [Figure 9](#) (TSQ Vantage, TSQ Quantum Access, or TSQ Quantum Access MAX mass spectrometer) and [Figure 10](#) (TSQ Quantum Ultra mass spectrometer).

Figure 9. Front panel LEDs of the TSQ Vantage, TSQ Quantum Access, or TSQ Quantum Access MAX mass spectrometer


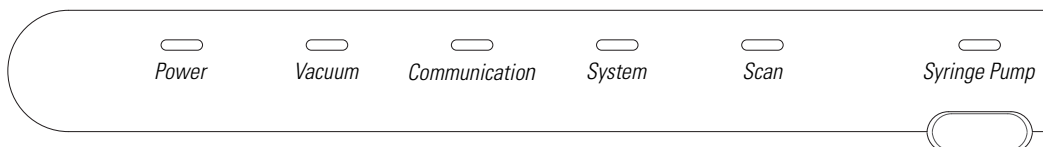


Figure 10. Front panel LEDs of the TSQ Quantum Ultra mass spectrometer

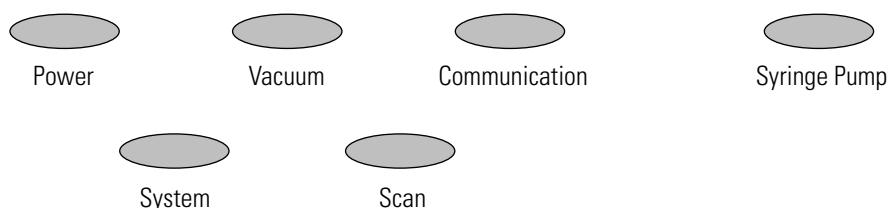

The Power LED illuminates green whenever power is supplied to the vacuum system and electronic assemblies of the mass spectrometer.

The Vacuum LED illuminates yellow when the turbomolecular pump is nearly at speed (80 percent of its operating speed of 750 MHz) and it is safe to turn on the ion gauge. The Vacuum LED is off if the turbomolecular pump is not at speed. The Vacuum LED illuminates green whenever the pressure in the analyzer chamber is at or below the value required to enable high voltages to the mass analyzer (7×10^{-4} Torr).

The Communication LED illuminates yellow when the mass spectrometer and the data system are trying to establish a communication link. The Communication LED illuminates green when the Ethernet communication link between the mass spectrometer and the data system has been made.


The System LED illuminates yellow whenever the mass spectrometer is in standby—that is, high voltage is not supplied to the ion source, mass analyzer, or ion detection system, but the mass spectrometer power is on. The System LED illuminates green whenever the high voltage is enabled and the system is in the On state. High voltage is enabled if the analyzer chamber is below 7×10^{-4} Torr.

The Scan LED flashes blue whenever the mass spectrometer is on and is scanning ions.

The syringe pump LED illuminates green whenever the syringe pump is pumping. The LED illuminates yellow if the syringe pump is at the end of its travel.

Two additional LEDs and a push-button switch are located on the front panel above the divert/inject valve. See [Figure 11](#). When the divert/inject valve is set up for loop injections, the divert/inject valve button switches the valve between load and inject modes and the labels *Load* and *Inject* apply.

Figure 11. Divert/inject valve button and LEDs

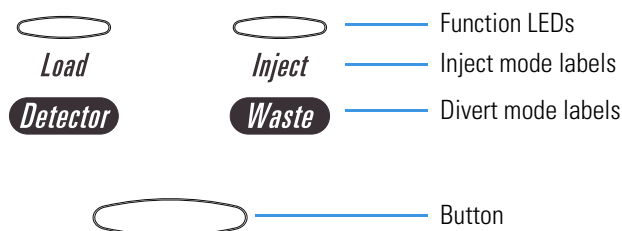

When the divert/inject valve is set up for divert valve operation, the divert/inject valve button switches the LC flow between the mass spectrometer and the waste container and the labels *Detector* and *Waste* apply.

The main power circuit breaker switch (labeled Main Power) is located on the power panel at the lower right corner of the right side panel of the mass spectrometer. See [Figure 12](#) and [Figure 13](#). In the Off (O) position, the circuit breaker removes all power to the mass spectrometer, including the vacuum pumps. In the On (I) position, power is supplied to the mass spectrometer. In the standard operational mode, the circuit breaker is kept in the On (I) position.

Figure 12. Power panel of the TSQ Vantage or TSQ Quantum Ultra mass spectrometer


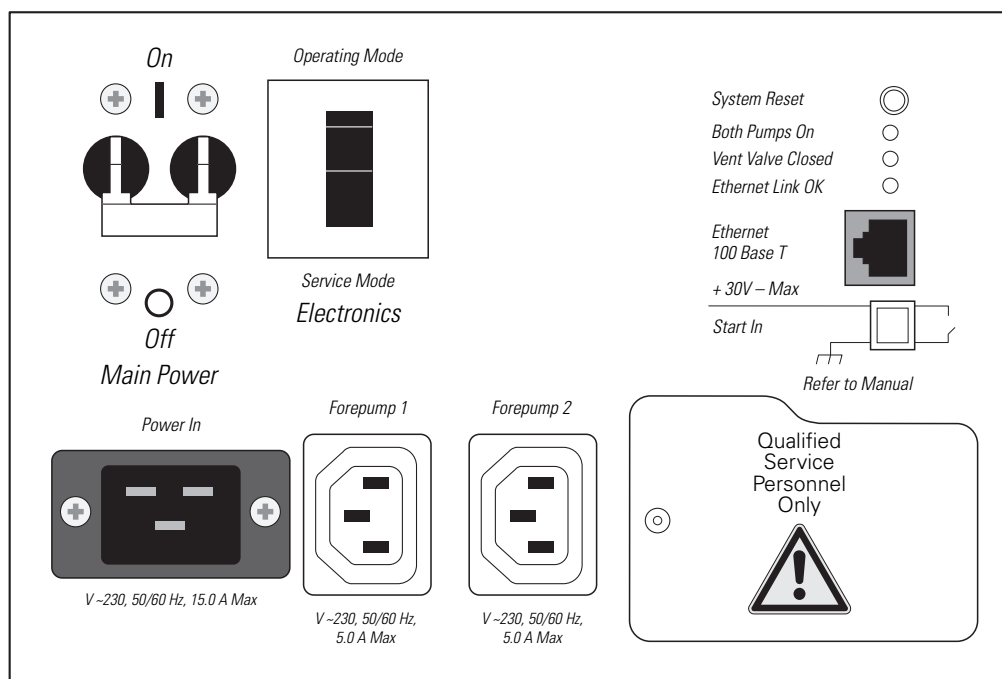


Figure 13. Power panel of the TSQ Quantum Access or TSQ Quantum Access MAX mass spectrometer, showing switches and LEDs

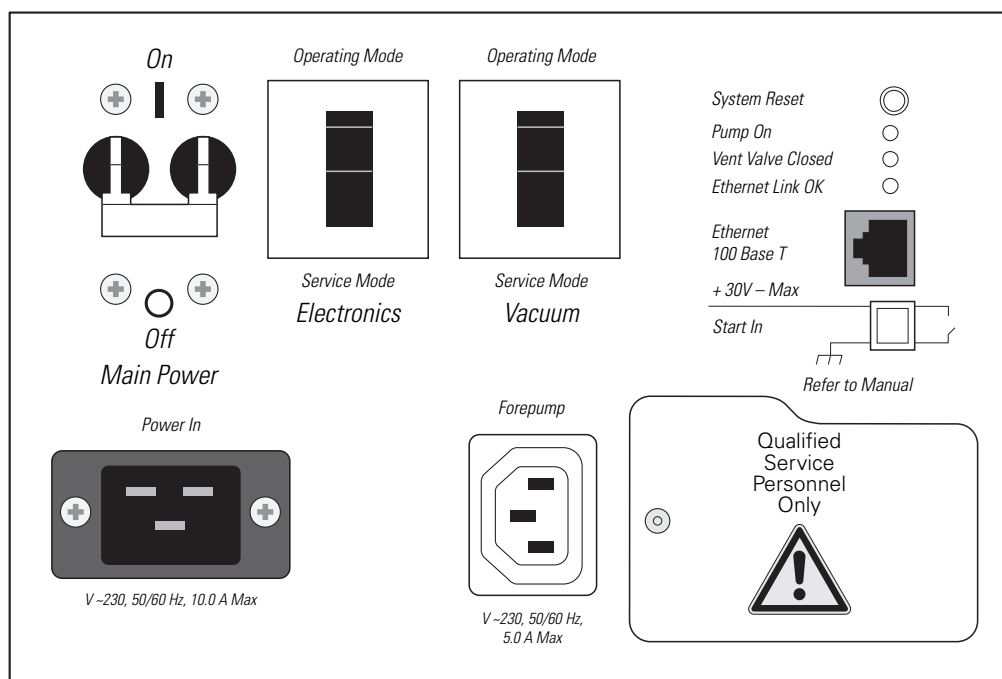

The electronics service switch (labeled Electronics) is located next to the main power circuit breaker on the mass spectrometer power panel (Figure 12 and Figure 13). In the Service Mode position the switch removes power to all components of the mass spectrometer other than the vacuum system. The Operating Mode position supplies power to all non-vacuum system components of the mass spectrometer.

On the TSQ Quantum Access and TSQ Quantum Access MAX mass spectrometers, a vacuum service switch (labeled Vacuum) is located next to the Electronics service switch on the mass spectrometer power panel (Figure 13). In the Service Mode position the switch removes power to all components of the vacuum system, including the forepump, turbomolecular pump, and turbomolecular pump controller. The switch in the Operating Mode position supplies power to all vacuum system components of the mass spectrometer.

The System Reset button is also located on the mass spectrometer power panel. Pressing the System Reset button causes the embedded computer on the System Control PCB to reboot. The TSQ mass spectrometer software is then reloaded from the data system. See “[Resetting the Mass Spectrometer](#)” on [page 57](#) for information on how to reset the mass spectrometer.

Three LEDs are located on the power panel:

- The Pump On LED illuminates green when the forepump current sensor detects current to the forepump. The LED is off when the forepump current sensor does not detect current to the forepump. If the current sensor detects a loss of current when the TSQ mass spectrometer is on, high voltage turns off and the vacuum system vents.

- The Vent Valve Closed LED illuminates green whenever the vent valve current sensor detects current through the vent valve and the vent valve is closed. The LED is off when the vent valve is open.
- The Ethernet Link OK LED illuminates green when the System Control PCB is communicating with the data system PC. The LED is off when there is no communication between the System Control PCB and the data system PC.



CAUTION In an emergency, to shut off all power to the mass spectrometer, place the Main Power circuit breaker switch (labeled Main Power) in the Off (O) position.


API Source

The atmospheric pressure ionization (API) source forms gas phase sample ions from sample molecules that elute from the LC or are introduced by the syringe pump. You can operate the API source in the electrospray ionization (ESI), heated electrospray ionization (H-ESI), nanospray ionization (NSI), atmospheric pressure photo ionization (APPI), or atmospheric pressure chemical ionization (APCI) mode. For more information regarding the API source, refer to the *Ion Max and Ion Max-S API Source Hardware Manual*, *HESI-II Probe User Guide*, *H-ESI Probe User Guide*, *Ion Max APPI Source Operator's Manual*, or *Nanospray Ion Source Operator's Manual*.

Ion Source Interface

The ion source interface consists of the components of the API source that are held under vacuum (except for the atmospheric pressure side of the ion sweep cone). The ion source interface includes an ion transfer tube, two cartridge heaters, a heater block, a platinum probe sensor, a vent prevent ball, and an ion sweep cone. See [Figure 14](#).

Figure 14. Cross-sectional view of the ion source interface, Q00 ion optics, and Q0 ion optics of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra mass spectrometer

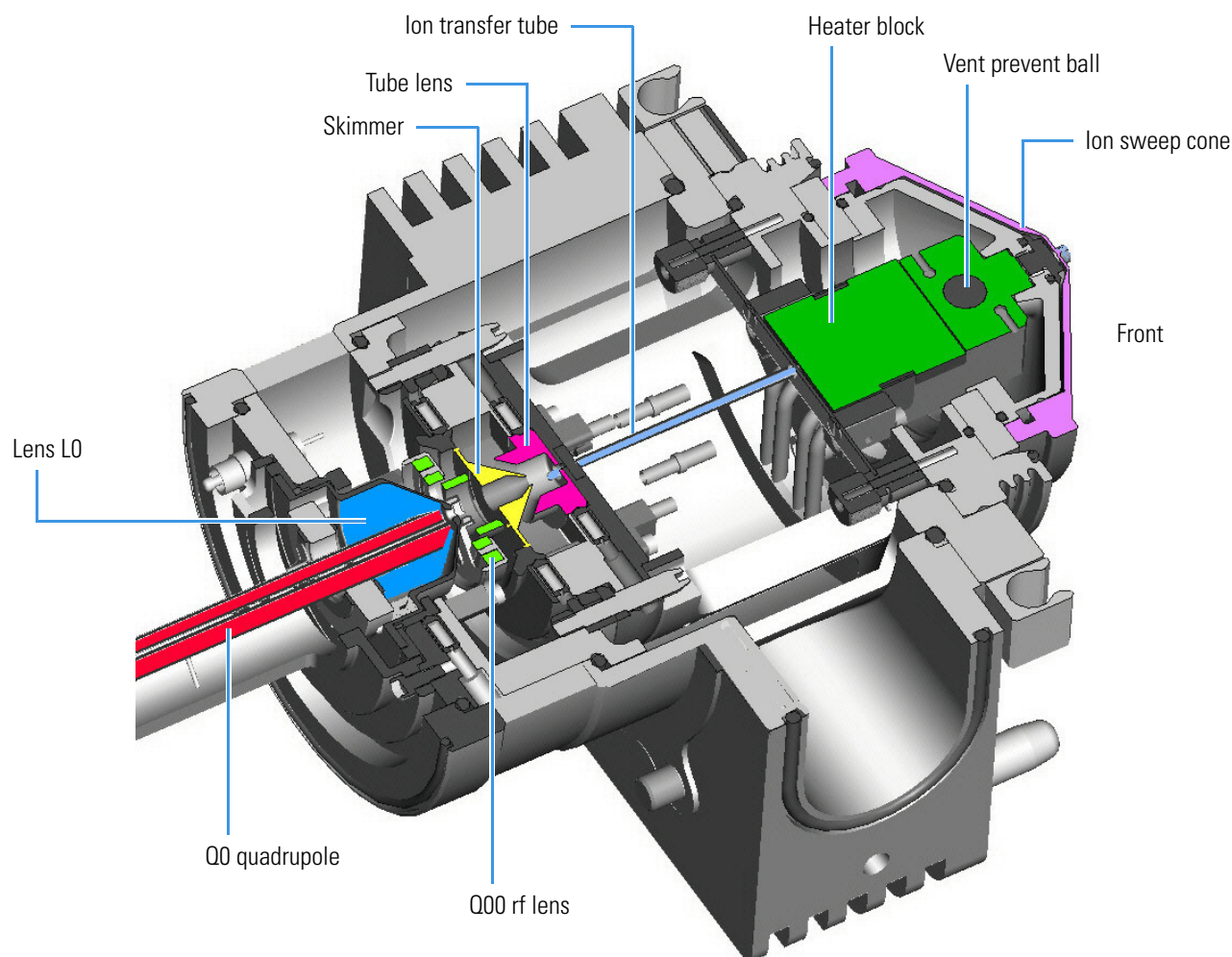


The ion transfer tube assists in desolvating ions that are produced by the API probe. The tube is an elongated, 4 in. (on the TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Quantum Ultra mass spectrometers) or 2.3 in. (on the TSQ Vantage mass spectrometer) cylindrical metal tube. See [Figure 15](#). Two heater cartridges are embedded in the heater block. The heater block surrounds the ion transfer tube and heats it to temperatures up to 400 °C. A platinum probe sensor measures the temperature of the heater block. Typical temperatures of the ion transfer tube are 270 °C for ESI and 250 °C for APCI, but these temperatures vary with flow rate and mobile phase composition. A decreasing pressure gradient draws ions into the ion transfer tube in the atmospheric pressure region and transports them to the ion transfer tube-skimmer region of the vacuum manifold. A potential of typically ± 35 V (positive for positive ions and negative for negative ions) assists in repelling ions from the ion transfer tube to the skimmer. The vent prevent ball falls into the space occupied by the ion transfer tube when the tube is removed, thus preventing air from entering the vacuum manifold. The vent prevent ball allows you to remove the ion transfer tube for cleaning without venting the system.

The ion sweep cone is a metallic cone over the ion transfer tube. The ion sweep cone channels the sweep gas towards the entrance of the tube.

The ion source interface is enclosed in a vacuum chamber that is evacuated by the forepump to a pressure of approximately 1.5 Torr.

Figure 15. Ion transfer tube

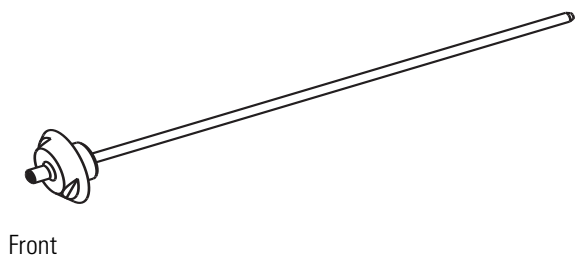

Ion Optics

The ion optics focus the ions produced in the API source and transmit them to the mass analyzer. The TSQ mass spectrometer uses two ion optics elements:


- Q00 ion optics
- Q0 ion optics

Q00 Ion Optics

The Q00 ion optics are the ion optics that are located closest to the API source. Q00 ion optics include either the tube lens and skimmer or the S-lens and exit lens, the Q00 rf lens, the ion source interface cage, and the lens L0. See [Figure 14](#).


On the TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Quantum Ultra mass spectrometers, ions from the ion transfer tube enter the tube lens. See [Figure 16](#). The tube lens has a mass-dependent potential applied to it to focus the ions toward the opening of the skimmer. An additional potential of between 0 and ± 250 V (positive for positive ions and negative for negative ions), called the tube lens offset voltage, can be applied to the tube lens. This accelerates the ions into background gas that is present in the ion transfer tube-skimmer region. Collisions with the background gas aid in the desolvation of the ions and increase sensitivity. If the tube lens offset voltage is too high, however, collisions with the background gas can be energetic enough to cause ion fragmentation. This fragmentation, called ion source collision-induced dissociation (CID), decreases sensitivity. When you tune the TSQ mass spectrometer, you adjust the tube lens offset voltage to maximize sensitivity by balancing desolvation with fragmentation.

Figure 16. Tube lens of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra mass spectrometer

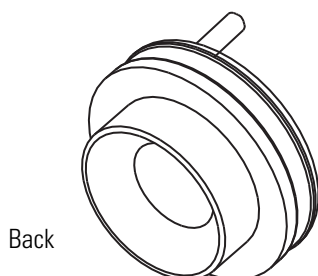

The TSQ Vantage mass spectrometer uses an S-lens and exit lens rather than a tube lens and skimmer. The S-lens is an ion transmission device consisting of progressively-spaced, stainless-steel electrodes. See [Figure 17](#). A 650 kHz rf voltage is applied to the electrodes, and adjacent electrodes have voltages of opposite phase. As the rf amplitude increases, ions of progressively higher mass-to-charge ratios pass through to the exit lens. During the tune procedure, the TSQ Vantage mass spectrometer determines the mass-dependent rf amplitudes for optimum transmission of ions through the lens. The maximum amplitude is 300 V peak-to-peak.

Figure 17. S-lens of the TSQ Vantage mass spectrometer

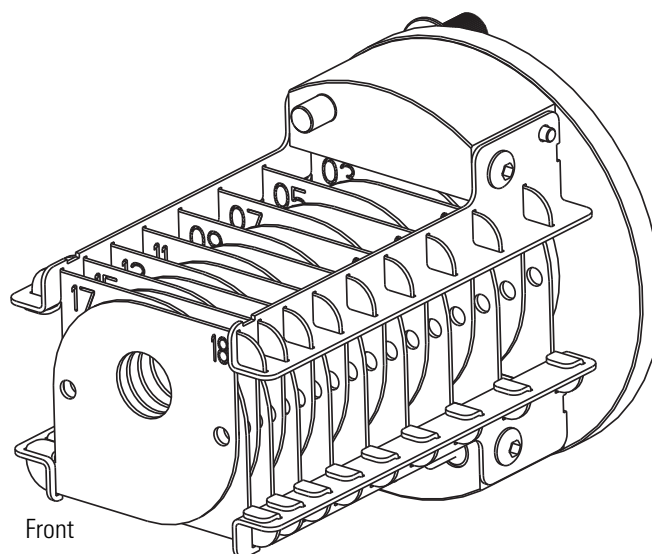

Ions from the tube lens or S-lens pass through the skimmer ([Figure 18](#)) or exit lens ([Figure 19](#)) and move toward the Q00 rf lens. The skimmer and exit lens act as a vacuum baffle between the higher pressure ion source interface region (at 1.5 Torr) and the lower pressure Q00 ion optics region (at 50 mTorr) of the vacuum manifold. The apertures of the skimmer and exit lens are offset with respect to the bore of the ion transfer tube. This arrangement reduces the number of large, charged particles that pass through the skimmer or exit lens into the mass analyzer, which reduces detector noise. The tube lens and skimmer, or the S-lens and exit lens, mount to the ion source interface cage.

Figure 18. Skimmer of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra mass spectrometer



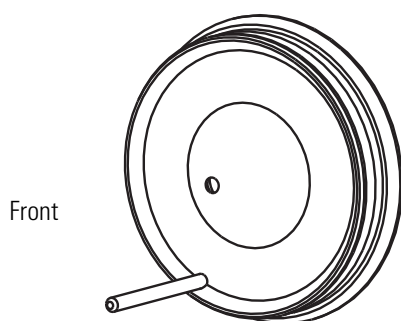


Figure 19. Exit lens of the TSQ Vantage mass spectrometer

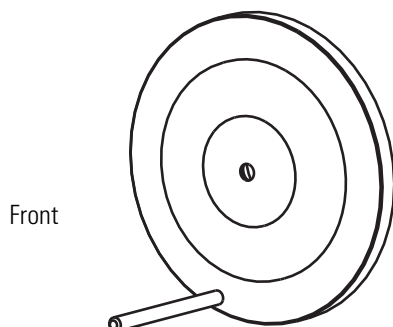

The Q00 rf lens is a square array of thin metal elements that act as an ion focusing device. See [Figure 20](#). An rf voltage that is applied to the elements gives rise to an electric field that focuses the ions along the axis of the lens. A dc voltage offset from ground applied to Q00 (called the Q00 offset voltage) increases the translational kinetic energy of ions emerging from the skimmer. During ion focusing, the offset voltage is negative for positive ions and positive for negative ions. Increasing the offset voltage increases the translational kinetic energy of the ions. Typical values of the Q00 offset voltage are 0 and ± 4 V (negative for positive ions and positive for negative ions).

Figure 20. Q00 rf lens

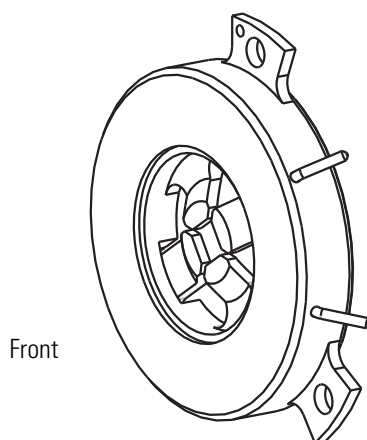

Source CID can also be accomplished by increasing the dc voltage offset between the skimmer, which is at ground potential, and the rest of the mass spectrometer beginning with Q00. You use the Source CID parameter in the Tune Master software to adjust this potential.


The lens L0 is a metal cylinder with a small hole in one end through which the ion beam can pass. A potential of between 0 and ± 3 V (negative for positive ions and positive for negative ions) is applied to lens L0 to aid in ion transmission. Lens L0 also acts as a vacuum baffle between the Q00 and Q0 ion optics chambers.

Q0 Ion Optics

The Q0 ion optics transmits ions from the Q00 ion optics to the mass analyzer. The Q0 ion optics includes the Q0 quadrupole and lenses L11 and L12.

The Q0 quadrupole is a square array of square-profile rods that act as an ion transmission device. See [Figure 21](#). An rf voltage applied to the rods gives rise to an electric field that guides the ions along the axis of the quadrupole. The Q0 offset voltage increases the translational kinetic energy of ions emerging from Q00.


Figure 21. Q0 quadrupole

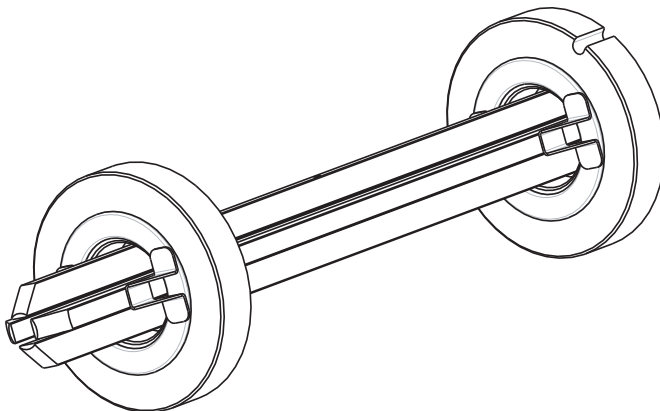

The Q0 offset voltage can also be used to fragment ions. In ion source collision-induced dissociation, the offset voltage (typically -30 V to $+30$ V) imparts enough kinetic energy to the ions so that, when they collide with solvent or air molecules, the ions dissociate to form product ions. The translational kinetic energy imparted to the ions determines the degree of dissociation. At low energies, adduct ions convert to sample ions by collision-induced dissociation without fragmenting. At higher energies, molecular ions can fragment and thereby enable primitive MS/MS/MS analysis.

The L11 and L12 lenses are metal disks with a circular hole in the center through which the ion beam can pass. See [Figure 22](#). Together they act as a two-element cone lens. An electrical potential applied to the lens accelerates (or decelerates) ions as they approach the lens to focus the ion beam as it passes through the lens. The value ranges between 0 and ± 300 V. Lenses L11 and L12 also act as a vacuum baffle between the Q0 ion optics chamber and the mass analyzer chamber.


Figure 22. Lenses L11 (left) and L12 (right)

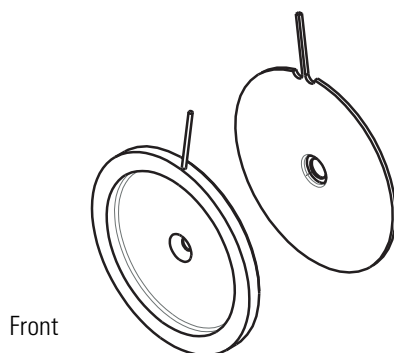

Mass Analyzer

The mass analyzer separates ions according to their mass-to-charge ratio and then passes them to the ion detection system. The mass analyzer on the TSQ mass spectrometer consists of three quadrupole rod assemblies (Q1, Q2, and Q3) and three lens sets. See [Figure 23](#).

Figure 23. Mass analyzer, ion detection system, and ion optics

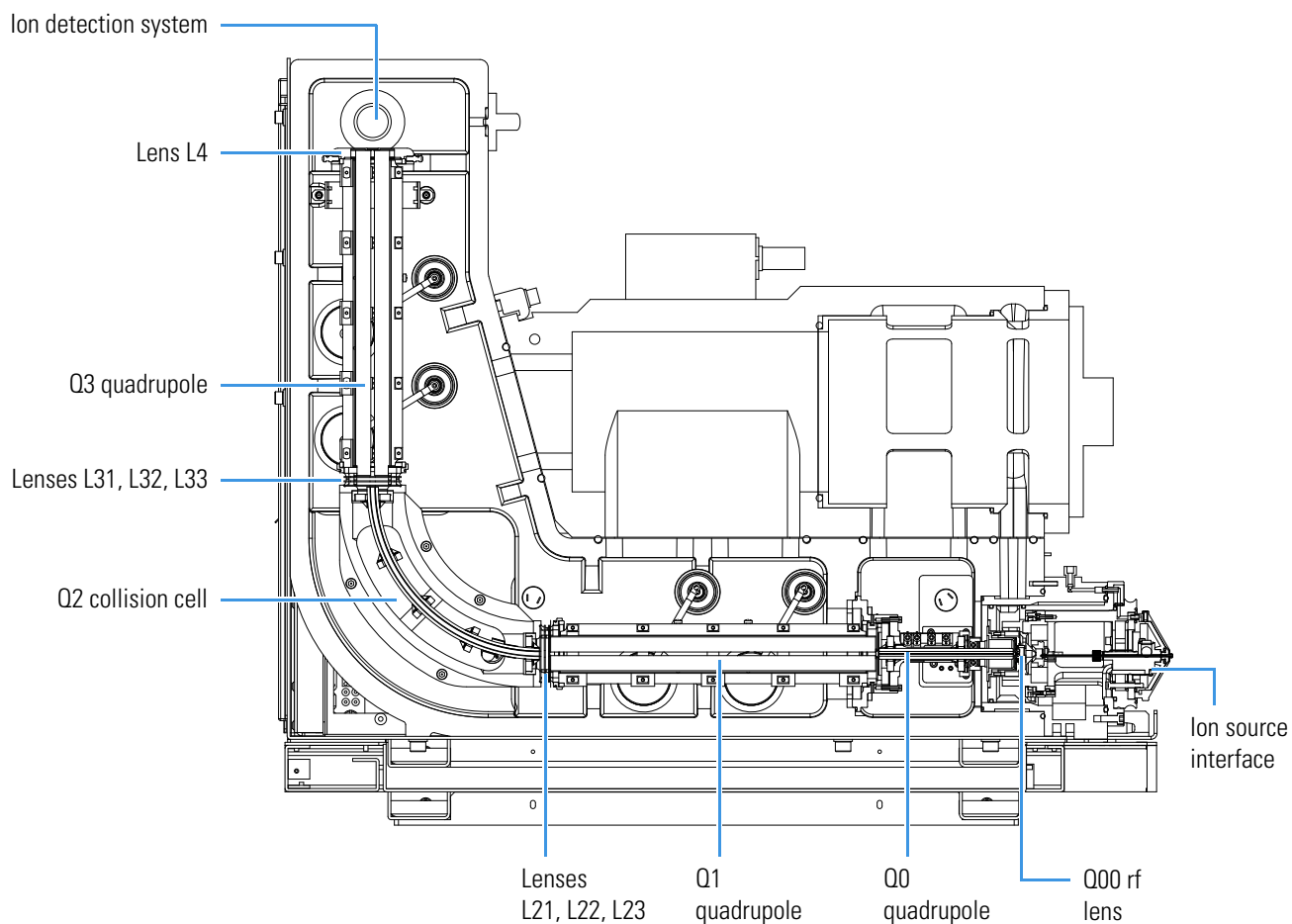

The following subtopics discuss the mass analyzer in detail:

- [Quadrupole Rod Assemblies](#)
- [RF and DC Fields Applied to the Quadrupoles](#)
- [Mass Analysis](#)
- [Collision Cell and CID Efficiency](#)
- [Quadrupole Offset Voltage](#)
- [Mass Analyzer Lenses](#)

Quadrupole Rod Assemblies

The three rod assemblies used in the TSQ mass spectrometer are numbered from the ion source end of the manifold and are designated Q1, Q2, and Q3. Q1 and Q3 are quadrupoles that enable high-resolution scans without signal loss.

Q1 and Q3 are square arrays of precision-machined and precision-aligned, true hyperbolic quadrupoles, called hyperquads (TSQ Vantage and TSQ Quantum Ultra mass spectrometers) or hyperbolic-profile round rods (TSQ Quantum Access and TSQ Quantum Access MAX mass spectrometers). See [Figure 24](#) and [Figure 25](#). Quartz spacers act as electrical insulators between adjacent rods.

Figure 24. Q1 or Q3 hyperquad of the TSQ Vantage or TSQ Quantum Ultra mass spectrometer



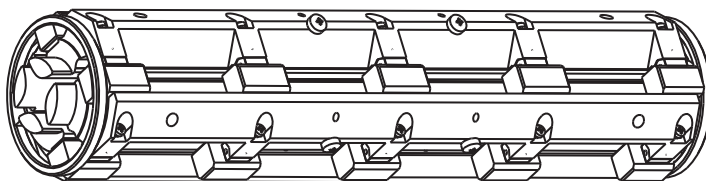


Figure 25. Q1 or Q3 quadrupole of the TSQ Quantum Access, or TSQ Quantum Access MAX mass spectrometer

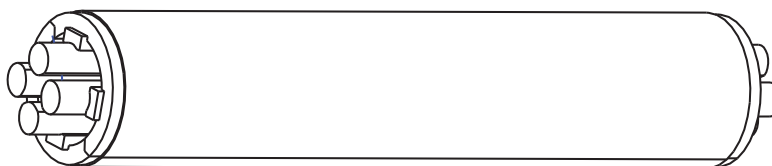

Q2 is a square-profile quadrupole rod assembly. Q2 always acts as an ion transmission device. The Q2 quadrupole rods are bent through a 90-degree arc. In addition to reducing the footprint of the instrument, this bending of the rods prevents the transmission of unwanted neutral species to the detector and dramatically lowers the noise level in the data. Q2 has become synonymous with the term *collision cell*. Technically, the collision cell is the chamber that encloses Q2 where collision-induced dissociation can take place if the argon collision gas is present. See [Figure 26](#).


Figure 26. Collision cell (top), Q2 quadrupole (bottom), and lenses

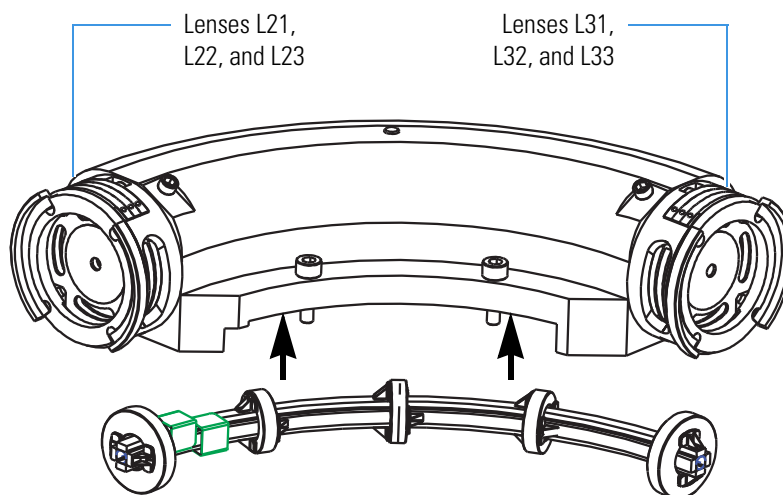

RF and DC Fields Applied to the Quadrupoles

In a quadrupole rod assembly, because rods opposite each other in the array connect electrically, the four rods can be considered as two pairs of two rods each. Ac and dc voltages are applied to the rods and these voltages are ramped during the scan. Voltages of the same amplitude and sign are applied to the rods of each pair. However, the voltages applied to the different rod pairs are equal in amplitude but opposite in sign. See [Figure 27](#).

Figure 27. Polarity of the rf and dc voltages applied to the rods of the Q1 and Q3 mass analyzers

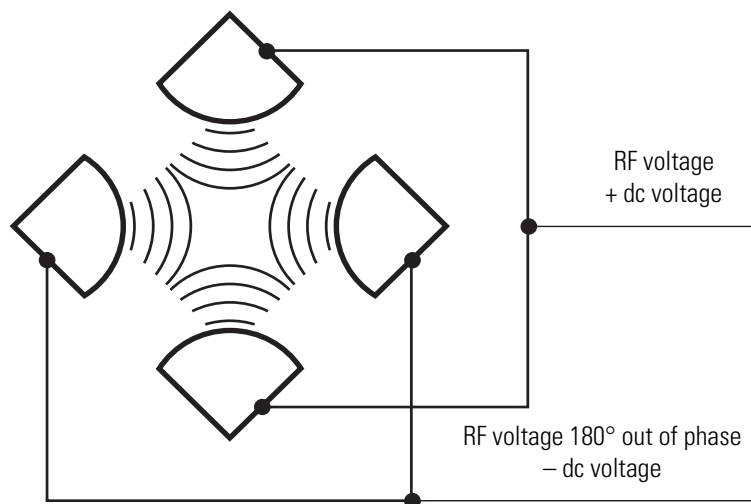


The ac voltage applied to the quadrupole rods is of constant frequency (1.123 MHz). The rf voltage applied to the rods varies from 0 to 10000 V peak-to-peak, and the dc voltage varies from 0 to ± 840 V. Voltages of the same amplitude and sign are applied to each rod pair. However, the voltages applied to the other rod pair are equal in amplitude but opposite in sign.

Because the frequency of this ac voltage is in the radio frequency range, it is referred to as rf voltage. In Figure 28, the solid line represents the combined rf and dc voltage applied to one rod pair, and the dashed line represents the combined rf and dc voltage applied to the other rod pair. The ratio of rf voltage to dc voltage determines the ability of the mass spectrometer to separate ions of different mass-to-charge ratios.

The first and third quadrupole rod assemblies (Q1 and Q3 quadrupoles) can act as mass analyzers or as ion transmission devices. When both rf and dc voltages are applied, Q1 and Q3 function as mass analyzers. When only rf voltage is applied, they act as ion transmission devices. In the ion transmission mode, the quadrupole rod assemblies allow ions in a wide window of mass-to-charge ratios to pass.

The square quadrupole rod assembly (Q2) operates in the ion transmission mode only. Surrounding Q2 is a collision cell where collision-induced dissociation (CID) can take place if the argon collision gas is present in the cell.

Figure 28. Magnitude of the rf and dc voltages applied to the rods of the Q1 and Q3 mass analyzers

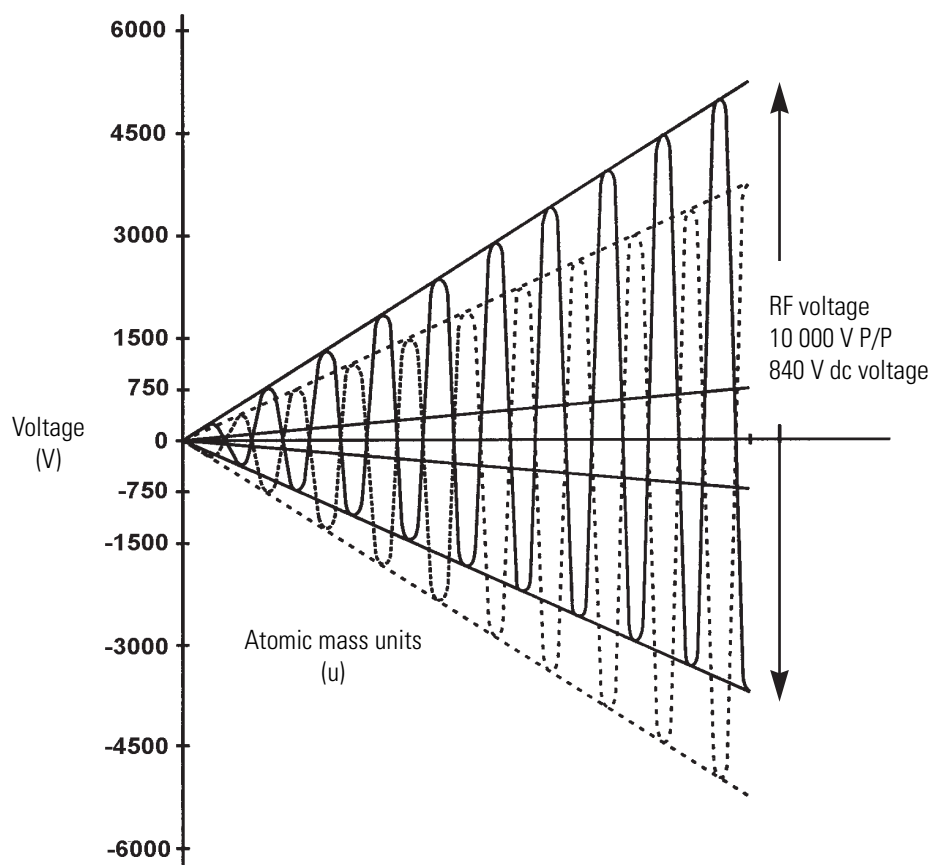

Mass Analysis

The mass analyzers (Q1 and Q3) are square arrays of precision-machined and precision-aligned hyperbolic or round-profile rods. The rods are charged with a variable ratio of rf voltage and dc voltage (Figure 28). These potentials give rise to an electrostatic field that gives stable oscillations to ions with a specific mass-to-charge ratio and unstable oscillations to all others.

At any given instant, one particular set of rf and dc voltage values is being applied to the mass analyzer rods. Under these conditions, only ions of one mass-to-charge ratio (for example, m/z 180) are maintained within bounded oscillations as their velocity carries them through the mass analyzer. At the same time, all other ions undergo unbounded oscillations. These ions strike one of the rod surfaces, become neutralized, and are pumped away, or they are ejected from the rod assembly.

Then, at a later time, both rf and dc voltages change, and ions of the next mass-to-charge ratio (for example, m/z 181) are allowed to pass, while all other ions (including m/z 180) become unstable and undergo unbounded oscillations. This process continues, with ions of one mass-to-charge ratio after another being transmitted, as the rf and dc voltages change in value. At the end of the scan, the rf and dc voltages are discharged to zero, and the process repeats.

The TSQ system can rapidly and precisely change the potentials on the quadrupole rods. The rf and dc voltages in the TSQ mass spectrometer can be scanned over the full mass range of the system, m/z 10 to 3000, in 0.85 seconds.

The more closely the electrostatic field generated by a set of quadrupole rods approximates a hyperbolic geometry, the better their operating characteristics are. As a result, the precision quadrupole rods of the TSQ mass spectrometer provide excellent sensitivity, peak shape, resolution, and high mass transmission.

Collision Cell and CID Efficiency

In the MS/MS scan modes, the TSQ mass spectrometer applies a large voltage of opposite polarity to the rod pairs between scans, which empties the collision cell. This process ensures that no ions remain in the collision cell from scan to scan.

The collision cell quadrupole rod assembly (Q2), which always acts as an ion transmission device, is a quadrupole array of square-profile rods. A variable rf voltage charges the rods, which creates an electrostatic field that gives stable oscillations to ions in a wide window of mass-to-charge ratios.

The collision cell surrounds Q2 and is usually pressurized from about 1×10^{-3} to 4×10^{-3} Torr with argon collision gas. The collision cell is where collision-induced dissociation takes place.

CID is a process in which an ion collides with a neutral atom or molecule and then, because of the collision, dissociates into smaller fragments. The mechanism of dissociation involves converting some of the translational kinetic energy (TKE) of the ion into internal energy. This collision places the ion in an excited state. If the internal energy is sufficient, the ion fragments.

Three expressions convey the efficiency of the CID process:

- Collection efficiency
- Fragmentation efficiency
- Overall CID efficiency

Collection efficiency: The ion flux ratio measured at the exit of the collision cell and at its entrance. With no collision gas present, the TSQ mass spectrometer obtains virtually 100 percent collection efficiency. Collection efficiency is a mass-dependent parameter. For example, with mid-range collision gas pressure, the collection efficiency might vary from about 50 percent for comparatively less massive ions (which are more prone to scatter) up to 75 percent for comparatively more massive ions (which are less prone to scatter).

Fragmentation efficiency: The fraction of the ion flux at the exit of the collision cell that results from fragmented ions. Fragmentation efficiency depends directly on the stability of the ion and indirectly on the mass of the ion. The more stable the ion, the less likely a given collision will fragment the ion. The more massive the ion, the greater its ability to distribute the vibrational energy imparted by a collision. As a result, ion fragmentation might decrease.

With mid-range collision gas pressure, fragmentation efficiency might vary from 15 percent to 65 percent for various compounds. As the collision gas pressure increases, the fragmentation efficiency for all compounds approaches 100 percent due to multiple collisions. The collection efficiency decreases, however, due to scattering.

Overall CID efficiency: The product of the collection efficiency and the fragmentation efficiency. The overall CID efficiency exhibits a maximum with intermediate pressure. As the pressure increases beyond the optimum value, more and more collisions take place, the probability of scattering increases, and fewer and fewer ions pass through the collision cell. Scattering results in the collection efficiency decreasing. The fragmentation efficiency also decreases as the pressure is decreased from its optimum value, because fewer and fewer collisions take place.

Quadrupole Offset Voltage

The quadrupole offset voltage is a dc potential applied to the quadrupole rods in addition to the ramping dc voltage. The offset voltage applied to the two rod pairs of the assemblies is equal in amplitude and equal in sign. The quadrupole offset voltage accelerates or decelerates ions and, therefore, sets the TKE of the ions as they enter the quadrupole rod assembly.

In general, for a given experiment, the TSQ mass spectrometer has fixed offset voltages for Q1 and Q2. However, in MS/MS experiments, the quadrupole offset voltage applied to Q3 usually varies as a scan proceeds. The TSQ mass spectrometer automatically computes the Q3 quadrupole offset voltage necessary and then varies the voltage, as appropriate, as each scan proceeds.

The offset voltage applied to Q2 (which contains the collision cell) is responsible for the collision energy. The collision energy is the difference in potential between the ion source (where parent ions are formed) and Q2 (where they collide with collision gas). As the offset voltage on Q2 increases, the TKE of the parent ions also increases. As a result, increases in the Q2 offset voltage increase the energy of ion/Ar collisions. The collision energy is generally set to one value for an entire scan and can be set from 0 to ± 200 V.

Before obtaining any mass spectra, the TSQ mass spectrometer tunes Q1 in the Q1MS scan mode (Q2 and Q3 rf voltage only), and tunes Q3 in the Q3MS scan mode (Q1 and Q2 rf voltage only). During tuning, the TSQ mass spectrometer determines the optimum quadrupole offset voltage for Q1 and for Q3.

Mass Analyzer Lenses

The TSQ system mass analyzer has three lens sets. See [Figure 26 on page 30](#). Those between Q1 and Q2 are designated L21, L22, L23; those between Q2 and Q3 are designated L31, L32, L33; and the lens between Q3 and the ion detection system is designated as L4 (or the exit lens). All of the lenses have circular holes in their centers through which the ion beam passes.

The lens assemblies also retain the three rod assemblies to ensure accurate and automatic axial alignment of the rod assemblies.

The L2x lens set (between Q1 and Q2) and the L3x lens set (between Q2 and Q3) serve these functions:

- To minimize the amount of collision gas that enters the mass analyzers (Q1 and Q3) from the collision cell (Q2). (For high-mass transmission, it is important to maintain a low pressure in the mass analyzers.)
- To retain the collision gas. Lenses L23 and L31 form two of the walls of the collision cell, so they tend to hold the collision gas in the collision cell. The collision gas escapes, however, through the same lens holes through which the ion beam passes.
- To prevent gas from entering the mass analyzers. Lenses L22 and L21 on one side of Q2 and lenses L32 and L33 on the other side of Q2 act as baffles to help prevent the gas that escapes from the collision cell from entering the mass analyzers.
- To shield Q1 from the rf voltage applied to Q2 and vice versa (L2x lens set) and to shield Q3 from the rf voltage applied to Q2 and vice versa (L3x lens set).
- To focus the ion beam. The three lenses between Q1 and Q2 (and those between Q2 and Q3) together form a three-element aperture lens. The first and third lenses are generally set to similar or identical values and the central lens is set to a value different (either higher or lower) from the other two.

The voltage applied to each of the lenses can vary from about -300 to $+300$ V. Typically, however, the voltage applied to the first and third elements of the L2x lens set is somewhat greater than the quadrupole offset voltage applied to Q1. Because the Q1 quadrupole offset voltage is generally set to about ± 5 V (depending on the charge of the ions of interest), the voltage applied to lenses L21 and L23 is typically about -10 V for positive ions and $+10$ V for negative ions. The voltage applied to the central lens of the L2x lens set is typically about ± 225 V.

In the Q3MS scan mode, the voltage applied to the lenses of the L3x lens set is about the same as that applied to the corresponding lens in the L2x lens set. Note, however, that in the MS/MS scan modes, the voltage applied to the L3x lens set automatically varies with the quadrupole offset voltage applied to Q3. As the Q3 quadrupole offset voltage ramps, the voltages applied to the lenses ramp correspondingly.

Lens L4 is located between Q3 and the ion detection system. L4 is held at ground potential. Its purpose is to shield Q3 from the high voltage applied to the ion detection system and to shield the ion detection system from the high rf voltages applied to Q3.

Ion Detection System

The TSQ mass spectrometer is equipped with a high-sensitivity, off-axis ion detection system. This system produces a high signal-to-noise ratio and allows for voltage polarity switching between positive ion and negative ion modes of operation. The ion detection system includes a 15 kV conversion dynode and a channel electron multiplier. See [Figure 29](#). The ion detection system is located at the back of the vacuum manifold behind the mass analyzer.

The conversion dynode is a concave metal surface located at a right angle to the ion beam. The TSQ mass spectrometer applies to the conversion dynode a potential of $+15$ kV for negative ion detection or -15 kV for positive ion detection. When an ion strikes the surface of the conversion dynode, one or more secondary particles are produced. These secondary particles can include positive ions, negative ions, electrons, and neutrals. When positive ions strike a negatively charged conversion dynode, the secondary particles of interest are negative ions and electrons. When negative ions strike a positively charged conversion dynode, the secondary particles of interest are positive ions. The curved surface of the conversion dynode focuses these secondary particles and a voltage gradient accelerates the particles into the electron multiplier.


The electron multiplier includes a cathode and an anode. The cathode of the electron multiplier is a lead-oxide, funnel-like resistor. The high voltage ring applies a potential of up to -2.5 kV to the cathode. The exit end of the cathode (at the anode) is near ground potential.

The anode of the electron multiplier is a small cup located at the exit end of the cathode. The anode collects the electrons produced by the cathode. The anode screws into the anode feedthrough in the base plate.

Secondary particles from the conversion dynode strike the inner walls of the electron multiplier cathode with sufficient energy to eject electrons. The ejected electrons are

accelerated farther into the cathode, drawn by the increasingly positive potential gradient. The funnel shape of the cathode causes the ejected electrons not to travel far before they again strike the inner surface of the cathode, which causes the emission of more electrons. A cascade of electrons is then created that finally results in a measurable current at the end of the cathode where the anode collects the electrons. The current collected by the anode is proportional to the number of secondary particles striking the cathode.

Figure 29. Ion detection system, showing the electron multiplier (top) and conversion dynode (bottom)

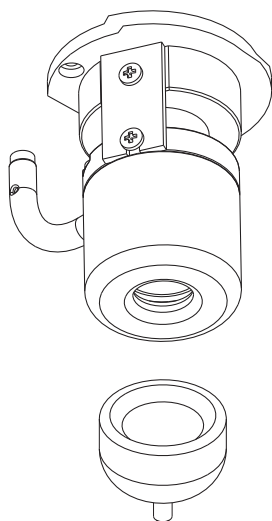

Typically, the electron multiplier is set to a gain of about 3×10^5 (that is, for each ion or electron that enters, 3×10^5 electrons exit) in MS mode and 2×10^6 in MS/MS mode. The electrometer circuit converts the current that leaves the electron multiplier via the anode to a voltage and the data system records the voltage.

The ion detection system of the TSQ mass spectrometer increases signal and decreases noise. The high voltage applied to the conversion dynode results in a high conversion efficiency and increased signal. That is, for each ion striking the conversion dynode, many secondary particles are produced. The increase in conversion efficiency is more pronounced for more massive ions than it is for less massive ions.

Because of the off-axis orientation of the ion detection system relative to the mass analyzer, neutral molecules from the mass analyzer tend not to strike the conversion dynode or electron multiplier. As a result, noise from neutral molecules is reduced.

Vacuum System and Inlet Gases Hardware

The vacuum system evacuates the region around the ion source interface, ion optics, mass analyzer, and ion detection system. The principal components of the vacuum system include the following:


- Vacuum manifold
- Turbomolecular pump
- Forepump
- Convectron™ gauges
- Ion gauges

The inlet gases hardware controls the flow of collision gas, sheath gas, auxiliary gas, and air (during venting) into the mass spectrometer. The inlet gasses hardware includes the following components:

- Vent valve
- Collision gas flow control valves
- Sheath gas valve
- Auxiliary gas valve

A functional block diagram of the vacuum system and inlet gasses hardware is shown in [Figure 30](#).

Figure 30. Functional block diagram of the vacuum system and inlet gasses hardware

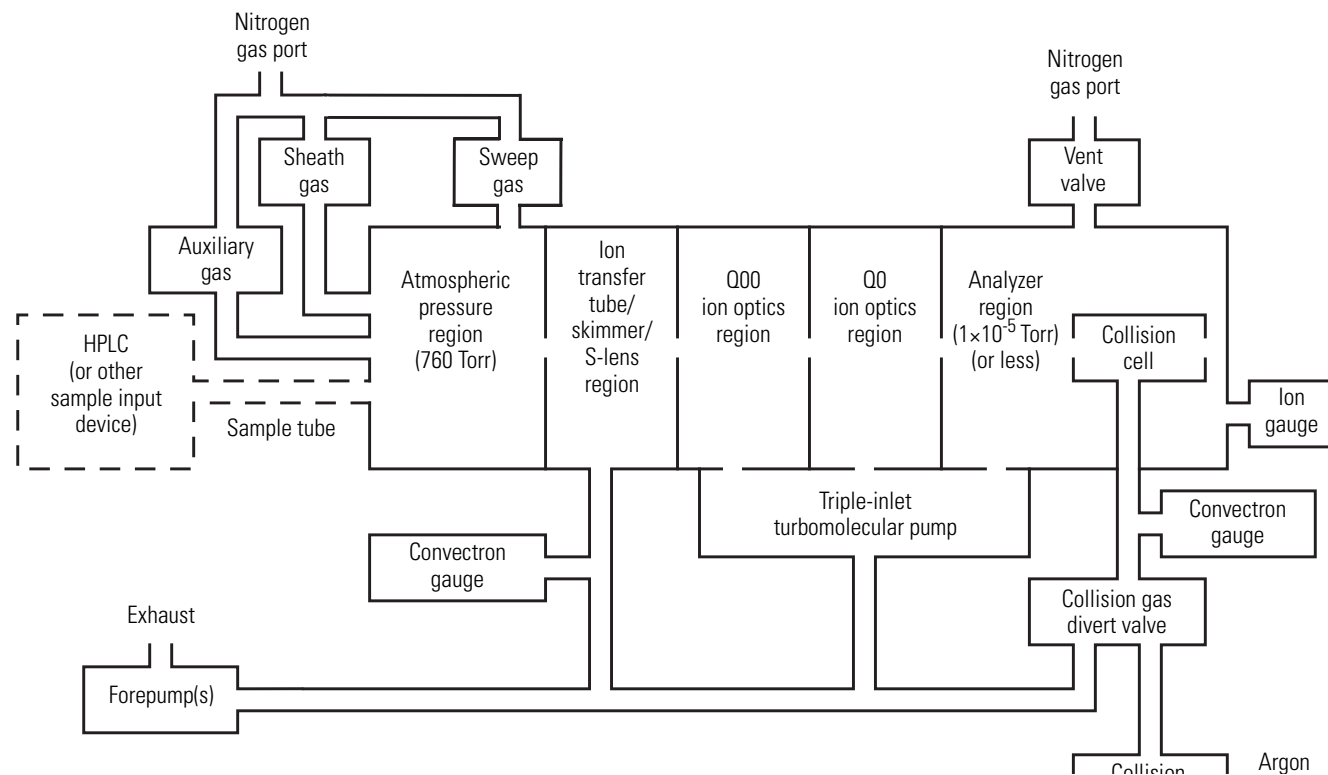


Vacuum Manifold

The vacuum manifold encloses the ion source interface, ion optics, mass analyzer, and ion detection system assemblies. The vacuum manifold is a thick-walled, aluminum chamber with two removable side cover plates, with openings on the front, sides, and top, and various electrical feedthroughs and gas inlets.

The vacuum manifold is divided into four chambers by three baffles. See [Figure 31](#). The forepump evacuates the region inside the first chamber, called the ion transfer tube/skimmer region. The third inlet in the molecular drag section of the triple-inlet turbomolecular vacuum pump evacuates the region inside the second chamber, called the Q00 ion optics region. The interstage port of the turbomolecular vacuum pump evacuates the region inside the third chamber, called the Q0 ion optics region. The high vacuum port of the turbomolecular pump evacuates the region inside the fourth chamber, called the analyzer region, to less than 10^{-5} Torr. The turbomolecular pump in turn discharges into the forepump through the foreline.

Figure 31. Vacuum manifold (interior)

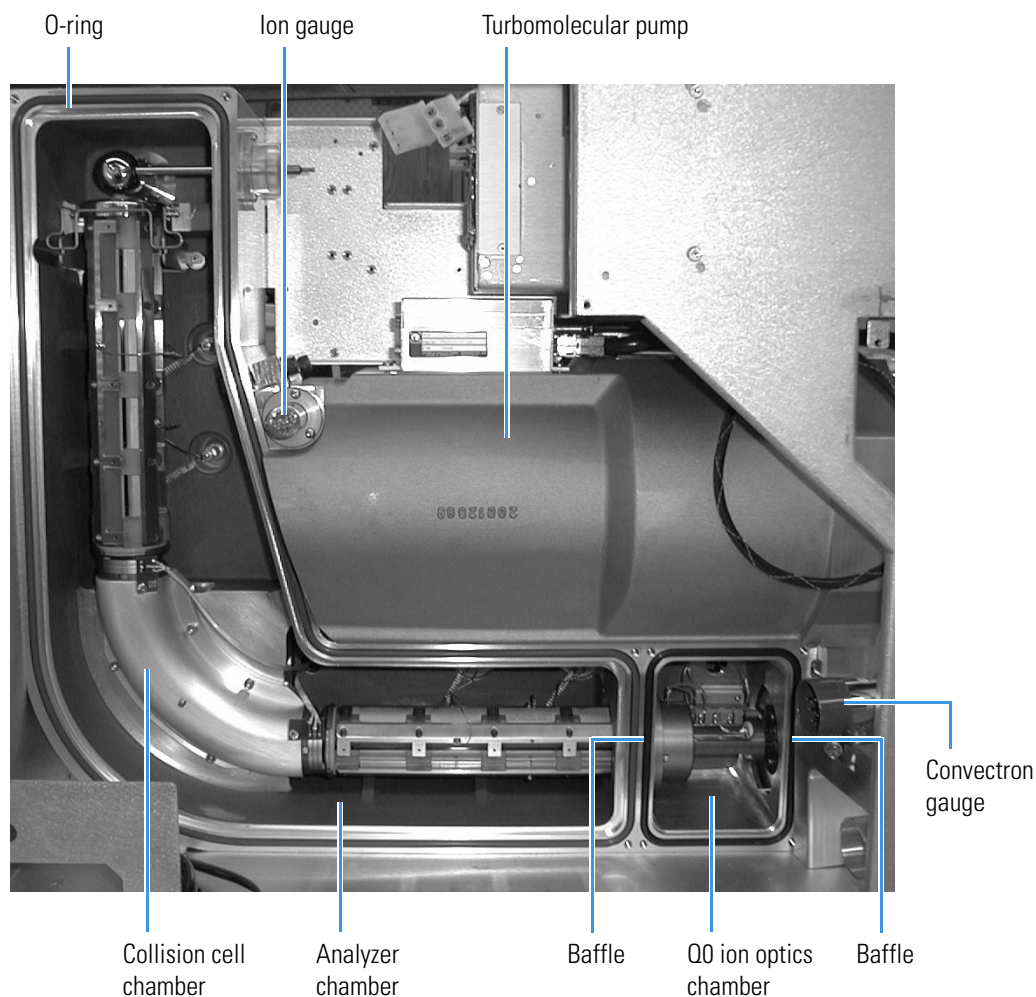

The collision cell chamber, inside the analyzer chamber, has a user controlled argon pressure of between one and four mTorr when CID is turned on. Argon in the collision cell is evacuated by the forepump when CID is turned off.

The vacuum manifold has the following feedthroughs and inlets:

- A feedthrough for the high voltage for the conversion dynode
- A feedthrough for the high voltage for the electron multiplier
- A feedthrough for the ion current signal from the anode of the electron multiplier
- Two feedthroughs for the Q1 quadrupole rf voltage
- Two feedthroughs for the Q3 quadrupole rf voltage
- A feedthrough for the Q2 quadrupole rf voltage
- A feedthrough for lens L21, L22, L23, L31, L32, and L33 voltages

- A feedthrough for lens L0, L11, and L12 voltages
- A feedthrough for the Q00 and Q0 ion optics rf voltages
- A feedthrough for the tube lens offset voltage, ion transfer tube offset voltage, and ion transfer tube heater voltages
- A vacuum connection for measuring the pressure in the analyzer region with the ion gauge
- An argon gas inlet into the collision cell
- A vent gas inlet

Two removable side cover plates on the left side of the vacuum manifold allow access to the Q0 ion optics, mass analyzer, and ion detection system. Two electrically conductive O-rings provide a vacuum-tight seal between the side cover plates and the vacuum manifold.

Turbomolecular Pump

A Leybold TW220/150/15S triple-inlet turbomolecular pump provides the vacuum for the Q00 ion optics region, Q0 ion optics region, and analyzer region of the vacuum manifold. The turbomolecular pump mounts onto the top of the vacuum manifold ([Figure 31](#)).

The turbomolecular pump has three pumping inlets:

- A high-vacuum inlet at the top of the rotor stack, which evacuates the analyzer chamber
- An interstage inlet about halfway down the rotor stack, which evacuates the Q0 ion optics chamber
- A third inlet in the molecular drag section of the pump, which evacuates the Q00 ion optics chamber

The turbomolecular pump is controlled by a Leybold TDS controller and powered by a +24 V dc (250 W) power supply. The Main Power circuit breaker switch and the Vacuum service switch, not the Electronics service switch, turn power for the turbomolecular pump off and on. The pump is air cooled by a fan that draws air in from the front of the instrument.

Power to the turbomolecular pump is shut off if the foreline pressure, as measured by the Convection gauge, is too high, or if the turbomolecular temperature is too high.

Forepumps

One (TSQ Quantum Access and TSQ Quantum Access MAX mass spectrometers) or two (TSQ Vantage and TSQ Quantum Ultra mass spectrometers) Edwards™ E2M30 forepumps (or roughing pumps) establish the vacuum necessary for the proper operation of the turbomolecular pump. The forepump also evacuates the inlet valve and the collision cell. The pump has a maximum displacement of 650 L/min and maintains a minimum pressure of approximately 1 Pa (0.01 Torr).

The forepump connects to the turbomolecular pump by a section of reinforced PVC tubing. The power cord of the forepump plugs into the outlet labeled *Forepump* on the power panel (see [Figure 13](#) on [page 21](#)). This outlet supplies power to the pump and is controlled by the Main Power circuit breaker switch and Vacuum service switch, not by the Electronics service switch.



CAUTION Always plug the forepump power cord into the outlets labeled *Forepump* on the right-side power panel of the mass spectrometer. Never plug it directly into a wall outlet.

Convectron Gauges

A Convectron gauge measures the pressure in the inlet valve and the foreline, which connects the turbomolecular pump and the forepump. The Convectron gauge uses a Wheatstone bridge with a temperature dependent resistor to measure pressure down to a fraction of a milliTor. A second Convectron gauge measures the pressure of argon collision gas in the collision cell.

Ion Gauges

A Granville-Phillips™ 342™ mini ion gauge measures the pressures in the analyzer region of the vacuum manifold. The ion gauge produces energetic electrons that cause the ionization of molecules in the ion gauge. Positive ions formed in the ion gauge are attracted to a collector. The collector current is related to the pressure in the vacuum manifold. The ion gauge is also involved in vacuum protection.

Vent Valve

The vent valve, located in the area of Q2, allows the vacuum manifold to be vented to dry nitrogen or filtered air. The vent valve is a solenoid-operated valve, which is controlled by the Vent Delay PCB. When the solenoid is energized, the vent valve closes. If the power fails or the main power circuit breaker is placed in the Off (O) position, a 4-farad capacitor, which is located in the Power Entry Module, keeps the solenoid in the closed position for several minutes. If power is not restored in this time, then the solenoid opens and the system is vented with dry nitrogen or filtered air. Dry nitrogen (35 kPa [5 psi] maximum, 99% minimum purity) enters the mass spectrometer through a 1/4 in. port on the left side of the mass spectrometer. See [Figure 32](#). The vent valve closes after power is restored to the mass spectrometer.

Collision Gas Flow Control Valves

The collision gas flow control valves control the flow of argon collision gas into and out of the Q2 collision cell. A solenoid valve acts to shut off argon gas flow to the cell. The collision gas pressure is regulated by a proportional valve that is controlled by the data system. You can set the collision gas pressure (0 to 5 mTor) in the Define Scan dialog box of the EZ Tune window.

Ions enter the Q2 collision cell, collide with the argon collision gas, and then, because of the collision, dissociate into smaller fragments. See “Collision Cell and CID Efficiency” on page 32.

Argon (135 ± 70 kPa [20 ± 10 psi], 99.995% minimum purity) enters the mass spectrometer through a 1/8-in. port on the left side of the mass spectrometer. See Figure 32.

A second proportional valve allows the forepump to evacuate the Q2 collision cell of argon and waste gases when CID is turned off. The TSQ mass spectrometer automatically opens and closes the collision gas evacuation valve, depending on whether Q2 is acting as an ion transmission device or a collision cell.


Sheath Gas Valve

The sheath gas valve controls the flow of sheath gas (nitrogen) into the API source. Dry nitrogen (690 ± 140 kPa [100 ± 20 psi], 99% purity) enters the mass spectrometer through a 1/4 in. port on the left side of the mass spectrometer. See Figure 32. A solenoid shutoff valve starts and stops the flow of nitrogen to the API source. The data system controls a valve that regulates the sheath gas pressure. You can set the sheath gas flow rate (0 to 100 in arbitrary units) in the Ion Source dialog box from the EZ Tune window. The sheath gas enters the API source through 1/8 in. ID tubing.

Auxiliary Gas Valve

The auxiliary gas valve controls the flow of auxiliary gas (nitrogen) into the API source. Dry nitrogen (690 ± 140 kPa [100 ± 20 psi], 99% purity) enters the mass spectrometer through a 1/4 in. port on the left side of the mass spectrometer. A solenoid shutoff valve starts and stops the flow of nitrogen to the API source. The auxiliary gas pressure is regulated by a flow valve that is controlled by the data system. You can set the auxiliary gas flow rate (0 to 60 in arbitrary units) in the Ion Source dialog box. The auxiliary gas enters the API source through 1/8 in. ID tubing.

Figure 32. Nitrogen and argon gas inlets

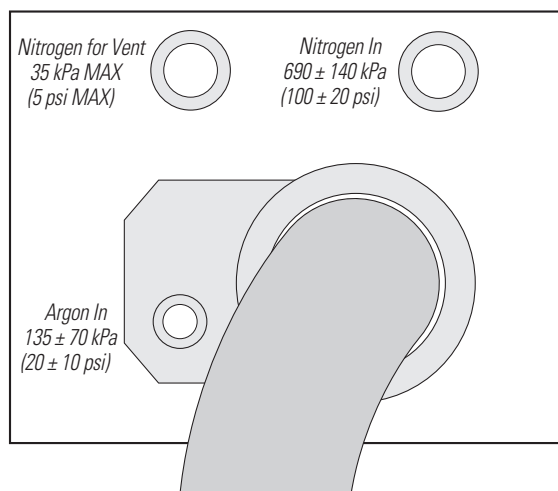

Electronic Assemblies

The electronic assemblies that control the operation of the mass spectrometer are distributed among various PCBs and other modules located in the tower, in the embedded computer, and on or around the vacuum manifold of the mass spectrometer.

The Power Entry Module provides mass spectrometer power control, a contact closure interface, vent valve control, an Ethernet 100 base-T connection from the System Control PCB to the data system PC, a mechanical pump failure protection circuit (linked to vent valve control), a system reset button, status LEDs, and service ports. The right-side power panel, shown in [Figure 13](#) on [page 21](#), is part of the Power Entry Module.

The Power Entry Module accepts line power, filters it, and provides it to various components of the mass spectrometer. The Power Module includes the following components:

- Main power circuit breaker switch
- Surge suppressor
- Line filter
- Electronics service switch
- Vacuum service switch

The “brains” of the TSQ mass spectrometer is the System Control PCB. The System Control PCB and embedded computer include the following:

- PowerPC processor
- Serial Peripheral Interconnect (SPI) bus
- I/O coprocessor
- Super Harvard Architecture Computer (Sharc) bus
- Scan generator DSP
- Acquisition processor DSP
- Interbus bridge
- 100 base-T Ethernet port

The rf voltage generation electronic assemblies produce the S-lens (TSQ Vantage mass spectrometer only), Q00, Q0, Q1, Q2, and Q3 rf voltages that enable ion transmission and mass analysis. The Analyzer Control PCB and the System Control PCB control all rf voltages.

The Q1 and Q3 rf voltage amplifier circuits are identical, and the circuits for Q0 and Q2 are similar.

The rf voltage generation electronic assemblies include the following components:

- RF oscillator
- RF Voltage Amplifier PCB
- Low Pass Filter PCB
- RF voltage coil
- RF voltage detector
- Mass DAC
- Integrating amplifier

The ion detection system electronic assemblies provide high voltage to the electron multiplier and conversion dynode of the ion detection system. They also receive the electron multiplier output current signal, convert it to a voltage (by the electrometer circuit), and pass it to the embedded computer.

The ion detection system electronic assemblies include the following:

- Electron multiplier power supply
- Conversion dynode power supply
- Electrometer PCB
- Acquisition DSP

The Analyzer Control PCB contains circuitry for controlling and monitoring the operation of the ion source, ion optics, mass analyzer, and ion detection system. These circuits are in turn monitored by the PowerPC processor of the System Control PCB via the SPI bus.

The Analyzer Control PCB controls and monitors the rf voltages for Q0, Q1, Q2, and Q3 quadrupoles. It also has lens voltage drivers that convert ± 330 V dc power from the DC Rod Driver PCB to dc voltages that are applied to the lenses.

Data System

The data system controls and monitors the TSQ mass spectrometer. The data system also processes data that the TSQ mass spectrometer acquires. The data system is composed of the following:

- [Computer Hardware](#)
- [Data System / Mass Spectrometer / LC Communication, Contact Closure, and Line Power Connections](#)
- [Data System / Local Area Network Interface](#)

Computer Hardware

The data system computer has the following major features:

- Intel™ Pentium™ IV processor
- High capacity hard disk drive
- Recordable/rewriteable CD drive
- Primary Ethernet port (data system to mass spectrometer)
- Secondary Ethernet port (data system to local area network)
- High performance video graphics card
- DVD drive
- 1280 × 1024 resolution color monitor
- Keyboard and mouse
- Laser printer

For more information about the computer, refer to the appropriate manuals.

Data System / Mass Spectrometer / LC Communication, Contact Closure, and Line Power Connections

Your Thermo Scientific mass spectrometry application supports LC instruments supplied by Thermo Fisher Scientific and several other manufacturers. The instrument control software for these instruments is provided on the LC Devices software CD.

[Figure 33](#) and [Figure 34](#) illustrate the communication, contact closure, and line power connections for two Thermo Scientific Accela™ LC/TSQ Series MS systems.

[Figure 33](#) shows the communication, contact closure, and line power connections for an LC/MS system with a TSQ Series mass spectrometer, Accela Open Autosampler, and Accela 600 or 1250 Pump. The TSQ Series mass spectrometer communicates with the data system computer through an Ethernet link by way of an Ethernet switch. The Accela Pump communicates with the data system computer through a USB link and the Accela Open Autosampler communicates with the data system computer through an RS232 link. Two contact closure cables, the Accela Open Autosampler interconnect cable and the Detector adapter cable, synchronize the system run signals.

[Figure 34](#) shows the communication, contact closure, and line power connections for an LC/MS system with a TSQ Series mass spectrometer, Accela Autosampler, Accela Pump, and Accela Detector. The TSQ Series mass spectrometer, Accela Autosampler, and Accela Detector (PDA or UV/Vis) communicate with the data system computer through an Ethernet link by

way of an Ethernet switch. The Accela pump (Accela Pump, Accela 600 Pump, or Accela 1250 Pump) communicates with the data system computer through a USB link. Two contact closure cables, the Accela interconnect cable and the TSQ adapter cable, synchronize the system run signals.

For information about connecting the communication cables to the Accela LC system devices and making the contact closure connections for your Accela LC /TSQ Series MS system, refer to the *Accela Getting Connected Guide*.

Data System / Local Area Network Interface

The data system computer contains a secondary Ethernet adapter. This secondary Ethernet adapter is not involved in data system/mass spectrometer or LC communications. You can use this secondary Ethernet adapter to access your local area network.

Printer

A high-resolution laser printer communicates with the data system via the local area network. Refer to the manual supplied by the manufacturer for details about the printer.

You set up the printer from the Print Setup dialog box. To open the Print Setup dialog box, choose **File > Print Setup** in any window.

Figure 33. LC/MS system with a TSQ Series mass spectrometer, an Accela Open Autosampler, and an Accela 600/1250 Pump


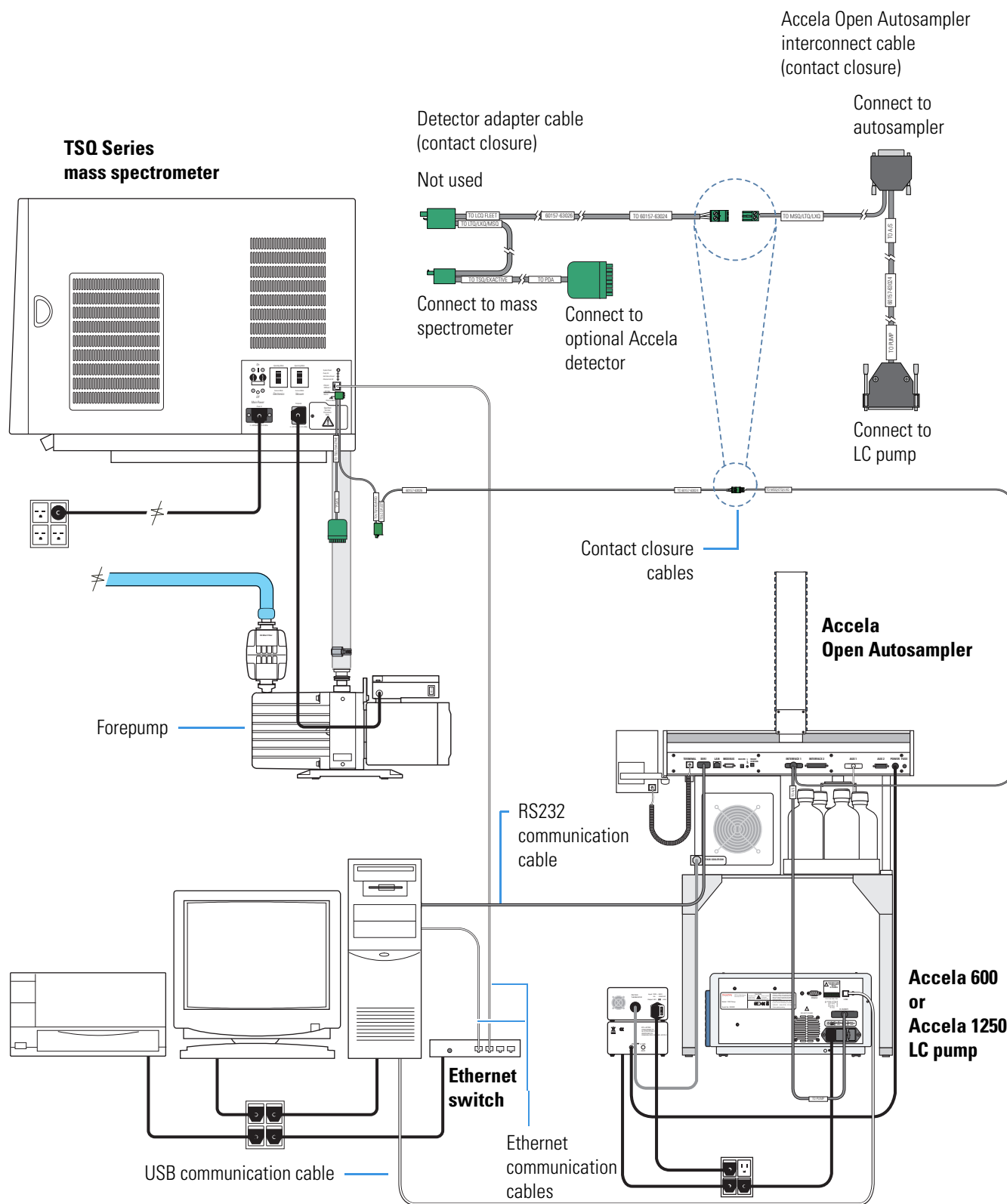


Figure 34. LC/MS system with a TSQ Series mass spectrometer, an Accela Autosampler, an Accela LC Pump, and an Accela PDA Detector

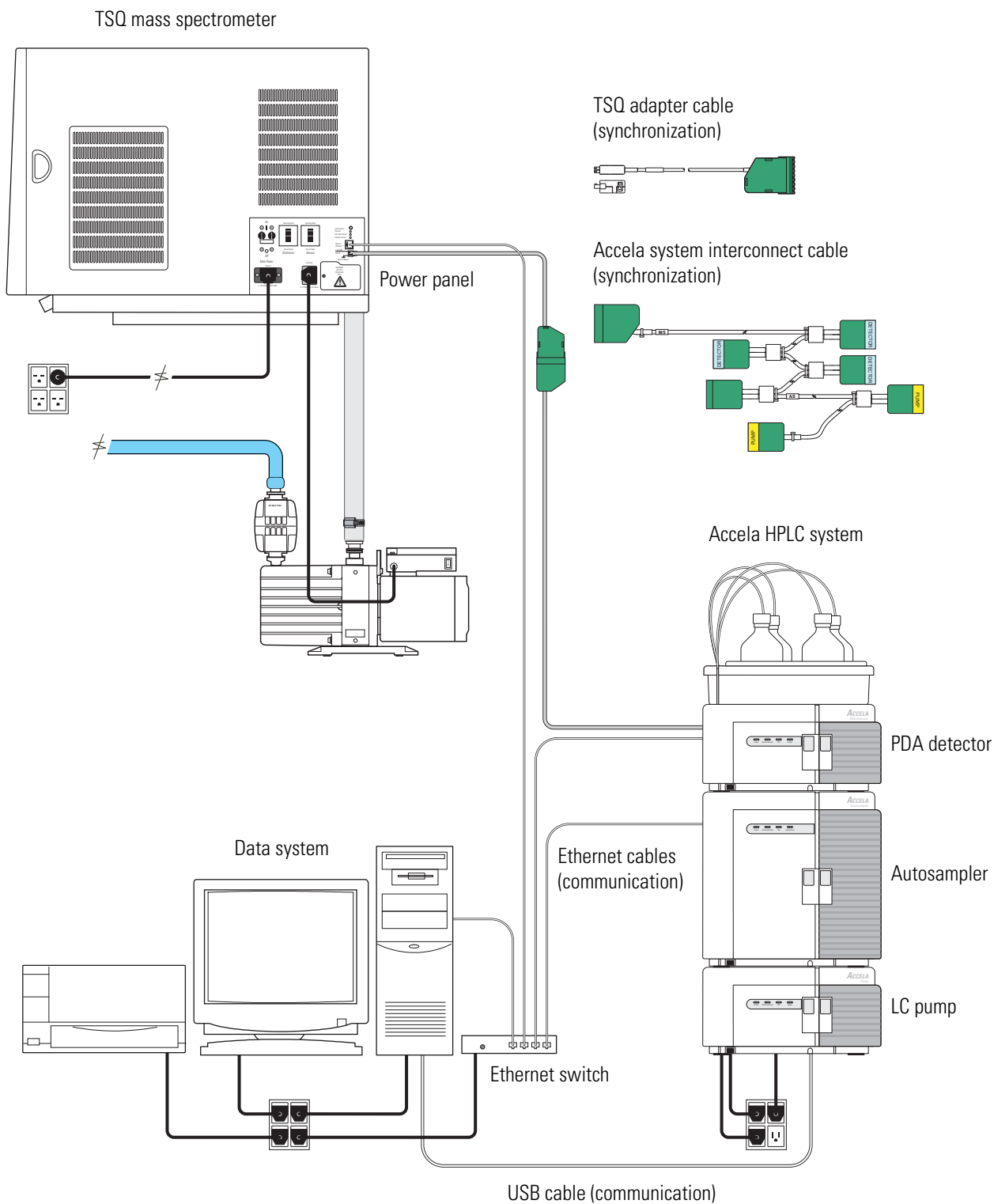

System Shutdown, Startup, and Reset

Many maintenance procedures for the TSQ system require that the mass spectrometer be shut down completely. In addition, you can place the TSQ mass spectrometer in standby mode if the system is not to be used for 12 hours or more.

Contents

- [Shutting Down the System in an Emergency](#)
- [Placing the System in Standby Condition](#)
- [Shutting Down the System Completely](#)
- [Starting Up the System after a Complete Shutdown](#)
- [Resetting the Mass Spectrometer](#)
- [Resetting the Data System](#)
- [Turning Off Selected Mass Spectrometer Components](#)

Shutting Down the System in an Emergency

If you need to turn off the mass spectrometer in an emergency, use the following process.



CAUTION Place the Main Power circuit breaker switch, located on the power panel on the right-side panel of the mass spectrometer, in the Off (O) position. See [Figure 35](#) and [Figure 36](#). This turns off all power to the mass spectrometer, including the vacuum pumps. Although removing power abruptly will not harm any component within the system, under normal conditions do not shut down the system with the main power circuit breaker switch. For the recommended procedure, see “[Shutting Down the System Completely](#)” on [page 52](#).

To turn off the LC, autosampler, and computer in an emergency, use the on/off switches on the LC, autosampler, and computer, respectively.

Figure 35. Power panel of the TSQ Vantage or TSQ Quantum Ultra mass spectrometer, showing the Main Power circuit breaker switch, Electronics service switch, and System Reset button


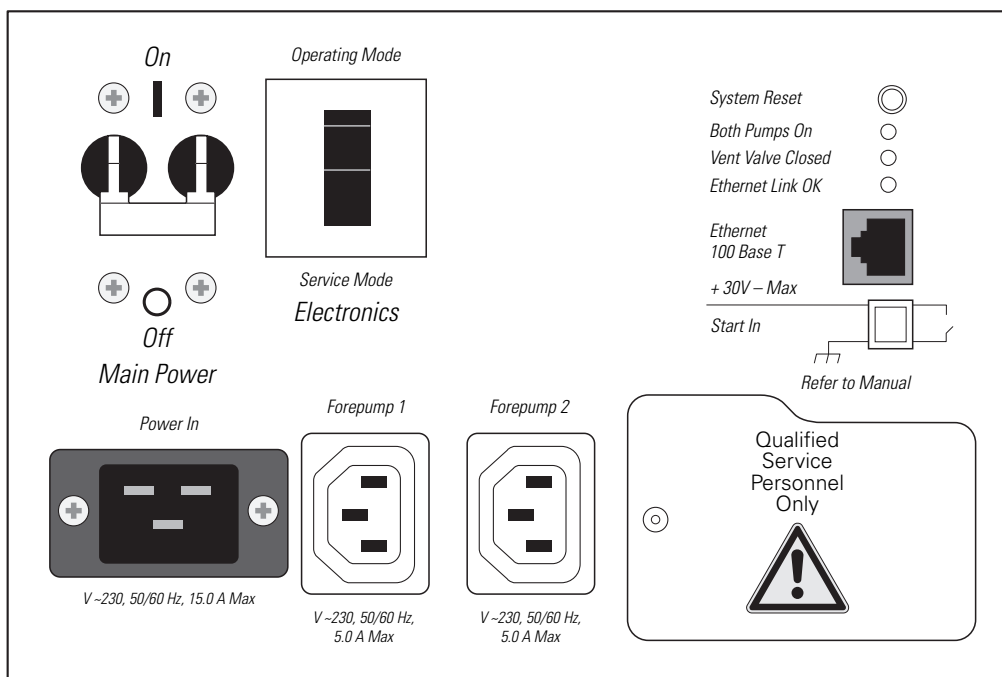



Figure 36. Power panel of the TSQ Quantum Access or TSQ Quantum Access MAX mass spectrometer, showing the Main Power circuit breaker switch, Electronics service switch, Vacuum service switch, and System Reset button

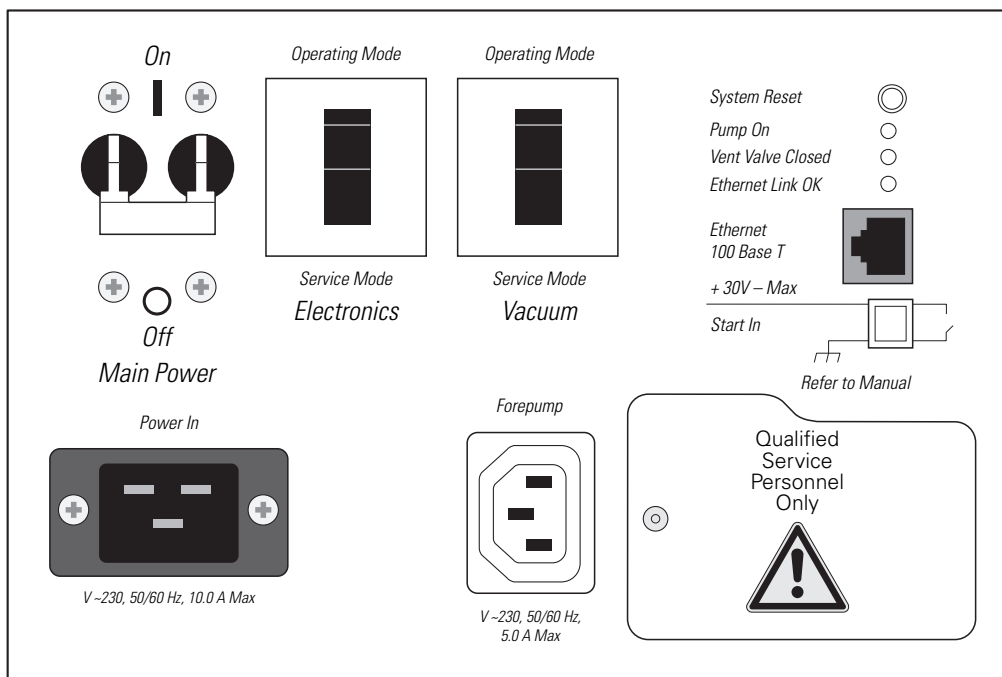

Placing the System in Standby Condition

If you are not going to use the TSQ system for a short period of time, such as overnight or over weekends, it does not need to be shut down completely. When you are not going to operate the system for 12 hours or more, you can leave the system in standby mode.

❖ To place the TSQ system in standby mode

1. Wait until data acquisition, if any, is complete.
2. Turn off the flow of liquid from the LC (or other sample introduction device) to the API source:
 - a. Choose **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune** to open the EZ Tune window.
 - b. In the EZ Tune window, choose **Setup > Inlet Direct Control**. The Inlet Direct Control view appears.
 - c. Click the LC's tab and click  (Stop) to stop the LC pump.
3. From the EZ Tune window, choose **Control > Standby** (or click the **On/Standby** button) to put the mass spectrometer in standby.



On



Off



Standby


When you choose Control > Standby, the TSQ system turns off the electron multiplier, conversion dynode, 8 kV power to the API source, and the mass analyzer and ion optics rf voltages. The TSQ system also turns off the auxiliary gas and sets the sheath gas flow to zero units. See [Table 2 on page 60](#) for the On/Off status of mass spectrometer components when the mass spectrometer is in standby mode. The System LED on the front panel of the mass spectrometer illuminates yellow when the system is in standby.

4. Flush the ion sweep cone and ion transfer tube of the API source as described in the topic [“Cleaning the Ion Sweep Cone and Ion Transfer Tube” on page 68](#).
5. Leave the mass spectrometer power on.
6. Leave the LC power on.
7. Leave the autosampler power on.
8. Leave the data system power on.

Shutting Down the System Completely

Shut down the TSQ system completely if not in use for an extended period only, or if it must be shut down for a maintenance or service procedure. For a short period of time, such as overnight or over weekends, you can place the system in standby. See [“Placing the System in Standby Condition.”](#)

❖ To shut down the TSQ system completely

1. Turn off the flow of liquid from the LC (or other sample introduction device) to the API source. To turn off the flow of liquid from the LC to the API source, do the following:
 - a. Choose **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune** to open the EZ Tune window.
 - b. In the EZ Tune window, choose **Setup > Inlet Direct Control**. The Inlet Direct Control view appears.
 - c. Click the LC's tab and click  (Stop) to stop the LC pump.
2. From the EZ Tune window, choose **Control > Off** (or click the **On/Standby/Off** button) to turn off the mass spectrometer.



On



Off



Standby

3. Place the Electronics service switch, located on the power panel (see [Figure 35](#) and [Figure 36](#)), in the Service Mode position.
4. On the TSQ Quantum Access or TSQ Quantum Access MAX, place the Vacuum service switch, located on the power panel (see [Figure 36](#) on [page 50](#)), in the Service Mode position.
5. Place the Main Power circuit breaker switch, located on the power panel (see [Figure 35](#) and [Figure 36](#)) in the Off (O) position. When you place the Main Power circuit breaker switch in the Off (O) position, the following occurs:
 - All power to the mass spectrometer is turned off. (All LEDs on the front panel of the mass spectrometer are off.)
 - A capacitor on the Vent Delay PCB provides power to the vent valve for two to four minutes (to allow the turbomolecular pump to spin down). After the capacitor discharges, power to the vent valve solenoid shuts off. When power to the vent valve solenoid shuts off, the vent valve opens and the vacuum manifold is vented to filtered air or nitrogen.
 - After about two minutes, the pressure of the vacuum manifold reaches atmospheric pressure.

6. Unplug the power cord for the mass spectrometer.



CAUTION Allow heated components to cool before servicing them.

Note If you are planning to perform routine or preventive system maintenance on the mass spectrometer only, you do not need to turn off the LC, argon and nitrogen gases, data system, or autosampler. In this case, the shutdown procedure is completed. However, if you do not plan to operate your system for an extended period of time, Thermo Fisher Scientific recommends that you turn off the LC, data system, autosampler, and gases as described in the following steps, 7 through 12.

7. Turn off the (optional) LC. Follow the procedure described in the manual that came with the LC.
8. Turn off the argon collision gas supply at the tank.
9. Turn off the nitrogen supply at the tank.
10. Turn off the data system:
 - a. Choose **Start > Shut Down** from the Windows™ taskbar. The Shut Down Windows dialog box appears.
 - b. To start the Windows shutdown procedure, select **Shut down** and click **OK**.
11. Turn off the laser printer by using the on/off switch.
12. Turn off the autosampler by using the Main Power on/off switch.

Starting Up the System after a Complete Shutdown

To start up the TSQ system after it has been shut down completely, do the following procedures:

- [Starting Up the LC](#) (optional)
- [Starting Up the Data System](#)
- [Starting Up the Mass Spectrometer](#)
- [Starting Up the Autosampler](#) (optional)
- [Setting Up Conditions for Operation](#)

Starting Up the LC

To start up the LC, follow the startup procedure described in the manual that came with the LC. If necessary, set up the LC as described in its *Getting Connected Guide*. If necessary, configure the LC in the Instrument Configuration window. To open the Instrument Configuration window, choose **Start > All Programs > Thermo Foundation > Instrument Configuration**.

Do not turn on the liquid flow to the mass spectrometer.

Starting Up the Data System

❖ To start up the data system

1. Turn on the monitor, computer, and printer.
2. Observe the Windows XP startup procedure on the monitor and press CTRL+ALT+DEL when you are prompted to do so.
3. In the Logon Information dialog box, click **OK** or enter your password (if you have one).

Starting Up the Mass Spectrometer

❖ To start up the mass spectrometer

Note The data system must be running before you start up the mass spectrometer. The mass spectrometer will not operate until software is received from the data system.

1. Turn on the flows of argon and nitrogen at the tanks if they are off.
2. Make sure that the Main Power circuit breaker switch is in the Off (O) position and that the Electronics service switch and the Vacuum service switch (TSQ Quantum Access or TSQ Quantum Access MAX) are both in the Service Mode position.
3. Plug in the power cord for the mass spectrometer.

4. Place the Main Power circuit breaker switch in the On (I) position. When you place the Main Power circuit breaker switch in the On (I) position, power is supplied to those mass spectrometer components that are not affected by the Vacuum service switch and the Electronics service switch.
5. On the TSQ Quantum Access or TSQ Quantum Access MAX, place the Vacuum service switch in the Operational position.
6. Place the Electronics service switch in the Operational position. When you place the Electronics service switch in the Operational position, the following occurs:
 - The Power LED on the mass spectrometer front panel illuminates green to indicate that power is provided to the mass spectrometer electronics. (The electron multiplier, conversion dynode, ion source, and mass analyzer and ion optics rf voltages remain off.)
 - The embedded computer reboots. After several seconds the Communication LED on the front panel illuminates yellow to indicate that the data system and the mass spectrometer have started to establish a communication link.
 - After several more seconds, the Communication LED illuminates green to indicate that the data system and the mass spectrometer have established a communication link. Software for the operation of the mass spectrometer is then transferred from the data system to the mass spectrometer.
 - After three minutes, the System LED illuminates yellow to indicate that the software transfer from the data system to the mass spectrometer is complete and that the instrument is in standby.
 - The Vacuum LED on the front panel of the mass spectrometer remains off until the turbomolecular pump reaches 80 percent of its operational speed of 750 Hz. At this time the ion gauge is turned on and the Vacuum LED illuminates yellow. The Vacuum LED illuminates green, and the high voltage can be turned on, only if the pressure in the mass analyzer region, as measured by the ion gauge, is below the maximum allowable pressure of 7×10^{-4} Torr.

If you have an autosampler, go to the next topic, [“Starting Up the Autosampler.”](#) If you do not have an autosampler, go to the topic [“Setting Up Conditions for Operation.”](#)

Starting Up the Autosampler

❖ To start up the autosampler

Place the power switch on the autosampler in the On position. If necessary, configure the autosampler. For procedures for placing sample vials, preparing solvent and waste bottles, installing syringes, and so on, refer to the manual that came with the autosampler. Refer also to its *Getting Connected Guide*.

Setting Up Conditions for Operation

❖ To set up your TSQ mass spectrometer for operation

1. Before you begin data acquisition with your TSQ system, allow the system to pump down for at least one hour. Operation of the system with excessive air and water in the vacuum manifold can cause reduced sensitivity, tuning problems, and reduced lifetime of the electron multiplier.
2. Ensure that the argon pressure and nitrogen pressure are within the operational limits [argon: 135 ± 70 kPa (20 ± 10 psig), nitrogen: 690 ± 140 kPa (100 ± 20 psig)].

Note Air in the argon line must be purged or given sufficient time to be purged for normal TSQ mass spectrometer performance.

3. In the EZ Tune window, choose **View > Instrument Status** and look at the Instrument Status view. Check that the ion gauge pressure is below about 5×10^{-6} Torr, and that the forepump pressure is below about 50 mTorr (with the ion transfer tube sealed with a septum cap).
4. Remove the septum cap from the orifice of the ion transfer tube.
5. Set up the API source for operation as described in the *Ion Max and Ion Max-S API Source Hardware Manual*, *HESI-II Probe User Guide*, *H-ESI Probe User Guide*, *Nanospray Ion Source Operator's Manual*, and *TSQ Series Getting Started Guide*.

Note You do not need to calibrate and tune the TSQ mass spectrometer each time you restart the TSQ system.

Calibration parameters are instrument parameters that affect the mass accuracy and resolution. Tune parameters are instrument parameters that affect the intensity of the ion signal. You must tune and calibrate the TSQ mass spectrometer (that is, optimize the tune and calibration parameters) perhaps once a quarter. Refer to “Automatic Tuning and Calibrating in the ESI/MS Mode” in the *TSQ Series Getting Started Guide* for a procedure for tuning and calibrating the TSQ mass spectrometer.

You must optimize the tune parameters (or change the Tune Method) whenever you change the type of experiment. For procedures on how to optimize the tune parameters for an ESI or APCI experiment, refer to “Optimizing the Mass Spectrometer with Your Compound in ESI/MS/MS Mode” or “Optimizing the Mass Spectrometer with Your Compound in APCI/MS/MS Mode” in the *TSQ Series Getting Started Guide*.

Resetting the Mass Spectrometer

If communication between the mass spectrometer and data system computer is lost, it might be necessary to reset the mass spectrometer using the System Reset button on the right-side power panel. Pressing the System Reset button creates an interrupt in the embedded computer. This causes the embedded computer to restart in a known (default) state. See [Figure 35](#) on [page 50](#) for the location of the System Reset button.

The procedure given here assumes that the mass spectrometer and data system computer are both powered on and operational. If the mass spectrometer, data system computer, or both are off, go to [“Starting Up the System after a Complete Shutdown”](#) on [page 54](#).

❖ To reset the mass spectrometer

Press the System Reset button located on the right-side power panel. Make sure the Communication LED is extinguished before releasing the System Reset button. When you press the Reset button, the following occurs:

- An interrupt on the embedded computer causes the CPU to reboot. All LEDs on the front panel of the mass spectrometer are off except the Power LED.
- After several seconds, the Communication LED illuminates yellow to indicate that the data system and the mass spectrometer are starting to establish a communication link.
- After several more seconds, the Communication LED illuminates green to indicate that the data system and the mass spectrometer have established a communication link. Software for the operation of the mass spectrometer is then transferred from the data system to the mass spectrometer.
- After three minutes the software transfer is complete. The System LED illuminates either green to indicate that the instrument is functional and the high voltages are on, or yellow to indicate that the instrument is functional and is in standby.

Resetting the Data System

There are two ways to reset the data system:

- [Resetting the Data System by Using the Windows Shutdown and Restart Procedure](#)
- [Resetting the Data System by Turning the Personal Computer Off Then On](#)

Resetting the Data System by Using the Windows Shutdown and Restart Procedure

If possible, use the Windows shutdown and restart procedure to shut down and restart the data system so that Windows can properly close programs and save changes to files.

❖ To reset the data system by using the Windows shutdown and restart procedure

1. Choose **Start > Shut Down** from the Windows taskbar. The Shut Down Windows dialog box appears.
2. To start the Windows shutdown and restart procedure, select **Restart** and click **OK**.
3. Observe the Windows shutdown and restart procedure on the monitor. Press CTRL+ALT+DELETE when you are prompted to do so.
4. To complete the shutdown and restart procedure, click **OK** or enter your password (if you have one) in the Logon Information dialog box.

Note The communications link between the data system and the mass spectrometer should automatically reestablish after you reset the data system. When this occurs the Communication LED on the front panel of the mass spectrometer illuminates yellow and then green. If the system is unable to reestablish the communications link, press the System Reset button on the power panel of the mass spectrometer.

Resetting the Data System by Turning the Personal Computer Off Then On

If you are unable to reset the data system by using the Windows shutdown and restart procedure, proceed as follows.

❖ To reset the data system by turning the PC off and on

1. Press the Power button on the personal computer to turn the personal computer off.
2. After several seconds, press the Power button on the personal computer to turn the personal computer on.
3. Observe the Windows XP startup procedure on the monitor and press CTRL+ALT+DELETE when you are prompted to do so.
4. To complete the shutdown and restart procedure, click **OK** or enter your password in the Logon Information dialog box.
5. When the shutdown and restart procedure has completed, choose **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune** to display the EZ Tune window.

Note The communications link between the data system and the mass spectrometer should automatically reestablish after you reset the data system. When this occurs the Communication LED on the front panel of the mass spectrometer illuminates yellow and then green. If the system is unable to reestablish the communications link, press the System Reset button on the right-side power panel of the Power Entry Module of the mass spectrometer. See [Figure 35](#) on [page 50](#).

Turning Off Selected Mass Spectrometer Components

There are different ways that you can turn off some or all of the mass spectrometer components:

- Turn off individual mass spectrometer components from the EZ Tune window. Turning off individual mass spectrometer components might be necessary when you are troubleshooting or when you are running certain diagnostic procedures.
- Place the mass spectrometer in standby mode. Standby is the normal condition to leave the mass spectrometer in when it is not in use. Choose **Control > Standby** (or turn the On/Standby button on or off) from the EZ Tune window to place the mass spectrometer in standby.
- Place the mass spectrometer in the off condition. The off condition is similar to standby, except all high voltage components of the mass spectrometer are turned off. Choose **Control > Off** from the EZ Tune window to place the mass spectrometer in the off condition.
- Place the Electronics service switch in the Service Mode position. The Electronics service switch turns off all components in the mass spectrometer other than the +24 V power supply, forepump, turbomolecular pump, Vent Delay PCB, and fans.
- Place the Vacuum service switch on the TSQ Quantum Access or TSQ Quantum Access MAX in the Service Mode position. The Vacuum service switch turns off all vacuum system components, including the +24 V power supply, forepump, turbomolecular pump, Vent Delay PCB, and fans.
- Place the Main Power circuit breaker switch in the Off (O) position. Placing the Main Power circuit breaker switch in the Off (O) position removes all power to the mass spectrometer, including the vacuum system.

3 System Shutdown, Startup, and Reset

Turning Off Selected Mass Spectrometer Components

Table 2 summarizes the on/off status of mass spectrometer components, voltages, and gas flows.

Table 2. On/Off status of mass spectrometer components, voltages, and gas flows (Sheet 1 of 2)





Mass spectrometer component	Standby 	Off 	Electronics service switch in Service Mode position	Vacuum service switch (TSQ Quantum Access or Access MAX) in Service Mode position	MS Main Power circuit breaker switch in Off position
Electron multiplier	Off	Off	Off	Off	Off
Conversion dynode	Off	Off	Off	Off	Off
Mass analyzer rf voltage	Off	Off	Off	Off	Off
Mass analyzer dc offset voltage	Off	Off	Off	Off	Off
Q00 and Q0 ion optics rf voltage	Off	Off	Off	Off	Off
Q00 and Q0 ion optics dc offset voltage	Off	Off	Off	Off	Off
Tube lens or S-lens	On	Off	Off	Off	Off
Ion transfer tube heater	On	Off	Off	Off	Off
Ion transfer tube dc offset	On	Off	Off	Off	Off
APCI corona discharge needle	Off	Off	Off	Off	Off
APCI vaporizer	Off	Off	Off	Off	Off
ESI needle	Off	Off	Off	Off	Off
Sheath gas	Off	Off	Off	Off	Off
Auxiliary gas	Off	Off	Off	Off	Off
Argon collision gas	Off	Off	Off	Off	Off
Vent valve	Closed	Closed	Closed	Open (after 2 to 4 min)	Open (after 2 to 4 min)
Turbomolecular pump	On	On	On	Off	Off
Forepump	On	On	On	Off	Off
Vent Delay PCB	On	On	On	Off	Off
Embedded computer	On	On	Off	On	Off
Turbomolecular pump controller	On	On	On	Off	Off
Power supply, electron multiplier, and conversion dynode	Off	Off	Off	Off	Off
Power supply, 8 kV	Off	Off	Off	Off	Off
PS1 power supply, +24 V	On	On	On	Off	Off

Table 2. On/Off status of mass spectrometer components, voltages, and gas flows (Sheet 2 of 2)

Mass spectrometer component	Standby 	Off 	Electronics service switch in Service Mode position	Vacuum service switch (TSQ Quantum Access or Access MAX) in Service Mode position	MS Main Power circuit breaker switch in Off position
PS2 power supply, +5, ± 15 , ± 24 V dc	On	On	Off	Off	Off
PS3 power supply, +36, -28 V dc	On	On	Off	Off	Off
Fan, turbomolecular pump	On	On	On	Off	Off
Fan, above manifold	On	On	On	Off	Off
Fan, center wall	On	On	On	Off	Off
Convectron gauge, foreline	On	On	On	On	Off
Convectron gauge, collision cell	On	On	Off	On	Off
Ion gauge	On	On	Off	Off	Off

Daily Operation

To ensure the proper operation of your system, Thermo Fisher Scientific recommends daily performance of these checks and cleaning procedures for the TSQ system.

Contents

- [Things to Do Before Operating the TSQ System](#)
- [Things to Do After Operating the TSQ System](#)

Note You do not need to tune and calibrate the mass spectrometer as part of your daily routine.

Tune parameters are instrument parameters that affect the intensity of the ion signal. Calibration parameters are instrument parameters that affect the mass accuracy and resolution of the mass spectrum. You must tune and calibrate the mass spectrometer (that is, optimize the tune and calibration parameters) perhaps once a quarter. For procedures on how to tune and calibrate your mass spectrometer, refer to “Automatic Tuning and Calibrating in the ESI/MS Mode” in the *TSQ Series Getting Started Guide*.

You must optimize the tune parameters (or change the Tune Method) whenever you change the type of experiment. For procedures on how to optimize the tune parameters for your ESI or APCI experiment, refer to “Optimizing the Mass Spectrometer with Your Compound in ESI/MS/MS Mode” or “Optimizing the Mass Spectrometer with Your Compound in APCI/MS/MS Mode” in the *TSQ Series Getting Started Guide*.

Things to Do Before Operating the TSQ System

Every day before starting analyses, verify that the instrument is ready for operation by doing the following:

- Check the argon and nitrogen supplies.
- Check the ESI fused-silica sample tube for elongation.
- Check the system vacuum levels.

Checking the Argon and Nitrogen Supplies

Check the argon supply on the regulator of the gas tank. Make sure that you have sufficient gas for your analysis. If necessary, install a new tank of argon. Verify that the pressure of argon reaching the mass spectrometer is between 135 ± 70 kPa (20 ± 10 psig). If necessary, adjust the pressure with the tank pressure regulator.

Check the nitrogen supply on the regulator of the nitrogen gas tank or liquid nitrogen boil-off tank. Make sure that you have sufficient gas for your analysis. Based on 24-hour per day operation, typical nitrogen consumption is 2800 liters per day (100 cubic feet per day). If necessary, replace the tank. Verify that the pressure of nitrogen reaching the mass spectrometer is between 690 ± 140 kPa (100 ± 20 psig). If necessary, adjust the pressure with the tank pressure regulator.



CAUTION Before you begin normal operation each day, ensure that you have sufficient nitrogen for your API source. The presence of oxygen in the ion source when the mass spectrometer is on could be unsafe. The TSQ system displays a pop-up message if the nitrogen pressure is too low.

Checking the ESI Fused-Silica Sample Tube for Elongation

Using acetonitrile in the mobile phase can cause elongation of the polyimide coating on the fused-silica sample tube. Elongation of the polyimide coating can degrade signal intensity and signal stability over time.

If you are running in ESI mode with a fused-silica sample tube, verify the sample tube is not elongated past the tip of the ESI spray needle. (This elongation does not occur with a metal needle.)

To cut and reposition the sample tube approximately 0.5 mm (between 0 and 1 mm) inside the end of the ESI needle, refer to the *Ion Max and Ion Max-S API Source Hardware Manual* for more information.

Checking the System Vacuum Levels

For proper performance, your TSQ system must operate at the proper vacuum levels. Operation of the system with poor vacuum levels can cause reduced sensitivity, tuning problems, and reduced lifetime of the electron multiplier. You should check your system for air leaks by checking the system vacuum levels before you begin your first acquisition.

You can check the current values of the pressures in the ion transfer tube-skimmer region and foreline (labeled *Forepump Pressure*) and in the analyzer region (labeled *Ion Gauge Pressure*) in the Instrument Status view of the EZ Tune window. To display the EZ Tune window, from the Windows taskbar choose **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune**. Choose **View > Instrument Status** to display the Instrument Status view.

Compare the current values of the pressures in the vacuum manifold with the values listed in [Table 3](#). If the current values are higher than normal, you might have an air leak.

Table 3. Typical pressure readings for TSQ Vantage and TSQ Quantum Ultra

Conditions	Convectron gauge reading (foreline, ion transfer tube- skimmer region)	Ion gauge reading (analyzer region)
Collision gas off, ion transfer tube orifice sealed	less than 0.05 Torr	Less than 3×10^{-6} Torr (after pumping for at least one hour)
Collision gas off, ion transfer tube orifice open	0.9 to 1.2 Torr (110 V at 60 Hz) 1.5 to 2.2 Torr (220 V at 50 Hz)	Less than 5×10^{-6} Torr
Collision gas set to 2 mTorr, ion transfer tube orifice open	0.9 to 1.2 Torr (110 V at 60 Hz) 1.5 to 2.2 Torr (220 V at 50 Hz)	Less than 3×10^{-5} Torr

Table 4. Typical pressure readings for TSQ Quantum Access and TSQ Quantum Access MAX

Conditions	Convectron gauge reading (foreline, ion transfer tube- skimmer region)	Ion gauge reading (analyzer region)
Collision gas off, ion transfer tube orifice sealed	less than 0.05 Torr	Less than 3×10^{-6} Torr (after pumping for at least one hour)
Collision gas off, ion transfer tube orifice open	1.2 to 1.8 Torr (110 V at 60 Hz) 1.5 to 2.2 Torr (220 V at 50 Hz)	Less than 5×10^{-6} Torr
Collision gas set to 2 mTorr, ion transfer tube orifice open	1.2 to 1.8 Torr (110 V at 60 Hz) 1.5 to 2.2 Torr (220 V at 50 Hz)	Less than 3×10^{-5} Torr

If the pressure is high (above 5×10^{-5} Torr in the analyzer region), and you have restarted the system within the last 30 to 60 minutes, wait an additional 30 minutes and recheck the pressure. If the pressure is decreasing with time, check the pressure periodically until it is within the typical pressure range of the mass spectrometer.

If the pressure remains high, your system might have an air leak. If you suspect an air leak, shut down the system as described in the topic [“Shutting Down the System Completely”](#) on [page 52](#). Make a visual inspection of the vacuum system and vacuum lines for leaks. Check each fitting and flange on the system for tightness, and tighten the fittings or flanges that are loose. Do not tighten fittings indiscriminately. Pay particular attention to fittings that have been changed recently or to fittings that have been subjected to heating and cooling. Make sure that the cover plates of the vacuum manifold are properly seated.

Things to Do After Operating the TSQ System

After operating the TSQ system, verify the instrument is ready for a period of inactivity by doing the following:



- [Flushing the Sample Transfer Line, Sample Tube, and API Probe](#)
- [Placing the System in Standby Condition](#)
- [Cleaning the Ion Sweep Cone and Ion Transfer Tube](#)
- [Emptying the Solvent Waste Bottle](#)

Flushing the Sample Transfer Line, Sample Tube, and API Probe

Thermo Fisher Scientific recommends flushing the sample transfer line, sample tube, and API probe at the end of each working day (or more often if you suspect they are contaminated) by flowing a 50:50 methanol/distilled water solution from the LC through the API source.

❖ To flush the sample transfer line, sample tube, and API probe

1. Wait until data acquisition, if any, is complete.
2. To display the EZ Tune window, from the Windows taskbar choose **Start > All Programs > Thermo Instruments > TSQ > TSQ Tune**.
3. From the EZ Tune window, choose **Control > On** (or click the **On/Standby** button) to turn on the voltages and gas flows to the API source.
 - If you are operating in APCI mode, go to step 4.
 - If you are operating in ESI mode, go to step 5.
4. Set up the APCI source:
 - a. In the EZ Tune window, click on the Ion Source button to display the Ion Source Devices dialog box.
 - b. Set the APCI vaporizer temperature to 500 °C:
Select **Vaporizer Temperature** in the device list, and then type **500** in the box and click **Apply**.
 - c. Set the sheath gas flow rate to 30:
Select **Sheath Gas Pressure** in the device list, and then type **30** in the box and click **Apply**.
 - d. Set the auxiliary gas flow rate to 5:
Select **Aux Valve Flow** in the device list, and then type **5** in the box and click **Apply**.
 - e. Set the spray current to 0:
Select **Spray Current** in the device list, and then type **0** in the box and click **Apply**.Go to step 6.

5. Set up the ESI source as follows:
 - a. In the EZ Tune window, click on the Ion Source button to display the Ion Source Devices dialog box.
 - b. Set the sheath gas flow rate to 30:
Select **Sheath Gas Pressure** in the device list, and then type **30** in the box and click **Apply**.
 - c. Set the auxiliary gas flow rate to 5:
Select **Aux Valve Flow** in the device list, and then type **5** in the box and click **Apply**.
 - d. Set the ESI spray voltage to 0:
Select **Spray Voltage** in the device list, and then type **0** in the box and click **Apply**.
6. Set up and start a flow of 50:50 methanol/water solution from the LC to the API source as follows:
 - a. In the EZ Tune window, choose **Setup > Inlet Direct Control**. The Inlet Direct Control view appears.
 - b. Click the LC's tab.
 - c. Set the flow rate to a value that is typical for your experiments.
 - d. Set the solvent proportions to 50% methanol and water.
 - e. Click  (Start) to start the LC pump.
7. Let the solution flow through the sample transfer line, sample tube, and API probe for 15 minutes. After 15 minutes, turn off the flow of liquid from the LC to the API source, as follows: Leave the API source (including the APCI vaporizer, sheath gas, and auxiliary gas) on for an additional 5 minutes. Click  (Stop) to stop the LC pump.
8. After 5 minutes, turn off the API source by placing the mass spectrometer in standby: From the EZ Tune window, choose **Control > Standby** (or click the **On/Standby** button) to put the mass spectrometer in standby.

Go to the next topic, [“Placing the System in Standby Condition.”](#)

Placing the System in Standby Condition

❖ To place the TSQ system in standby condition

1. From the EZ Tune window, choose **Control > Standby** (or click the **On/Standby** button) to put the mass spectrometer in standby. The System LED on the front panel of the mass spectrometer is illuminated yellow when the system is in standby.
2. Leave the mass spectrometer power on.
3. Leave the LC power on.
4. Leave the autosampler power on.
5. Leave the data system power on.

Go to the next topic, [“Cleaning the Ion Sweep Cone and Ion Transfer Tube.”](#)

Cleaning the Ion Sweep Cone and Ion Transfer Tube

You must clean the ion sweep cone and the ion transfer tube on a regular basis to prevent corrosion and to maintain optimum performance of your API source. If you use nonvolatile buffers in your solvent system or high concentrations of sample, you might need to clean the ion sweep cone and ion transfer tube more often. See the topic, [“Maintaining the Ion Sweep Cone and Ion Transfer Tube”](#) on page 79.

Go to the next topic, [“Emptying the Solvent Waste Bottle.”](#)

Emptying the Solvent Waste Bottle

Thermo Fisher Scientific recommends checking the solvent level in the solvent waste bottle on a daily basis. If necessary, empty the solvent waste bottle. Dispose of the solvent waste in accordance with local and federal regulations.

Removing and Reinstalling the Ion Source Housing

You must remove the Ion Max™ (on the TSQ Quantum Access MAX, TSQ Vantage, or TSQ Quantum Ultra) or Ion Max-S (on the TSQ Quantum Access) ion source housing before performing maintenance on the ion optics or ion source interface, or to access the ion sweep cone.

Contents

- [Removing the Ion Max or Ion Max-S Ion Source Housing](#)
- [Installing the Ion Max or Ion Max-S Ion Source Housing](#)
- [Ion Max and Ion Max-S Housing Maintenance](#)

Removing the Ion Max or Ion Max-S Ion Source Housing

Note If an ion source probe is still installed in the ion source housing, disconnect the external liquid lines to the probe before removing the ion source housing.

❖ To remove the ion source housing

1. Remove the drain tube from the ion source housing drain. See [Figure 37](#) and [Figure 38](#).
2. Rotate the ion source housing locking levers 90 degrees to release the ion source housing from the ion source mount assembly.
3. Remove the ion source housing by pulling it straight off the ion source mount assembly, and place the housing in a safe location for temporary storage.

5 Removing and Reinstalling the Ion Source Housing

Removing the Ion Max or Ion Max-S Ion Source Housing

Figure 37. Ion Max source housing (front)


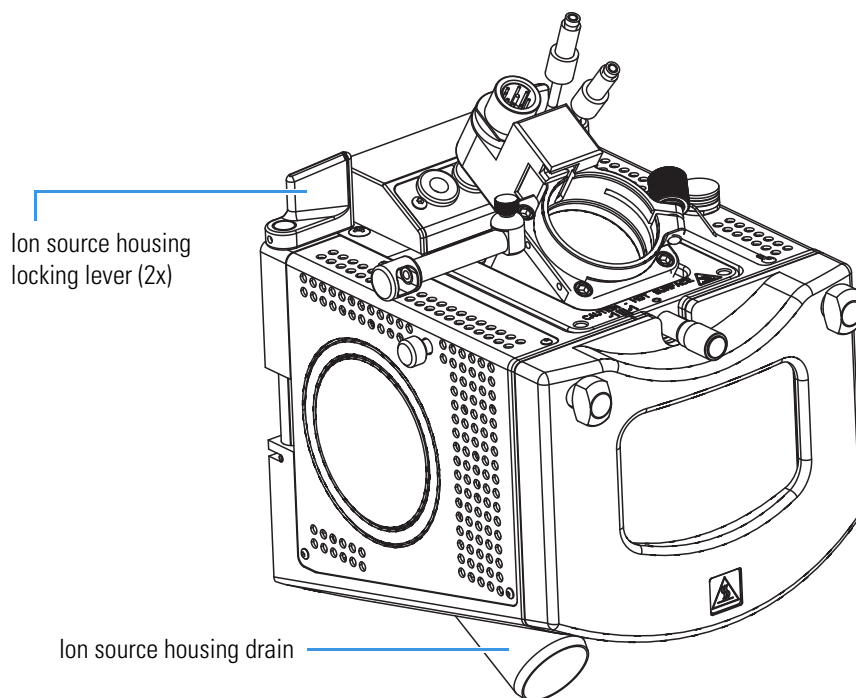


Figure 38. Ion Max-S source housing (front)

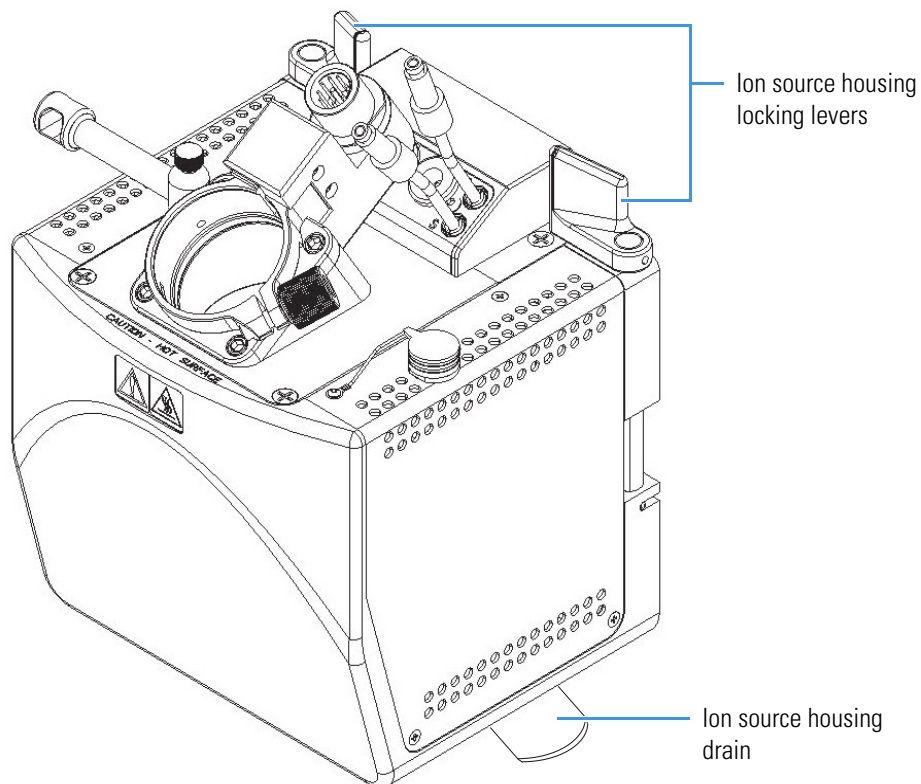


Installing the Ion Max or Ion Max-S Ion Source Housing

❖ **To reinstall the Ion Max or Ion Max-S ion source housing**

1. Align the two guide pin holes on the of the ion source housing with the ion source housing guide pins on the mass spectrometer. Carefully press the ion source housing onto the ion source mount. See [Figure 39](#) and [Figure 40](#).


Figure 39. Ion source housing

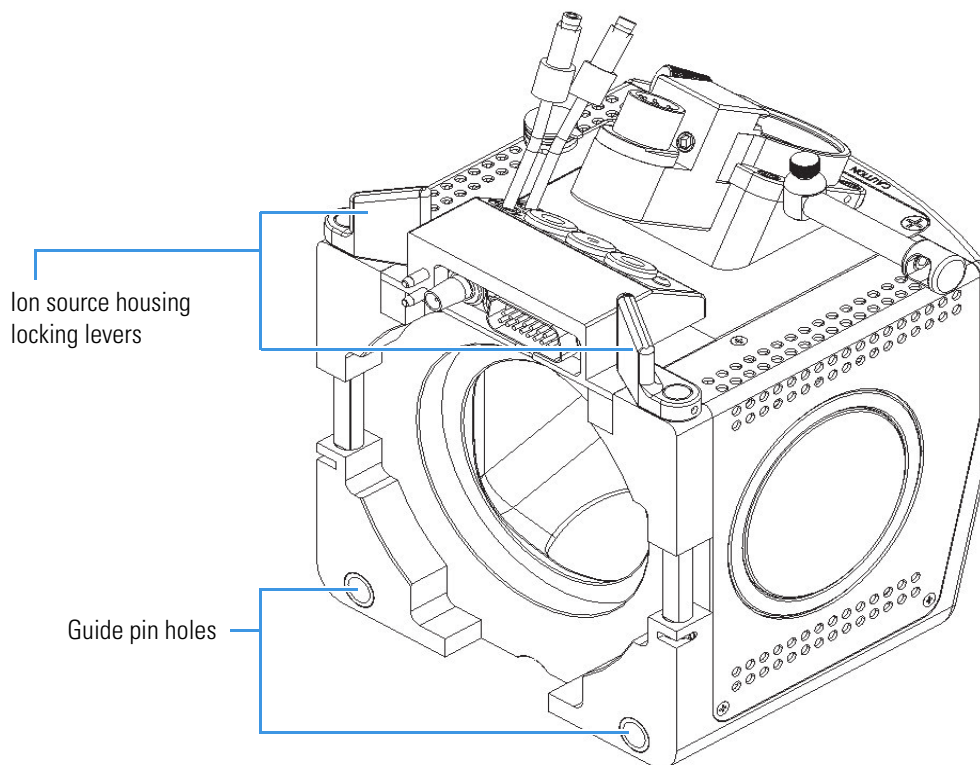

5 Removing and Reinstalling the Ion Source Housing

Installing the Ion Max or Ion Max-S Ion Source Housing

Figure 40. Ion source mount, showing the ion source housing guide pins and ion source interface release lever

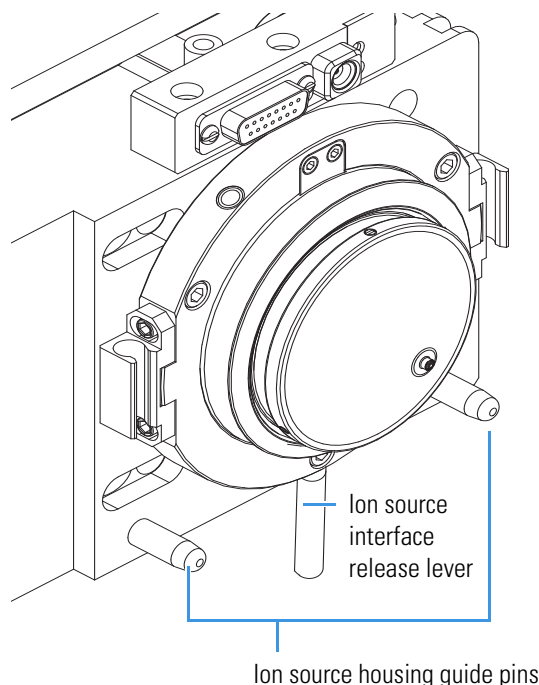

2. Rotate the ion source housing locking levers 90 degrees to lock the ion source housing onto the ion source mount assembly.



CAUTION Prevent solvent waste from backing up into the ion source and mass spectrometer. Always ensure that liquid in the drain tube is able to drain to a waste container.

3. Reinstall the ion source housing drain tube:



CAUTION Do not vent the ion source drain tube (or any vent tubing connected to the waste container) to the same fume exhaust system to which you have connected the forepump. The ion source chamber and analyzer optics can become contaminated with mechanical pump oil if the ion source drain tube and the (blue) forepump exhaust tubing are connected to the same fume exhaust system.

Your laboratory must be equipped with at least two fume exhaust systems. Route the (blue) forepump exhaust tubing to a dedicated fume exhaust system. Route the drain tube from the ion source to a waste container. Vent the waste container to a dedicated fume exhaust system.

- a. Connect the 1 in. ID Tygon™ tubing (P/N 00301-22922) to the ion source housing drain fitting.
- b. Attach the free end of the hose to a dedicated drain system. Vent the drain system to a fume exhaust system.

The Ion Max or Ion Max-S ion source is now properly installed on the mass spectrometer.

Ion Max and Ion Max-S Housing Maintenance

The Ion Max and Ion Max-S ion source housings are designed to be serviced by Thermo Fisher Scientific customer support engineers. User maintenance is limited to cleaning the ion source housing as necessary. To clean the housing, remove the housing from the instrument, and in an appropriate fume hood rinse the interior of the housing with methanol. Allow the housing to dry before you install it on the mass spectrometer. Be sure to follow all safety precautions in the sections regarding the installation and removal of the housing. For any additional service that might be required, contact your local Thermo Fisher Scientific service representative.

Maintenance

This chapter describes routine maintenance procedures that you must perform to ensure optimum performance of the instrument. Optimum performance of the TSQ mass spectrometer depends on the maintenance of all parts of the instrument. **You are responsible for maintaining your system properly by performing the system maintenance procedures on a regular basis.**

Note When you perform maintenance procedures, be methodical; always wear clean, lint-free gloves when handling the components of the ion source; always place the components on a clean, lint-free surface; and never overtighten a screw or use excessive force.

Contents

- [Frequency of Cleaning](#)
- [Maintaining the Ion Sweep Cone and Ion Transfer Tube](#)
- [Maintaining the Ion Source Interface Assembly](#)
- [Cleaning the Q00 Ion Optics](#)
- [Cleaning the Fan Filters](#)

[Table 5](#) lists routine and infrequent mass spectrometer maintenance procedures. Most of the procedures involve cleaning. For example, procedures are provided for cleaning the ion transfer tube, the ion sweep cone, the skimmer and tube lens of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra, the exit lens and S-lens of the TSQ Vantage, the Q00 rf lens, the lens L0, and the fan filters. Procedures for replacing the ion transfer tube and the capillary heater assembly are also provided. For instructions on maintaining the LC or autosampler, refer to the manual that comes with the LC or autosampler.

Table 5. Mass spectrometer maintenance procedures (Sheet 1 of 2)

Mass spectrometer component	Procedure	Frequency	Procedure location
API source	Flush (clean) sample transfer line, sample tube, and API probe	Daily	page 66
	Remove and clean ion transfer tube	Weekly, or if ion transfer tube bore is contaminated or obstructed	page 79
	Remove and clean ion sweep cone	As needed ^a	page 79
	Clean Ion Max or Ion Max-S housing	As needed ^a	page 73
	Replace ion transfer tube	If ion transfer tube bore is corroded	page 79
Q00 ion optics	Clean S-lens and exit lens (TSQ Vantage)	As needed ^a	page 87
	Clean tube lens and skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Quantum Ultra)	As needed ^a	page 88
	Clean Q00 rf lens and lens L0	Every 3 to 4 months ^a	page 97
APCI, ESI, or H-ESI probe	Replace fused-silica sample tube	If sample tube is broken or obstructed	<i>Ion Max and Ion Max-S API Source Hardware Manual or H-ESI Probe User Guide</i>
ESI or H-ESI probe	Trim fused-silica sample tube	If polyimide coating on the end of the sample tube has elongated	<i>Ion Max and Ion Max-S API Source Hardware Manual or H-ESI Probe User Guide</i>
HESI-II probe	Replace needle insert	If metal needle is plugged	<i>HESI-II Probe User Guide</i>
Forepump	Purge (decontaminate) oil	If the oil is cloudy or discolored	Manufacturer's documentation
	Add oil	If the oil level is below the MIN mark	Manufacturer's documentation
	Change oil	Every four months, or if oil remains cloudy or discolored after purging	Manufacturer's documentation

Table 5. Mass spectrometer maintenance procedures (Sheet 2 of 2)

Mass spectrometer component	Procedure	Frequency	Procedure location
Q0 ion optics	Clean Q0 quadrupole and lenses L11 and L12 ^b	As needed ^a	N/A
Mass analyzer	Clean Q1, Q2, and Q3 quadrupoles and lenses ^b	As needed ^a	N/A
Ion detection system	Clean ion detection system (electron multiplier and conversion dynode) ^b	Whenever the side cover plate of the vacuum manifold is removed	N/A
Cooling fans	Clean fan filters	Every 4 months	page 101
Ion detection system	Replace electron multiplier assembly ^b	If noise in spectrum is excessive or proper electron multiplier gain can not be achieved	N/A
Electronic modules	Replace electronic module ^b	If electronic module fails	N/A
PCBs	Replace PCB ^b	If PCB fails	N/A

^a The frequency of cleaning the components of the mass spectrometer depends on the types and amounts of samples and solvents that are introduced into the instrument. Cleaning of the Q0, Q1, Q2, and Q3 quadrupoles is rarely (if ever) required.

^b A Thermo Fisher Scientific service engineer should perform this maintenance procedure.

Frequency of Cleaning

The frequency of cleaning the components of the mass spectrometer depends on the types and amounts of samples and solvents that are introduced into the instrument. In general, for a given sample and ionization technique, the closer a mass spectrometer component is to the source of the ions, the more rapidly it becomes dirty.

- Clean the sample transfer line, sample tube, and API probe at the end of each operating day to remove any residual salts from buffered mobile phases or other contamination that might accumulate during normal operation. See [“Flushing the Sample Transfer Line, Sample Tube, and API Probe”](#) on [page 66](#).
- Periodically remove and clean the ion transfer tube and ion sweep cone of the API source. See [“Maintaining the Ion Sweep Cone and Ion Transfer Tube”](#) on [page 79](#).
- The S-lens and exit lens (on the TSQ Vantage) or the tube lens and skimmer (on the TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Quantum Ultra) of the Q00 ion optics become dirty at a slower rate than the API probe, ion sweep cone, and ion transfer tube. See [“Cleaning the S-Lens and Exit Lens \(TSQ Vantage\)”](#) on [page 87](#).
- The rf lens of the Q00 ion optics becomes dirty at a rate significantly slower than the API source, tube lens and skimmer, or S-lens and exit lens. See [“Cleaning the Q00 Ion Optics”](#) on [page 91](#).

When the performance of your system decreases significantly due to contamination, clean the components of the mass spectrometer in the following order:

1. Clean the API probe, ion sweep cone, and ion transfer tube.
2. Clean the S-lens and exit lens (on the TSQ Vantage) or the tube lens and skimmer (on the TSQ Quantum Access, TSQ Quantum Access MAX, and TSQ Quantum Ultra).
3. Clean the Q00 rf lens and lens L0.

Generally, do not open the vacuum manifold. Cleaning the Q0, Q1, Q2, and Q3 quadrupoles is rarely (if ever) required. (Q0 can be removed and cleaned along with Q00.) If you believe that the Q1, Q2, or Q3 quadrupoles need cleaning, contact Thermo Fisher Scientific and schedule a visit by a Thermo Fisher Scientific service engineer.

If you open the vacuum manifold, clean the electron multiplier and conversion dynode by blowing them with a clean, dry gas. If you do not open the vacuum manifold, the electron multiplier and conversion dynode do not normally require cleaning, although rarely the electron multiplier might require replacement.



CAUTION As with all chemicals, store and handle solvents and reagents according to standard safety procedures. Dispose of them according to local and federal regulations.

Maintaining the Ion Sweep Cone and Ion Transfer Tube

To prevent corrosion and to maintain optimum performance of the API source, remove and clean the ion sweep cone and the ion transfer tube on a regular basis.

Removing and Cleaning the Ion Transfer Tube

Buffer salts or high concentrations of sample can block the bore of the ion transfer tube. You can easily remove the ion transfer tube for cleaning, however, as it does not require venting the system before removal.

If the pressure in the ion transfer tube-skimmer region (as measured by the Convectron gauge) drops considerably below 1 Torr, a blocked ion transfer tube could be the cause. Check the Convectron gauge pressure by choosing **View > Instrument Status** from the EZ Tune window. The pressure is listed under Forepump Pressure.

❖ To remove and clean the ion transfer tube

1. Place the instrument in standby:
 - a. Choose **Start > Programs > Thermo Instruments > TSQ > TSQ Tune** to open the EZ Tune window.
 - b. From the EZ Tune window, choose **Control > Standby** (or click the **On/Standby** button) to put the mass spectrometer in Standby mode.




On



Off



Standby

2. Turn off the flow of liquid from the LC (or other sample introduction device) to the API source:
 - a. In the EZ Tune window, choose **Setup > Inlet Direct Control**. The Inlet Direct Control dialog box appears.
 - b. Click the **LC** tab and click  (Stop) to stop the LC pump.




CAUTION Lower the temperature of the ion transfer tube to 50 °C or less before you remove it. Handling a hot ion transfer tube or ion sweep cone can cause burns.

Attempting to remove a hot ion transfer tube might cause it to bind and break in the heater assembly or strip the threads of the ion source interface.

3. Lower the temperature of the ion transfer tube (capillary):

- a. In the EZ Tune window, click

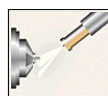

to open the Ion Source Devices dialog box.

See [Figure 41](#).

6 Maintenance

Maintaining the Ion Sweep Cone and Ion Transfer Tube

Figure 41. Ion Source Devices dialog box showing the ion transfer tube (capillary) at 270 °C

The dialog box titled "Ion Source Devices" contains a table of device parameters. The "Capillary Temperature" row is highlighted. Below the table, the "Tube Lens Offset" is set to 270.

Device	Value	Readback
✓ Spray Voltage	3500	3502
✓ Sheath Gas Pressure	0	0
✓ Ion Sweep Gas Pressure	0.0	0.1
✓ Aux Gas Pressure	0	0
✓ Capillary Temperature	270	270
✓ Capillary Offset	35	35
✓ Tube Lens Offset	101	102
✓ Skimmer Offset	-0	-0

LC Flow Rate (ul/min): 1.00 [Apply]

Tube Lens Offset: 270 [Apply] [Close] [Help]

- Click **Capillary Temperature** in the Device list.
- Set the ion transfer tube (capillary) temperature to 50 °C or lower in the Capillary Temperature box and click **Apply**.
- Allow the ion transfer tube (capillary) to cool to 50 °C or lower. See [Figure 42](#).

Figure 42. Ion Source Devices dialog box showing the ion transfer tube at 50 °C

The dialog box titled "Ion Source Devices" shows the "Capillary Temperature" row highlighted. The temperature has been changed to 50 °C. The "Capillary Offset" is now 35, and the "Tube Lens Offset" is 103.

Device	Value	Readback
✓ Spray Voltage	3500	3501
✓ Sheath Gas Pressure	0	0
✓ Ion Sweep Gas Pressure	0.0	0.1
✓ Aux Gas Pressure	0	0
✓ Capillary Temperature	50	50
✓ Capillary Offset	35	35
✓ Tube Lens Offset	101	103
✓ Skimmer Offset	-0	-0

LC Flow Rate (ul/min): 1.00 [Apply]

Capillary Temperature: 50 [Apply] [Close] [Help]

4. Remove the Ion Max or Ion Max-S ion source housing from the front of the mass spectrometer. See [“Removing the Ion Max or Ion Max-S Ion Source Housing”](#) on page 69.
5. Remove the ion sweep cone by pulling it off the ion source interface.
6. With the custom removal tool from the accessory kit (P/N 70111-20258), remove the ion transfer tube by turning it counterclockwise until you can pull it free from the ion source interface. See [Figure 43](#).

Figure 43. Removal tool attached to the ion transfer tube


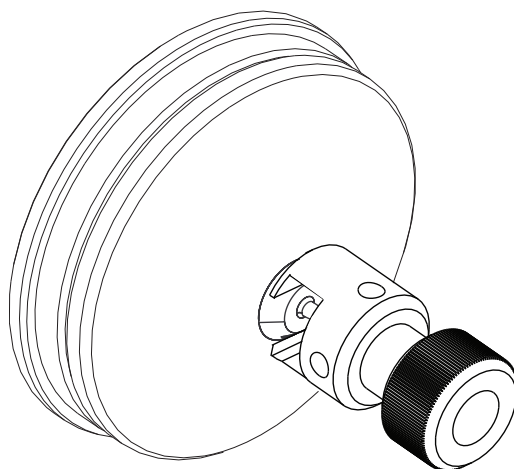


Figure 44. Spray cone, ion transfer tube, O-ring, and ion sweep cone of the ion source interface

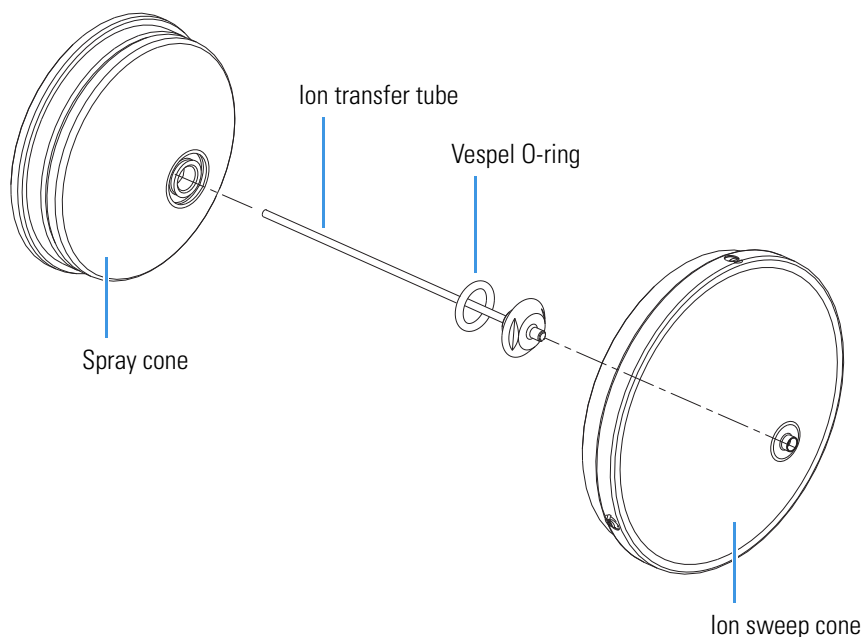

7. Soak the ion transfer tube in a dilute solution of nitric acid to remove contaminants.

6 Maintenance

Maintaining the Ion Sweep Cone and Ion Transfer Tube

8. Sonicate the ion transfer tube in distilled water.
9. Clean the ion sweep cone by wiping the inside and outside with methanol and a Kimwipe.
10. Remove, clean with methanol, and inspect the Vespel™ O-ring (P/N 97055-20442) that is seated in the spray cone under the entrance end of the ion transfer tube. Replace it if necessary.
11. Reseat the O-ring in the spray cone.



CAUTION When you reinstall the ion transfer tube, take these precautions:

- Ensure that everything is properly aligned to prevent stripping the threads on the ion transfer tube.
- For mass spectrometers other than the TSQ Vantage, be careful not to bend the ion transfer tube. Rotate the tube as you insert it. (The ion transfer tube of the TSQ Vantage will bend slightly when you insert it.)

12. Insert the ion transfer tube into the heater block. Rotate the tube as you insert it. After you insert it, turn the tube clockwise until it is fingertight.
13. Align the gas inlet on the ion sweep cone with the sweep gas supply port on the ion source mount. Firmly press the ion sweep cone into the ion source mount. See [Figure 45](#) and [Figure 46](#).
14. Reinstall the Ion Max or Ion Max-S source housing on the mass spectrometer as described in the topic “[Installing the Ion Max or Ion Max-S Ion Source Housing](#)” on [page 71](#).

Figure 45. Ion source interface, showing the sweep gas supply port in the API cone seal


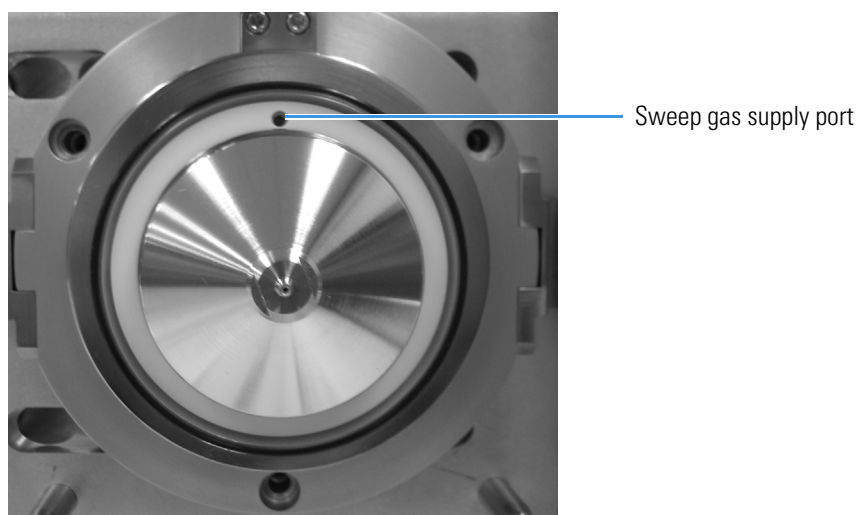


Figure 46. Ion sweep cones for FAIMS and APPI (left) and H-ESI, ESI, and APCI (right), showing the gas inlets

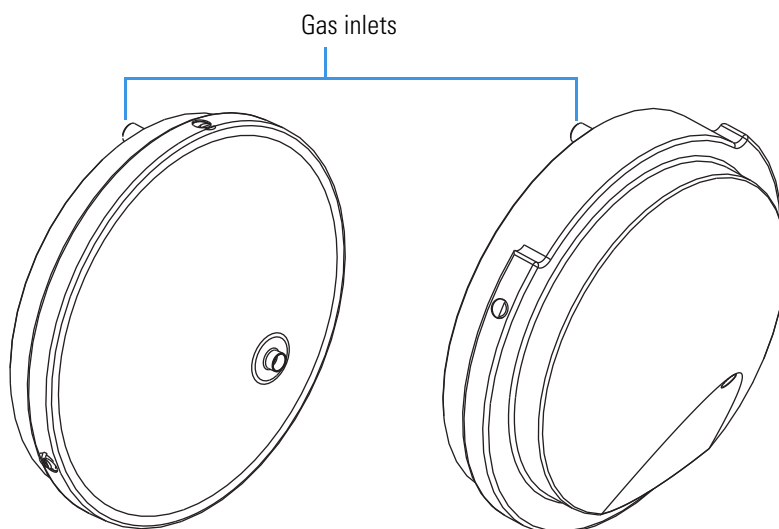

Note Unblocking the ion transfer tube should increase the Convectron gauge pressure to a normal value (approximately 1.5 Torr). If this method fails to clear the ion transfer tube, replace it. See [Table 6](#) for the correct ion transfer tube part number for your instrument.

Table 6. Ion transfer tube part numbers

Instrument	Ion transfer tube part number
TSQ Vantage series	70005-20423
TSQ Quantum Access series	70111-20972
TSQ Quantum Ultra series	97055-20199
TSQ Quantum Discovery Max	70111-20396
TSQ Quantum Discovery	70111-20100
TSQ Quantum AM	70111-20100
TSQ Quantum	70111-20100

Maintaining the Ion Source Interface Assembly

The ion source interface assembly includes the ion sweep cone, ion transfer tube, capillary heater assembly, S-lens or tube lens, and exit lens or skimmer. The ion transfer tube has a finite lifetime. You must replace the ion transfer tube if the ion transfer tube bore becomes corroded.

You can easily remove the ion sweep cone and ion transfer tube to clean or replace them without venting the system. See [“Maintaining the Ion Sweep Cone and Ion Transfer Tube”](#) on [page 79](#).

To clean the ion source interface assembly components, other than the ion sweep cone and ion transfer tube, follow these procedures:

1. [Shutting Down and Venting the System](#)
2. [Removing the Ion Max or Ion Max-S Ion Source Housing](#)
3. [Removing the Ion Source Interface](#)
4. [Removing the S-Lens and Exit Lens \(TSQ Vantage\)](#)
5. [Cleaning the S-Lens and Exit Lens \(TSQ Vantage\)](#)
6. [Removing the Tube Lens and Skimmer \(TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra\)](#)
7. [Cleaning the Tube Lens and Skimmer \(TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra\)](#)
8. [Reinstalling the S-Lens and Exit Lens \(TSQ Vantage\)](#)
9. [Reinstalling the Tube Lens and Skimmer \(TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra\)](#)
10. [Reinstalling the Ion Source Interface Assembly](#)
11. [Reinstalling the Ion Max or Ion Max-S Ion Source Housing](#)
12. [Starting Up the System](#)

Shutting Down and Venting the System

Shut down and vent the system. See [“Shutting Down the System Completely”](#) on [page 52](#).

Removing the Ion Max or Ion Max-S Ion Source Housing

You must remove the Ion Max or Ion Max-S ion source housing to access the ion source interface assembly. See [“Removing the Ion Max or Ion Max-S Ion Source Housing”](#) on [page 69](#).

Removing the Ion Source Interface

You must remove the ion source interface to access the exit lens or skimmer, the S-lens or tube lens, and the capillary heater assembly.




CAUTION Wait for the ion source interface to cool to ambient temperature before you remove it.

❖ To remove the ion source interface

1. Wear clean, talc-free gloves.
2. Using a right-to-left motion, move the ion source interface release lever several times to loosen the ion source interface from the ion optics cage assembly. See [Figure 47](#).
3. Grab the exposed outer grooves of the ion source interface and firmly pull the assembly straight out.
4. Place the assembly on a clean, lint-free surface.

Figure 47. Removing the ion source interface

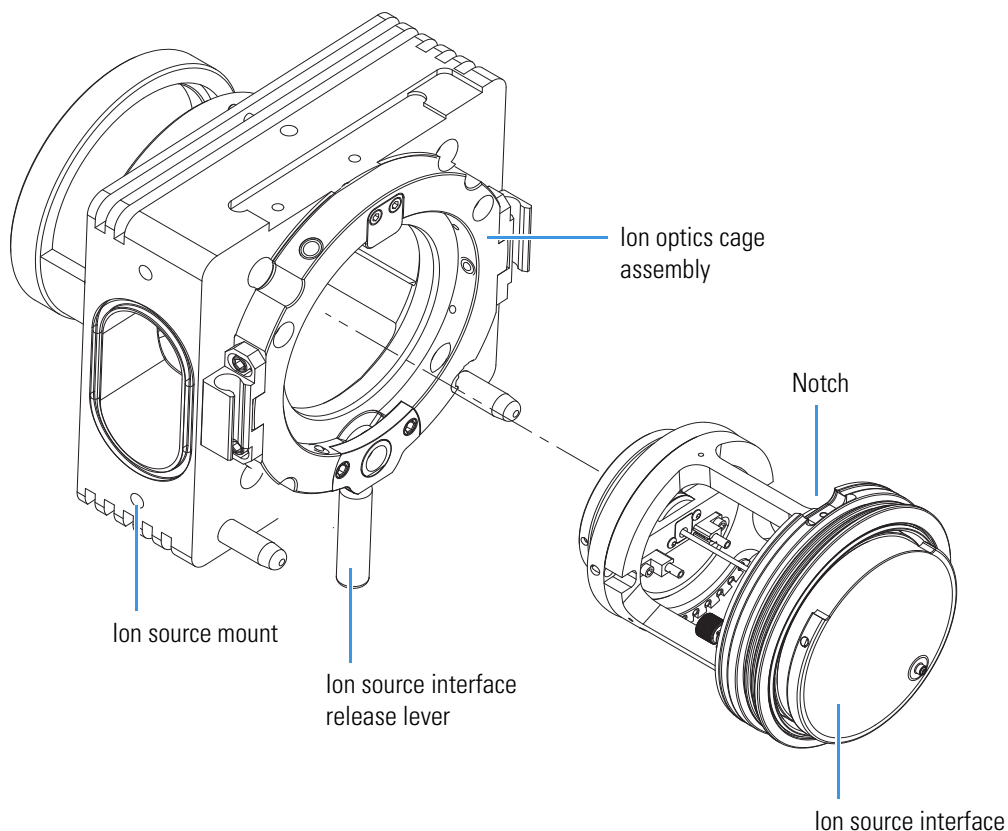


Removing the S-Lens and Exit Lens (TSQ Vantage)

You must remove the S-lens and exit lens from the ion source interface cage before they can be cleaned.

❖ To remove the S-lens and exit lens from the ion source interface assembly

1. Wearing clean, lint- and powder- free gloves, loosen and extend the two thumbscrews that secure the S-lens to the ion source interface cage and the exit lens to the S-lens. See [Figure 48](#).
2. Grasp the two thumbscrews and carefully pull the S-lens straight out of the ion source interface cage. Place it on a clean, lint-free surface.
3. Remove the exit lens from the S-lens and place it on a clean, lint-free surface.

Figure 48. Removing the exit lens and S-lens from the ion source interface cage of the TSQ Vantage

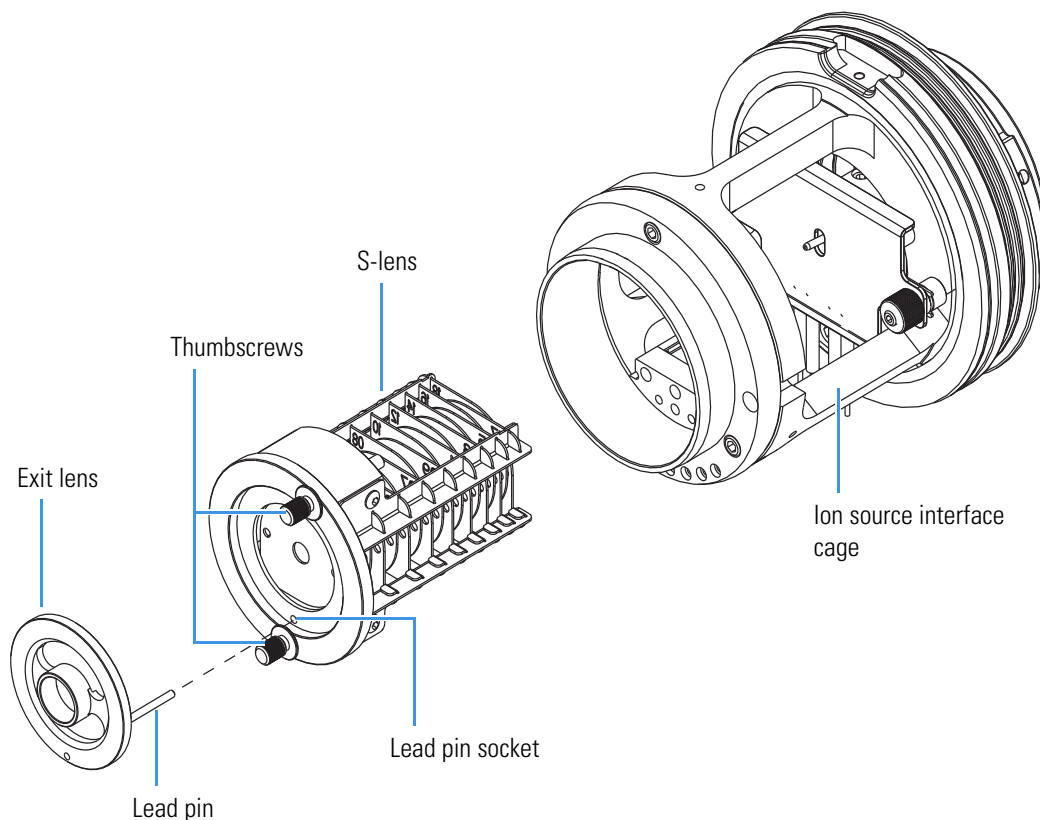

Cleaning the S-Lens and Exit Lens (TSQ Vantage)

❖ To clean the S-lens and exit lens

Note Wear clean gloves when you handle the S-lens and exit lens.



CAUTION Do not clean the S-lens or exit lens with detergents, acidic or caustic substances, or abrasives.

1. Inspect the S-lens and exit lens under magnification for any lint, particulates, and sample buildup or coatings.
2. Sonicate the S-lens and exit lens for 10–15 min. in a 50:50 methanol/water solution or a 1% solution of Liquinox in water. If a sonicator is not available, do the following:
 - a. To clean the exit lens, use a soft toothbrush with a 1% solution of Liquinox in water.
 - b. To clean the S-lens, use a chamois-tip swab with a 1% solution of Liquinox in water, and use a 6000 grit MICRO-MESH polishing swab (P/N 00301-01911) to clean the areas that are inaccessible to the chamois-tip swab.
3. For the exit lens, clean the bore by using the 6000 grit MICRO-MESH polishing swab.
4. Rinse the S-lens and exit lens thoroughly with water.
5. Rinse the S-lens and exit lens with fresh methanol.
6. Air dry the S-lens and exit lens or blow dry them with nitrogen gas. Ensure that all solvent has evaporated from the components before reassembly.
7. Inspect the S-lens and exit lens under magnification for any lint or particulates.


Removing the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)

You must remove the tube lens and skimmer from the ion source interface assembly before they can be cleaned.

❖ To remove the tube lens and skimmer from the ion source interface assembly

1. Wearing clean, lint- and powder- free gloves, reach behind the skimmer with your fingers and gently press the skimmer out of the contact ring support. If necessary, loosen the set screws. See [Figure 49](#).
2. Place the skimmer on a clean, lint-free surface.
3. Push the tube lens, from the back, out of the contact ring support.
4. Place the tube lens on a clean, lint-free surface.

Figure 49. Removing the skimmer and tube lens from the ion source interface cage of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra

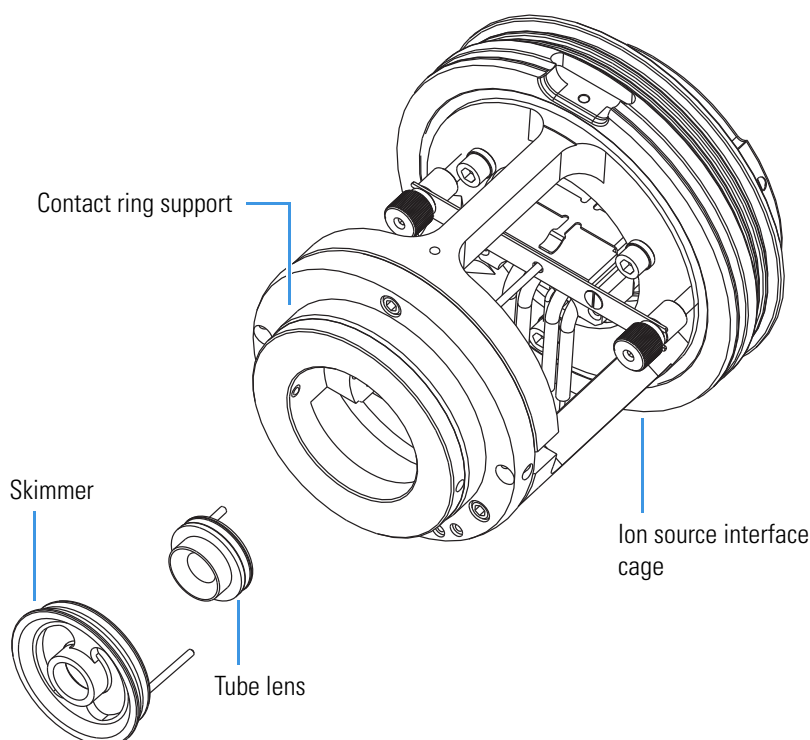

Cleaning the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)

❖ To clean the tube lens and skimmer

Note Wear clean gloves when you handle the tube lens and skimmer.



CAUTION Do not clean the tube lens or skimmer with detergents, acidic or caustic substances, or abrasives.

The sharp edge of the skimmer cone is fragile and any deformation might result in a loss of efficient ion transmission. Use care not to drop the skimmer when cleaning it. Sonicate the skimmer alone with the cone side up.

Sonicate the tube lens and skimmer alone in organic or aqueous solution. For most cleaning applications, HPLC-grade methanol is the solvent of choice. However, use of buffers or salt solutions might require that you use an aqueous solution. If you need to use a solvent other than methanol, after cleaning the component, flush the component with distilled water and then flush it with methanol as a final wash. You can air dry the skimmer and tube lens or blow dry them with nitrogen gas. In all cases, ensure that all solvent has evaporated from the components before reassembly.

Reinstalling the S-Lens and Exit Lens (TSQ Vantage)

❖ To reinstall the S-lens and exit lens

1. Reinstall the exit lens onto the S-lens:
 - a. Wearing clean, lint- and powder- free gloves, extend the thumbscrews on the S-lens. See [Figure 48](#) on [page 86](#).
 - b. Insert the exit lens lead pin into the lead pin socket and seat the exit lens onto the S-lens.
2. Reinstall the exit lens and S-lens into the ion source interface cage:
 - a. Hold the S-lens by the thumbscrews and orient the thumbscrews toward the thumbscrew holes on the ion source interface. Only one orientation will work.
 - b. Carefully insert the S-lens into the ion source interface cage.
 - c. Tighten the thumbscrews until fingertight.

Reinstalling the Tube Lens and Skimmer (TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra)

❖ To reinstall the tube lens and skimmer


1. Reinstall the tube lens into the ion source interface cage (see [Figure 50](#)):
 - a. Orient the tube lens so that the lead pin points toward the socket on the tube lens connection wire in the contact ring support.
 - b. Insert the lead pin into the socket and firmly press the tube lens into the contact ring support.



CAUTION Ensure that you do not scratch or nick the skimmer cone.

2. Reinstall the skimmer into the ion source interface assembly:
 - a. Orient the skimmer so that the lead pin points toward the socket on the tube lens connection wire in the contact ring support.
 - b. Insert the lead pin into the socket and firmly press the skimmer into the contact ring support.

Figure 50. Ion source interface cage of the TSQ Quantum Access, TSQ Quantum Access MAX, or TSQ Quantum Ultra, showing the skimmer and tube lens lead pin sockets

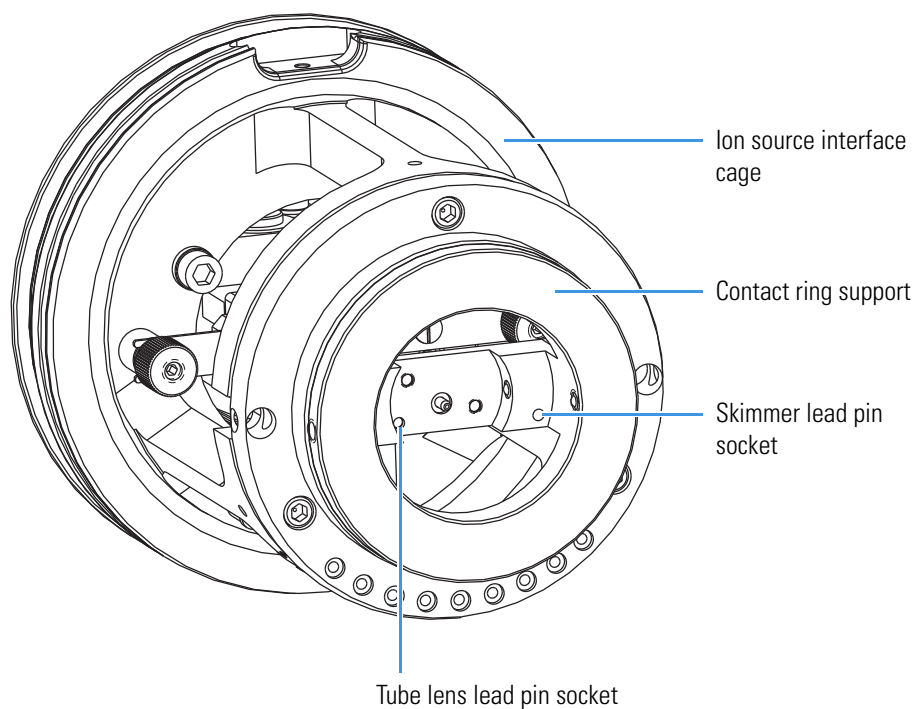

Reinstalling the Ion Source Interface Assembly

❖ To reinstall the ion source interface assembly

1. Orient the ion source interface assembly so that the sweep gas inlet notch on the ion source interface is aligned with the sweep gas supply block on the ion source mount.
2. Carefully slide the ion source interface assembly into the ion optics cage assembly.

Reinstalling the Ion Max or Ion Max-S Ion Source Housing

Reinstall the Ion Max or Ion Max-S ion source housing as described in the topic [“Installing the Ion Max or Ion Max-S Ion Source Housing”](#) on [page 71](#).

You might also need to reinstall an ion source probe and LC liquid lines, as appropriate, as described in the corresponding topics in the *TSQ Series Getting Connected Guide*.

Starting Up the System

Start up the system as described in the topic [“Starting Up the System after a Complete Shutdown”](#) on [page 54](#).

Cleaning the Q00 Ion Optics

Because an accumulation of chemicals on the surfaces of the tube lens, skimmer, and Q00 rf lens forms an insulating layer that can modify the electrical fields that control ion transmission, clean ion optics components are essential for the proper operation of the instrument. The tube lens and skimmer require cleaning less often than the API source. The Q00 rf lens requires cleaning less often than the tube lens and skimmer. The frequency of cleaning depends on the type and quantity of the compounds that you analyze.

Clean or replace the Q00 ion optics components by doing the following:

1. [Shutting Down and Venting the System](#)
2. [Removing the Ion Max or Ion Max-S Ion Source Housing](#)
3. [Removing the Ion Source Interface](#)
4. [Removing the Ion Optics Cage Assembly](#)
5. [Disassembling the Ion Optics Cage Assembly](#)
6. [Cleaning Q00 and L0](#)
7. [Reassembling the Ion Optics Cage Assembly](#)
8. [Reinstalling the Ion Optics Cage Assembly](#)
9. [Reinstalling the Ion Source Interface Assembly](#)
10. [Reinstalling the Ion Max or Ion Max-S Ion Source Housing](#)
11. [Starting Up the System](#)

Shutting Down and Venting the System

Shut down and vent the system as described in the topic [“Shutting Down the System Completely”](#) on [page 52](#).

Removing the Ion Max or Ion Max-S Ion Source Housing

Remove the Ion Max or Ion Max-S ion source housing to access the ion source interface assembly. Remove the Ion Max or Ion Max-S ion source housing as described in the topic [“Removing the Ion Max or Ion Max-S Ion Source Housing”](#) on [page 69](#).

Removing the Ion Source Interface Assembly


Remove the ion source interface assembly to access the Q00 ion optics assembly. Remove the ion source interface assembly as described in the topic [“Removing the Ion Source Interface”](#) on [page 85](#).

Removing the Ion Optics Cage Assembly

❖ To remove the ion optics cage assembly


1. Using a Phillips screwdriver, remove the three screws fastening the ion source interface cover to the mass spectrometer, and place the cover in a safe place. See [Figure 51](#).

Figure 51. Removing the ion source interface cover screws

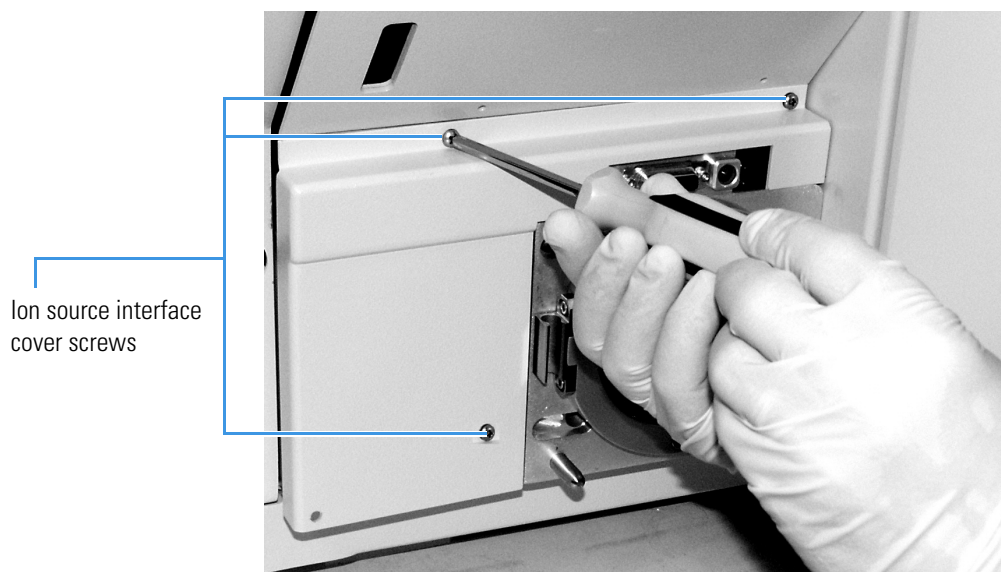

2. Using a hex ball driver, remove the roughing pump adapter access plate. See [Figure 52](#).

Figure 52. Removing the roughing pump adapter access plate screws

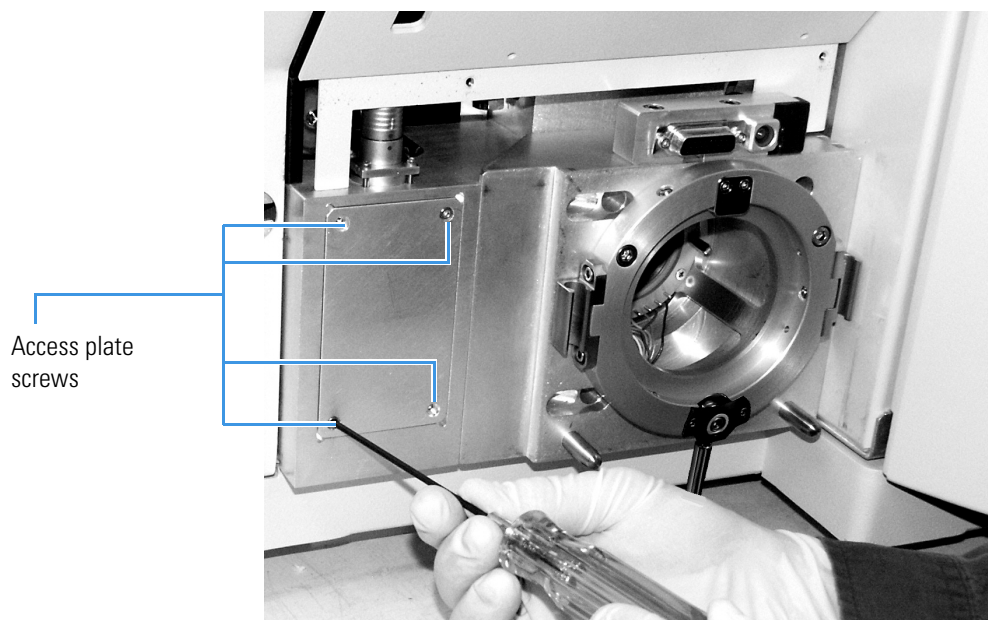


3. Unplug the ion source cable harness from the connector. See [Figure 53](#).

Figure 53. Unplugging the ion source cable harness connector



4. Remove the ion optics cage assembly:
 - a. Using a hex ball driver, remove the three screws fastening the ion optics cage assembly to the ion source mount assembly. See [Figure 54](#).

Figure 54. Removing the ion optics cage assembly retaining screws

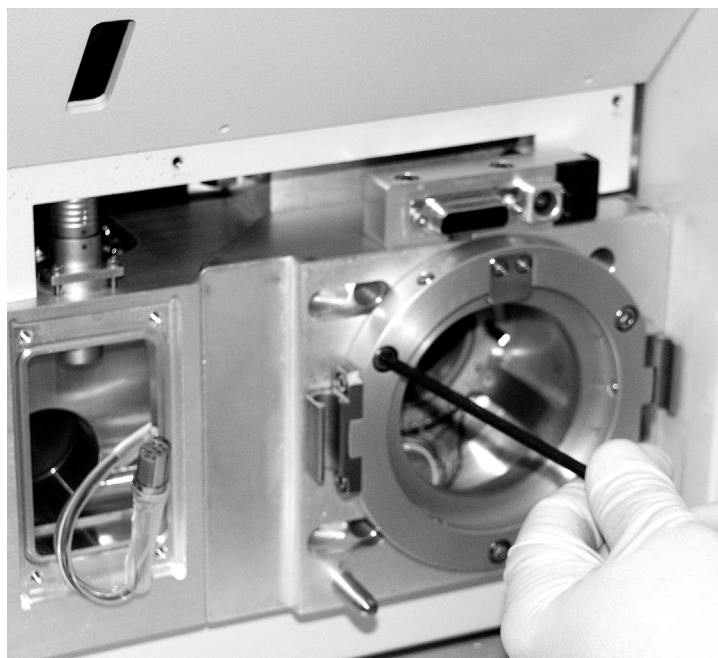

- b. Carefully feed the cable harness back through the roughing pump adapter and out of the front of the cage assembly to ensure that it does not get trapped between the cage and the ion source mount assembly while you are removing the cage assembly. See [Figure 55](#).

Figure 55. Pulling the ion source cable harness out of the ion optics cage assembly



- c. Carefully slide the ion optics cage assembly out of the ion source mount assembly. See [Figure 56](#).

Figure 56. Removing the ion optics cage assembly



- d. Place the ion optics cage assembly on a clean, lint-free surface.

Disassembling the Ion Optics Cage Assembly

The Q00 rf lens and L0 lens must be removed from the ion optics cage assembly before they can be cleaned.


Note Make careful note of the proper orientation of all wiring before removing any electrical lead from its contact.

❖ To disassemble the ion optics cage assembly

Note Wear clean gloves when you handle the ion optics components.


1. Using a hex ball driver, remove the three screws that fasten the Q0 mount assembly to the ion optics cage. See [Figure 57](#).

Figure 57. Ion optics cage assembly with Q0 mount assembly and Q00/L0 mount assembly removed

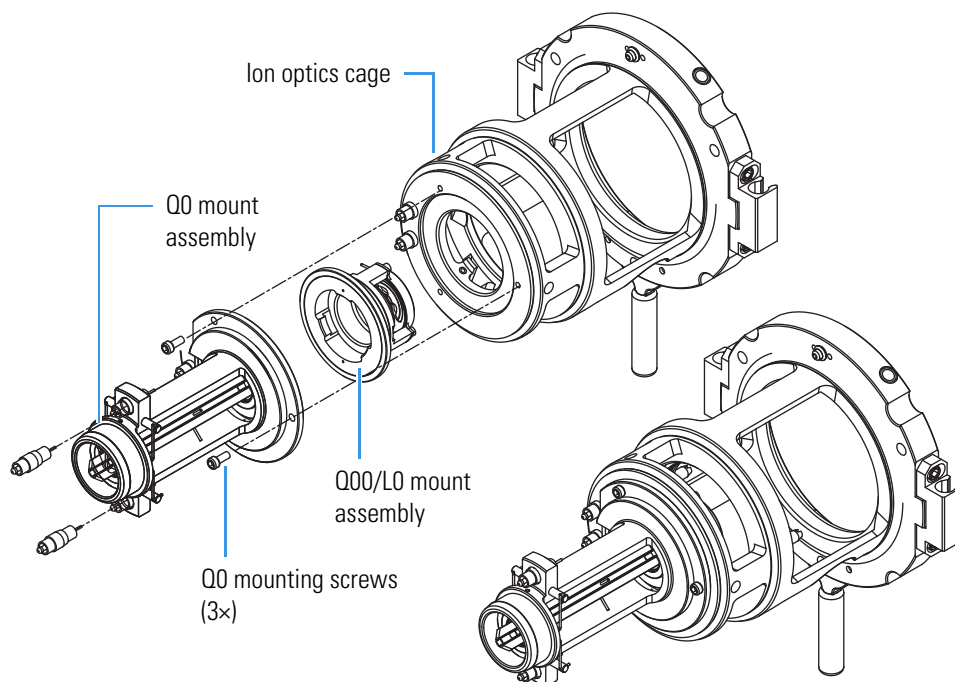

2. Remove the Q0 mount assembly. Place it on a clean, lint-free surface.
3. Using needle nose pliers, disconnect the two rf lead wires from the Q00 lead pins.
4. Carefully push the Q00/L0 mount assembly out of the ion optics cage by pressing on the mount assembly with your fingers.
5. Using a hex ball driver, remove the three Q00 mounting screws that fasten the Q00 rf lens assembly to the Q00/L0 mount. Leave the lens L0 mounted in the PEEK™ Q00/L0 mount.

Figure 58. Exploded diagram of the Q00/L0 mount assembly

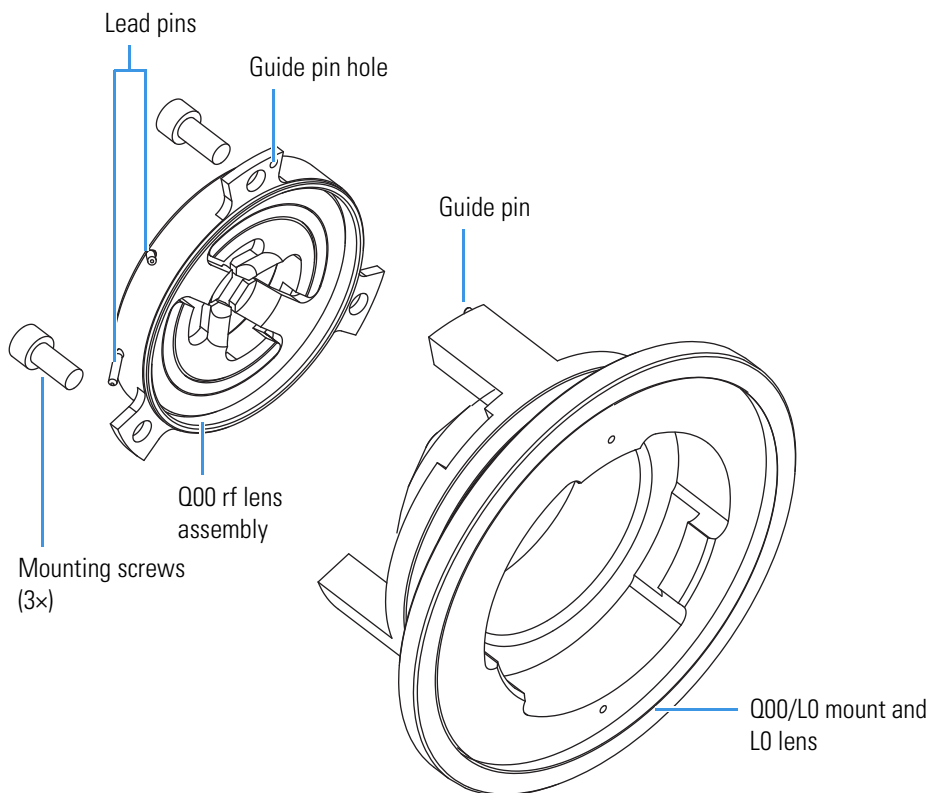

6. Place all components onto a clean, lint-free surface.



CAUTION Do not disassemble the Q00 rf lens assembly any further.

Cleaning Q00 and L0

❖ To clean the Q00 rf lens and L0 lens

1. Clean the Q00 rf lens and L0 lens with a Kimwipe™ soaked in methanol.



CAUTION Do not use abrasives to clean the Q00 rf lens. Do not soak the Q00 rf lens and L0 lens in solvent.

Note Wear clean gloves to handle the parts after you clean them.

2. Dry the parts with a rapid stream of nitrogen gas.
3. Inspect each part for contamination and dust. If necessary, repeat the cleaning procedure.

Reassembling the Ion Optics Cage Assembly

❖ To reassemble and mount the Q00/L0 assembly


1. Reinstall the Q00 assembly on the Q00/L0 mount:
 - a. Orient the Q00 assembly so that the guide pin on the Q00/L0 mount inserts into the guide pin hole in the Q00 assembly.
 - b. Secure the Q00 assembly with the 3 Q00 mounting screws.
2. Reinstall the Q00/L0 mount assembly in the ion optics cage:
 - a. Orient the Q00/L0 mount assembly so that the lead pins are closest to the lead wires in the ion optics cage.
 - b. Press the Q00/L0 assembly into the ion optics cage. See [Figure 57](#) on [page 96](#).
3. Use needle nose pliers to carefully reconnect the top rf lead wire to the top Q00 lead pin and the bottom rf lead wire to the bottom Q00 lead pin.
4. Reinstall the Q0 mount assembly onto the ion optics cage:
 - a. Orient the Q0 mount assembly so that the flat side is on the same side as the two contacts.
 - b. Secure the Q0 mount assembly on the ion optics cage with the 3 Q0 mounting screws. See [Figure 57](#) on [page 96](#).

Reinstalling the Ion Optics Cage Assembly

❖ To reinstall the ion optics cage assembly


1. Reinstall the 3.25 in. ID Viton™ O-ring into the groove on the ion source mount assembly. See [Figure 59](#).

Figure 59. Ion source mount assembly, showing groove for O-ring

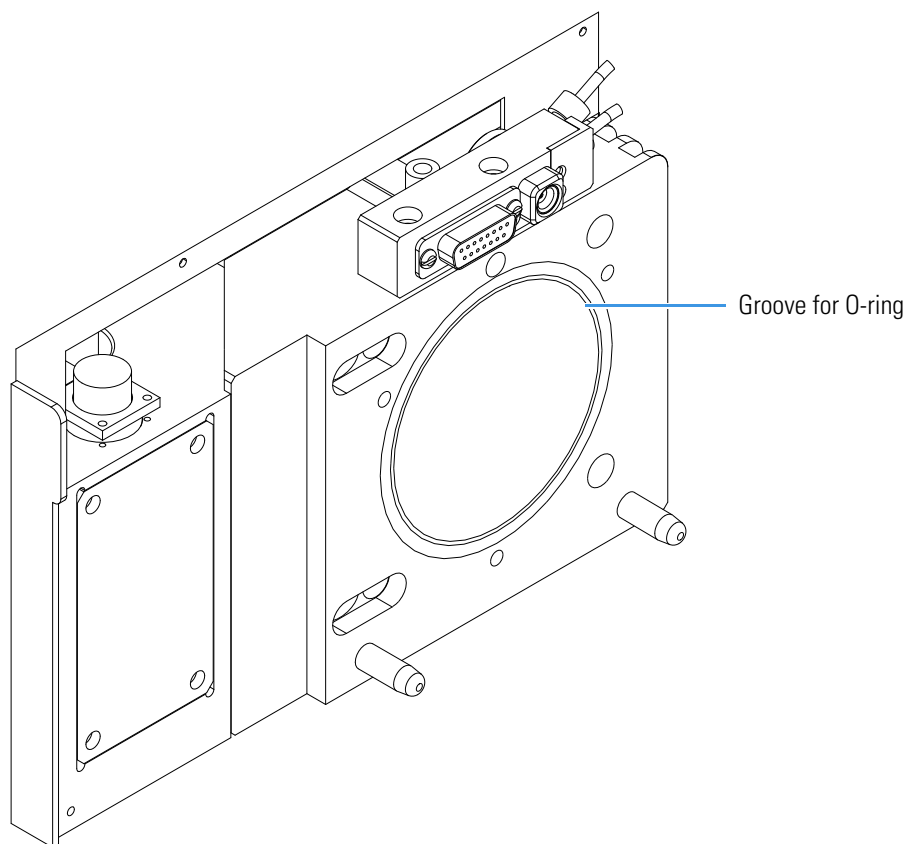

2. Orient the ion optics cage assembly with the contact pins down, and then partially insert it into the ion source mount assembly.
3. Feed the cable harness into the roughing pump adapter and reconnect the cable. Rotate the ion optics cage assembly if necessary.
4. Insert the ion optics cage assembly completely so that the two interstage contact pins are inserted into the sockets in the ion source mount assembly.
5. Ensure that the 3.25 in. ID Viton O-ring is properly seated between the ion optics cage assembly and the ion source mount assembly.
6. Carefully press the wiring harness down into the recess in the bottom of the ion source mount assembly so that the wires are out of the path of the ion source interface assembly that you will install in the next procedure. See [Figure 60](#).

Figure 60. Ion optics cage assembly, installed

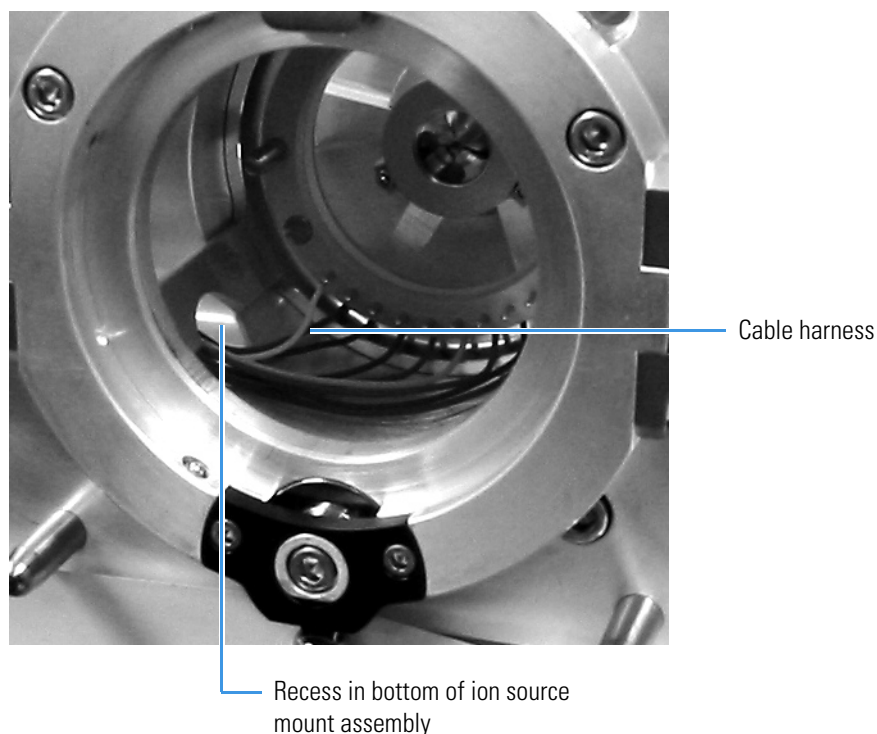

7. Reinstall the three ion optics cage assembly screws that fasten the cage assembly to the ion source mount assembly. See [Figure 54](#) on [page 93](#).
8. Plug in the ion source cable harness connector. Secure the connector. See [Figure 53](#) on [page 93](#).
9. Reinstall the roughing pump adapter access plate and screws. See [Figure 52](#) on [page 92](#).
10. Reinstall the ion source interface cover and screws. See [Figure 51](#) on [page 92](#).

Reinstalling the Ion Source Interface Assembly

Reinstall the ion source interface assembly as described in the topic “[Reinstalling the Ion Source Interface Assembly](#)” on [page 90](#).

Reinstalling the Ion Max or Ion Max-S Ion Source Housing

Reinstall the Ion Max or Ion Max-S ion source housing as described in the topic “[Installing the Ion Max or Ion Max-S Ion Source Housing](#)” on [page 71](#).

Starting Up the System

Start up the system as described in the topic “[Starting Up the System after a Complete Shutdown](#)” on [page 54](#).

Maintaining the Forepump

The forepump is located under the TSQ mass spectrometer work table. Inspecting, adding, purging, and changing the oil are all that is required to maintain the forepump.

Forepump oil is a translucent, light amber color, which you must check often. During normal operation, oil must always be visible in the oil level sight glass between the MIN and MAX marks. If the oil level is below the MIN mark, add oil. If the oil is cloudy or discolored, purge the oil to decontaminate dissolved solvents. If the pump oil is still discolored, change it. Plan to change the pump oil after every 3000 hours (about four months) of operation. Refer to the manufacturer's documentation for procedures for purging, adding, and changing the forepump oil.

Cleaning the Fan Filters

You must clean the two fan filters every four months. The fan filters are located on the of the front cover and on the right side cover of the TSQ Series mass spectrometer.

❖ To clean the fan filters

1. Remove the fan filter from the right side cover:
 - a. On the TSQ Vantage, using a Phillips screwdriver, remove the six screws that hold the blue appliqué to the right-side cover.
 - b. Remove the appliqué.
 - c. Loosen and remove the four screws that hold the filter bracket to the right-side cover. Note that from the outside of the mass spectrometer the filter bracket appears to be a grille.
 - d. Remove the filter and filter bracket.
2. Remove the fan filter from the front cover:
 - a. Remove the front cover of the TSQ mass spectrometer:
 - i. Disconnect all tubing to the divert/inject valve and syringe pump.
 - ii. Depress the two (blue) spring catches that are located on either side of front cover.
 - iii. Pull the front cover straight off by about 6 in. (15 cm) to access the Front Panel PCB cable and the two grounding wires.
 - iv. Disconnect the Front Panel PCB cable from the TSQ mass spectrometer by pushing the connector retainer up and then pulling the cable free.
 - v. Pull the two grounding wires free from their respective spade connectors.
 - vi. Remove the front cover.

- b. Loosen and remove the four screws that hold the filter bracket to the front cover.
 - c. Remove the filter and filter bracket.
3. Wash the fan filters in a solution of soap and water.
4. Rinse the fan filters with tap water.
5. Squeeze the water from the fan filters and allow them to air dry.
6. Reinstall the fan filter on the right-side cover:
 - a. Place the filter in the filter bracket and position the filter bracket and filter on the right-side cover.
 - b. Install and tighten the four screws that hold the filter bracket to the right-side cover.
 - c. On the TSQ Vantage, position the blue appliqué on the right-side cover.
 - d. Install and tighten the six screws that hold the appliqué to the right-side cover.
7. Reinstall the fan filter on the front cover:
 - a. Place the filter in the filter bracket and position the filter bracket and filter on the of the front cover.
 - b. Install and tighten the four screws that hold the filter bracket to the front cover.
 - c. Reinstall the front cover of the TSQ mass spectrometer.
 - i. Position the front cover about 6 in. (15 cm) from the TSQ mass spectrometer.
 - ii. Reconnect the grounding wire to the space connector.
 - iii. Reconnect the Front Panel PCB cable to the TSQ mass spectrometer by inserting the cable into the connector. Push the connector retainer down to secure the cable.
 - iv. Align the front cover with the two guide pins on the TSQ mass spectrometer.
 - v. Push the front cover against the TSQ mass spectrometer until the two spring catches lock.
 - vi. Reconnect all tubing to the divert/inject valve and the syringe pump.

Diagnostics and PCB and Assembly Replacement

The TSQ system diagnostics can test many components of the TSQ mass spectrometer. If there is a problem with the instrument electronics, the diagnostics can often locate the problem. Replacing a faulty PCB or assembly can usually correct the problem. After the PCB or assembly is replaced, the diagnostic tests are rerun to verify that the instrument is functioning properly.

A subset of the diagnostic tools that are part of the Tune application targets specific functions of the entire mass spectrometer. In general, these system evaluation tools do not contain pass or fail criteria nor do they rate the performance. Instead, they provide data that can be interpreted by a Thermo Fisher Scientific service engineer. Because the system evaluations examine complex interactions, the service engineer uses the data in conjunction with other tools and tests to form a diagnosis.

Note Three levels of protection are possible:

- No protection—All operators can access all workspaces.
- Automatic protection—Tune Master uses the default password, lctsq, to protect the secure workspaces.
- Custom password protection—The Key Operator (or Laboratory Administrator or Manager) can select a password to protect the secure workspaces.

If your TSQ system has been password protected, you must obtain the password before you can access the secure workspaces (including the System Tune and Calibration workspace). If the instrument password is lost, you must reinstall the TSQ software to reset the default password (lctsq).

Contents

- [Running the TSQ System Diagnostics](#)
- [Replacing a Fuse](#)
- [Replacing PCBs and Power Supplies](#)

Running the TSQ System Diagnostics

The TSQ system diagnostics are used to test the major electronic circuits within the instrument and indicate whether the circuits pass or fail the tests. If there is a problem with the instrument electronics, the TSQ system diagnostics can often locate the problem.



CAUTION Non-expert users should run only the following diagnostics because certain diagnostics can overwrite system parameters.

- Ion Source Board
- System Control
- Analyzer Control Board

The TSQ system diagnostics do not diagnose problems that are not electrical in nature. For example, they do not diagnose poor sensitivity due to misaligned or dirty components or to improper tuning. Therefore, it is important that the person running the diagnostics be familiar with system operation and basic hardware theory as well as the details of the diagnostics.

Typically, only a Thermo Fisher Scientific service engineer runs diagnostic tests because certain tests can overwrite system parameters. However, before calling a Thermo Fisher Scientific service engineer to run diagnostics, consider the following:

- Did the system fail when you were running samples?
- Did problems occur after you performed maintenance on the instrument, data system, or peripherals?
- Did you change the configuration of the system, cables, or peripherals just before the problem occurred?

If the answer is yes to the first item above, there is the possibility of a hardware failure, and running the diagnostics is appropriate.

If the answer is yes to either of the last two questions above, the problem is probably mechanical, not electrical. Verify again that alignment, configurations, and cable connections are correct before you call the Thermo Fisher Scientific service engineer. Keep careful notes documenting the nature of the problem and the corrective steps you have taken. If you are not successful in correcting the problem, you can email this information to your service engineer. The Field Service department can then do a preliminary evaluation of the problem before an engineer arrives at your site.

❖ To run diagnostics



CAUTION Non-expert users should run only the following diagnostics:

- Ion Source Board
- System Control
- Analyzer Control Board

1. Open TSQ Tune:
 - a. Right-click the **Start** button and choose **Explore**.
 - b. Browse to the **C:\Thermo\Instruments\TSQ\system\programs** folder.
 - c. Double-click **TSQTune**.
2. Choose **Workspace > Full Instrument Control** to display the Full Instrument Control Workspace.
3. Choose **Workspace > Diagnostics** to display the Diagnostics Workspace.
4. Select Test Name: **Ion Source Board**, **System Control**, or **Analyzer Control Board**.
5. Select how many times you want to run the tests.
6. Click **Start** to start the diagnostics.

Replacing a Fuse

Fuses protect the various circuits by opening the circuits whenever overcurrent occurs. On the TSQ mass spectrometer, a failed fuse indicates a failed board or electronic module that must be replaced by a Thermo Fisher Scientific service engineer.



CAUTION Due to the complexity of accessing the TSQ mass spectrometer electronic assemblies, only a Thermo Fisher Scientific service engineer can replace fuses.

Replacing PCBs and Power Supplies

The electronic assemblies that control the operation of the mass spectrometer are distributed among various PCBs and other modules located in the tower, in the embedded computer, and on or around the vacuum manifold of the mass spectrometer.



CAUTION The TSQ mass spectrometer electronic assemblies are close-packed to minimize the size of the system. Due to the complexity of removing and reinstalling the TSQ mass spectrometer electronic assemblies, only a Thermo Fisher Scientific service engineer can replace electronic assemblies.

Replaceable Parts

This chapter contains part numbers for replaceable and consumable parts for the mass spectrometer and kits. To ensure proper results in servicing the TSQ system, order only the parts listed or their equivalent.

Contents

- [MS Accessory Kit](#)
- [TSQ Vantage Instrument Specific Accessory Kit](#)
- [TSQ Quantum Ultra, Ultra AM, and Ultra EMR Instrument Specific Accessory Kit](#)
- [HESI-II Probe Kit](#)
- [TSQ Quantum Access Instrument Specific Accessory Kit](#)
- [TSQ Accurate Mass Calibration Compounds Kit](#)
- [Fittings, Ferrules, and Sample Loops](#)

MS Accessory Kit

Accessory Kit	70111-62034
Fuse, 0.25 A, 250 V, 5×20	00006-01700
Fuse, 0.25 A, 250 V, 5×20	00006-05080
Fuse, 0.50 A, 250 V, 5×20	00006-07608
Fuse, 1.0 A, 250 V, 5×20	00006-07610
Fuse, 2.0 A, 250 V	00006-09102
Fuse, 1.0 A, 250 V, 5×20	00006-10510
Fuse, 4.0 A, 250 V	00006-10705
Fuse, 0.16 A, 250 V, 5×20	00006-11204
Fuse, 4 A, 250 V	00006-11420
Fuse, 1.0 A, 250 V	00006-14015
Fuse, 6.3 A, 250 V, 5×20	00006-11450
Fitting, ferrule, Swagelok™, Back, 1/8 in.	00101-02500
Fitting, ferrule, Swagelok, Back, 1/4 in.	00101-04000
Fitting, ferrule, Swagelok, Front, 1/8 in.	00101-08500
Fitting, ferrule, Swagelok, Front, 1/4 in.	00101-10000

8 Replaceable Parts

TSQ Vantage Instrument Specific Accessory Kit

Fitting, Swagelok, Nut, 1/4 in., brass	00101-12500
Fitting, Swagelok, Nut, 1/8 in., brass	00101-15500
Fitting, HPLC adaptor, 10-32 × 1/4-28	00101-18080
Ferrule, 0.008 in. ID, KEL-F™, HPLC	00101-18114
Fitting, Fingertight 2, Upchurch™	00101-18195
Ferrule, Fingertight 2, Upchurch	00101-18196
Fitting, HPLC, Tee, 0.020 orifice, PEEK	00101-18204
Tubing, fused silica, 0.150 mm ID × 0.390 mm OD	00106-10498
Tubing, fused silica, 0.10 mm ID × 0.19 mm OD	00106-10499
Tubing, fused silica, 0.05 mm ID × 0.19 mm OD, deactivated	00106-10502
Tube, hypodermic 28-gauge × 10 in. length	00106-20000
Fitting, HPLC, 10-32, long, one piece	00109-99-00016
Septa	00301-16999
Syringe, 500 µL, Gastight™, Removable needle	00301-19016
Tubing, PEEK, 0.005 ID × 1/16 in. OD, Red, 5 ft. (1.5 m long)	00301-22912
Tubing, Teflon™, 0.030 in. ID × 1/16 in. OD	00301-22915
Syringe Adapter Kit	70005-62011
Needle, Corona, APCI	70005-98033
Tool, ion transfer tube removal	70111-20258
O-ring, graphite Vespel (for ion transfer tube)	97055-20442

TSQ Vantage Instrument Specific Accessory Kit

Accessory Kit	70111-62081
5/32 in. Ball driver	00025-10020
Reserpine	00301-12901
Ultramark 1621 (EMR only)	00301-12200
Ion transfer tube	70005-20423
Polytyrosine, liquid	00301-22924
Polytyrosine, solid	00301-22925
Cable, LC/MS	60053-63035

TSQ Quantum Ultra, Ultra AM, and Ultra EMR Instrument Specific Accessory Kit

Accessory Kit	70111-62057
5/32 in. Ball driver	00025-10020
Ultramark 1621 (EMR only)	00301-12200
Reserpine	00301-12901
Needle seal	07055-20271
ESI needle	07055-20273
Kit, metal needle, H-ESI	97055-62026
Ion transfer tube	97055-20199
Polytyrosine, liquid	00301-22924
Polytyrosine, solid	00301-22925
Cable, LC/MS	60053-63035

HESI-II Probe Kit

HESI-II Probe Kit	OPTON -20037
High-flow needle insert assembly	OPTON-53010
Low-flow needle insert assembly	OPTON-20037

The HESI-II Probe Kit contains the HESI-II probe with a factory-installed high-flow needle insert. The high-flow metal needle insert is suitable for flow rates equal to or greater than 5 µL/min. For low-flow rate applications, order the low-flow needle insert assembly.

TSQ Quantum Access Instrument Specific Accessory Kit

Accessory Kit	70111-62059
Reserpine	00301-12901
Ultramark 1621	00301-12200
Needle seal	00950-00952
ESI needle	00950-00990
Ion transfer tube	70111-20972
Polytyrosine, liquid	00301-22924
Polytyrosine, solid	00301-22925
Kit, metal needle, high flow	70005-62013
Kit, metal needle, low flow	97144-62080
Cable, LC/MS	60053-63035
CsI	HAZMAT-01-00004

TSQ Accurate Mass Calibration Compounds Kit

Calibration Compounds Kit	70111-62029
Polyethylene glycol, 1000 molecular weight, 5 g	00301-07710
Polyethylene glycol, 200 molecular weight, 5 g	00301-07712
Polyethylene glycol, 400 molecular weight, 5 g	00301-07714
Polyethylene glycol, 600 molecular weight, 5 g	00301-07716

Fittings, Ferrules, and Sample Loops

Ferrule, Fingertight 2, Upchurch. (for: PEEK tubing and Teflon tube)	00101-18196
Ferrule, 0.016 in. ID, PEEK, Upchurch (for: Fused-silica capillary infusion line)	00101-18120
Ferrule, 0.012 in. ID, Kel-F, Upchurch (for: High flow metal needle and low flow metal needle)	00101-18114
Ferrule, 0.008 in. ID, Kel-F, Upchurch (for: Fused-silica capillary sample line)	00101-18116
Ferrule, LC, 1/16 in., stainless steel, Rheodyne	2522-3830
Fitting, Fingertight, Upchurch.	00101-18195
Fitting, Fingertight, Upchurch.	00101-18081

8 Replaceable Parts

Fittings, Ferrules, and Sample Loops

Fitting, adapter, Kel-F, 10-32 × 1/4-28, Upchurch (connects directly to ESI probe inlet).....	00101-18080
Nut, LC, 1/16 in., stainless steel, Rheodyne.....	2522-0066
Fitting, LC union, 0.010 in. orifice, PEEK.....	00101-18202
Fitting, LC TEE union, 0.020 in. orifice, PEEK.....	00101-18204
Fitting, adapter union, PEEK, Upchurch.....	00101-18206
5 µL sample loop, stainless steel.	00110-22010

Index

A

Accela System, interconnect diagram 48
accessory kit 107
air leaks, checking for 64
analyzer assembly, lenses, voltages applied to 35
analyzer chamber, location (figure) 39
Analyzer Control PCB, discussion 44
analyzer region, vacuum manifold, description 38
anode, electron multiplier, description 35
API probe, flushing 66
argon inlet, location (figure) 42
argon supply
 checking 64
 pressure 64
atmospheric pressure ionization (API) source
 description 71
 ion source interface 22
autosampler
 about 14
 configuring 14
 interconnect diagram (Accela System) 48
 startup 55
auxiliary gas pressure 42
auxiliary gas valve, description 42
auxiliary gas, on/off status 60

C

calibration
 and H-SRM (note) 63
 discussed 63
 frequency (note) 56, 63
capillary, heated. *See* ion transfer tube
cathode, electron multiplier, description 35
Cautions
 capillary heater cage assembly temperature 85
 emergency shutdown 49
 forepump and line power 41

 lower ion transfer tube temperature before
 removing 79
 nitrogen supply and API source operation 64
 replacing fuses 105
 replacing PCBs and power supplies 105
 scratching the skimmer cone 89
 solvent waste backing into ion source housing 72
centroid scan, defined 12
CID. *See* collision-induced dissociation
circuit breaker 20
cleaning procedures
 API probe, flushing 66
 ESI probe, flushing 66
 exit lens 87
 frequency 77
 ion sweep cone 79
 ion transfer tube 68, 79
 Q00 ion optics 91
 Q00 rf lens 91, 97
 sample transfer line 66
 sample tube 66
 skimmer 87–88
 S-lens 87
 sweep cone 68
 tube lens 87–88
collision cell
 figure 30
 location (figure) 39
collision energy (Q2 offset voltage) 34
collision gas valve, description 41
collision gas, description 32
collision-induced dissociation (CID), discussion 32
Communication LED
 description 19
 mass spectrometer reset 57
 mass spectrometer startup 55
compliance
 EMC iii
 FCC v

- regulatory compliance [iii](#)
- WEEE [vii](#)
- computer, features [45](#)
- contacting us [xxi](#)
- Convection gauge
 - description [41](#)
 - location (figure) [39](#)
 - pressure readings (table) [65, 83](#)
- conversion dynode
 - description [35](#)
 - figure [36](#)
 - on/off status [60](#)

D

- daily maintenance
 - API probe, flushing [66](#)
 - emptying solvent waste bottle [68](#)
 - ion sweep cone, flushing [68, 79](#)
 - ion transfer tube, flushing [68, 79](#)
 - sample tube, flushing [66](#)
 - transfer line, flushing [66](#)
- daily system checks
 - air leaks [64](#)
 - argon supply [64](#)
 - check list [66](#)
 - list [63](#)
 - nitrogen supply [64](#)
 - overview [63](#)
- data system
 - description [44](#)
 - LAN interface [46](#)
 - personal computer [45](#)
 - resetting [57](#)
 - shutdown [53](#)
- data types, discussed [11](#)
- data-dependent scan mode, discussed [9](#)
- diagnostics
 - Caution [104](#)
 - running [104](#)
- divert/inject valve
 - description [16](#)
 - location (figure) [17](#)
 - loop injection [16](#)
 - plumbing (figure) [17](#)
- dynode [35](#)

E

- electron multiplier
 - anode, description [35](#)
 - description [35](#)
 - on/off status [60](#)

- electronic assemblies
 - description [43, 105](#)
 - ion detection system [44](#)
 - Power Entry Module, description [43](#)
 - rf voltage generation, description [43](#)
- electronics service switch
 - and complete shutdown [52](#)
 - location [50](#)
 - mass spectrometer components on/off status [60](#)
- embedded computer
 - description [43](#)
 - resetting [57](#)
- EMC compliance [iii](#)
- emergency shutdown procedure [49](#)
- ESI probe, flushing [66](#)
- Ethernet Link OK LED [22](#)
- exit lens
 - cleaning [87](#)
 - reinstalling [89](#)
 - removing [86](#)

F

- FCC compliance [v](#)
- filters [43](#)
- fittings, replaceable parts [109](#)
- forepump
 - description [40](#)
 - maintenance [101](#)
 - on/off status [60](#)
 - power cord (Caution) [41](#)
- front panel LEDs
 - Communication [19](#)
 - description [19](#)
 - figure [19](#)
 - Power [19](#)
 - Scan [19](#)
 - System [19](#)
 - Vacuum [19](#)
- full scan type [10](#)
- functional description
 - data system [44](#)
 - divert/inject valve [16](#)
 - liquid chromatograph [15](#)
 - mass analyzer [28](#)
 - mass spectrometer [18](#)
 - syringe pump [15](#)
 - TSQ mass spectrometer [13](#)
- fuses, mass spectrometer, replacing [105](#)

I

- inlet gasses hardware
 - auxiliary gas valve [42](#)
 - collision gas valve [41](#)
 - description [37](#)
 - sheath gas valve [42](#)
- Instrument Configuration window, opening [14](#)
- interconnect diagram, Accela System [48](#)
- ion detection system
 - conversion dynode [35](#)
 - electron multiplier [35](#)
 - electron multiplier gain [36](#)
 - electronic assemblies [44](#)
- ion gauge
 - description [41](#)
 - pressure readings (table) [65](#), [83](#)
- Ion Max API source
 - housing (figure) [70](#)
 - housing, rear (figure) [71](#)
 - installing [71](#)
 - maintenance [75](#)
- Ion Max-S API source
 - housing (figure) [70](#)
 - installing [71](#)
- ion optics
 - cage assembly location (figure) [85–86](#), [88](#)
 - description [24](#)
 - Q0 quadrupole, description [27](#)
 - Q00 quadrupole, description [24](#)
- ion optics cage assembly
 - disassembling [96](#)
 - exploded view (figure) [96](#)
 - reinstalling [99](#)
- ion polarity modes, discussed [3](#)
- ion source interface
 - cross-sectional view [23](#)
 - description [22](#)
 - figure [85–86](#), [88](#)
 - ion transfer tube [23](#)
 - location (figure) [82](#)
 - maintenance [84](#)
 - reinstalling [90](#)
 - release lever (location) [72](#)
 - removing [85](#)
- ion source mount
 - figure [72](#)
 - guide pins [72](#)
 - location (figure) [85–86](#), [88](#)
- ion source, description [3](#)
- ion sweep cone
 - cleaning [82](#)
 - description [24](#)

- figure [82](#)
 - flushing [68](#), [79](#)
 - location (figure) [23](#)
- ion transfer tube
 - description [23](#)
 - figure [24](#)
 - flushing [68](#), [79](#)
 - location (figure) [23](#)
 - lower temperature before removing (Caution) [79](#)
 - part numbers [83](#)
 - removing and cleaning [79](#)
 - setting temperature [79](#)
- ion transmission device
 - defined [4](#)
 - rod assembly [4](#)

L

- leaks, checking for [64](#)
- LEDs
 - Communication LED [19](#)
 - Communication LED, mass spectrometer reset [57](#)
 - Communication LED, mass spectrometer startup [55](#)
 - divert/inject valve LEDs [20](#)
 - Ethernet Link OK [22](#)
 - Power LED [19](#)
 - Power LED, mass spectrometer reset [57](#)
 - Power LED, mass spectrometer startup [55](#)
 - Pump On, location (figure) [22](#)
 - Scan LED [19](#)
 - syringe pump LED [16](#), [19](#)
 - System LED [19](#)
 - System LED, mass spectrometer reset [57](#)
 - System LED, mass spectrometer startup [55](#)
 - Vacuum LED [19](#)
 - Vent Valve Closed [21](#)
- lens L0, description [27](#)
- lenses
 - analyzer assembly, voltages applied to [35](#)
 - L21, L22, L23, figure [30](#)
 - L31, L32, L33, figure [30](#)
- liquid chromatograph (LC)
 - about [15](#)
 - configuring [15](#)
 - functional description [15](#)
 - startup [54](#)
- liquid chromatograph, interconnect diagram (Accela System) [48](#)
- loop injection, divert/inject valve [16](#)

M

main power circuit breaker, description 20

maintenance

- API probe, flushing 66
- exit lens, cleaning 87
- forepump 101
- frequency (table) 76
- Ion Max API source 75
- Ion Max-S API source 75
- ion source interface 84
- ion sweep cone, cleaning 79
- ion transfer tube, cleaning 68, 79
- ion transfer tube, removing and cleaning 79
- Q00 ion optics, cleaning 91
- Q00 rf lens, cleaning 91, 97
- sample transfer line, flushing 66
- sample tube, flushing 66
- skimmer, cleaning 87–88
- S-lens, cleaning 87
- sweep cone, cleaning 68
- tube lens, cleaning 87–88

mass analysis

- collision-induced dissociation (CID) 32
- discussion 32
- rf and dc fields (figure) 30

mass analyzer

- description 4, 28
- on/off status 60

mass ranges 12

mass spectrometer

- collision gas valve 41
- diagnostics 104
- electronic assemblies 43, 105
- emergency shutdown 49
- frequency of cleaning 77
- front panel LEDs 19
- functional description 18
- fuses, replacing 105
- inlet gasses hardware 37
- interconnect diagram (Accela System) 48
- ion detection system electronic assemblies 44
- ion gauge 41
- LEDs and system startup 57
- maintenance procedures (table) 76
- mass analyzer 28
- Off condition 60
- Power Entry Module, description 43
- pressure readings (table) 65, 83
- resetting 57
- RF voltage generation electronic assemblies 43
- shutdown 52
- Standby condition 51
- starting up 54

turbomolecular pump 40

vacuum manifold 38

vacuum system 37

vent valve 41

mechanical pump 40

MS main power circuit breaker, mass spectrometer components on/off status 60

MS/MS scan modes

- neutral loss 7
- parent 6
- product 5

N

neutral loss scan mode

- discussed 7
- example (figure) 9
- illustration (figure) 8

nitrogen inlet, location (figure) 42

nitrogen supply

- Caution 64
- checking 64
- pressure 42, 64
- rate of consumption 64

O

Off condition, on/off status of components 60

offset voltage, quadrupole 33

P

parent scan mode

- discussed 6
- illustration (figure) 6

part numbers of replaceable parts 107

password 103

personal computer, features 45

power cord, forepump (Caution) 41

Power Entry Module, description 43

Power LED

- description 19
- mass spectrometer reset 57
- mass spectrometer startup 55

power panel, figure 20–21

pressure levels, checking 64

printer, setting up 46

procedures

- cleaning the exit lens 87
- cleaning the skimmer 88
- cleaning the S-lens 87
- cleaning the tube lens 88
- complete system shutdown 52

- configuring the LC 15
- disassembling the ion optics cage assembly 96
- installing the Ion Max API source 71
- installing the Ion Max-S API source 71
- maintaining the forepump 101
- placing system in standby 51
- reinstalling the exit lens 89
- reinstalling the ion optics cage assembly 99
- reinstalling the ion source interface 90
- reinstalling the skimmer 89
- reinstalling the S-lens 89
- reinstalling the tube lens 89
- removing skimmer 87
- removing the exit lens 86
- removing the S-lens 86
- removing tube lens 87
- resetting mass spectrometer 57
- product scan mode
 - discussed 5
 - illustration (figure) 6
- profile scan, defined 12
- pumps
 - forepump 40
 - turbomolecular pump 40

Q

- Q0 ion optics chamber, location (figure) 39
- Q0 ion optics, discussion 27
- Q0 mount assembly, location (figure) 96
- Q0 quadrupole
 - description 27
 - figure 27
- Q00 ion optics
 - cleaning 91
 - cleaning components 97
 - cross-sectional view 23
 - reassembling 98
 - skimmer 25
- Q00 RF lens
 - description 26
 - figure 26
 - on/off status 60
- Q00 rf lens
 - cleaning 91, 97
 - location (figure) 23
- Q00/L0 mount assembly
 - exploded view (figure) 97
 - location (figure) 96
- Q1 quadrupole, figure 29
- Q1MS and Q3MS scan modes 5
- Q2 rod assembly, figure 30
- Q3 quadrupole, figure 29

- quadrupole mass analyzer
 - functional description 29, 32
 - quadrupole offset voltage 33
- quadrupoles
 - mass analysis 29, 32
 - Q0, description 27
 - Q1 and Q3 29
 - RF and dc fields (figure) 31
 - rf and dc fields (figure) 30

R

- regulatory compliance iii
- replaceable parts
 - accessory kit 107
 - fittings 109
 - part numbers 107
 - sample loops 109
- RF voltage generation, discussion 43
- right side power panel
 - discussion 21
 - figure 21
- rod assemblies
 - description 30
 - ion transmission 4
 - mass analysis 4
 - mass analyzer(s), quadrupole 29, 32
 - Q1, Q2, and Q3 (note) 5
- roughing pump 40

S

- safety precautions xviii
- sample loop, replaceable parts 109
- sample transfer line, flushing 66
- sample tube, flushing 66
- scan data types
 - centroid scan 12
 - profile scan 12
- Scan LED, description 19
- scan modes
 - data-dependent 9
 - discussed 3
 - mass spectrometer 3
 - neutral loss 7
 - neutral loss, example (figure) 9
 - neutral loss, illustrated (figure) 8
 - parent 6
 - parent, illustrated (figure) 7
 - product 5
 - Product, illustrated (figure) 6
 - Q1MS and Q3MS 5
 - summary (table) 4

scan types
 discussed 10
 full-scan 10
 selected ion monitoring (SIM) 10
 selected reaction monitoring (SRM) 11
 selected ion monitoring (SIM) scan type 10
 selected reaction monitoring (SRM) scan type 11
 sheath gas pressure 42
 sheath gas valve, description 42
 sheath gas, on/off status 60
 shutdown
 data system 53
 emergency procedure 49
 mass spectrometer 52
 nonemergency procedure 52
 side cover plate (vacuum manifold), description 40
 skimmer
 cleaning 88
 description 25
 figure 26
 location (figure) 23
 reinstalling 89
 removing 87
 S-lens
 cleaning 87
 on/off status 60
 reinstalling 89
 removing 86
 solvent waste bottle, emptying 68
 solvents
 cleaning 88
 purity requirements xix
 Standby condition
 on/off status of components 60
 placing system in 51
 startup
 autosampler 55
 liquid chromatograph 54
 mass spectrometer 54
 operating conditions, setting 56
 sweep gas supply port, location (figure) 82
 switches 21
 syringe pump
 description 15
 figure 16
 system checks
 argon and nitrogen supplies 64
 sample tube polyimide elongation 64
 vacuum levels 64
 System Control PCB, description 43

System LED
 mass spectrometer reset 57
 mass spectrometer startup 55
 system reset button
 description 21
 location (figure) 21
 mass spectrometer reset 57
 system shutdown
 emergency procedure 49
 nonemergency procedure 52

T

tables
 mass spectrometer components on/off status 60
 mass spectrometer maintenance procedures 76
 pressure readings 65, 83
 summary of scan modes 4
 troubleshooting 103
 TSQ system
 autosampler 14
 calibrating 63
 calibration, frequency 56
 complete shutdown 52
 daily system checks 63
 data system 44
 data types 11
 diagnostics 104
 divert/inject valve 16
 emergency shutdown 49
 front panel LEDs 19
 functional block diagram (figure) 13–14
 functional description 13
 ion polarity modes 3
 LEDs and system startup 57
 liquid chromatograph 15
 mass ranges 12
 mass spectrometer 18
 operating conditions, setting 56
 overview 1
 regulatory compliance iii
 replaceable parts 107
 scan modes 3
 scan types 10
 shutdown, emergency procedure 49
 shutdown, nonemergency procedure 52
 Standby condition 51
 starting up 54
 syringe pump 15
 tuning 63
 tuning, frequency 56

- tube lens
 - cleaning [88](#)
 - description [24](#)
 - figure [24](#)
 - location (figure) [23](#)
 - on/off status [60](#)
 - reinstalling [89](#)
 - removing [87](#)
 - voltages [24](#)
- Tune Master window, opening [16](#)
- tuning
 - discussed [63](#)
 - frequency (note) [56](#)
 - tube lens offset voltage [24](#)
- turbomolecular pump
 - description [40](#)
 - on/off status [60](#)

V

- Vacuum LED
 - and analyzer region pressure [19](#)
 - description [19](#)
- vacuum levels, checking [64](#)
- vacuum manifold
 - analyzer chamber, location (figure) [39](#)
 - collision cell chamber, location (figure) [39](#)
 - description [38](#)
 - feedthroughs [39](#)
 - interior (figure) [39](#)
 - pressure readings (table) [65](#), [83](#)
 - Q0 ion optics chamber, location (figure) [39](#)
 - side cover plate, description [40](#)
- vacuum pumps
 - forepump [40](#)
 - turbomolecular pump [40](#)
- vacuum service switch
 - and mass spectrometer shutdown [52](#)
 - location [50](#)
 - mass spectrometer components on/off status [60](#)
- vacuum system
 - auxiliary gas valve [42](#)
 - block diagram [38](#)
 - checking pressure levels [64](#)
 - collision gas valve [41](#)
 - description [37](#)
 - forepump [40](#)
 - ion gauge [41](#)
 - powering off [21](#)
 - pressure readings [65](#), [83](#)
 - sheath gas valve [42](#)
 - turbomolecular pump [40](#)
 - Vacuum LED [19](#)

- vacuum service switch [21](#)
- vent valve [41](#)
- vent valve
 - description [41](#)
 - LED [21](#)
- Vent Valve Closed LED [21](#)
- voltages
 - conversion dynode [35](#)
 - ion transfer tube [23](#)
 - Q0 ion optics [27](#)
 - Q0 offset [27](#)
 - Q00 offset [26](#)
 - quadrupole rf and dc [30](#)
 - tube lens offset [24](#)

W

- WEEE compliance [vii](#)

